

softWoRx[®]

Imaging Workstation

User's Manual

Revision C

built with precisionware[®]

AppliedPrecision[®]

Legal Notices

Revision C of the User's Manual for the *softWoRx* Imaging Workstation. Part number 04-720103-000 Rev C.

© 1999-2008 Applied Precision, Inc. All rights reserved. No part of this manual may be reproduced, transmitted, stored in a retrieval system, or translated into any language in any form by any means without the written permission of Applied Precision, Inc.

Information in this document is subject to change without notice.

DeltaVision, Applied Precision, and *softWoRx* are registered trademarks of Applied Precision, Inc. All other registered names and trademarks referred to in this manual are the property of their respective companies.

Applied Precision, Inc
1040 12th Ave. NW
Issaquah, WA 98027
(425) 557-1000 FAX: (425) 557-1055

Other Manuals and Guides

The following documents are provided for *softWoRx*.

Document	Purpose	Available for...
Online Help	Provides reference information for <i>softWoRx</i> and procedures that show how to use <i>softWoRx</i> tools	All <i>softWoRx</i> workstations
Product Notes	Provide examples and tips for using <i>softWoRx</i>	All <i>softWoRx</i> users (online at www.appliedprecision.com)
<i>The DeltaVision RT Restoration Microscopy System User's Manual</i>	Shows how to acquire data and how to maintain the data acquisition system	Acquisition workstations
<i>Getting Started with QLM</i>	Shows how to acquire photokinetic data with the QLM module	Acquisition workstations that have the optional QLM module
<i>RedHat Linux Bible</i>	Shows how to use Linux	All <i>softWoRx</i> Users. (This is a third party manual.)

Contents

Preface	i
About This Manual.....	i
Document Conventions	ii
Lists	ii
Notes, Warnings and Cautions	ii
User Interface Description Conventions.....	ii
Contacting Applied Precision, Inc.....	iii
Customer Service Hotline	iii
Corporate Office	iii
Acknowledgements	iii
Introduction	1
What is <i>softWoRx</i> ?	1
What Can You Use <i>softWoRx</i> for?.....	2
Acquiring Data	2
Processing Data	3
Visualizing and Presenting Data	3
Analyzing Results	4
Optional Components.....	5
The Quantitative Laser Module	5
<i>softWoRx</i> Explorer	5
Part One: Processing Data	7
1. Deconvolving Image Data	9
About Deconvolution Processing.....	10
Deconvolution Tools	10
Deconvolving an Image	11
Deconvolving Several Images.....	13
Common Deconvolution Options	14
More about the <i>softWoRx</i> Deconvolution Tools	16
2. Correcting Images	17
About Correcting Images	17
Correcting Z Section Image Data	18
Equalizing Intensities in Time-lapse Image Data.....	19
Calibrating Image Data.....	20
Aligning Adjacent Images	21
Correcting Chromatic Aberration	23

3. Stitching	25
Stitching Images That Have a Single Z Section	26
Stitching Images That Have Multiple Z Sections	27
4. Importing Data	31
Converting TIFF Images	32
Converting ISee Files.....	33
ISee Series Conversion.....	34
Converting Pic Files.....	35
Converting STK Files.....	36
5. Data and Task Manipulation	39
Selecting Data	40
Selecting Rectangular Data Regions	40
Selecting Irregular Data Regions.....	42
Cropping and Trimming Data	45
Cropping a Rectangular Region.....	45
Cropping an Irregular Data Region	47
Trimming Time Data	50
Combining Data of Two Images.....	52
Setting Up Process Chains with Task Builder	54
Using Ratio Imaging.....	57
Using the Multiplexed Wavelength Option.....	62
Setting Up the Multiplexed Wavelength Option.....	63
Designing a Multiplexed Wavelength Experiment	68
6. DMS Integration.....	73
What is DMS?	73
Connecting to a DMS Database	74
Uploading Images to a DMS Database	74
Uploading Images from the Image Window.....	74
Uploading Images Using Task Builder	77
Auto-uploading Images after Acquisition.....	80
Downloading Files from DMS	81
Browsing and Locating Images in a DMS Database	82
Browsing Image Files using P/D/I Hierarchy.....	82
Browsing Image Files using CG/C/I Hierarchy.....	83
Searching for Image Files Based on Annotation	84

Part Two: Visualizing & Presenting Data	85
7. Viewing Image Data.....	87
Opening an Image	88
The Image Window	89
Viewing 5D Images	90
Viewing Z Sections and Time Points.....	90
Viewing Areas	92
Displaying or Hiding Channels	94
Adjusting Brightness and Contrast.....	95
Assigning Colors or Grayscale to Channels	99
Grayscale Mode.....	100
Color Mode	101
Blended Color Mode.....	101
Controlling the Image Window Display	102
Hiding or Displaying Image Window Border Tools.....	103
The Image Window Scale Bar.....	103
Resizing or Reorienting an Image	105
Resizing an Image	105
Rotating an Image	107
Viewing Cross Sections.....	111
8. Viewing Movies	113
Viewing Volumetric or Time-Lapse Movies.....	113
Tracking Particle Movement with Trails Movies.....	114
9. Viewing Projections and Volumes	119
Creating 2D Projections	119
Creating Volume Projections	122
Volume Rendering.....	122
Axes of Rotation	123
Methods for Projecting Volumes	125
10. Filtering Image Data	137
About <i>softWoRx</i> Filters	138
Using Convolution Filters	138
Enhancing Object Boundaries.....	139
Using 2D Statistical Filters.....	141
Using Image Arithmetic	142
Scaling Pixel Intensity to Enhance Local Contrast.....	143
Setting an Intensity Threshold	144

11. Saving, Exporting, and Printing	147
Image Data Files and Image Graphic Files.....	148
Saving <i>DeltaVision</i> Files.....	148
Saving to the Data Management System (DMS).....	150
Exporting <i>DeltaVision</i> Files.....	151
Exporting to TIFF Files.....	151
Exporting to PhotoShop Files.....	153
Exporting to Movie Files.....	153
Exporting to JPEG Files.....	155
Capturing Screen Shots.....	156
Archiving Files to CD/DVD.....	157
Printing Images.....	159
Part Three: Analyzing Results	161
12. Examining Intensity Data	162
Examining Point Values.....	162
Examining Intensity Data with Data Inspector.....	163
Selecting a Region of Interest.....	165
Viewing Intensity Line Profiles.....	166
Viewing the Line Intensity of a Row or Column.....	167
Viewing the Line intensity in Any Direction.....	168
Calculating Statistics.....	169
Calculating Statistics for Selected Areas.....	169
Calculating Statistics for Irregular Areas.....	171
13. Measuring Distance and Velocity	175
Measuring Distances.....	175
Measuring Velocity.....	177
14. Volume Modeling	181
About Volume Modeling.....	182
Edit Polygon Dialog Box.....	182
2D Polygon Finder.....	185
3D Object Builder.....	187
Creating and Viewing the 3-D Object.....	188
Volume Modeling Example.....	190
15. Detecting and Analyzing Colocalization	193
Examining the Entire Image.....	194
Identifying Potential Colocalized Areas.....	197
Detecting Colocalization with ROIs.....	201

16. Other Applications.....	203
About Photokinetics.....	203
Analyzing Fluorescence Recovery After Photo-bleaching	204
About FRAP Experiments.....	205
Analyzing Fluorescence Resonance Energy Transfer.....	208
Using the FRET Analysis Tool	208
Appendix A: Image Quality.....	215
Using Deconvolution Residuals	216
Visually Evaluating Images	217
Index	221

Preface

This manual shows how to use *softWoRx* to process, visualize, and analyze image data.

- *About This Manual* describes the information in the manual.
- *Document Conventions* explains the typography, notes, and other conventions used in this manual.
- *Contacting Applied Precision, Inc.* provides information about how to contact customer support.

About This Manual

This manual is divided into three parts that contain the following information:

- Part One includes instructions for processing and importing data.
- Part Two shows how to visualize data and prepare it for presentations. It also shows how to save or export data in a variety of formats.
- Part Three shows how to use *softWoRx* tools to perform quantitative analysis.

The manual also includes an appendix that shows how to analyze image quality.

Document Conventions


In order to make the information in this manual as easy as possible for you to locate and use, the following conventions are observed.

Lists


- Round bullets indicate options in procedures.
- 1. Numbered items are sequential steps for completing a procedure.
- Square bullets indicate items in a list.
- ▶ Arrows indicate single step procedures.

Notes, Warnings and Cautions

 **Note** Indicates information about the previous paragraph or step in a procedure.


 **Important** Indicates important or critical information about the previous paragraph or step in a procedure.

 **Tip** Indicates helpful advice.

 **WARNING:** Indicates important information regarding potential injury.

 **WARNING:** Indicates risk of explosion.

 **WARNING:** Indicates risk of shock.

 **CAUTION:** Indicates important information regarding potential damage to equipment or software.

User Interface Description Conventions

Boldface indicates the names of buttons, menus, dialog box options, and fields.

Initial Capitals indicate the names of windows, dialog boxes, and tabs.

ALL CAPITALS SAN SERIF indicates the name of a key on your keyboard, such as ENTER or DELETE.

Uniform width font indicates text to enter on a command line or in the GUI.

Contacting Applied Precision, Inc

If you have questions about *DeltaVision*, first refer to this manual or consult the online Help system. If you don't find the information you need, contact us at one of the following addresses.

Customer Service Hotline

Phone: 800-862-5166

email: servicehotline@api.com

Hours: 8:00 AM – 5:00 PM, Pacific Time, Monday – Friday

Corporate Office

Applied Precision, Inc

1040 12th Avenue NW

Issaquah, WA 98027

USA

Phone: (425) 557-1000

Fax: (425) 557-1055

Internet address: www.appliedprecision.com

Acknowledgements

Applied Precision would like to thank Adrian Quintanilla and Dave McDonald of the Fred Hutchinson Cancer Research Center for providing a data acquisition workflow and other content that is used in this manual.

Introduction

This introduction provides an overview of how you can use *softWoRx* to process, visualize, and analyze multidimensional microscopy data. It also introduces optional *softWoRx* components.

In the Introduction

What is <i>softWoRx</i> ?	1
What Can You Use <i>softWoRx</i> for?	2
Optional Components	5

What is *softWoRx*?

softWoRx is a comprehensive software package designed for the analysis of multi-dimensional microscopy data. Although originally developed for use as a component of Applied Precision's *DeltaVision* Restoration Microscope System, *softWoRx* is now also available on a stand-alone analysis workstation, giving you a powerful yet friendly environment for exploring and refining your understanding of specimen structure. The flexibility of the software makes it ideal for the study of

images from fluorescence, brightfield, Differential Interference Contrast (DIC), and electron microscopy.

softWoRx is available on two types of Linux workstations:

- The Acquisition workstation is part of the *DeltaVision* data acquisition system and is used to control the system. You can also use it to process, analyze, and visualize data. (Refer to the *DeltaVision Core and personalDV Restoration Microscopy System User's Manual* for details on image acquisition.)
- The Analysis workstation is a stand-alone workstation. You can use it only to process, visualize, and analyze data.



The *DeltaVision Core* data acquisition system

What Can You Use *softWoRx* for?

You can use *softWoRx* to acquire, process, visualize, and analyze multi-dimensional renderings of a fluorescent specimen. You can also use it to save data in a variety of formats.

Acquiring Data

If you are using an acquisition workstation, you can acquire images with the *DeltaVision* Restoration Microscope System. The *softWoRx* Resolve3D module provides various options for acquiring time-lapse data, data with multiple Z sections, and data from multiple channels. (Refer to the *DeltaVision Core and*

personalDV Restoration Microscopy System User's Manual for details on image acquisition.)

If your system has the QLM module, you can use it to acquire photokinetic data for a variety of experiments. (*softWoRx* photokinetics data includes photo-bleaching or photo-activation that results from a laser pulse. See the *QLM Getting Started Guide*.)

Processing Data

Process image data to prepare it for visual examination and analysis. *softWoRx* provides several types of modules for processing image data.

Deconvolving Image Data

Deconvolve image data acquired with the *DeltaVision* system. Deconvolving improves contrast by relocating signal scatter and out-of-focus data.

Correcting Images

Correct image data for chromatic aberration (color shift) that results from oil matching and other environmental conditions. You can also correct data collection errors and equalize intensity values across Z sections.

Stitching

Stitch "panel" images collected with *DeltaVision* to generate a larger overall field of view. Stitched images are organized as either a series of time points or Z sections.

Importing

Import data from the TIFF format (16-bit grayscale), BioRad's Pic format, InoVision's ISee format, or MetaMorph's STK format.

Selecting, Cropping, and Combining data

Select data to crop it or to combine it with other data.

Visualizing and Presenting Data

After processing data, you can view and present it in a variety of ways. *softWoRx* provides several tools that you can use to visualize data and prepare it for presentations.

Viewing Image Data

Open data files in *softWoRx* and adjust the way that the image is displayed (e.g., display a scale bar, set grayscale or color modes, or adjust brightness and contrast). You can also rotate or resize image data or view data cross-sections.

Viewing Projections and Volumes

Render volumes and create 2D projections to visualize and explore three-dimensional data. Several methods for rendering volumes that you can

interactively rotate are available. The 2D projections quickly combine information from multiple Z Sections into a single section.

Viewing Movies

Create movies of volume rendered data or time-lapse data. You can also create movies to trace particle movement.

Filtering

Choose from several filters to improve the visual presentation of data, prepare data for modeling, or for other types of analysis. You can use statistical filters that are useful for removing noise from the image, threshold filters, and convolution filters.

Saving, Exporting, and Printing

Save and present data in a variety of formats. Export images to PhotoShop or JPEG formats or save image data in a *DeltaVision* file, a TIFF file, or a tabular format that can be opened in a spreadsheet. You can also save time-lapse or volume-rendered data as MPEG movies. All files can be archived to CD or DVD. If your system is configured with a printer, you can print *DeltaVision* files from *softWoRx*. If you have *softWoRx Explorer*, you can print *DeltaVision* files from a Mac or PC computer.

Analyzing Results

You can use measuring and modeling tools to perform quantitative analysis.

Examining Intensity Data

Study area and line profiles, calculate statistics, and display single point values.

Measuring Distance and Velocity

Measure features on an XY plane or across Z sections. You can also measure the velocity of particle movement.

Modeling

Use tools to create line models or volume models.

Detecting Colocalization

Use Colocalization modules to create a scatter plots and measure the Pearson Coefficient of Correlation to help determine whether colocalization is occurring.

Analyzing Fluorescence Resonance Energy Transfer Data

Use the Fluorescence Resonance Energy Transfer (FRET) module to analyze FRET data.

Analyzing Fluorescence Recovery After Photo-bleaching Data

Use the Analyze Fluorescence Recovery After Photo-bleaching (FRAP) module to analyze FRAP data.

Optional Components

You can purchase two optional components for *softWoRx*:

The Quantitative Laser Module

The Quantitative Laser Module (QLM) is a *DeltaVision* component that adds a laser beam into the back aperture of the microscope objective to provide a focused illumination spot in the center of the optical field. This optional component mounts to the Fiber Optic Module of a *DeltaVision Core* microscope, a *DeltaVision RT* microscope, or a microscope that is upgraded to the *DeltaVision 3.9* level.

If your system has the QLM hardware module, you can use *softWoRx* to analyze Photokinetic (photo-bleaching and photo-activation) experiments. The *softWoRx* FRAP analysis module discussed in Chapter 16 is only available for systems that have QLM hardware.

softWoRx Explorer

softWoRx[®] Explorer is a cross-platform image viewer available for many commonly used operating systems.

softWoRx Explorer allows you to view and explore *DeltaVision* images and images from other sources that contain spatial, temporal, and spectral ranges. In addition to displaying data in the X and Y plane, you can scroll through Z sections and time-lapse data. Individual spectrum (i.e., channels or fluorescent wavelengths) can be hidden or displayed in a variety of colors.

Part One

PROCESSING DATA

You will typically need to process data before you view or analyze it. Part one includes instructions for processing and importing data.

In Part One

Chapter 1. Deconvolving Image Data	9
Chapter 2. Correcting Images	17
Chapter 3. Stitching	25
Chapter 4. Importing Data	31
Chapter 5. Data and Task Manipulation	39
Chapter 6. DMS Integration	39

1. Deconvolving Image Data

This chapter shows how to use the *softWoRx* Deconvolve tool to remove blur in fluorescence optical sections.


You can deconvolve and view a single image or you can set up a queue to deconvolve several images. You can also set options to deconvolve a region of an image, select which wavelengths to include, or select the deconvolution method.

In This Chapter

Converting TIFF Images.....	32
Converting ISee Files	33
Converting Pic Files	35
Converting STK Files	36

About Deconvolution Processing

After you acquire images, you'll need to process them to remove blurred data. The process of relocating signal scatter and out-of-focus information present in digital images is known as Deconvolution. Applied Precision's proprietary deconvolution algorithms preserve the amount of light throughout the entire Z-stack. Blurred data is not dropped out. It is reassigned to its original location. This process increases contrast and comparative intensities within each Z-stack image, making for extremely sharp 3D reconstructions.

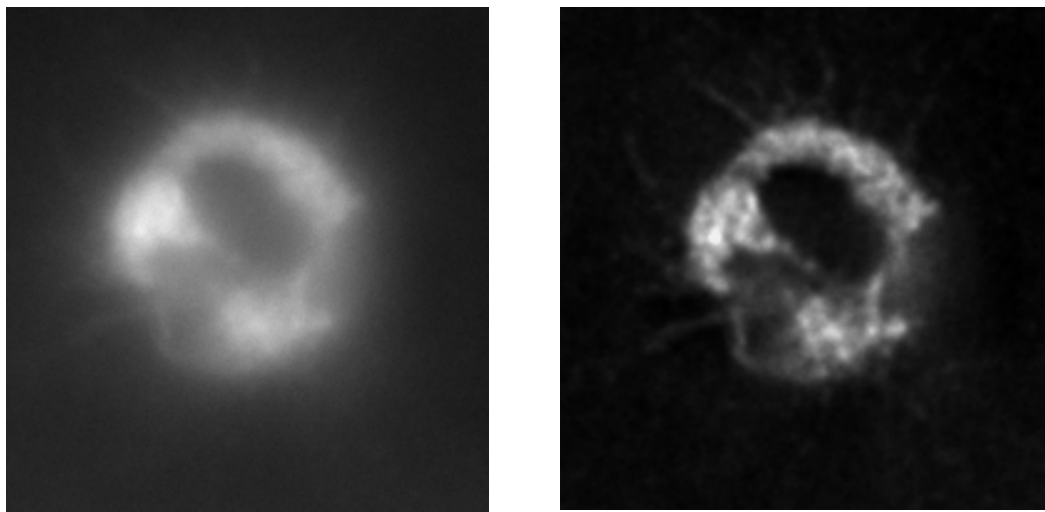

Unprocessed data on the left is deconvolved to create the image on the right

Deconvolution Tools

softWoRx provides two tools for deconvolving images: Deconvolve and Nearest Neighbor. The Deconvolve tool (described in this chapter) provides the best results for most applications. This method uses the iterative-constrained algorithm described by Agard¹. This tool should be used for experiments where the quantification of intensities is required.


The Nearest Neighbor tool (described only in the online Help) uses an approximate deconvolution approach, commonly referred to as *deblurring*, for removing blur from optical sections. The Nearest Neighbor tool should not be used in experiments where quantification of intensities is required.

¹ See: Agard, D.A. (1984) Optical Sectioning Microscopy: Cellular Architecture in Three Dimensions. *Ann. Rev. Biophys. Bioeng.* 13:191-219.

Deconvolving an Image

To deconvolve an image:

1. On the *softWoRx* main menu, choose **Process | Deconvolve** to open the Deconvolve dialog box.

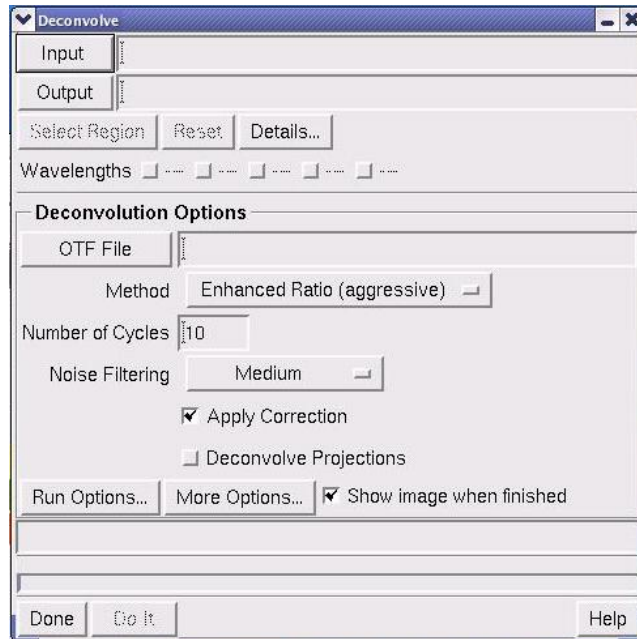

2. Enter the original `_R3D.dv` image file (for example, `/usr/local/softWoRx/data/samples/oocyte_R3D.dv`) in the **Input** field. You can do this in the following ways:

- From the Linux file Manager, drag and drop the file in the **Input** field.
- Click **Input** and browse to the file.
- Type the path and file name into the **Input** field.


softWoRx creates an output file name by appending the `_D3D` extension to the input file name. The new name (for example, `oocyte_R3D_D3D.dv`) is displayed in the **Output** field.

3. Enter the `.otf` file (for example, the `oocyte.otf` file in the `usr/local/softWoRx/data/samples` directory) into the **OTF File** field. You can use the same methods (drag and drop, browse, or type) to enter this file as you used in Step 2. For *DeltaVision* files, this is done automatically.

Note The `.otf` file is an Optical Transfer Function (OTF) file. In many microscopy systems, there is only one OTF per objective lens and the correct OTF is simply the one that corresponds to the lens used for measuring the optical

sections. Refer to the lens identification number in the measured data and the OTF to verify that the correct OTF is being used.

4. Click **Do It**.

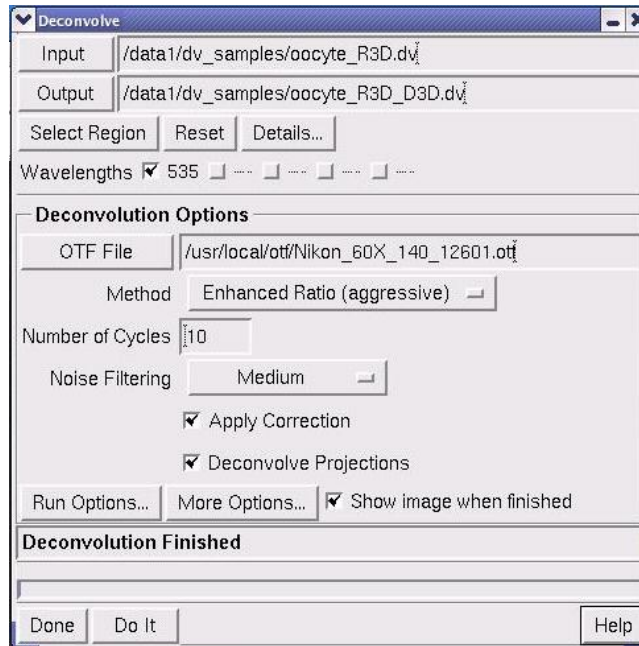


The deconvolution status is displayed in the Deconvolve Output window and on the bottom of the Deconvolve dialog box. When the deconvolution is finished, messages appear in each of these windows and the deconvolved image is displayed in the Image window.



The deconvolved image

- If you did not have the **Show image when finished** checkbox selected (default) on the Deconvolve dialog box, choose **File | Open** to open the View Image dialog box.
- Enter the file name (for example, `usr/local/softWoRx/data/samples/oocyte_R3D_D3D.dv`) into the **Input** field.

 **Note** You can use the same methods (drag and drop, browse, or type) to enter this file as you used to enter the `_R3D.dv` file in Step 2 on page 11.


- Click **Do It** to open the file in the Image window.

Deconvolving Several Images

You can create a queue to deconvolve several images and specify a time to start the deconvolution.

To deconvolve several images:

- On the *softWoRx* main menu bar, click **Process | Deconvolve** to open the Deconvolve dialog box.
- In the **Input** field, enter the `_R3D.dv` file (for example, `/usr/local/softWoRx/data/samples/oocyte_R3D.dv`). You can use the same methods (drag and drop, browse, or type) to enter this file as those used in Step 2 of the previous procedure.
- In the **OTF File** field, enter the `.otf` file (for example, `/usr/local/softWoRx/data/samples/Olympus_60X_142_10612.otf`).
- In the Deconvolve dialog box, click **Run Options** to open the Deconvolution Run Options dialog box.

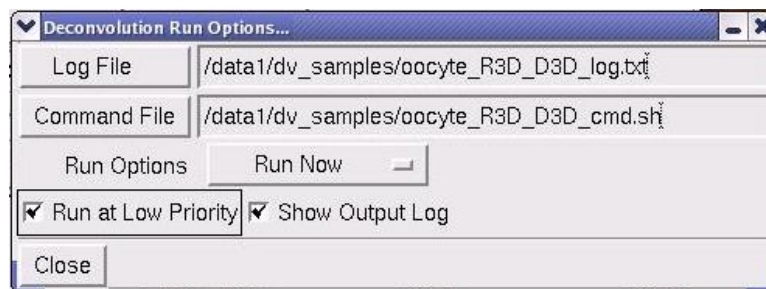


- In the **Run Options** pull-down list, select **Add to Queue**. Then click **Close**.

6. In the Deconvolve dialog box, click **Do It**. The file is added to the queue and displayed in the Queue Manager.



7. Repeat Steps 2 and 3 and click **Do It** in the Deconvolve dialog box for each of the remaining files.

 **Tip** If your file names are all the same except for the last digit (for example 040600aq01, 040600aq02, etc.), you can simply overwrite the last digit and press **Do It** for each file.

8. In the Queue Manager dialog box, choose one of the following options:

To perform the deconvolutions immediately, click **Start Now**.

or


To perform the deconvolutions later, click **Start Later** and select a time on the clock that appears in the dialog box.

As the files are deconvolved, the deconvolution status is displayed in the Queue Manager.

9. Click **Quit** to close the Queue Manager dialog box.

Common Deconvolution Options

You can set options to deconvolve a region of an image, deconvolve only data in specified wavelengths, or change the deconvolution method.

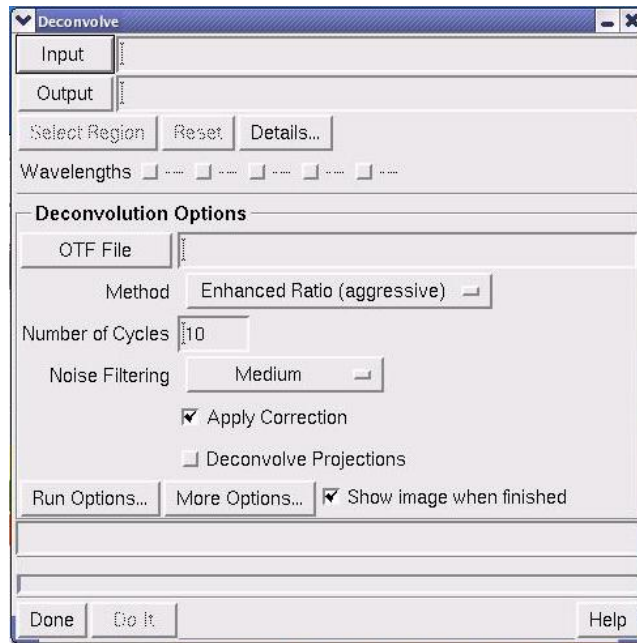

Use the Deconvolve dialog box to specify options for deconvolving

To	Do This
Deconvolve only part of the image	Click Select Region and use your mouse to define a specific region of the image to deconvolve. (This option is only available when the input data comes from a window.)
Select which wavelengths to include	Select the Wavelengths options.
Select a deconvolution method	<p>Choose one of the following deconvolution methods from the Method list:</p> <p>Ratio (conservative) method uses a more conservative algorithm that generally finds an accurate solution. Images with punctate fluorescence may deconvolve better using this method.</p> <p>Enhanced Ratio (the default method) is quicker because the residuals stabilize in fewer iterations, typically 10 or less.</p> <p>Additive uses a more conservative algorithm that generally finds an accurate solution.</p> <p>Enhanced Additive is faster than Additive because it requires fewer iterations (typically ten or fewer).</p> <p>The Additive and Enhanced Additive options are the preferred deconvolution methods for data acquired with the EMCCD electron multiplication camera. These methods are more tolerant of noisy data (images with higher noise levels).</p>
Display deconvolved images immediately after processing	Select the Show image when finished checkbox.

You can typically use the default settings for the rest of the options in the Deconvolve dialog box, including the options displayed when you click **More Options**.

More about the *softWoRx* Deconvolution Tools

softWoRx provides two tools for deconvolving images: the standard Deconvolve tool described in this section and the Nearest Neighbor Deconvolution tool. In most instances, the standard Deconvolve tool provides the best results. The Nearest Neighbor tool provides an approximate deconvolution approach for removing blur from optical sections. If you wish to learn more about Nearest Neighbor Deconvolution, refer to the online Help.

The standard Deconvolve tool uses the Constrained Iterative Deconvolution algorithm to remove the out-of-focus blur in fluorescence optical sections. This algorithm calculates a result using the following four steps:

1. The algorithm estimates what the object looks like.
2. The estimate is mathematically blurred to simulate the effects of the microscope's limited aperture.
3. The blurred estimate is compared to the actual image. The difference between the images is then used to modify the estimate.
4. The modified estimate is constrained to be non-negative, by setting pixels with negative intensity to 0.

The algorithm repeats this sequence of steps until the estimate, convolved with the point spread, closely approximates the actual image (see Agard²).

² See: Agard, D.A. (1984) Optical Sectioning Microscopy: Cellular Architecture in Three Dimensions. *Ann. Rev. Biophys. Bioeng.* 13:191-219.

2. Correcting Images

softWoRx provides several tools for correcting image data. You may need to correct data to prepare it for visualization and analysis.

In This Chapter

About Correcting Images	17
Correcting Z Section Image Data	18
Equalizing Intensities in Time-lapse Image Data	19
Calibrating Image Data	20
Aligning Adjacent Images	21
Correcting Chromatic Aberration.....	23

About Correcting Images

Correct Image is used to correct errors that are caused by photo-bleaching, inconsistent illumination intensity, or CCD defects. The Correct Image tool corrects the intensity values of sections within a Z series. With the exception of photo-bleaching, the tool can also correct intensity values between time points of a time-lapse experiment. By default, *softWoRx* automatically applies this tool during the deconvolution process. (See *Correcting Z Section Image Data* on Page 18.)

Equalize Time Points equalizes intensities of all time points to a reference time point that you select. The tool uses the mean image intensity to help generate a more uniform intensity display. You can use this tool to normalize time-lapse data for display purposes. In general, the data generated by this tool should be used for display only and should not be used for quantitative purposes. (See *Equalizing Intensities in Time-lapse Image Data* on Page 19.)

Calibrate calibrates a raw image when you have a calibration file and a bad pixel file that applies to the camera and conditions (array size, pixel size, and wavelength) that were used to collect the image. (See *Calibrating Image Data* on Page 20.)

Align Image corrects single wavelength images that have motion artifacts, problems with Z sectioning, or problems with time series. It allows you to align adjacent images by applying an XY shift with an optional rotation. (See *Aligning Adjacent Images* on Page 21.)

Chromatic Aberration Corrector allows you to adjust channels relative to each other to correct for shifts in color that result from oil matching, objective anomalies, and other environmental conditions that use X-Z and Y-Z image profiles. (See *Correcting Chromatic Aberration* on Page 23.)


Correcting Z Section Image Data

Correct Image options are used to correct systematic errors that occur during data collection. The three basic systematic errors are caused by photo-bleaching, inconsistent illumination intensity, and CCD defects.

By default, *softWoRx* automatically applies these options to images during the deconvolution process (These options are specified in the More Deconvolution Options dialog box).

If you are analyzing unprocessed images (images that are not deconvolved), use the Correct options tool to apply these options to the images before you process them. Applying correction options is especially important when you are performing quantitative analysis on images that contain multiple Z sections or Time series.

1. Open the Correct Image dialog box by choosing **Process | Correct** from the *softWoRx* main menu.

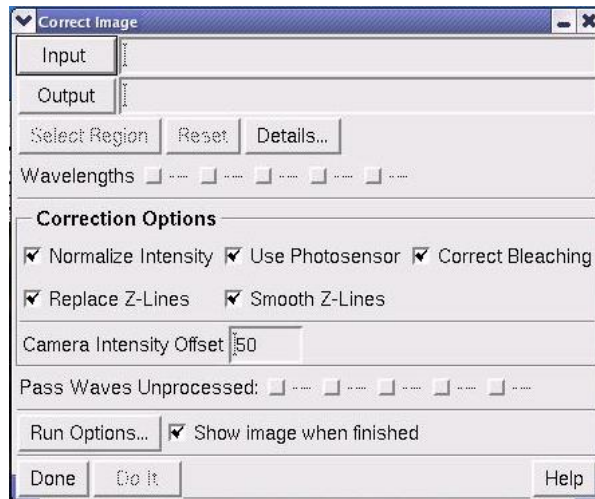

2. Enter the original `_R3D.dv` image file in the **Input** field.
3. Use the **Correction Option** toggles to select the desired image correction. See the *softWoRx* online help for a description of each of the **Correction** and **Run** options available from this dialog box.
4. When you are satisfied with your selections, select **Do It** to perform the correction process. Details of the process are displayed in a Correct Image Output window.

Equalizing Intensities in Time-lapse Image Data

Use Equalize Time Points to choose a reference time point and normalize the intensities of all other time points to it. The tool uses the mean image intensity to help generate a more uniform intensity display.


Note In general, the data generated by the Equalize Time Points tool should be used for display only and not for quantitative purposes.

The Correct tool and the Correction options of the Deconvolve tool make corrections to the intensity values of sections within a Z series and, with the exception of photo-bleaching inconsistencies, can also make corrections to intensity values between time points of a time-lapse experiment.

Note This tool is **not** intended to correct for photo-bleaching over time.

To equalize intensities to a time point:

1. Choose **Process | Equalize Time Points** from the main *softWoRx* menu to open the Equalize Time Points dialog box.

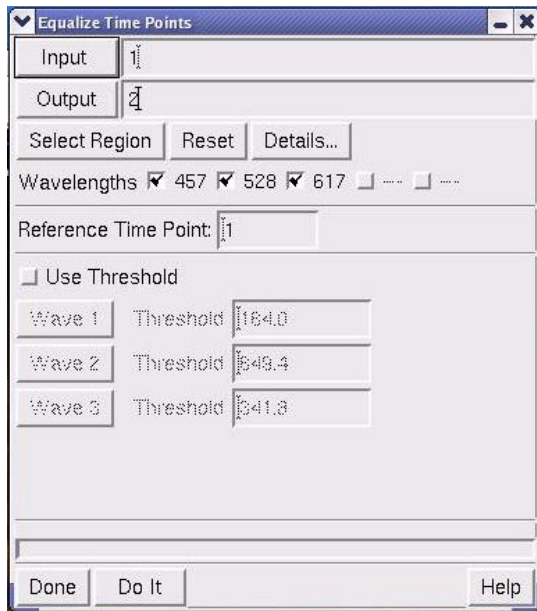

2. Enter the original `_R3D.dv` image file in the **Input** field.
3. Enter the time point to use as the standard for adjusting intensity values of other time points in the **Reference Time Point** field.
4. To perform equalization based on background and minimize the influence of the variation of higher intensities (signal), select **Use Threshold**. (This specifies to use only values less than the **Threshold** value for each wavelength when collecting **Min/Max/Mean** statistics for equalization.)
5. If you are using a threshold, specify the wavelengths in the **Wave** fields and specify the threshold for each wavelength.

Note The number of Wave fields in the Equalize Time Points dialog box adjusts automatically to the number of wavelengths in the selected image.


6. Click **Do It** to equalize the time points and then click **Done**.

Calibrating Image Data

Use the following instructions to calibrate an unprocessed image when you have a calibration file and a bad pixel file that applies to the camera and conditions (array size, pixel size, and wavelength) used to collect this image. If you do not have a calibration file, you must create one before you calibrate the image.

To calibrate an image:

1. Choose **Process | Calibrate** from the main *softWoRx* menu to open the Calibrate dialog box.

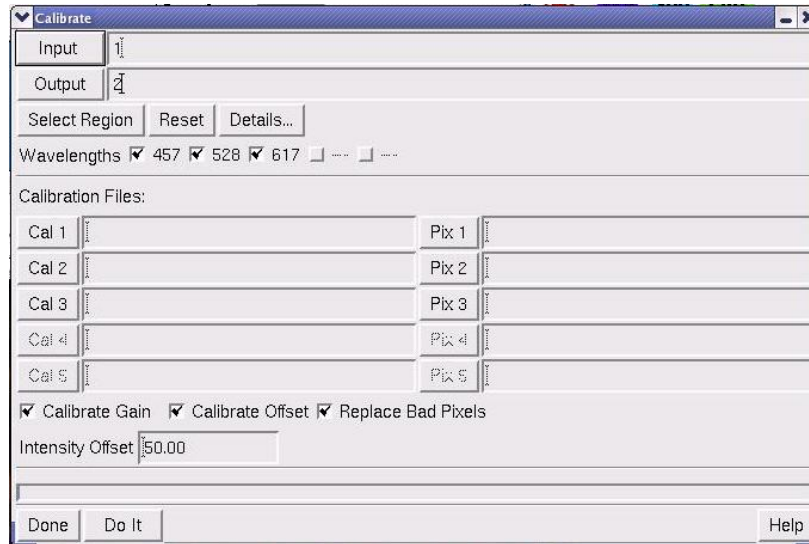


2. Click **Input** and browse to the file that you want to calibrate.
3. To select a region of the file, click **Details** and use the Region Details dialog box to select the region.
4. Click the **Cal** button for each channel and browse to the calibration file to use for this image.
5. Click the **Pix** button for each channel and browse to the bad pixel file to use for this image.
6. Select the types of calibration to perform (**Calibrate Gain**, **Calibrate Offset**, or **Replace Bad Pixels**).
7. Click **Do It** to calibrate the image.

Aligning Adjacent Images

Use Align Image to correct motion artifacts, problems with Z sectioning, or problems with time series. This tool allows you to align adjacent images by applying an XY shift with an optional rotation. Use Align Image only for images that have a single wavelength.


To align images:

1. Choose **Process** | **Align Image** on the *softWoRx* main menu to open the Align Image dialog box.

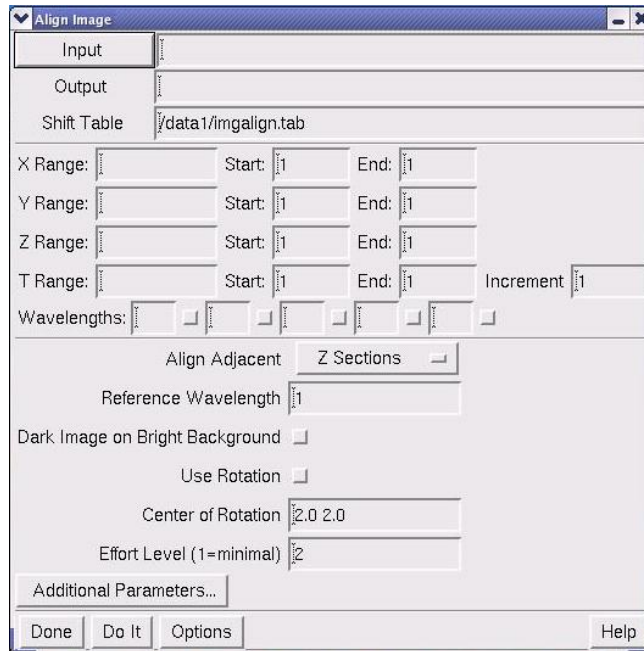

2. Click **Input** and browse to the appropriate file. The range fields and wavelengths are filled in automatically when a file is selected.

Note This dialog box requires an image name for processing. You cannot provide a window number for this field.

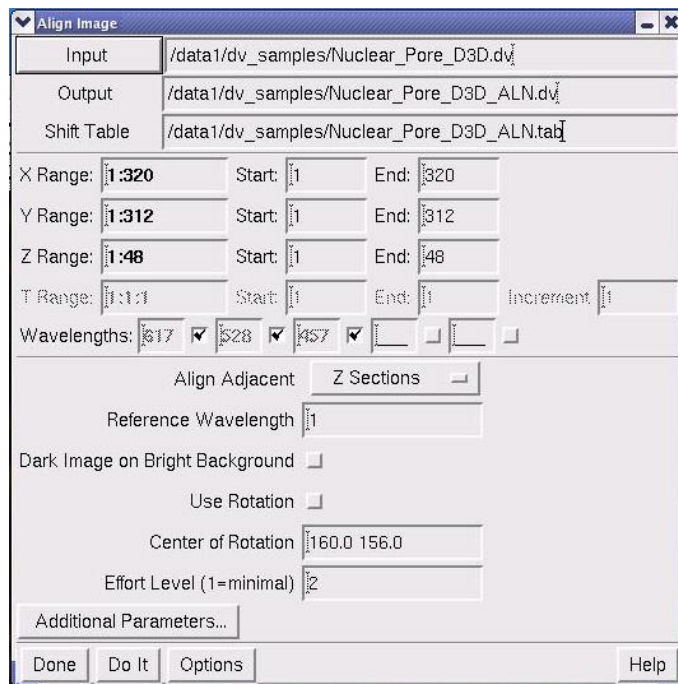


3. In the **Align Adjacent** list, select whether to align adjacent **Z Sections**, **Wavelengths**, or **Time Points**.
4. Click **Do It** to align the image.

Correcting Chromatic Aberration

Use the Chromatic Aberration Corrector to adjust channels relative to each other. This tool allows you to correct for shifts in color that result from oil matching and other environmental and optical conditions.

To correct Chromatic aberration:

1. Choose **Measure | Chromatic Correction** from the *softWoRx* main menu to open the Chromatic Aberration Corrector.

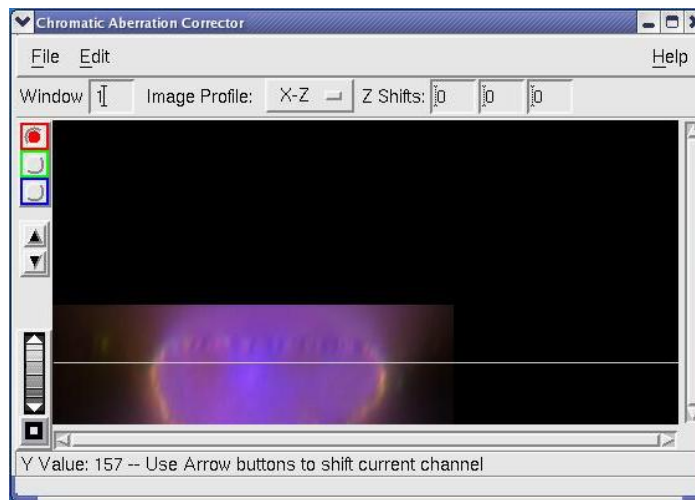


2. Select a multi-channel Image window to reference.
3. In the **Image Profile:** field, choose **X-Z** or **Y-Z** as the vertical profile to inspect.
4. Drag the yellow line in the Image window to adjust the X or Y position of the profile.
5. In the Chromatic Aberration Corrector, use the colored toggle buttons on the left to specify which channel you wish to adjust, relative to the others.
6. Click the up and down arrow buttons on the Chromatic Aberration Corrector to move the selected channel up or down relative to the others.
7. Select **File | Save Image with Corrections**.

Note For correcting chromatic aberration introduced by the optics, it is recommended that you make the measurement using a multi-colored bead so as not to bias the data. Measure the offset using the bead, then apply the same corrections to actual sample files.

3. Stitching

You can stitch images together to display images that are larger than a single field of view. This is especially useful when you want to collect data at a high magnification over a large area. You can also use it to display a sequence of time points in a time-lapse image.

If you are using a *DeltaVision* Acquisition workstation, you can create stitched images that are organized as either a series of time points or Z sections. Each time point, or Z section, is treated as a panel of the stitched image.


 **Note** Image stitching is only possible with certain types of *DeltaVision* image files. In particular, the image file must contain a series of images, along with a corresponding set of XY coordinates. To obtain images suitable for stitching, use the Panel Collection feature of Resolve3D.

In this Chapter

Stitching Images That Have a Single Z Section.....	26
Stitching Images That Have Multiple Z Sections	27

Stitching Images That Have a Single Z Section

You can use stitching for simple 2D images or for time-lapse images.

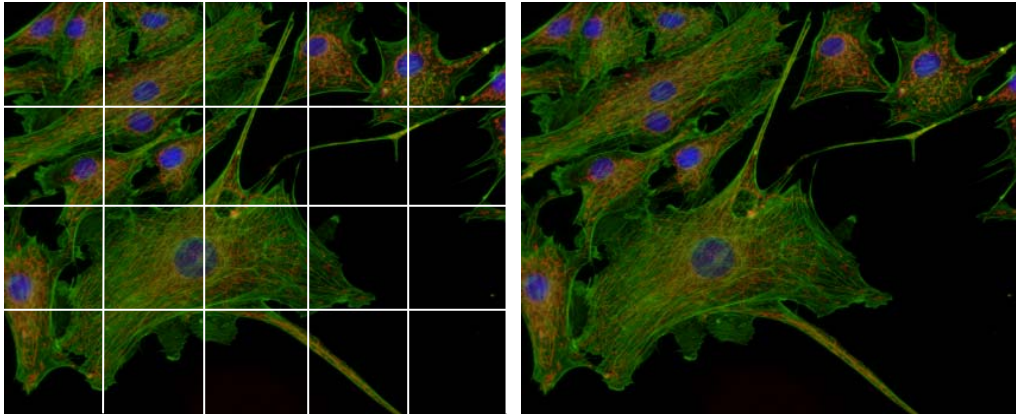


Individual panels were stitched to create the final image on the right

Before you start

Collect Panel data with the *DeltaVision* Acquisition workstation.


To stitch an image:

1. Click **View | Stitch** in the *softWoRx* main menu. The Stitch Image dialog box is displayed.

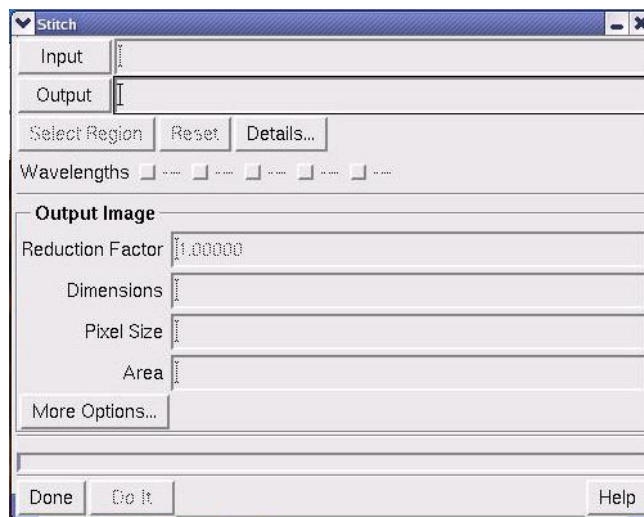

2. Click **Input** and browse to the file that you want to stitch. *softWoRx* reads the wavelengths from the file header and automatically selects the appropriate wavelengths. Change these settings only if your application does not require a stitched file for a particular wavelength, in which case you can de-select the appropriate wavelengths.
3. Click **Output** and browse to the same file previously selected. The file will appear in the Output field with “_STC” added to the file name. (With this

naming convention, the input source file is always associated with the output stitch file.)

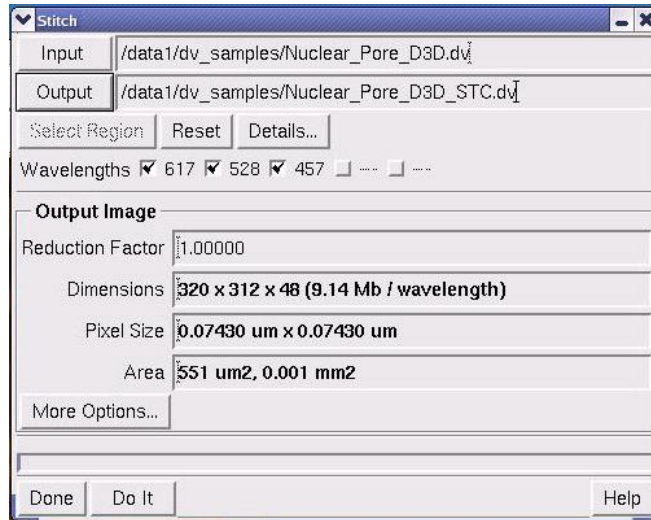

4. If desired, click **More Options** to display additional stitch options. For most applications, the default settings should work quite well. When finished with this dialog box, click **Done** to return to the main Stitch Image dialog box.
5. When all of the options have been specified, click **Do It**.

Stitching Images That Have Multiple Z Sections

For images that contain multiple Z sections, you will need to collect the images with the Panel tool and crop them before you stitch them.


Before you start

Collect Panel data with the *DeltaVision* Acquisition workstation.


Determine the width of the border rolloff (in voxels) for the images. (To minimize edge effects, the border rolloff is automatically set to about 1.5% of the image dimensions.)

To crop a multilayered image:

1. Collect 3D panel images.
2. Deconvolve the `_R3D.dv` file that you collected.
3. Open the deconvolved file in an Image window (this file has a `_R3D_D3D.dv` file extension).
4. Choose **Edit | Copy Region** on the main *softWorRx* menu. The Copy Region dialog box is displayed.

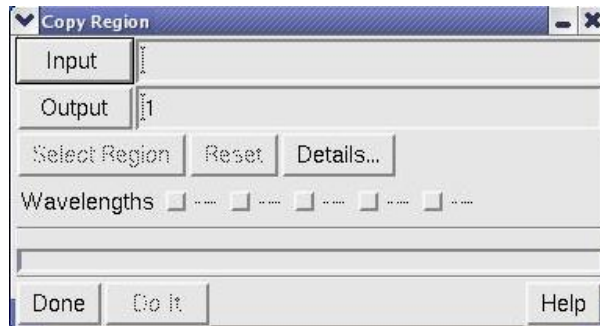

5. Drag the deconvolved file into the **Input** field and click **Details**. The Region Details dialog box is displayed.

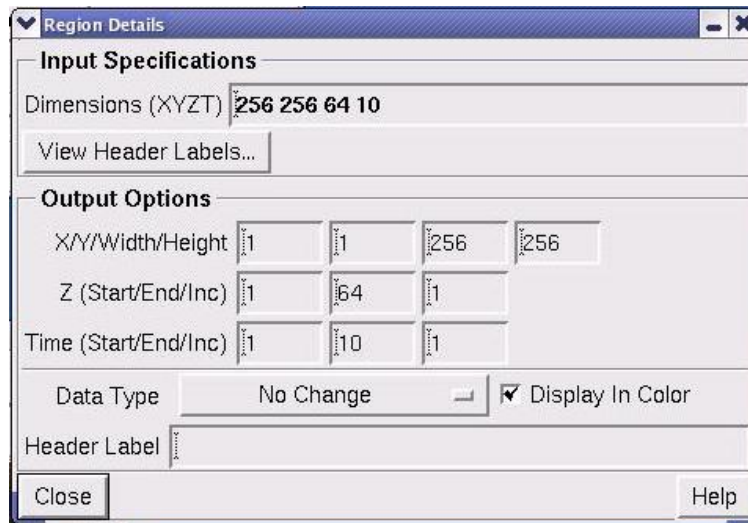

6. Enter values in the **X/Y/Width/Height**, **Z (Start, End, Inc)**, and **Time (Start/End/Inc)** fields under the **Output Options** section. (The **Dimensions (XYZT)** field displays the dimensions of the panel.) Use the following equation to determine the **Width** and **Height** values under **Output Options**:

$$\text{Width} = x - 2n$$

$$\text{Height} = y - 2n$$

Where x and y are dimensions x and y respectively and n is the number of border rolloff voxels.


7. Click **Close** to close the Region Details dialog box. Then click **Do It** in the Copy Region dialog box. (The cropped panel stack is displayed in a new Image window.)
8. Choose **File | Save** on the Image window menu to save the new cropped panel stack.

9. Stitch the cropped deconvolved file as shown in *Stitching Images That Have a Single Z Section* on Page 27.

4. Importing Data

You can convert the following image file formats to the *DeltaVision* file format.

- TIFF
- Inovision ISee™
- BioRad MRC-600 Pic
- UIC MetaMorph STK

 **Note** Since Applied Precision does not own the STK, PIC, ISEE, or TIFF formats, changes to those formats may occur that could make them incompatible with *softWoRx*. If this occurs, try saving the file as an older version of the format.

In This Chapter

Converting TIFF Images.....	32
Converting ISee Files	33
Converting Pic Files	35
Converting STK Files	36

Converting TIFF Images


You can convert a TIFF image file or series of TIFF files to a *DeltaVision* file format. When converting a series of TIFF files, the files are converted to a Z section stack.

DeltaVision expects each TIFF file to be 16 bits of grayscale data representing a single wavelength. If you have a multiple wavelength data set, you will need to create single-wavelength *DeltaVision* files and merge them using the Copy Region or Image Fusion tools.

When converting TIFF files to the DV format, you will need to provide pixel dimensions and a wavelength value for the output file. Other information may be added or modified using the Edit Image Header utility. You can also use this utility to reorganize the description of the data as a series of time points instead of Z sections.

To convert a TIFF image file to a DV image file:

1. Click **Conversions | Import from TIFF** in the *softWoRx* main menu. The Convert TIFF to DeltaVision dialog box is displayed.

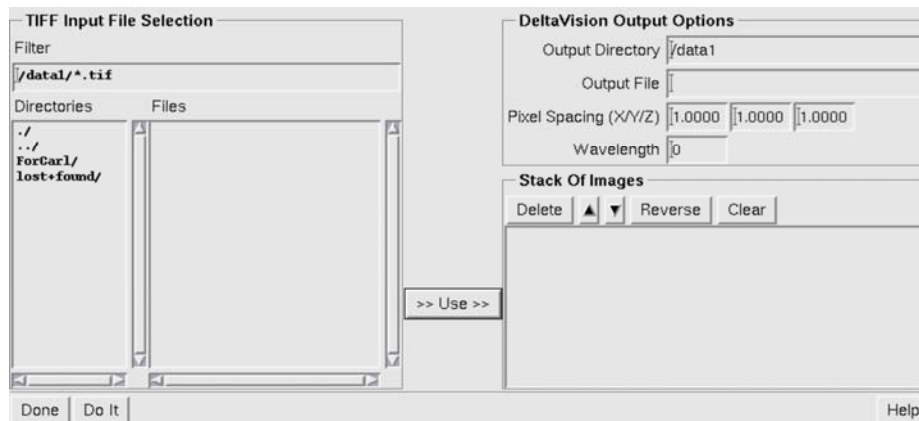

2. Select the file that you want to convert using the **TIFF Input File Selection** options.
3. Type the name of the directory in which to place the TIFF file in the **Output Directory** field.
4. Type a filename for the converted file into the **Output File** field.
5. Enter the X, Y, and Z spacing of the pixels in the data set (in microns) into the **Pixel Spacing (X/Y/Z)** fields.

Note The deconvolution program and other *softWoRx* software rely upon the presence of accurate wavelength and pixel spacing. *Not all TIFF files contain*

accurate pixel size and wavelength information, so it may be necessary to manually enter some of the fields in TIFF conversion.

6. Enter the wavelength of the data to be converted (in nanometers) into the **Wavelength** field.
7. Use the **Stack of Images** list and buttons to position the images in the desired order.
8. Click **Do It**.

Converting ISee Files


The ISee conversion tool is used to convert Inovision ISee™ images to the *DeltaVision* format.

The deconvolution program and other *softWoRx* software rely upon the presence of accurate wavelength and pixel spacing. Not all ISee files contain accurate pixel size and wavelength information. It may be necessary to manually enter values in the ISee Conversion fields.

Note The image wavelength for a *DeltaVision* file indicates the wavelength of the light imaged by the camera, rather than the illumination wavelength.

X/Y pixel spacing can be obtained in two ways: it can be measured with a test target or it can be approximated from the CCD detector element size and the total image magnification. For example, if the CCD detector has 6.7 μm pixels and the image was acquired with a 100X lens and a 1.5X optivar, then the pixel size is approximately $6.7 \mu\text{m}/(100 \times 1.5) = 0.045 \mu\text{m}$. The Z pixel spacing is the distance between adjacent optical sections.

Figure 1. ISee Converter

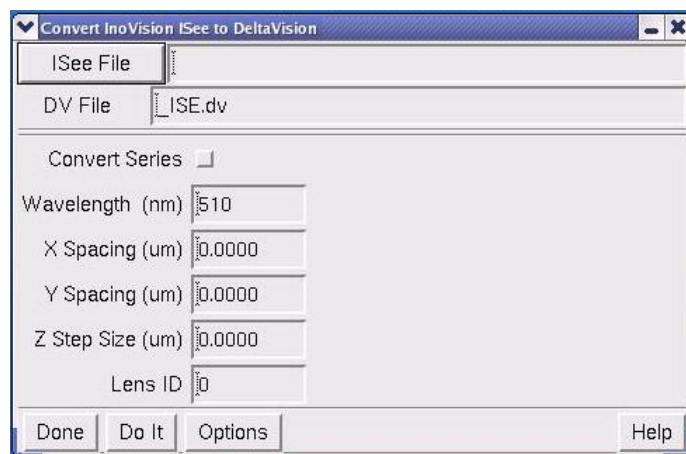

Options in this dialog box are described briefly in the following paragraphs. For additional information regarding ISee file conversion, refer to your online Help system.

ISee Series Conversion

The **Convert Series** option combines all files with a similar name into one *DeltaVision* file. The conversion program looks for files that have the same prefix as the input file. The program assumes that a series of Inovision files will be the same, except for the sequence numbers in the last 3 characters.

For example, the following files would be automatically combined into a single *DeltaVision* file:

my_file_name.001

my_file_name.002

my_file_name.003

my_file_name.004

Any of the above files could be entered into the **ISee File** option as the file to convert. The use of wildcards, such as "*", is not supported.

To convert an ISee image file to a DV image file:

1. Click **Conversions | Import from ISee** in the *softWoRx* main menu. The convert Inovision ISee to DeltaVision dialog box is displayed, as shown in Figure 1.
2. Select the file that you want to convert using the **ISee File** button and data-entry field.
3. Type a filename for the converted file into the **DV File** field.
4. Enter the wavelength of the data to be converted (usually in nanometers) into the **Wavelength** field.
5. Enter the X, Y, and Z spacing of the pixels in the data set (usually in microns) into the appropriate fields.
6. Enter the correct lens number in the **Lens ID** field.
7. Click **Do It**.

Converting Pic Files


The Pic conversion tool is used to convert BioRad MRC-600 Pic™ images to *DeltaVision* format.

The deconvolution program and other *softWoRx* software rely upon the presence of accurate wavelength and pixel spacing. Not all Pic files contain accurate pixel size and wavelength information, so it may be necessary to manually enter values in the Pic Conversion fields.

Note The image wavelength for a *DeltaVision* file indicates the wavelength of the light imaged by the camera, rather than the illumination wavelength.

X/Y pixel spacing can be obtained in two ways: it can be measured with a test target or it can be approximated from the CCD detector element size and the total image magnification. For example, if the CCD detector has 6.7 μm pixels and the image was acquired with a 100X lens and a 1.5X optivar, then the pixel size is approximately $6.7 \mu\text{m}/(100 \times 1.5) = 0.045 \mu\text{m}$. The Z pixel spacing is the distance between adjacent optical sections.

Figure 2

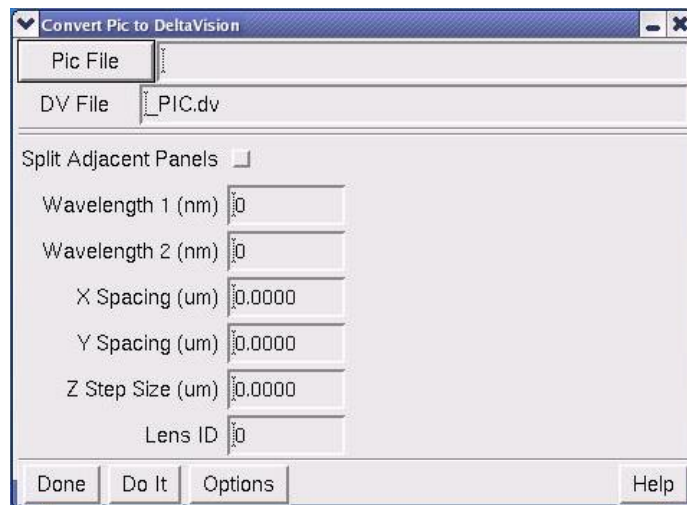

Options in this dialog box are described briefly in the following paragraphs. For additional information regarding Pic file conversion, refer to your online Help system.

To convert a Pic image file to a DV image file:

1. Click **Conversions | Import from Pic** in the *softWoRx* main menu. The Convert Pic to DeltaVision dialog box is displayed, as shown in Figure 2.

2. Select the file that you want to convert using the **Pic File** button and data-entry field.
3. Type a filename for the converted file into the **DV File** field.
4. If the image file consists of two wavelengths that are arranged in two adjacent panels, enable the **Split Adjacent Panels** option.
5. Enter the wavelength of the first image (or panel) into the **Wavelength 1 (nm)** field.
6. If necessary, enter the wavelength of the second image (or panel) into the **Wavelength 2 (nm)** field.
7. Enter the X, Y, and Z spacing of the pixels in the data set (usually in microns) into the appropriate fields.
8. Enter the correct lens number in the **Lens ID** field.
9. Click **Do It**.

Converting STK Files


The STK conversion tool is used to convert MetaMorph STK images to *DeltaVision* format. Unlike the ISee and Pic converters, the STK converter attempts to read the wavelength and pixel size values from the STK file's header immediately after you specify the name of the input file. After reading these values from the header, the converter enters this data into the fields in STK Conversion. If necessary, you may change this information manually before you click **Do It**. (The ISee and Pic converters do not read the input file until you click **Do It**.)

The deconvolution program and other *softWoRx* software rely upon the presence of accurate wavelength and pixel spacing. As with ISee and Pic files, not all STK files contain accurate pixel size and wavelength information, and it may be necessary to manually enter some of the fields in STK Conversion.

Note The image wavelength for a *DeltaVision* file indicates the wavelength of the light imaged by the camera, rather than the illumination wavelength.

X/Y pixel spacing can be obtained in two ways: it can be measured with a test target or it can be approximated from the CCD detector element size and the total image magnification. For example, if the CCD detector has 6.7 μm pixels and the image was acquired with a 100X lens and a 1.5X optivar, then the pixel size is approximately $6.7 \mu\text{m}/(100 \times 1.5) = 0.045 \mu\text{m}$. The Z pixel spacing is the distance between adjacent optical sections.

Figure 3

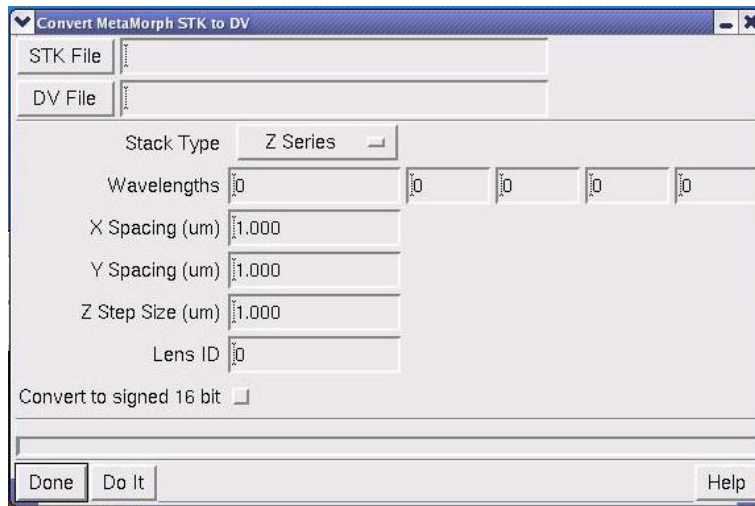

Options in this dialog box are described briefly in the following paragraphs. For additional information regarding STK file conversion, refer to your online Help system.

To convert an STK image file to a DV image file:

1. Click **Conversions | Import from MetaMorph STK** in the *softWoRx* main menu. The Convert MetaMorph STK to DV dialog box is displayed, as shown in Figure 3.
2. Select the file that you want to convert using the **STK File** button and data-entry field.
3. Type a filename for the converted file into the **DV File** field.
4. Enter the wavelengths (in nm) of the light collected by the camera for each channel into the **Wavelengths** fields.
5. If necessary, modify the X, Y, and Z spacing of the pixels in the data.
6. Enter the correct lens number in the **Lens ID** field.
7. Click **Do It**.
8. Proceed to enter the appropriate values into the rest of the dialog box and complete the desired manipulation.

5. Data and Task Manipulation

This chapter describes how to select data, crop data from images, and combine images.

In this Chapter

Selecting Data	40
Cropping and Trimming Data	45
Cropping an Irregular Data Region	47
Combining Data of Two Images.....	52
Setting Up Process Chains with Task Builder	54
Using Ratio Imaging.....	57
Using the Multiplexed Wavelength Option.....	62
Connecting to a DMS Database	74
Uploading Images to a DMS Database.....	74
Downloading Files from DMS	81
Browsing and Locating Images in a DMS Database.....	82

Selecting Data

You can select either rectangular or irregular data regions:


- Use the **Select Region** and **Details** buttons to select rectangular data regions. These buttons are included at the top of *softWoRx* dialog boxes that allow you to save, export, or select data.
- Use the Edit Polygon tool to select irregular data regions.

Selecting Rectangular Data Regions


Use **Select Region** to select data for volume rendering, Rotate3D tool applications, modeling, and other applications that require intensive processing. You can also use this tool to crop data and save it in a new file (see *Cropping a Rectangular Region* on Page 45).

To select a rectangular data region:


1. On any *softWoRx* process window that includes the **Select Region** button, click **Select Region** and drag the mouse across the image to select a rectangular area. The selected region is indicated by a dotted line.

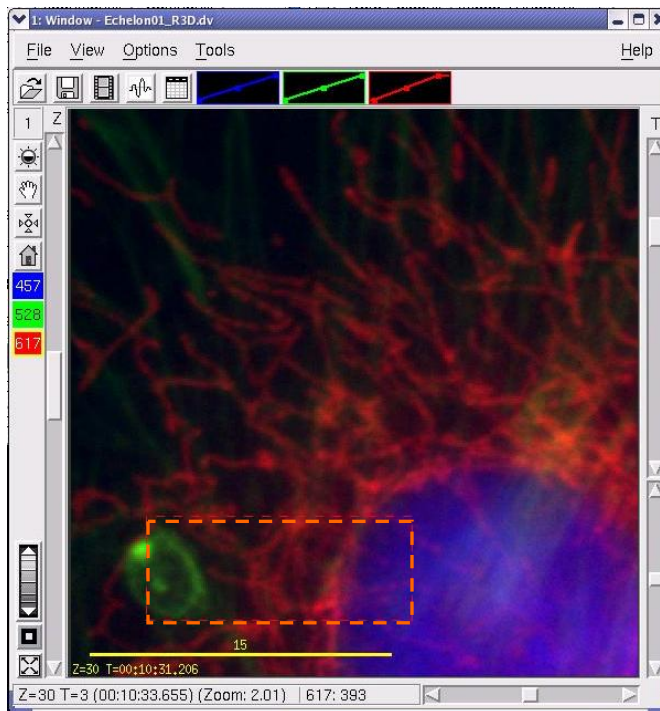

2. Click **Details** to open the Region Details dialog box.

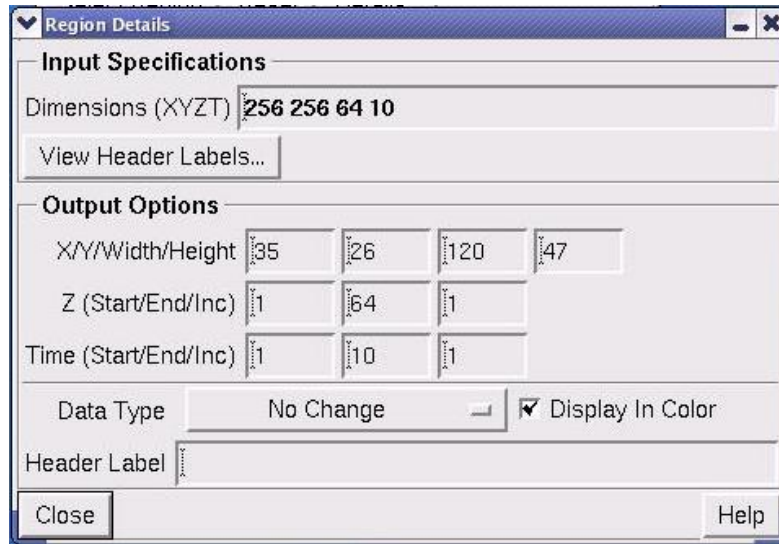

3. In the **Z (Start/End/Inc)** field, select a Z section range to include. **Start** and **End** are the beginning and end points. **Inc** (incremental) allows you to skip points (e.g., entering an **Inc** value of 2 skips every other point).
4. In the **Time(Start/End/Inc)** fields, select a time data range to include. Then click **Close**.
5. Select **Do It** on the process window you're using. The selected region is displayed in the chosen output window.

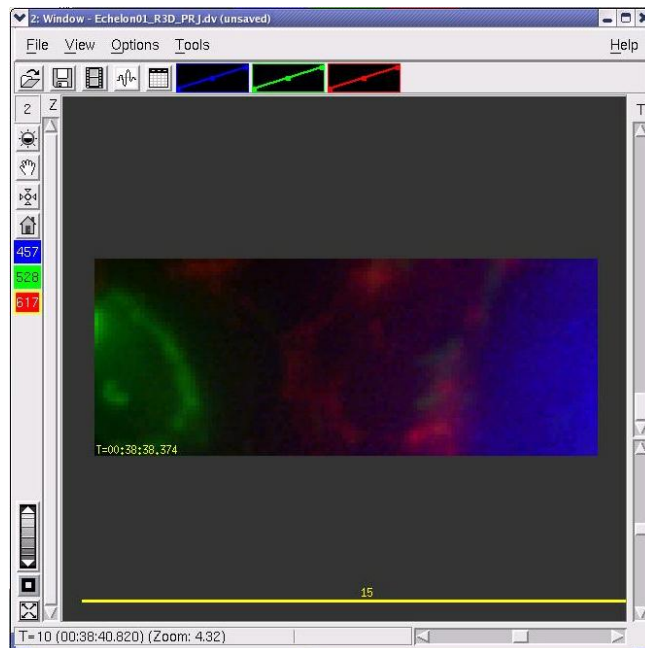


Selecting Irregular Data Regions


Use the Edit Polygon tool to select irregular data regions. This tool is applied differently than the tools for selecting rectangular regions. While selecting rectangular regions is usually used within other tools, the Polygon Editor stands alone. You must use the Polygon Editor tool with Cut Mask if you want to then apply another tool only to the selected region.


To select an irregular data region:


1. Open the image in the Image window.

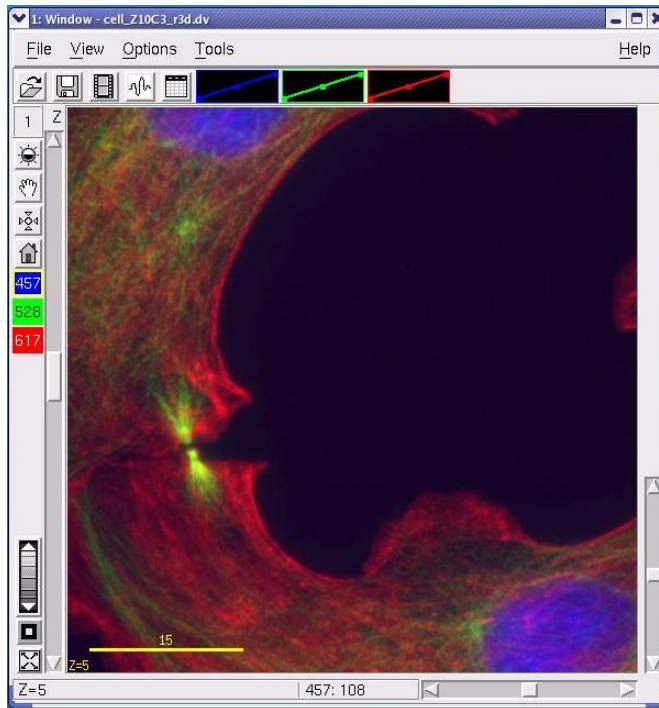


2. Choose **Model | Edit Polygon** to open the Edit Polygon menu.

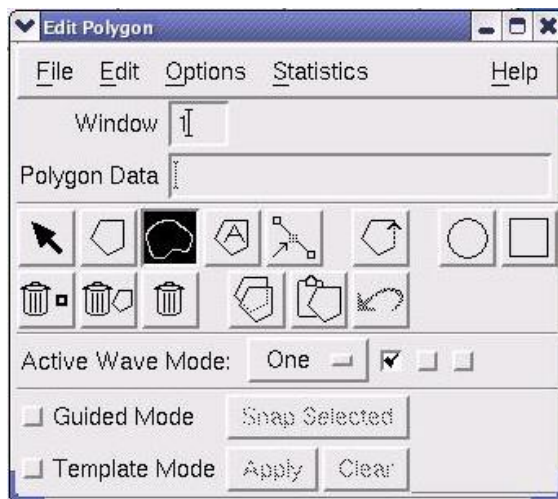


- Choose a selection tool (e.g., ) from the Edit Polygon menu. Then press and hold the left mouse button to draw a polygon around the area of interest within the Image window.

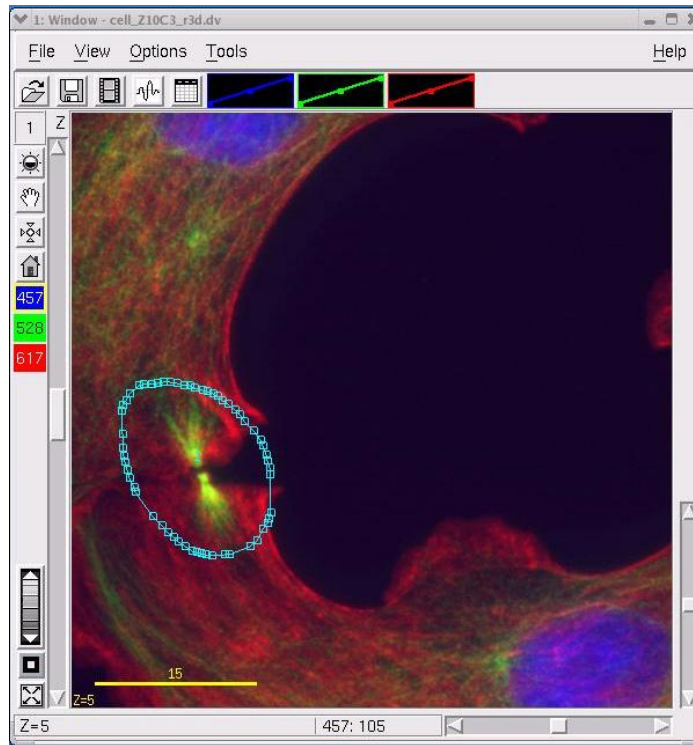


- To copy this region across wavelengths, time points, or through Z sections, choose **Edit | Propagate Polygons** from the Edit Polygon menu and enter the appropriate ranges.

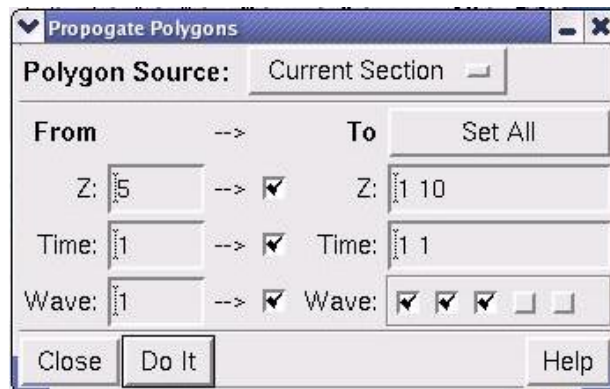



 **Tip** you can also select time points. You can select ranges or you can enter selected points (e.g., 1, 3-5, 20-25).

- Click **Set All** if you want to copy the polygons through all of the Z sections, time points, and wavelengths. Click **Do It** to copy the polygons. Then view the range of Z sections or time points to make sure that all of the data is included

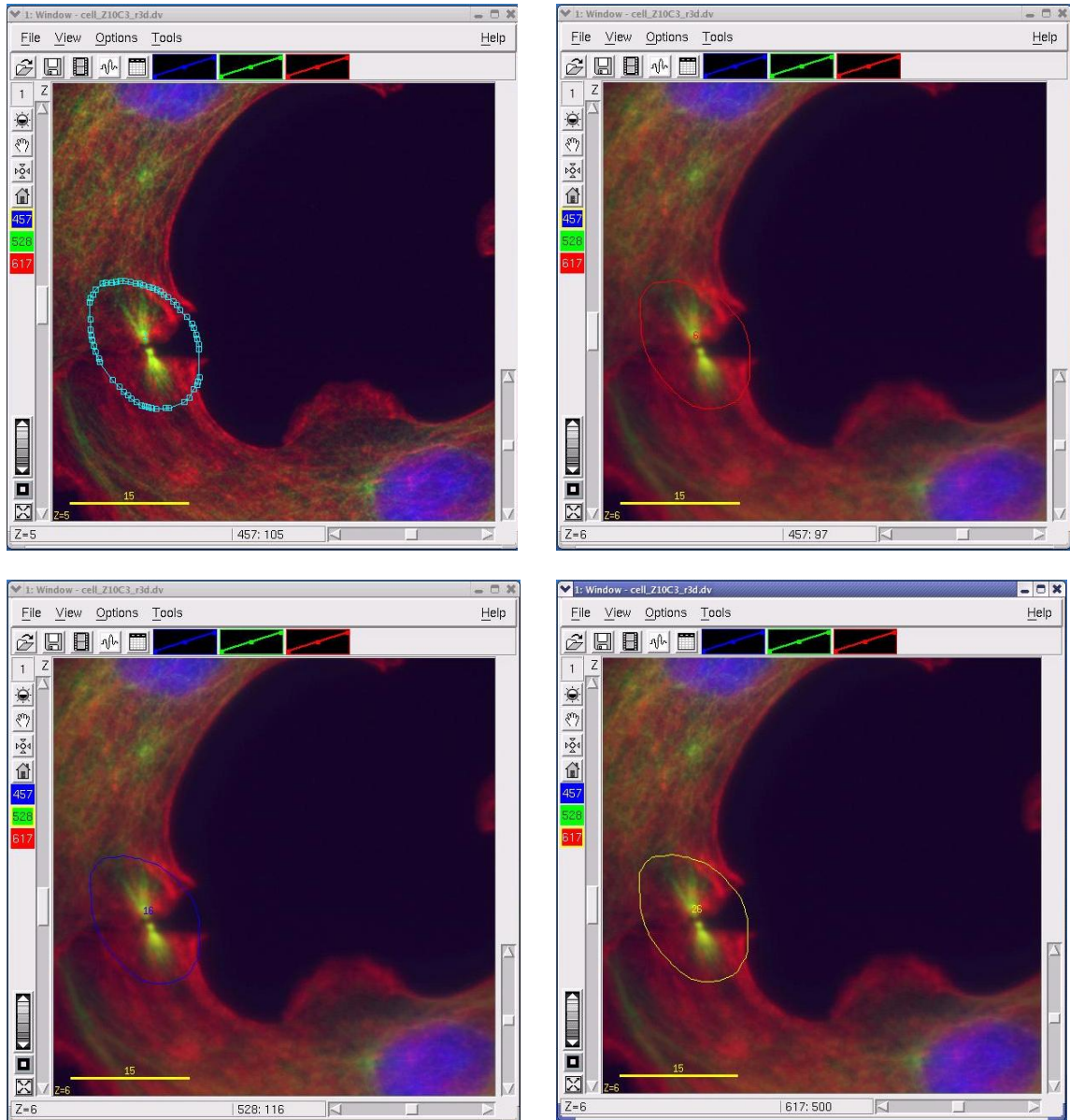in the polygon for each section. Use the channel selectors to view the polygons for each of the selected wavelengths.



In this example, all of the data for each of the selected wavelengths and between the selected cross sections is within the polygons for each Z section.

Tip You can change selected points on a polygon using the  button. You can move a selected polygon by selecting it with the  tool and dragging it to a new location.


Cropping and Trimming Data

You can selectively crop areas, Z sections, and channels from data files. You can also trim time points from time-lapse data. Cropping and trimming are useful for presenting data. It also helps prepare data for volume rendering, 3D rotation, modeling, and other types of visualization.


Cropping a Rectangular Region

To crop a rectangular region to a new file:


1. Open an Image window. From the Image window menu, choose **File | Save**.

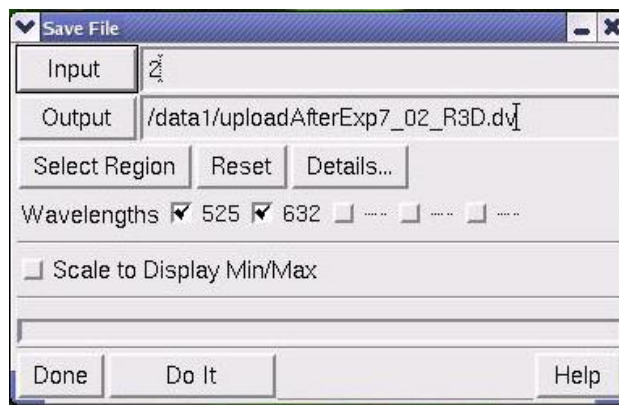

2. In the **Input** field, enter either a window number or an image file name. In the **Output** field, specify an image file or window as output.
3. If your input is a window, you can select a region. To do this, click **Select Region** and drag the mouse across the area to select it. Adjust the rectangle that you've created until it contains the desired area. Then click outside the Image window with the mouse.

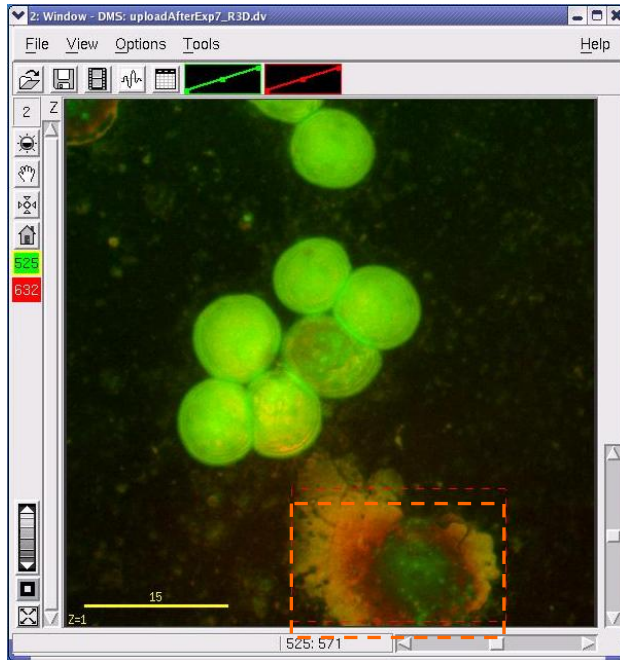


4. Click **Details** to open the Region Details dialog box.

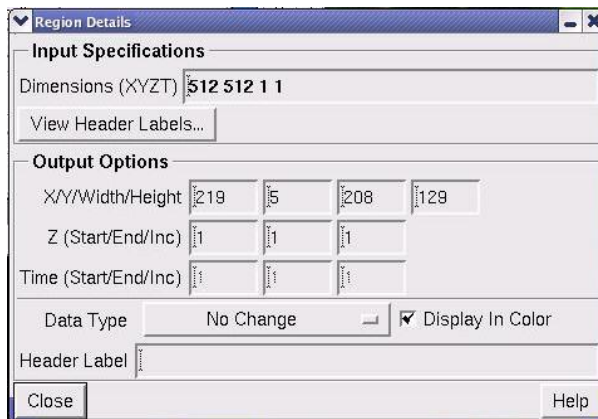

5. In the **Z (Start/End/Inc)** field, select a Z section range to include. **Start** and **End** are the beginning and end points. **Inc** allows you to skip points (e.g., entering an **Inc** value of 2 skips every other point).
6. In the **Time(Start/End/Inc)** fields, select a time data range to include. Then click **Close**.
7. In the Save File dialog box **Wavelengths** field, choose which wavelengths of the input data to process and include in the output data set. If you don't have the option of which wavelength to include, the toggle buttons are dimmed.


8. Click **Do It** to save the cropped image file. Then open the saved file in another Image window and view the results of the selections.

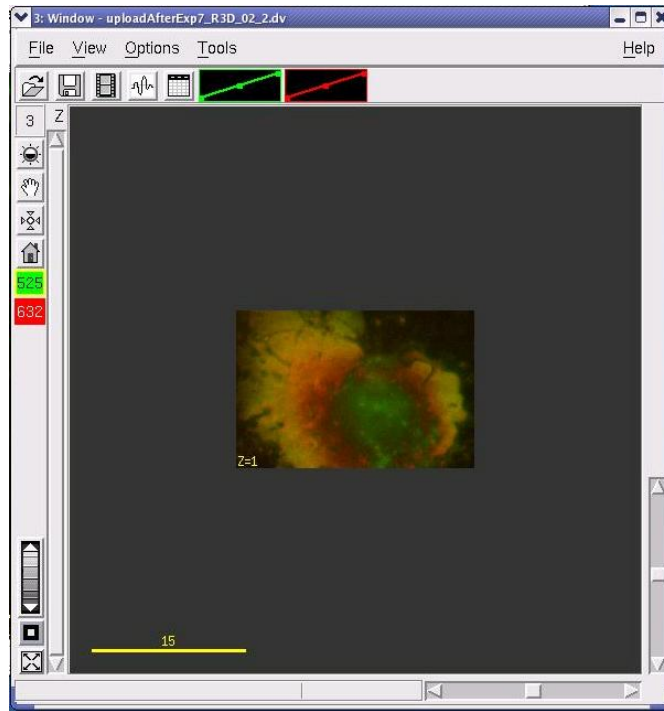


Cropping an Irregular Data Region


You can crop irregular data regions from image files.

To crop an irregular data region:

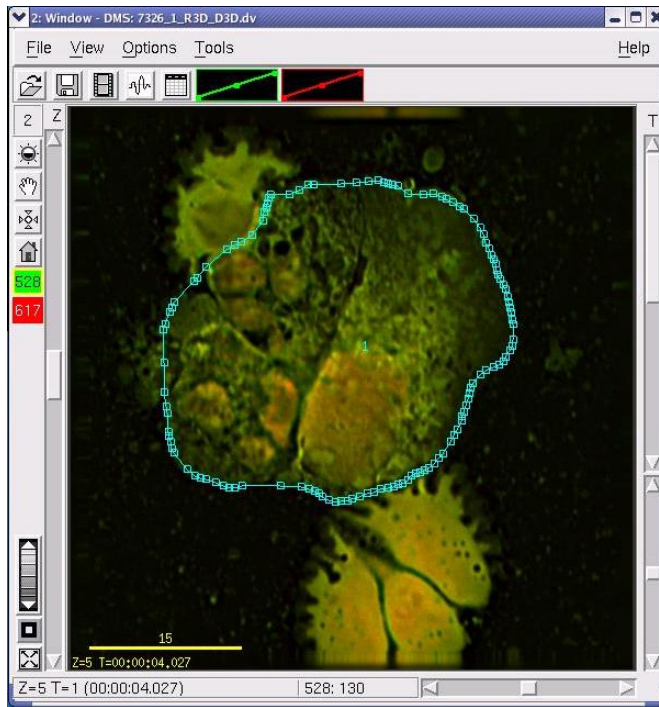


1. Open the image in the Image window.
2. Choose **Model | Edit Polygon** to open the Edit Polygon window.
3. Choose a selection tool (e.g. ) and draw a polygon around the area of interest.




- To copy this region across wavelengths, time points, or through Z sections, choose **Edit | Propagate Polygons** from the Edit Polygon menu and enter the appropriate ranges.

 **Tip** You can also select time points. You can select ranges or you can enter selected points (e.g., 1, 3-5, 20-25).


- Click **Do It** to copy the polygons. Then view the range of Z sections to make sure that all of the data is included in the polygon on each section.

 **Tip** You can change selected points on a polygon using the  button. You can move a selected polygon by selecting it with the  tool and dragging it to a new location.

- From the main *softWoRx* menu, choose **Edit | Cut Mask**.

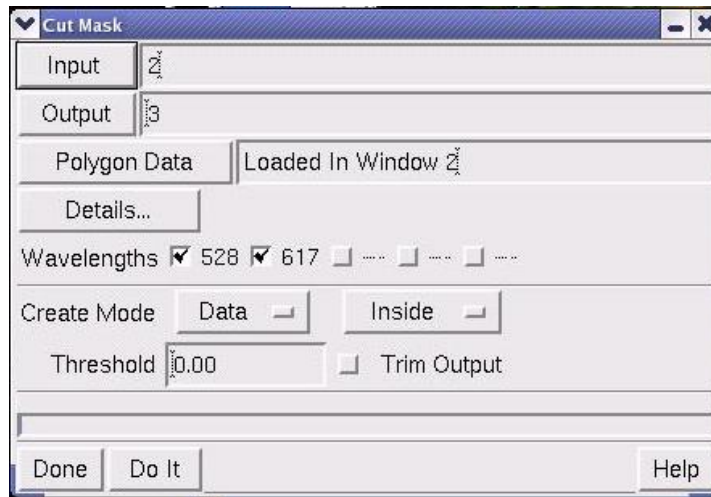


- In the **Input** field, enter the Image window number. Then click **Details** to open the Region Details dialog box and enter the Z or T sections to include. Click **Close**.

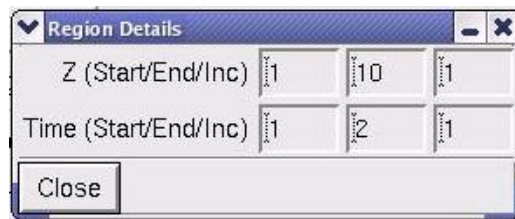

- In the Cut Mask window, choose which wavelengths to include. In **Create Mode**, choose one of the following modes and specify whether to act on the inside or the outside of the polygons.
 - Choosing **Data** cuts all of the data inside or outside of each polygon and copies it to the output destination.
 - Choosing **Mask** creates an output file with only 1's and 0's representing either the inside or the outside of the polygons.
- In the **Threshold** field, set a background intensity to remove from the selection. (For example, setting a threshold of 200 selects only data with an intensity value greater than 200.)


Note With Trim Output selected, the smallest area containing all the polygons is the area written to the output window. With Trim Output unselected, the size of the file has the same x-y dimensions as the original file, but only the part defined by polygons has intensity.

- Click **Do It** to crop the image.

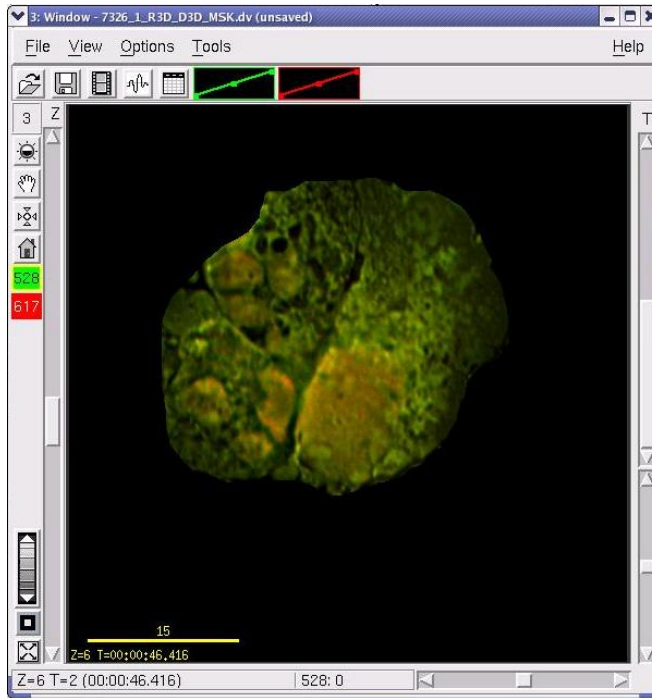

Trimming Time Data


You can trim time points from time-lapse images.

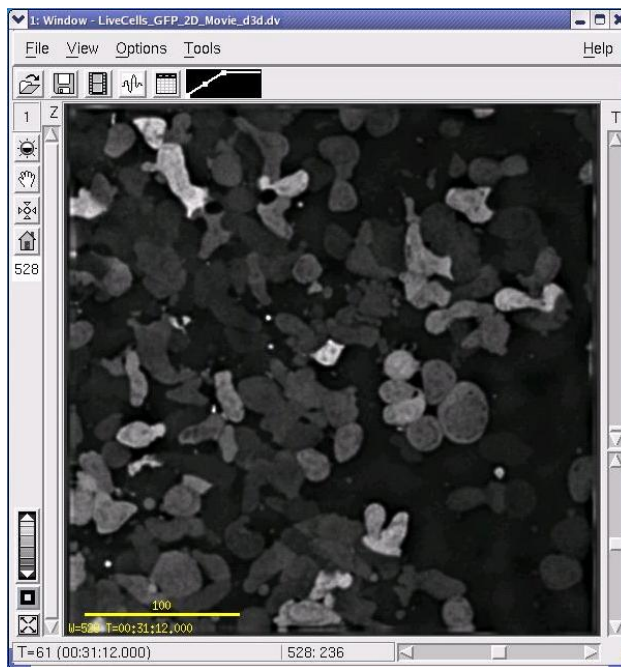


Starting image with 61 time points

To trim time points:


1. Open the image in the Image window.
2. From the main *softWoRx* menu, choose **Edit | Copy Region**.

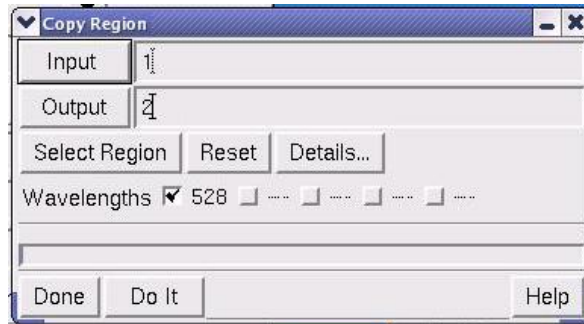

3. In the **Input** field, enter the Image window number.
4. Click **Details** to open the Region Details dialog box.

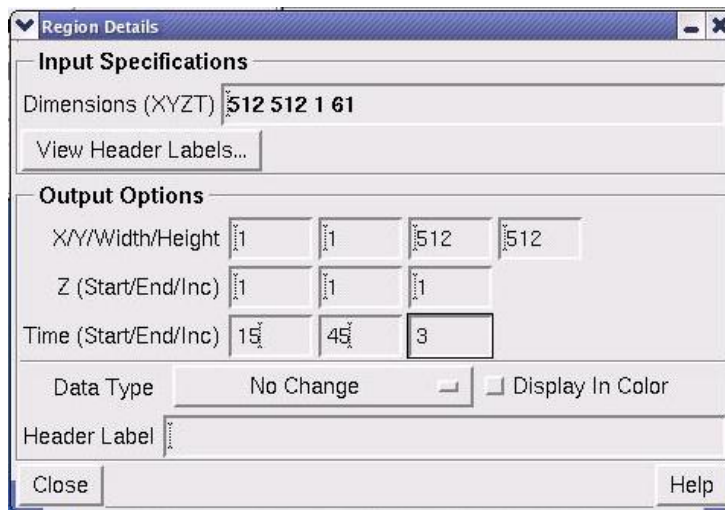

5. In the **Time (Start/End/Inc)** field, enter the first and last time points to include and the increment between points. For example, entering 15, 45, 1 includes all of the points between 15 and 45. You could skip every other point in this interval by entering 15, 45, 2. (In the example above, every third point is included.)
6. Click **Close** to quit Region Details.
7. In the Copy Region dialog box, click **Do It** to create the new image.

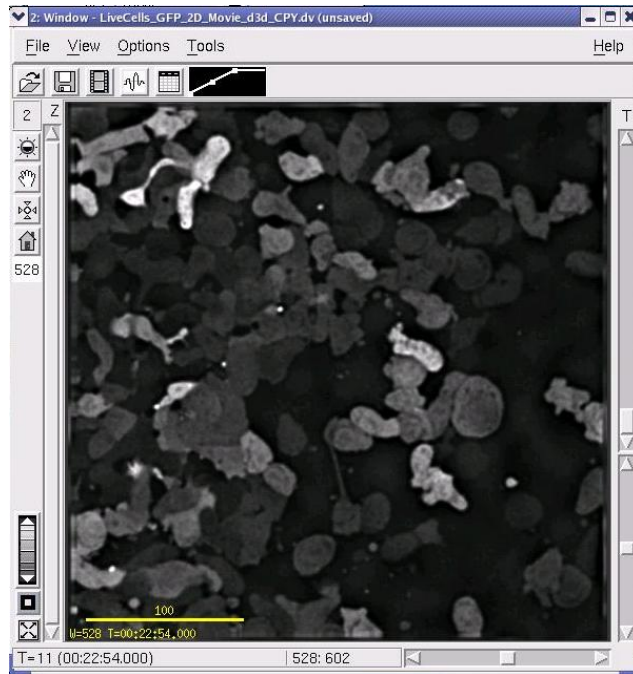

Ending image with 11 time points


Combining Data of Two Images

Use the Image Fusion dialog box to combine time points, Z sections, or wavelengths from two *DeltaVision* images into one output file. The input images may come from windows or files. After selecting the input images, you can specify exactly which wavelengths, time points, and Z sections you want to combine.

You can either append selected wavelengths to a single file or fuse time points or Z sections of the same wavelength, creating a single data set for each output wavelength.

To combine data of two image files:

1. Choose **Edit | Image Fusion** from the main *softWoRx* menu to open the Image Fusion dialog box.

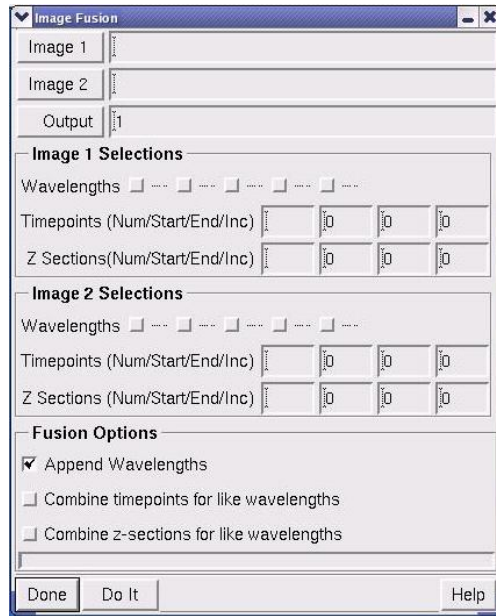

2. Enter image file names or window numbers for the two files that you want to combine in the **Image 1** and **Image 2** boxes.
3. Select which wavelengths, time points, and Z sections to combine from the first image in **Image 1 Selections**.
4. Select which wavelengths, time points, and Z sections to combine from the second image in **Image 2 Selections**.
5. Specify how to combine the data under **Fusion Options** as follows:
 - To append all selected wavelengths to the output data set, choose **Append Wavelengths**. (With this option, if Image 1 had wave 490 selected and Image 2 had wave 490 selected, the output data set would have two separate 490 wavelengths.)
 - To combine timepoints from all selected matching wavelengths into a single final wavelength data set, choose **Combine timepoints for like wavelengths**.
 - To combine Z sections from all selected matching wavelengths into a single final wavelength data set, choose **Combine z-sections for like wavelengths**.
6. Click **Do It**.

Notes

#1 If error messages are displayed referring to differences in file types that include image size, pixel size, lens info, data type, etc., use Copy Region and Edit Header to manipulate these items.


#2 If the number of sections varies from wavelength to wavelength during an operation, **Blank Z sections have been added to the image** is displayed. These blank sections are added to balance the number of sections between each wavelength of the output image. Each one is a zero intensity image added to the end of the appropriate wavelength.

Setting Up Process Chains with Task Builder


A *process chain* is a series of tasks that are predefined for a given collection of data. Task Builder is a unique feature of *softWoRx* that allows you to set up process chains for one or several tasks to be performed on a single set, or multiple sets of data. You can use the Task Builder dialog box to select files and define multiple operations to be performed. When you've finished providing the data information and the processes you want to accomplish, you can choose to either start the jobs immediately or start the jobs at a specified time.

To set up process chains with Task Builder:

1. Select **Process | Task Builder** from the main *softWoRx* menu to open the Task Builder dialog box.

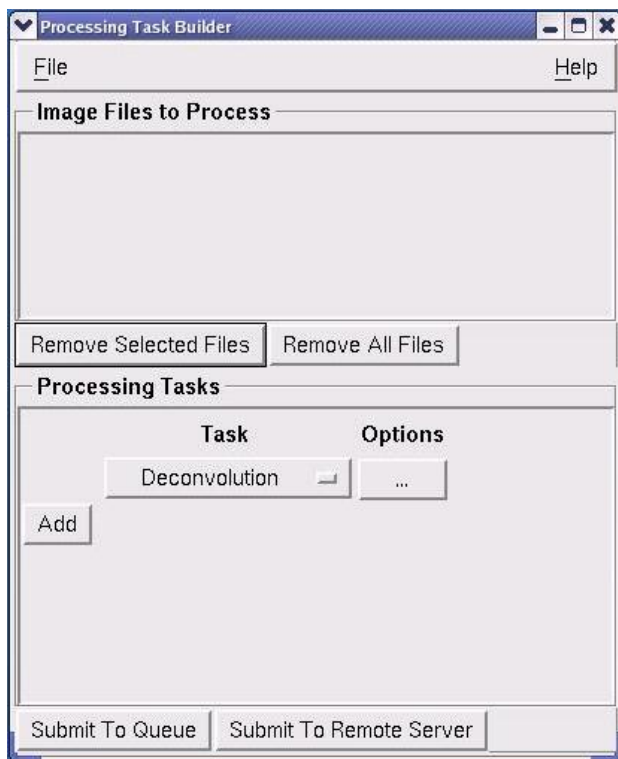


2. On the dialog box, select **File | Add Files**. You are presented with a list of files from which you can choose the specific files you want to add to your process chains.

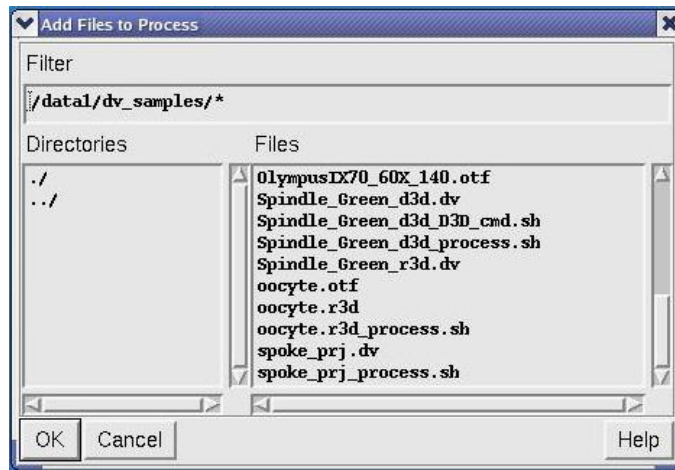

3. Select the file(s) you want to add to the Task Builder dialog box and click **OK**. The files are displayed in the Input Files section of the dialog box.


Note You can select files from this dialog box using the SHIFT key to select multiple contiguous files or the CTRL key to select multiple files from various parts of the list.

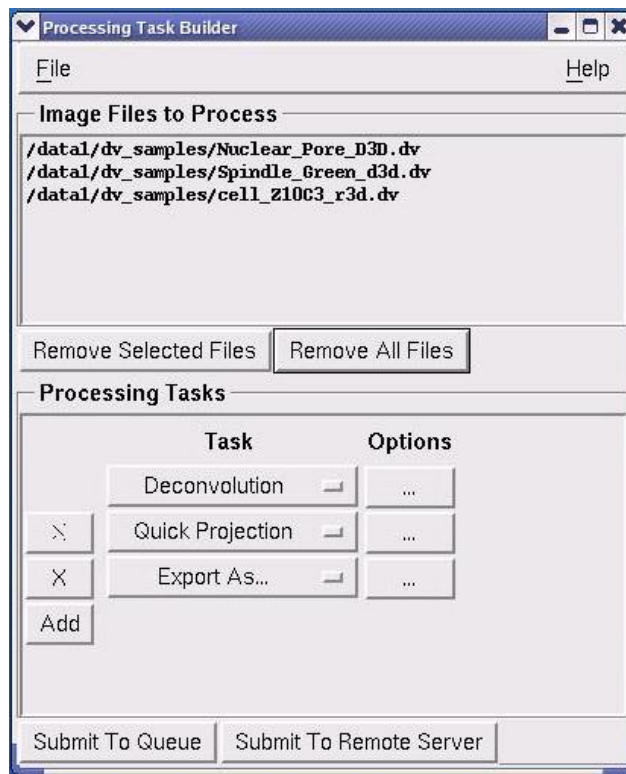


4. Next, use the **Task** options in the Processing Tasks section of the dialog box to select the tasks and the order in which you want these processes performed on the selected file(s). The Task options to choose from are **Deconvolution**, **Correction**, **Crop Image**, **Quick Projection**, **Volume Rendering**, and **Export As**. Use the **Add** button to include additional tasks and the **X** buttons to

remove tasks from the chain. Use the **Options** buttons next to each task to view a dialog box of available options for the specific task.


The task options you specify will be performed on each file in the exact order they appear in the Task Builder dialog box. Each of the selected files is run through the entire list of tasks before the Queue Manager moves on to the next file in the list.

 **Tip** You can use the left mouse button to drag image file icon, a group of file icons, or folder icons to the Task Builder or the Queue Manager.

- When you are satisfied with your selections and the order in which the tasks will occur, click **Submit Tasks to the Queue**. The *softWoRx* Queue Manager dialog box is displayed.

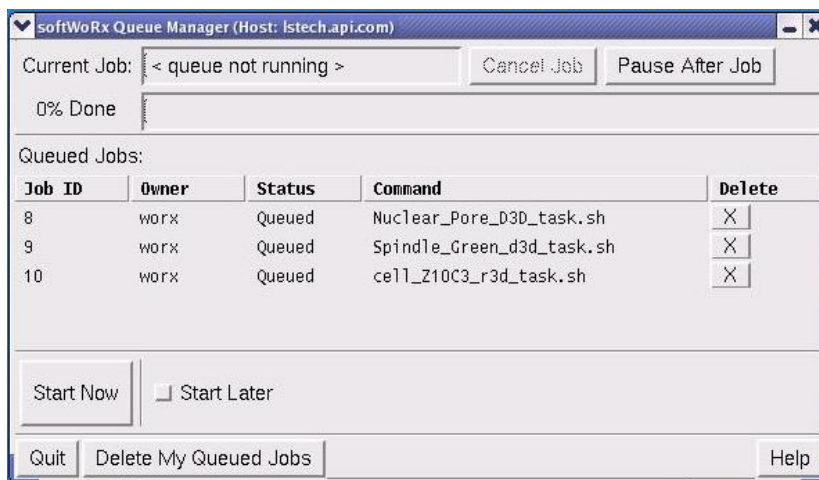


- To start the process chain,
 - Immediately*, click **Start Now**.
 - At a later time*, click **Start Later**. When you choose this option, a set of time option buttons is displayed as shown in the example below. Set the time at which you want the process chain to start. You can use the **Change It** button if you decide you want to change the time to begin the process chain.

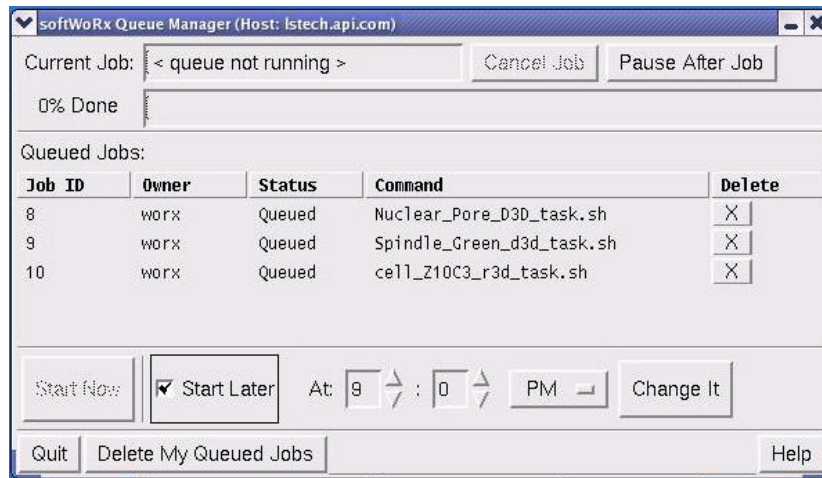

7. Select **Quit** to exit the Queue Manager dialog box.

Using Ratio Imaging


softWoRx provides a ratio imaging acquisition function that allows you to view a graphic representation of the ratio of two channels as the images are being collected. In addition, a ratio graph displays the mean value (of an area in the middle of the image) vs. time. Both the ratio image and the ratio graph are for monitoring purposes only. The ratio imaging experiment results in a two-channel time-lapse image.

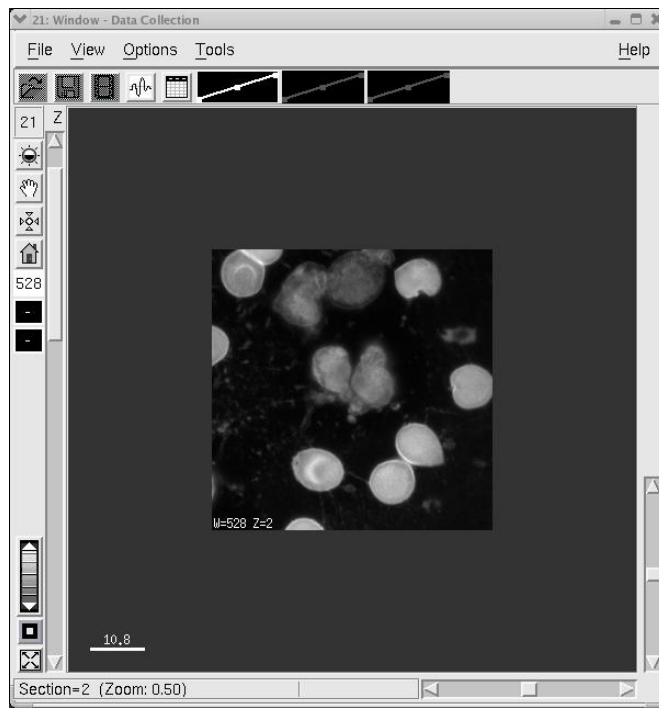

Sample image for ratio imaging experiment


To set up a ratio imaging experiment:

1. From the softWoRx main menu, select **File | Acquire (Resolve3D)** to open the Resolve3D window.

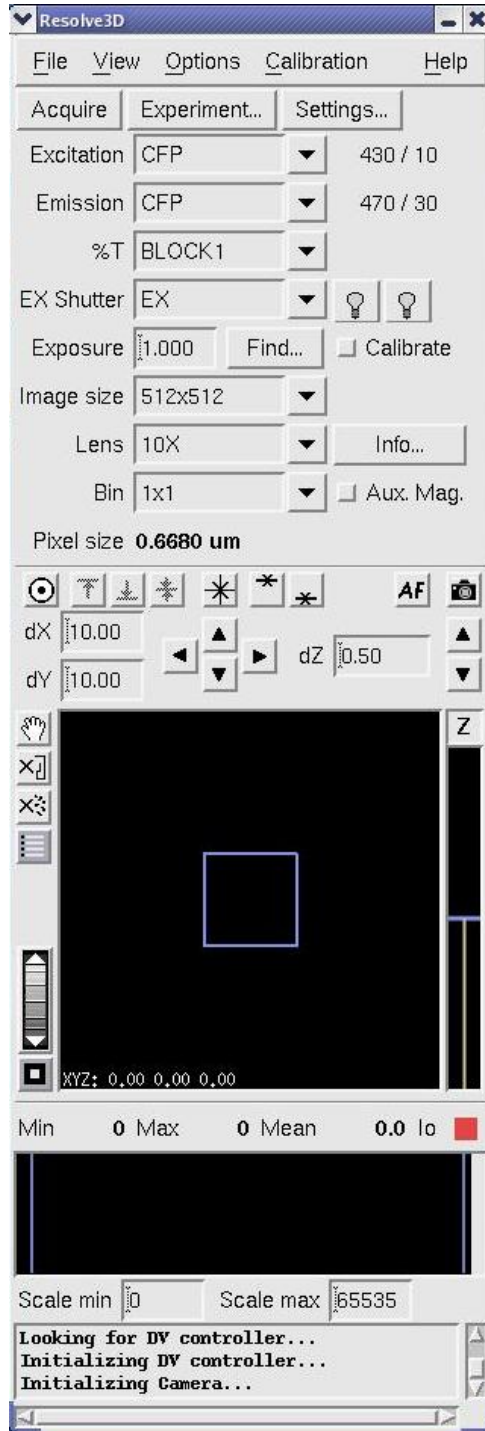


2. On the Resolve3D window, click the **Experiment** button to open the Design/Run Experiment dialog box.


3. Select the Sectioning tab and unselect the **Z Sectioning** toggle.

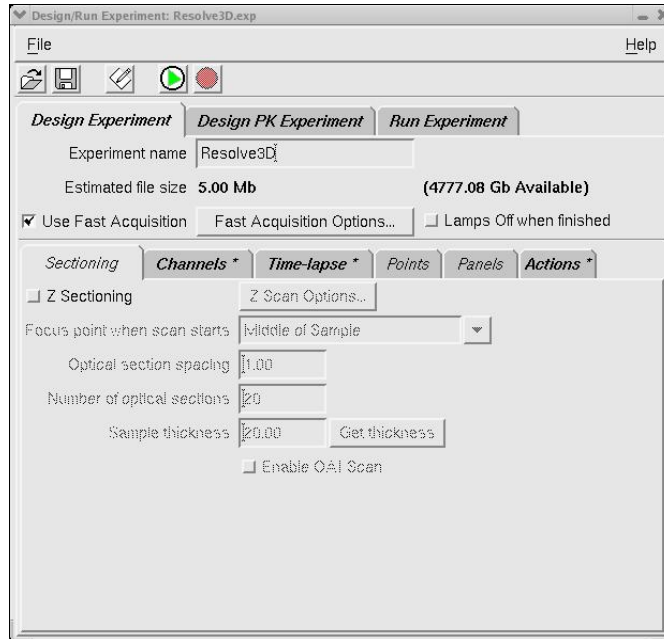


4. Select the Channels tab and specify the two channels you want to use for this experiment.

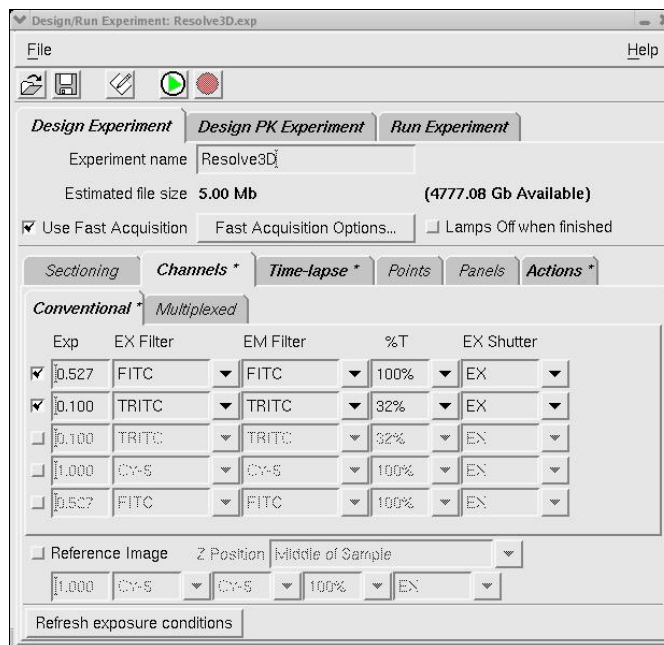


5. Select the Time-lapse tab and specify the time-lapse and total time for this experiment.

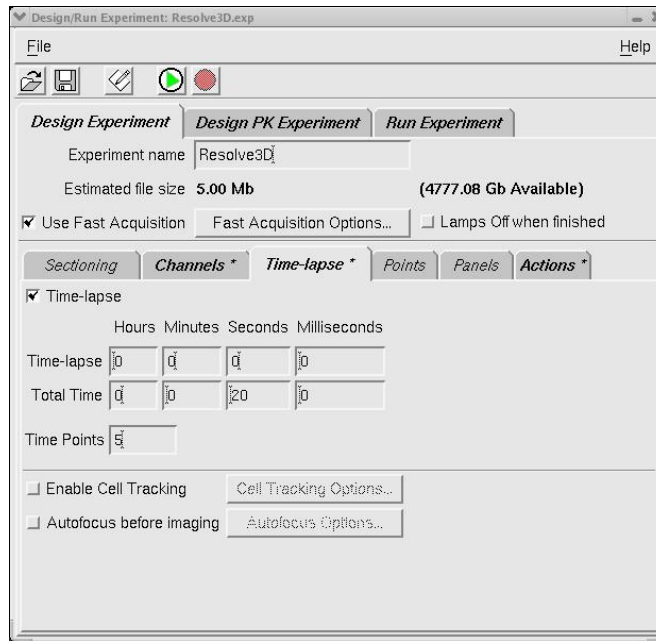


6. Select the **Actions** tab and select **Ratio Image** as the action for this experiment. The Time Points specification will default to **all** and the When control will be **After Imaging**.

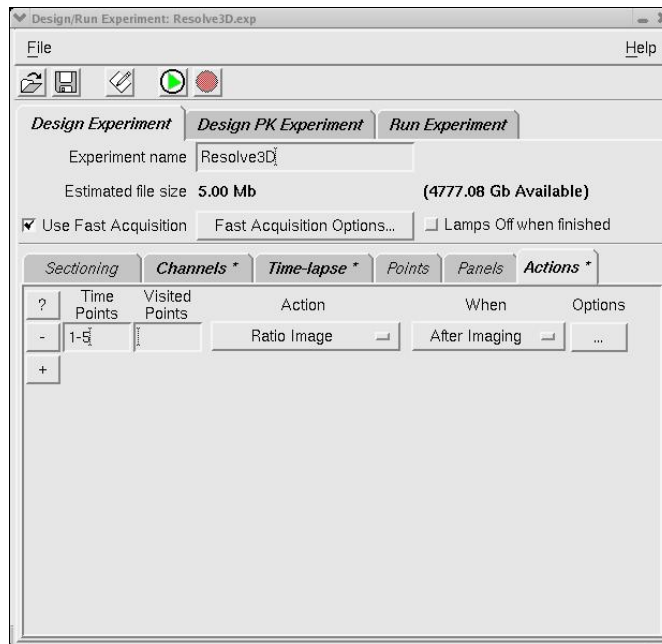

7. Select the **Run Experiment** tab and enter the image file name and a title for the ratio image. You can also enter text into the **Add note to log** field to include the text in your image log file.

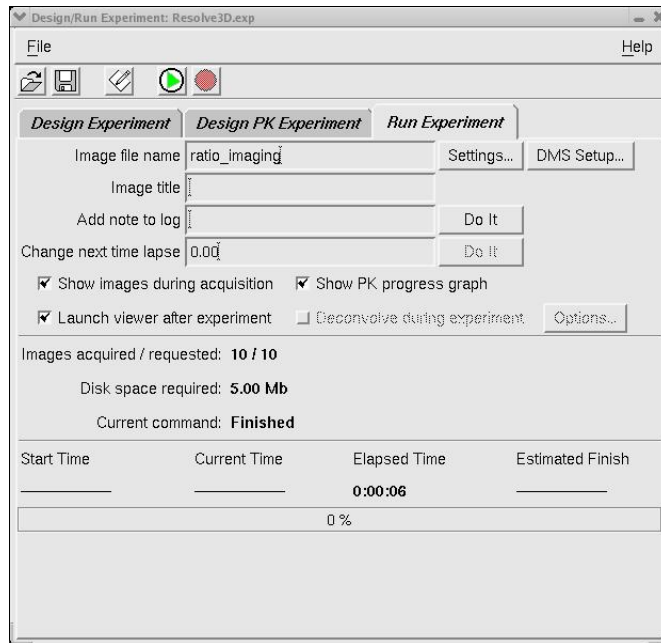


8. Click the **Start Scan** button to begin the imaging process. The ratio imaging process will occur in a separate Image viewer similar to the following:

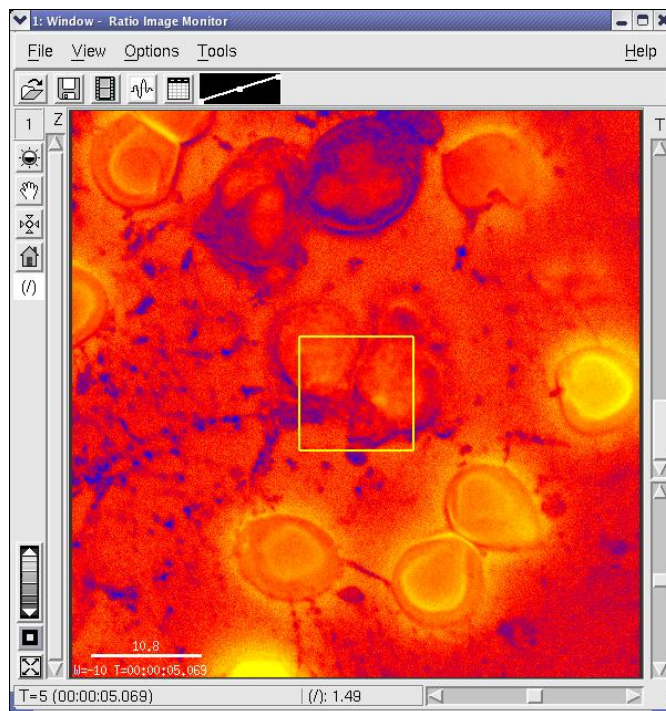

The square outline in the center of the image represents the mean value of the image data.


When the ratio imaging experiment is complete, a two-channel time-lapse image is displayed along with a ratio graph showing the mean value (the outlined area in the middle of the image) vs. time. The square outline

represents the portion of the image used to calculate the mean for a ratio graph.

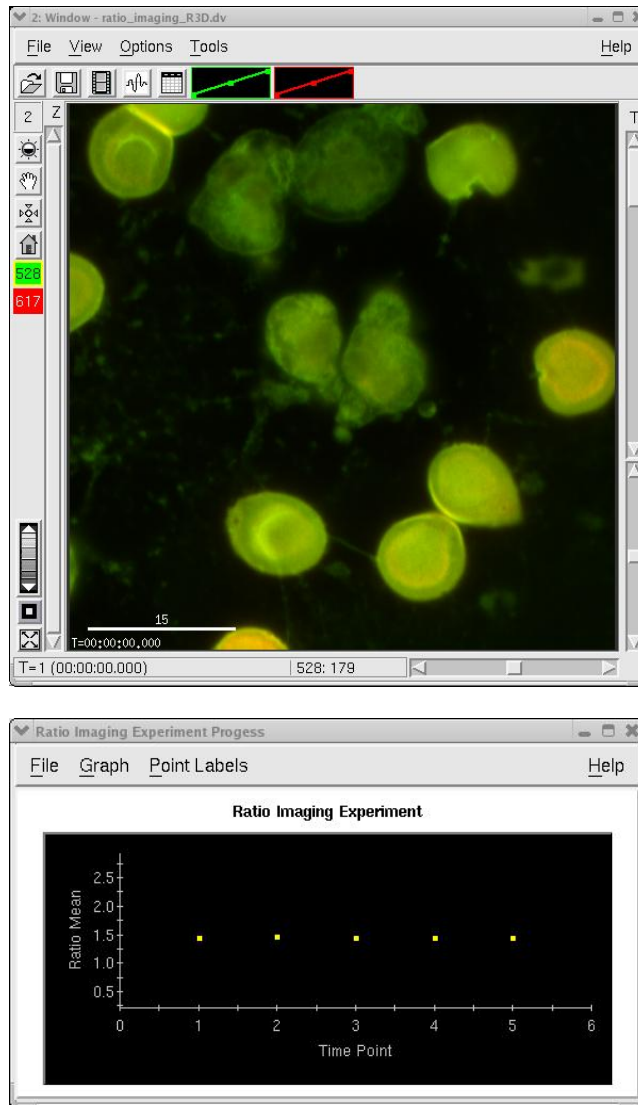

Using the Multiplexed Wavelength Option

The optional Multiplexed Wavelength module for the *DeltaVision* system allows you to perform nearly simultaneous two-channel imaging without the drawbacks associated with true simultaneous two-channel imaging. This option uses two shuttered illumination sources and a dual-band emission filter to eliminate filter wheel movement between channels, and therefore greatly reduces the time required for the *DeltaVision* system to acquire a set of two-channel images. The combined light path ensures no registration artifacts are introduced and independent excitation of probes helps to ensure minimal crosstalk.

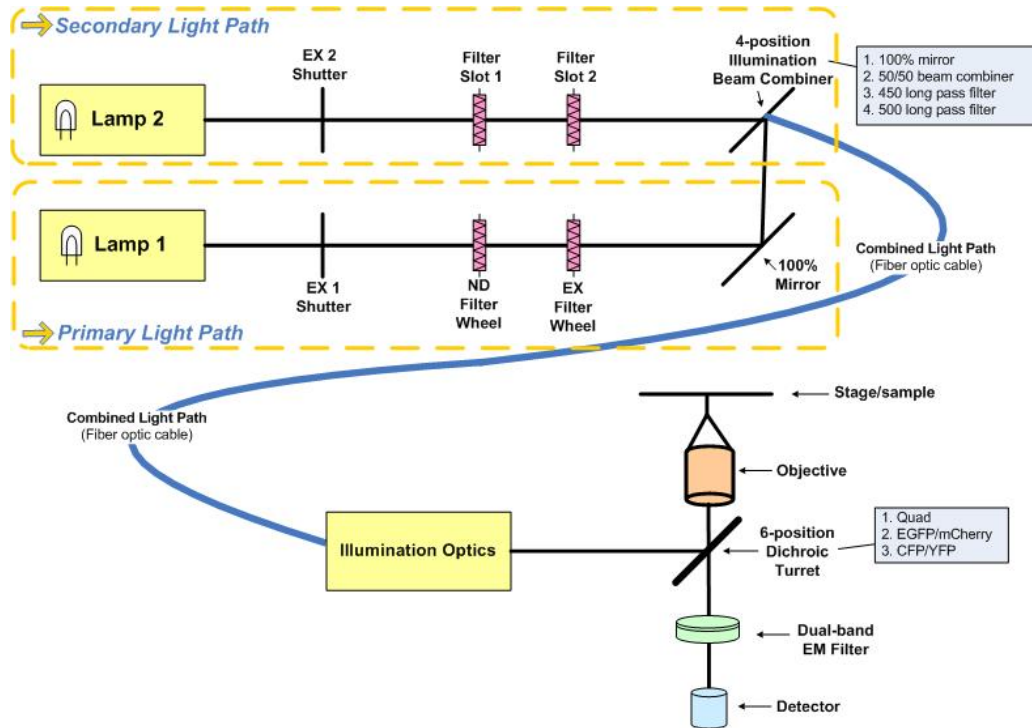

Conceptual view of Multiplexed Wavelength functionality

Before you use the Multiplexed Wavelength option, you must first have it installed and configured correctly. Your Applied Precision representative will assist you in setting up this option and help you ensure that all hardware and software to support Multiplexed Wavelength functionality is installed properly.


After the option has been installed and configured, the menus, tools, and other infrastructure necessary to use the feature will be available on your workstation.

Setting Up the Multiplexed Wavelength Option

Before you begin designing your Multiplexed Wavelength experiment, you'll need to perform the steps described in the following procedures to activate a Multiplexed Wavelength filter set and prepare the DeltaVision system for Multiplexed Wavelength operation.


To activate the Multiplexed Wavelength filter set:

1. To change the active filter set to a filter set that is Multiplex capable, select **Settings** in the Resolve3D main menu.

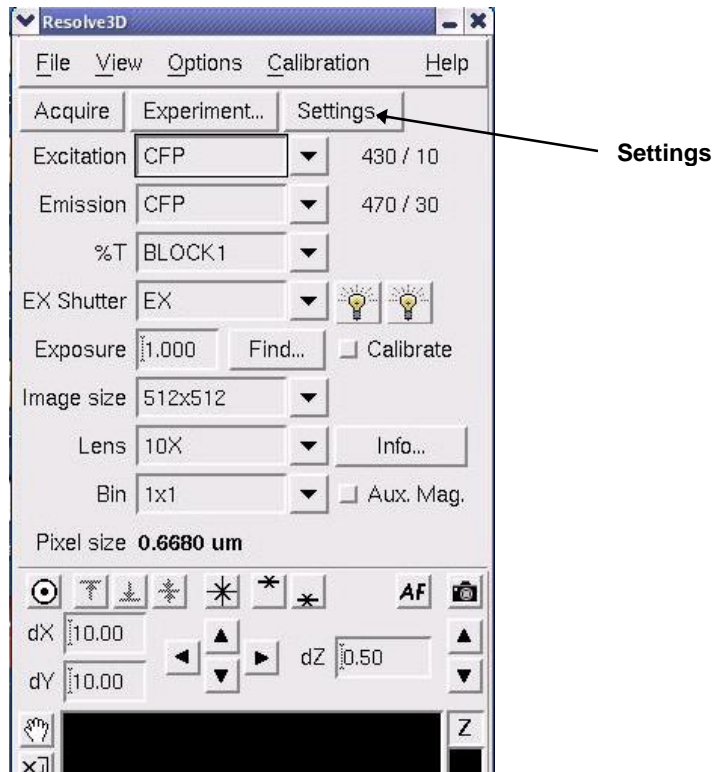


The Resolve3D Settings window is displayed.

2. In the Resolve3D Settings window, click on the Misc tab.

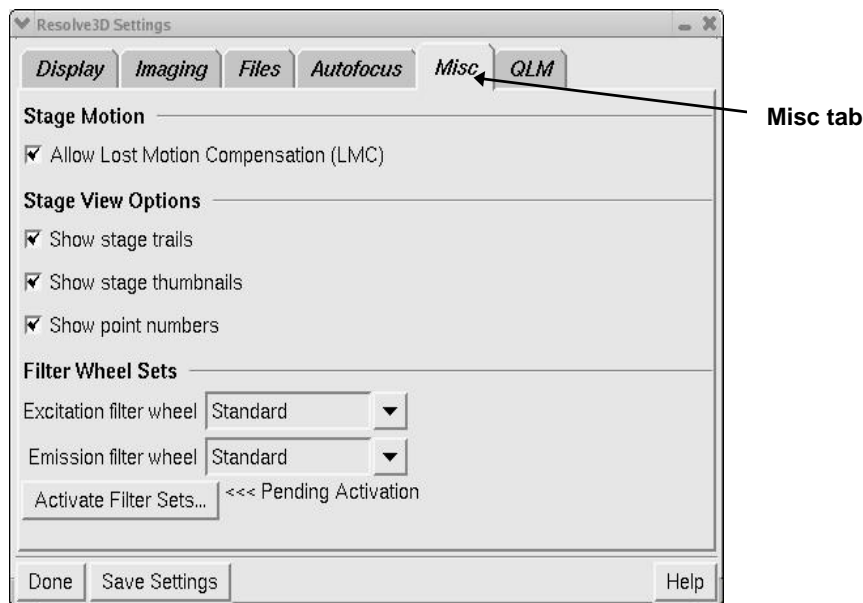


3. In the Resolve3D Settings window, select the EX and EM filter sets you want to use. When these fields are changed, the <<<Pending Activation message is displayed in the window (as shown).

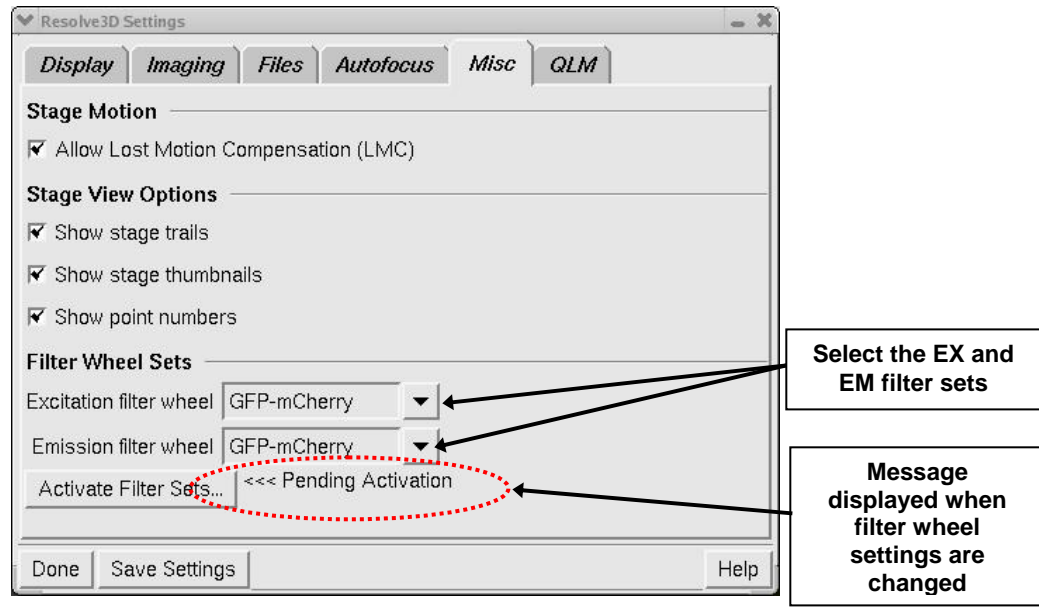

4. Click **Activate Filter Sets**. The following confirmation window is displayed.


Note If you select filter sets for the Excitation filter wheel and Emission filter wheel fields and then click **Done** in this window, your selections are retained until you either activate the filter sets or exit Resolve3D.

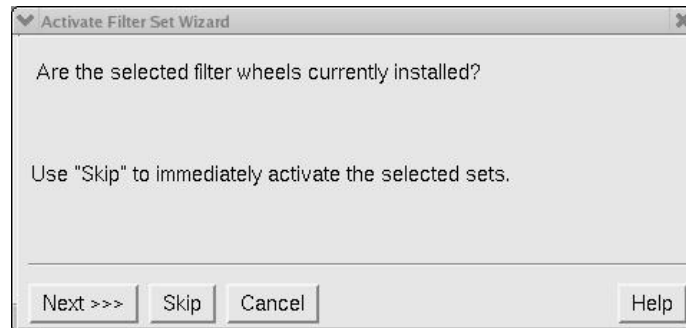


Note For the filter sets to be activated, the selected multiplexed filter set filters must exist in the currently installed excitation and emission filter wheels.

5. If the selected filter wheels are installed on your *DeltaVision*, click **Next** to continue.

Note If you click **Skip** from this window, the selected filter wheels are activated immediately and the remainder of the activation wizard is skipped.

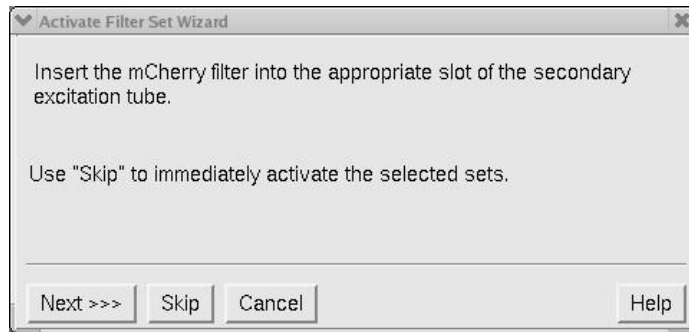


6. Insert the selected secondary filter insert (mCherry is used in this example) into Filter Slot 1 of the secondary light path and click **Next** to continue.

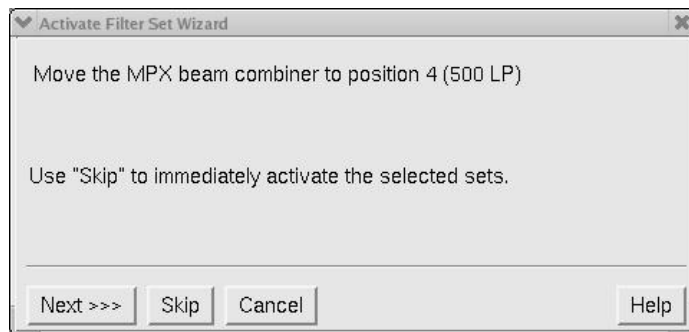


The system gathers the information for this window (in this case, "position 4 (500 LP)" from the `MXWSetup.ini` file, not from the Instrument Controller.

7. Move the beam combiner to the appropriate position and click **Next** to continue.

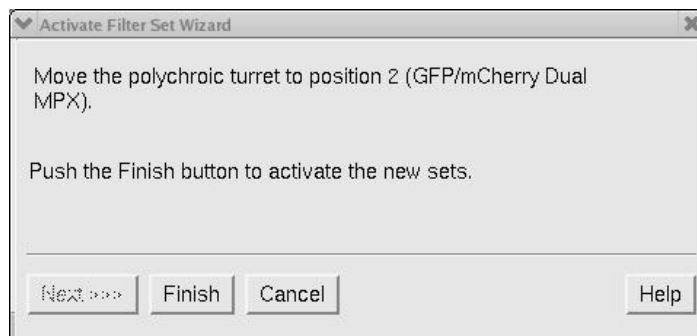

Again, *softWoRx* gets the information for this window (in this case, "position 2 (GFP/mCherry)") from the `MXWSetup.ini` file, not from the Instrument Controller.


8. Click **Finish** to complete the Multiplexed Wavelength filter activation process. After the selected filter set is activated, the Design/Run Experiment window will look similar to the following.

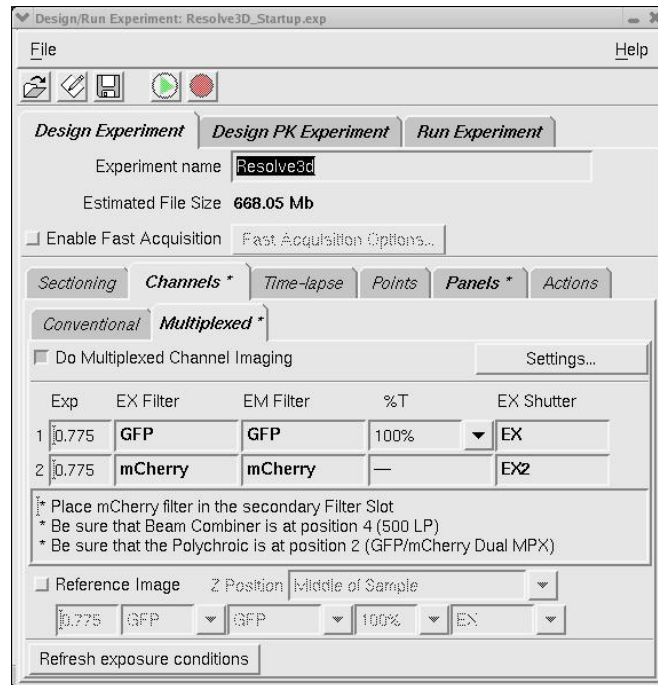


At this point, you have completed activating the Multiplex Wavelength filter set. You should now continue with the steps in the next procedure for viewing a sample with the Multiplexed Wavelength operation.

To view a sample using the Multiplexed Wavelength option:

1. Rotate the eyepiece filter wheel to the POL or BLANK position.
2. From the Resolve3D main menu, select the Excitation filter currently in the primary light path (CFP or GFP).

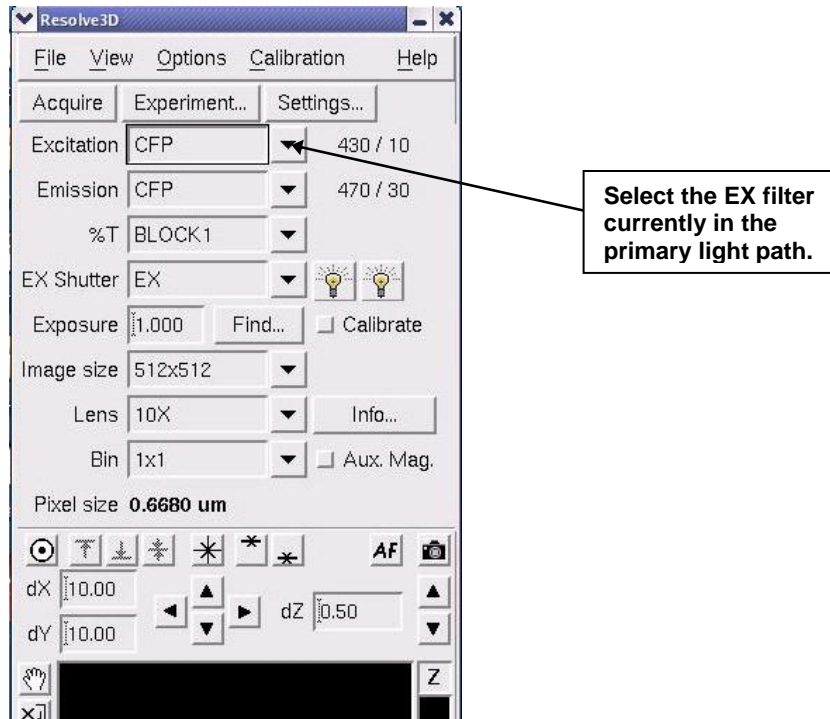

3. Use the EX button on the keypad to open the primary EX shutter and view the primary light path.
4. Use the EX2 (formerly CAMERA SHUTTER) button on the keypad to open the EX2 shutter and view the secondary light path.


 **Note** With the Multiplexed Wavelength option, you can view both selected wavelengths simultaneously by opening both EX shutters at the same time.

Designing a Multiplexed Wavelength Experiment


Use the following procedure to begin the design process for Multiplexed Wavelength experiments.

To design a Multiplexed Wavelength experiment:


1. From the Resolve3D main menu, click the **Experiment** button to open the Design/Run Experiment window.

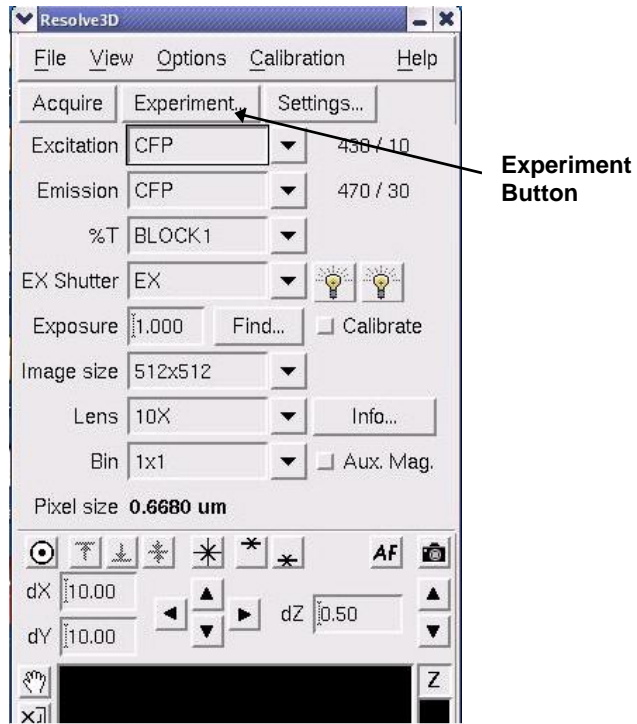

- If the Multiplexed Wavelength option is enabled on your *DeltaVision* system, you'll see the Multiplexed tab in the Design/Run Experiment window. Click on the Multiplexed tab to view the options for Multiplexed Wavelength experiments.

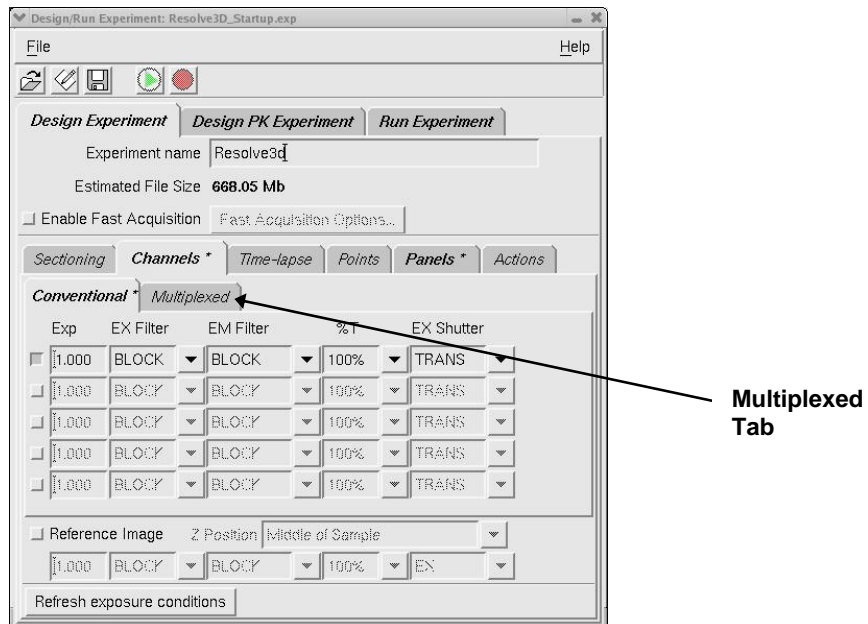


The Multiplexed tab of the Design/Run Experiment window is displayed as shown.

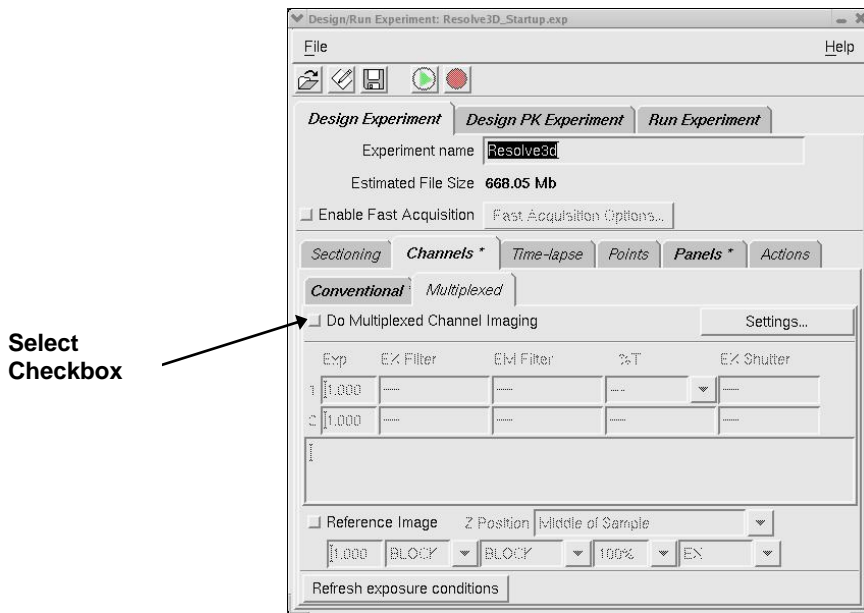


- From the Multiplexed tab, select the **Do Multiplexed Channel Imaging** checkbox.

If the currently active filter set is *not* Multiplex capable, the following window is displayed.

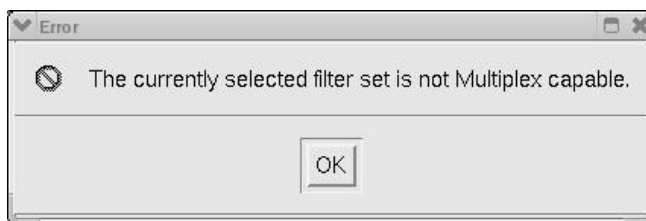

You will need to change the active filter set to continue. Press **OK** to return to the Design/Run Experiment window. To change the active filter set for Multiplexed Wavelength experiments, see the procedure for activating a multiplexed wavelength filter set in "Setting Up the Multiplexed Wavelength Option."

If the currently active filter set *is* Multiplex capable, the Design/Run Experiment window is displayed and will look similar to the following.

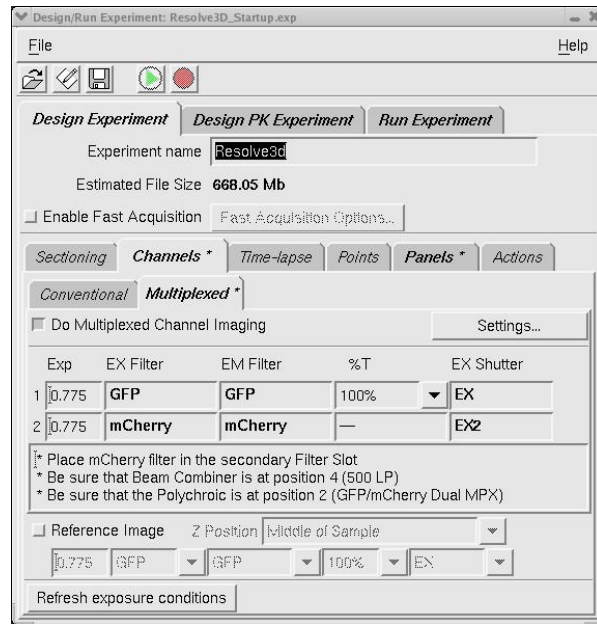

At this point, you have completed the initial portion of the Multiplexed Wavelength experiment design setup. You should now continue with the standard steps for the remainder of the experiment design, such as Sectioning, Timelapse, and so on.

Note As soon as you activate the **Do Multiplexed Channel Imaging** checkbox, all conventional imaging settings are disabled. This is also true of all multiplexed settings when you reactivate conventional imaging.

6. DMS Integration

This chapter describes how to use the DMS (Data Management System) with *softWoRx* to manage image files for your particular application.

What is DMS?

The DMS (Data Management Solution) product provides a functional infrastructure for the storage of biological images and their associated metadata. The DMS Server contains a data management system that centralizes all image data management. Once configured, the DMS Server becomes a repository for all data generated by a laboratory's image acquisition system(s). All visualization and analysis processes are performed on client workstations connected to the DMS Server.

In This Chapter


Connecting to a DMS Database	74
Uploading Images to a DMS Database	74
Downloading Files from DMS	81
Browsing and Locating Images in a DMS Database	82

Connecting to a DMS Database

Before you attempt to connect to a DMS database, you'll need to have a user name and password for the database you want to use. Check with your system administrator for this information.

To connect to a DMS database:

1. From the *softWoRx* main menu, select **File | Connect to DMS Database**. The Connect to DMS Database login box is displayed.

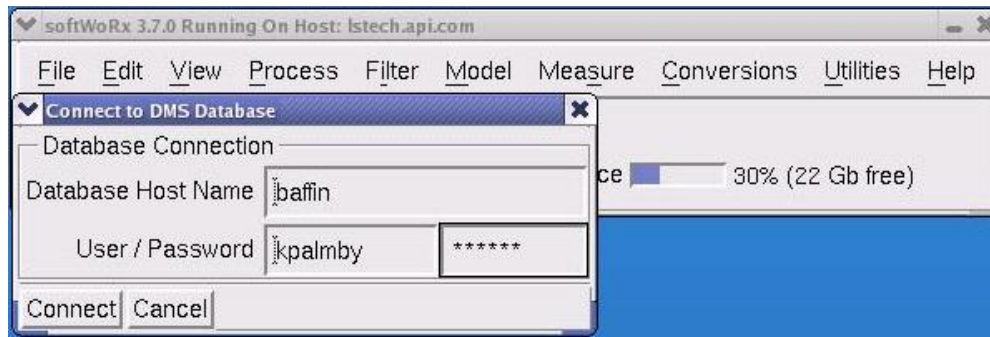

2. Enter the name of host computer in the Database Host Name field.
3. Enter your user name and password in the appropriate fields and click **Connect**. A pop-up message confirms that you are connected to the database.

If the connection fails, make sure you are using the correct user name and password combination for the specific database.

Uploading Images to a DMS Database

You can upload image data from your local file system to the DMS database in a number of ways. You can upload image data:

- Directly from an Image Window.
- As part of a Task Builder processing chain.
- As part of an experiment, in which the data is auto-uploaded at the conclusion of the experiment.


Each of these methods is discussed in the following sections.

Uploading Images from the Image Window


The simplest method for uploading image data to the DMS database from your local file system is to upload the data directly from the Image Window.

To upload image data from the Image Window:


1. From the *softWoRx* main menu, select **File | Open**. When the Open Image window is displayed, click the Files tab and select the appropriate directory.

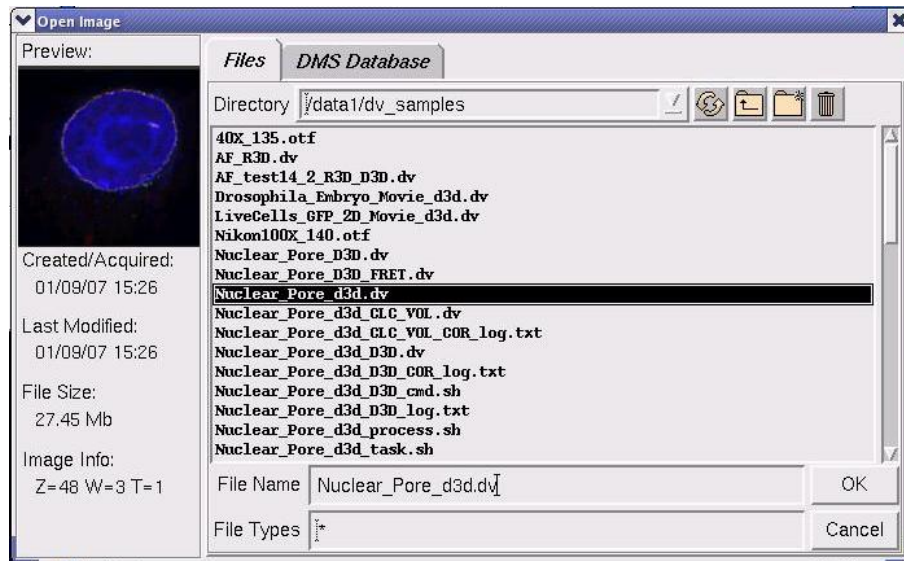

2. Select the file you want to upload and click **OK**. The Image Window is displayed with the file you selected.

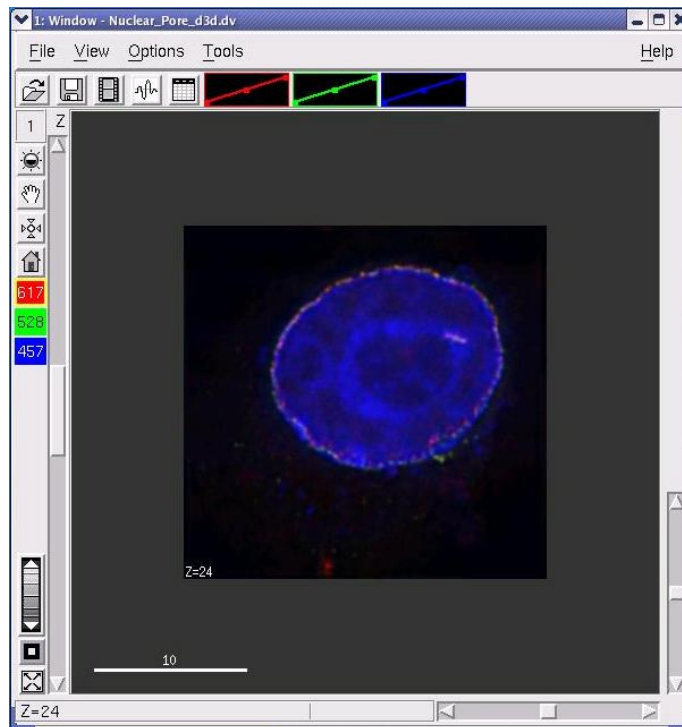


3. From the Image Window, select **File | Save to DMS**. The Save To Data Management System menu is displayed as shown.

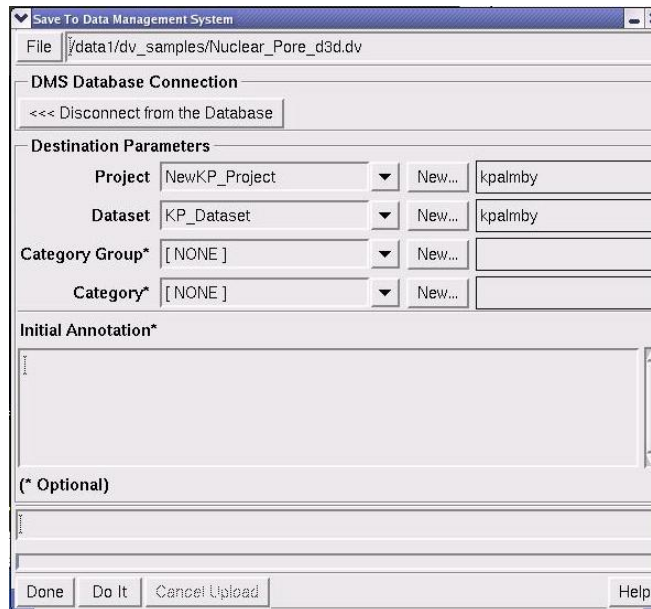


From this menu, select the Project and Dataset to which you want to save the data on the DMS database. You can optionally select a Category Group and Category for the saved image data and add initial annotation to the file for future reference. When you are satisfied with the image data to be uploaded, click **Do It**.

When the upload is complete, “Finished” is displayed on the status line as shown.

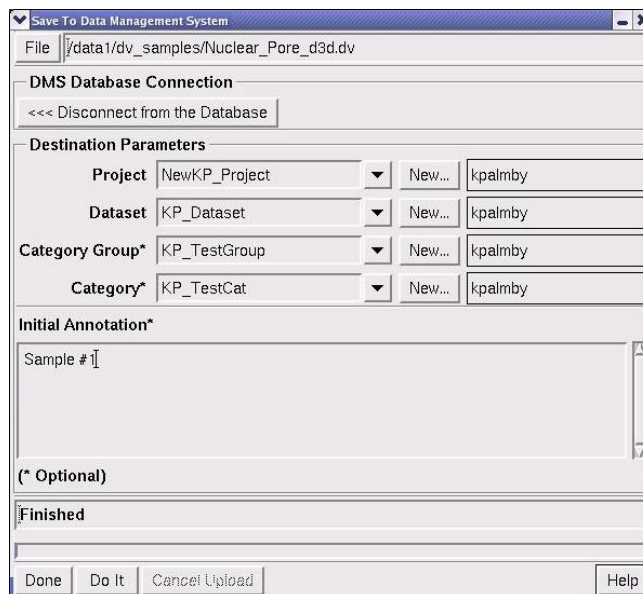

4. Click **Done** to exit the Save To Data Management System menu.
5. To confirm that the selected image file has been uploaded to the DMS database, select **File | Open** from the *softWoRx* main menu. When the Open Image window is displayed, click the DMS Database tab.

Note To activate the DMS Database tab, you must be currently connected to a DMS database. Refer to the previous section, “Connecting to a DMS Database” for details.

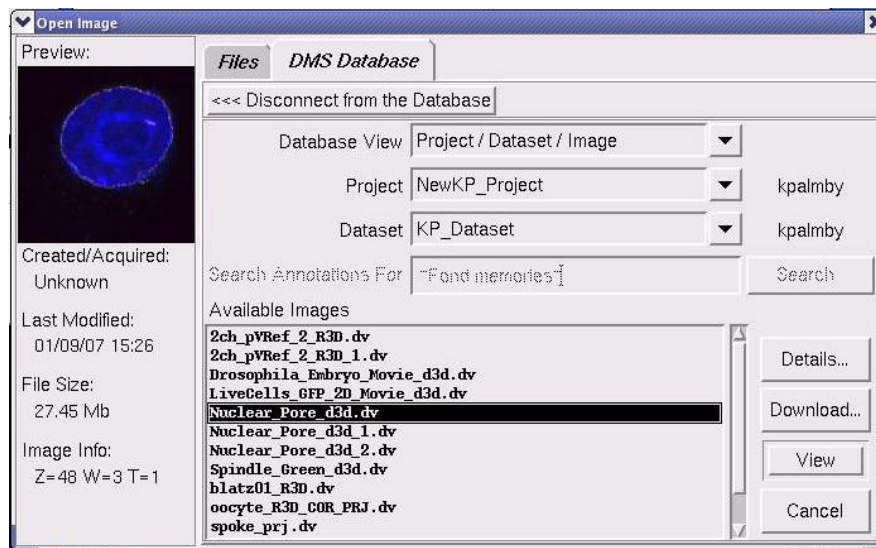


The file you uploaded should appear in the displayed list of files.

Uploading Images Using Task Builder


To upload image data from your local file system to the DMS database from within Task Builder, use the following procedure.

To upload image data to the DMS database using Task Builder:


1. From the *softWoRx* main menu, select **Process | Task Builder**. The Processing Task Builder menu is displayed.

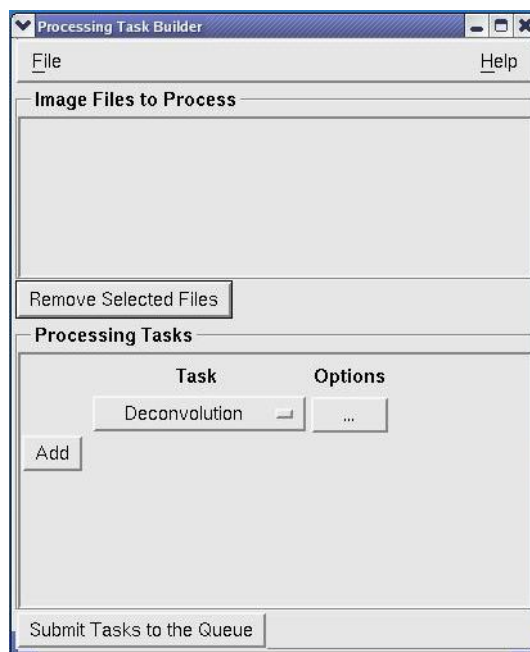


- In the Processing Task Builder menu, select the file(s) you want to process and upload to the DMS database. You can accomplish this by simply dragging and dropping the file or folder icon(s) into the **Image Files to Process** area of the menu. Alternatively, you can use **File | Add Files** from this menu to select the file(s) to include.

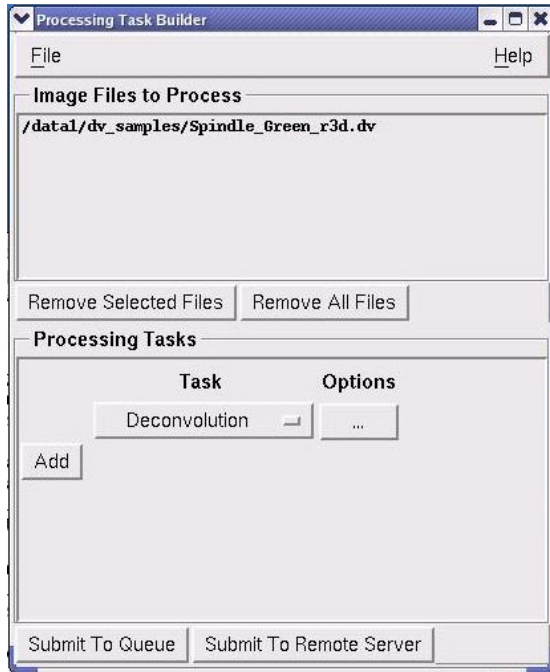


- Next, select the processes to perform on the image file(s). For the final task, select **Save to DMS** as shown.

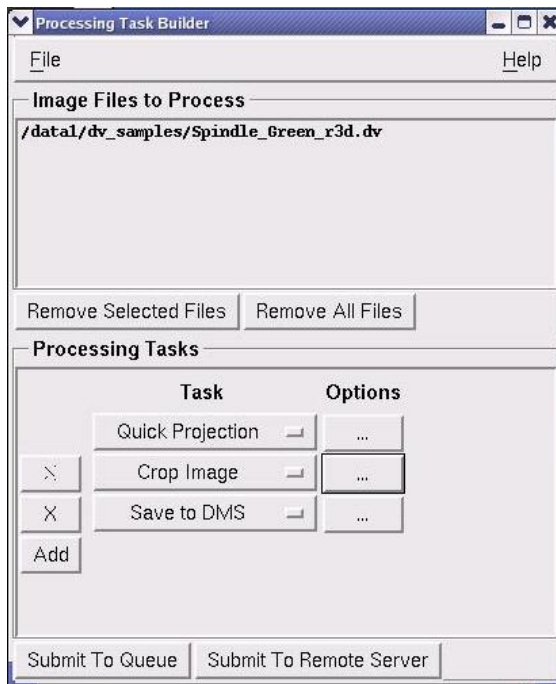

- Click the **Options** button next to the **Save to DMS** selection. The Save to DMS Options menu is displayed.

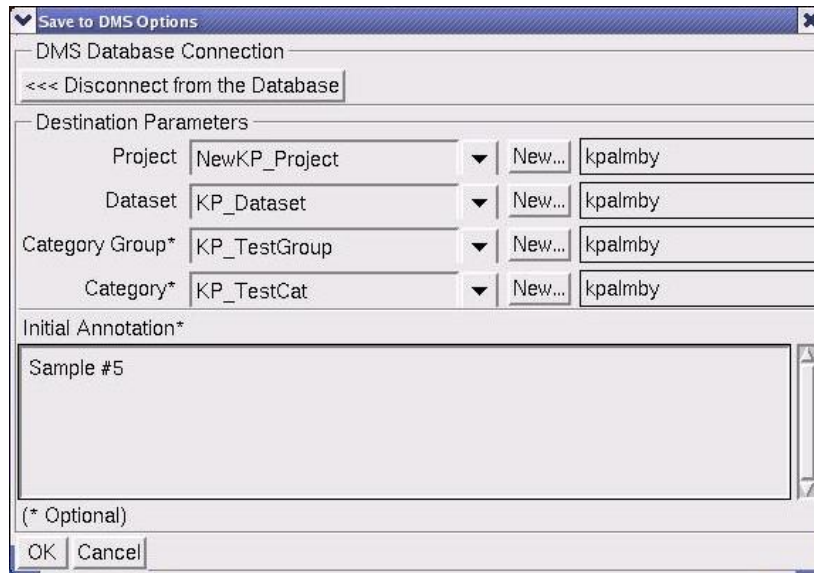

5. From this menu, select the Project and Dataset to which you want to save the data on the DMS database. You can optionally select the Category Groups and Category for the saved image data and add initial annotation to the file for future reference. When you are satisfied with the image data to be uploaded, click **OK**. You are returned to the Processing Task Builder menu.
6. From the Processing Task Builder menu, click **Submit to Queue** to open the *softWoRx* Queue Manager dialog box.

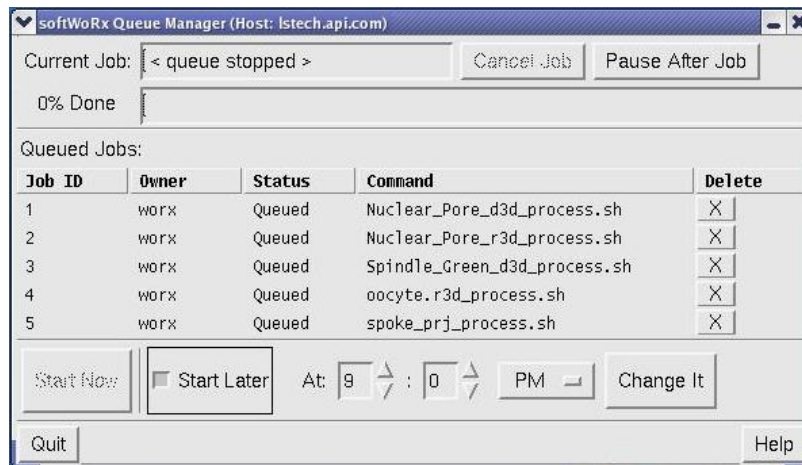


7. To start the process chain,
 - *Immediately*, click **Start Now**.
 - *At a later time*, click **Start Later**. When you choose this option, a set of time option buttons is displayed. Set the time at which you want the process chain to start. You can use the **Change It** button if you decide you want to change the time to begin the process chain.
8. Select **Quit** to exit the Queue Manager dialog box.

Auto-uploading Images after Acquisition


To auto-upload image data to the DMS database after an experiment is completed, use the following procedure.

To auto-upload image data immediately following an acquisition:

1. Set up your experiment as normal by clicking **Experiment** from the Resolve3D window to open the Design/Run Experiment window.
2. On the Design Experiment tab, enter the experiment name and configure the experiment. (For experiment setup details, see the "Setting Up and Running Experiments" Chapter in your *DeltaVision System User's Manual*.)
3. Select the Run Experiment tab and enter the image file name (or drag and drop from another location), the image title, and any annotation you want to add.

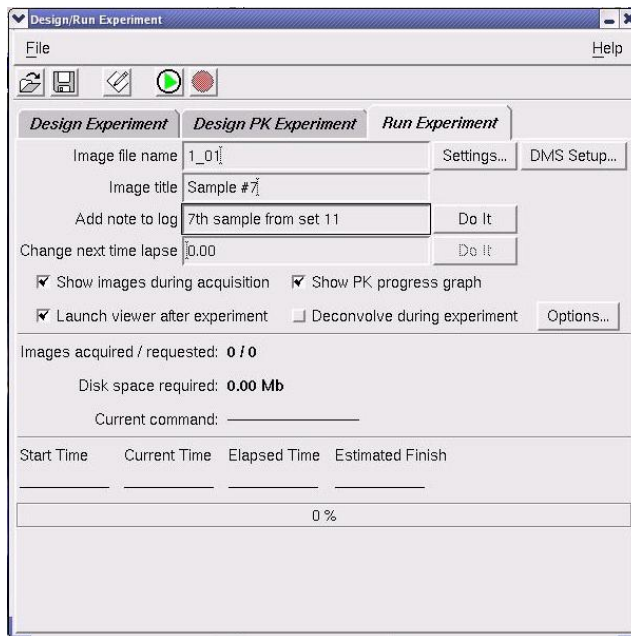

4. Click the **DMS Setup** button. The DMS Destination Setup window is displayed.

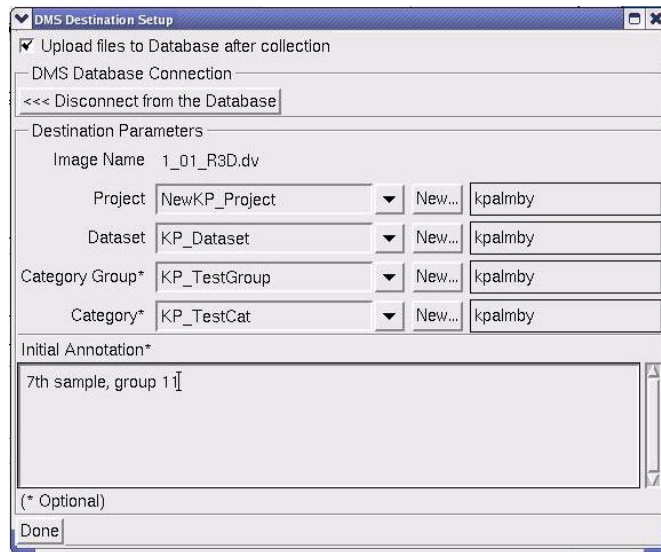


5. Be sure that the **Upload files to Database after collection** checkbox (at the top of this screen) is activated.
6. Enter the project and dataset where you want the acquired image data to upload.
7. Optionally, you can also enter the category group and the category for the uploaded data and include any annotation necessary.
8. When you are satisfied with the information entered in the DMS Destination Setup window, click **Done**.
9. Run the experiment macro as normal. The resulting image acquisition will be added to the DMS database as specified.

Downloading Files from DMS


You can use *softWoRx* to download image data from the DMS database to your local file system.

To download a file from the DMS database:

1. From the *softWoRx* main menu, select **File | Open**. When the Open Image window is displayed, click the DMS Database tab and select the file you want to download.

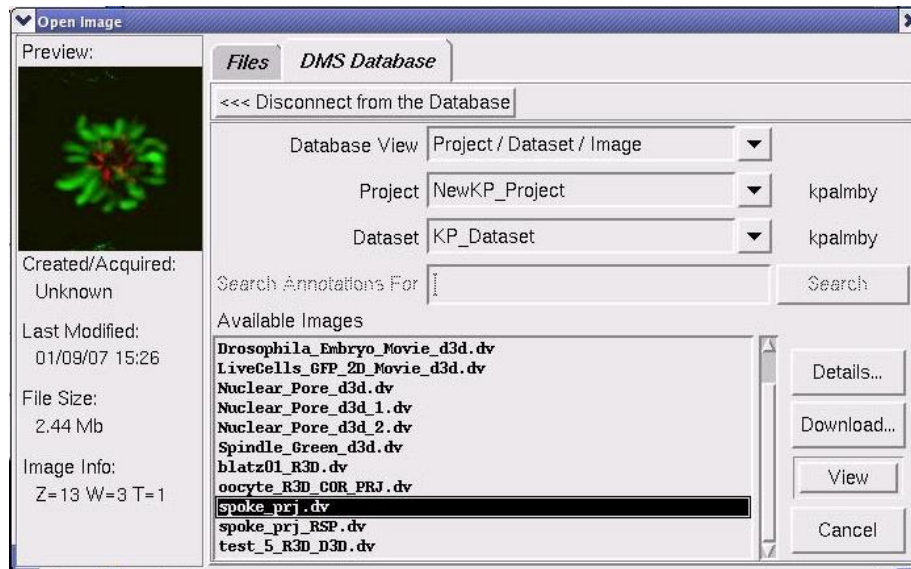

2. Click the **Download** button. The Choose Destination Folder dialog box is displayed.

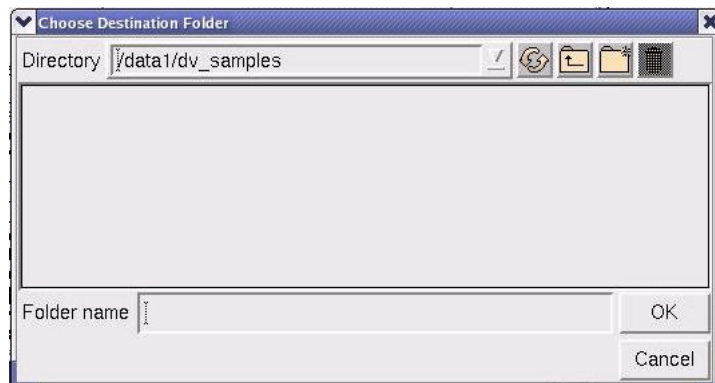

3. Select the directory on your local file system where you want the selected image file downloaded and click **OK**. The image file is downloaded from the DMS database to the specified location in your local file system.


Browsing and Locating Images in a DMS Database

The *softWoRx* software provides the ability to browse images within a DMS database by either the Project/Dataset/Image or the Category Group/Category/Image hierarchies. You can also search for specific images based on the image file's annotation.

Browsing Image Files using P/D/I Hierarchy

To browse through image files by Project/Dataset/Image:

1. From the *softWoRx* main menu, select **File | Open**. When the Open Image window is displayed, click the DMS Database tab.

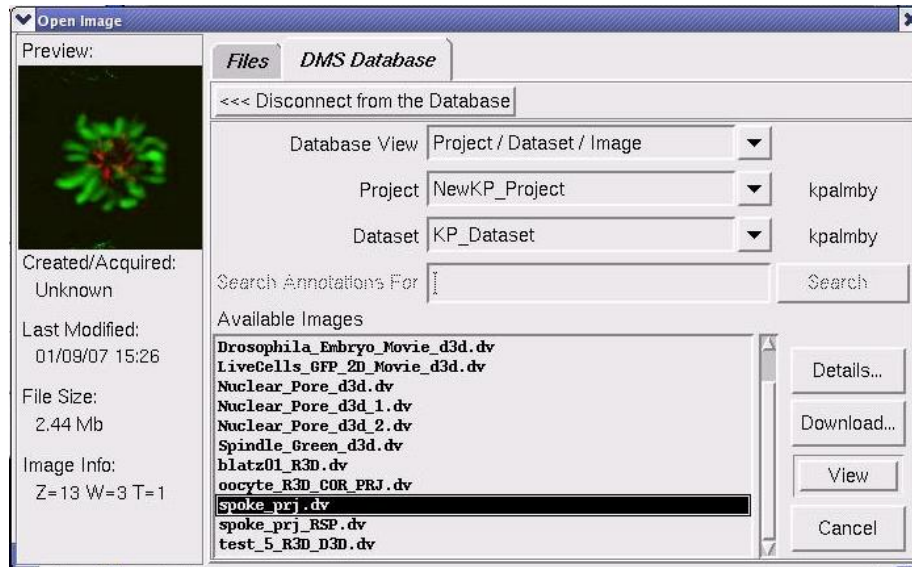


2. Use the drop-down menu in the Database View field to select **Project/Dataset/Image**.
3. Select the specific projects and datasets to browse using the dropdown menus in the Project and Dataset fields.

Browsing Image Files using CG/C/I Hierarchy

To browse through image files by Category-Group/Category/Image:

1. From the *softWoRx* main menu, select **File | Open**. When the Open Image window is displayed, click the DMS Database tab.

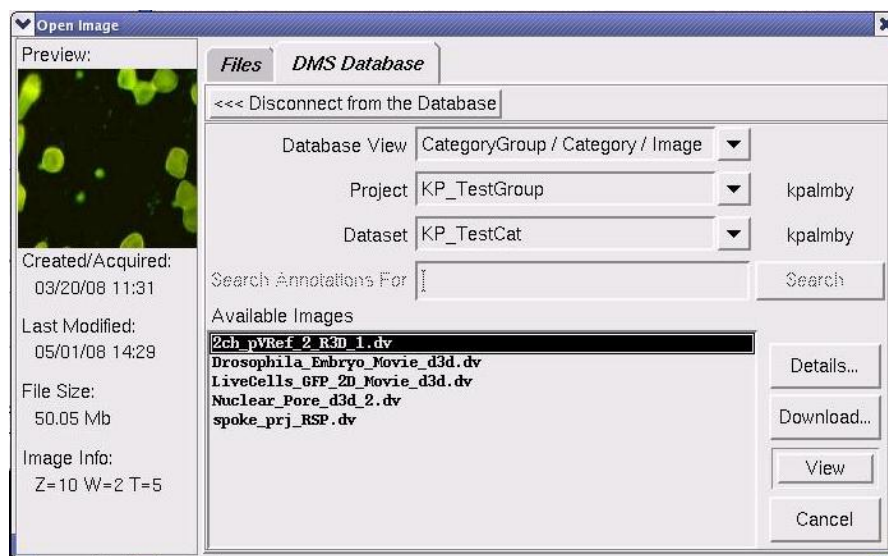

2. Use the drop-down menu in the Database View field to select **CategoryGroup/Category/Image**.


3. Select the specific category groups and categories to browse using the drop-down menus in the Category Group and Category fields.

Searching for Image Files Based on Annotation

You can locate images in a DMS database by performing a search based on the image file's annotation.

To search for an image file based on its annotation:

1. From the *softWoRx* main menu, select **File | Open**. When the Open Image window is displayed, click the DMS Database tab.

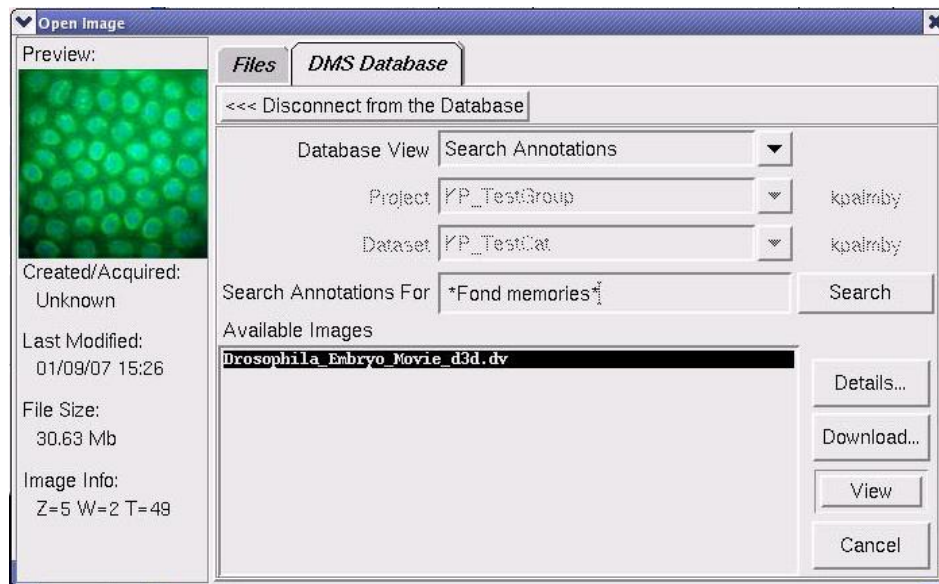

2. Use the drop-down menu in the Database View field to select **Search Annotations**.
3. In the Search Annotations For field, enter the annotation for the desired image file and click the **Search** button.

Note Holding your mouse cursor over the Search Annotations For field displays a tooltip window containing specific wildcard use information.

Part Two

VISUALIZING & PRESENTING DATA

softWoRx provides several tools that you can use to visualize data and prepare it for presentations. You can also save or export data in a variety of formats.

In Part Two

Chapter 7. Viewing Image Data	87
Chapter 8. Viewing Movies.....	113
Chapter 9. Viewing Projections and Volumes.....	119
Chapter	

10. Filtering Image Data.....137

Chapter

11. Saving, Exporting, and Printing147

7. Viewing Image Data

softWoRx provides several options for viewing data. You can open *DeltaVision* or TIFF files in the Image window and use slide bars and controls to view different Z sections and time points. You can also adjust the brightness and contrast of each channel in an image and assign a color or grayscale to each channel. To prepare your data for presentations, you can display a scale bar on the image and hide the Image window controls. You can also resample image data to change the size or orientation of the display and view cross sections of the data.


In This Chapter


Opening an Image.....	88
The Image Window	89
Viewing 5D Images.....	90
Adjusting Brightness and Contrast	95
Assigning Colors or Grayscale to Channels.....	99
Controlling the Image Window Display	102
Resizing or Reorienting an Image	105
Viewing Cross Sections.....	111


Opening an Image


You can open image data files in *DeltaVision* or TIFF image file formats.

To open an image data file:


1. Choose **Applied Precision | Start softWoRx** from the CentOS  menu. The softWoRx main menu is displayed.

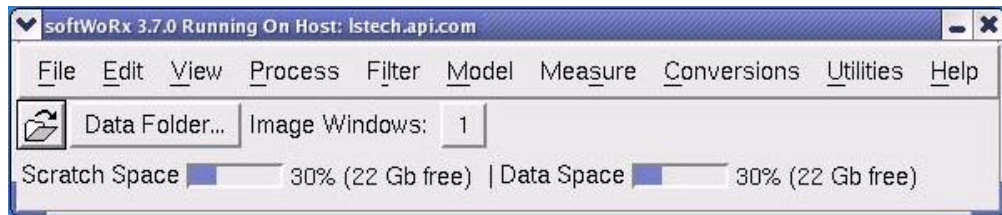


2. Click the  icon on the *softWoRx* toolbar, or **File | Open** from the menu. The Open Image dialog box is displayed. Click the Files tab.

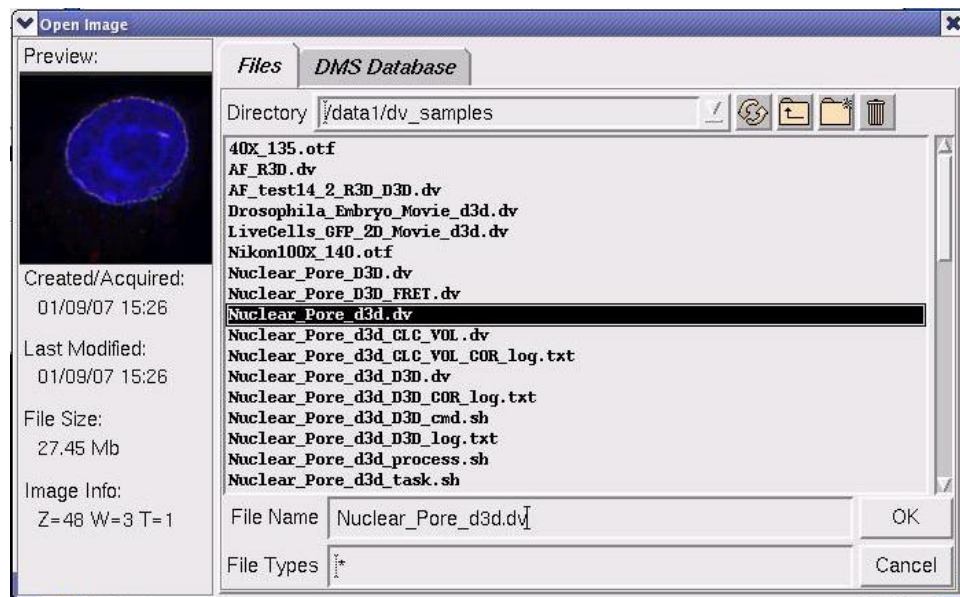

Select and preview *DeltaVision* images in the Open Image dialog box.

 **Tip** For *DeltaVision* images, you can adjust the thumbnail image in the Preview area. Change the brightness by dragging the left mouse across the preview image. View different Z sections or time points by pressing the right and middle mouse buttons.

3. Select an image file (for example, `Nuclear_PoreD3D.dv`) and click **OK** to open the image in the Image window.

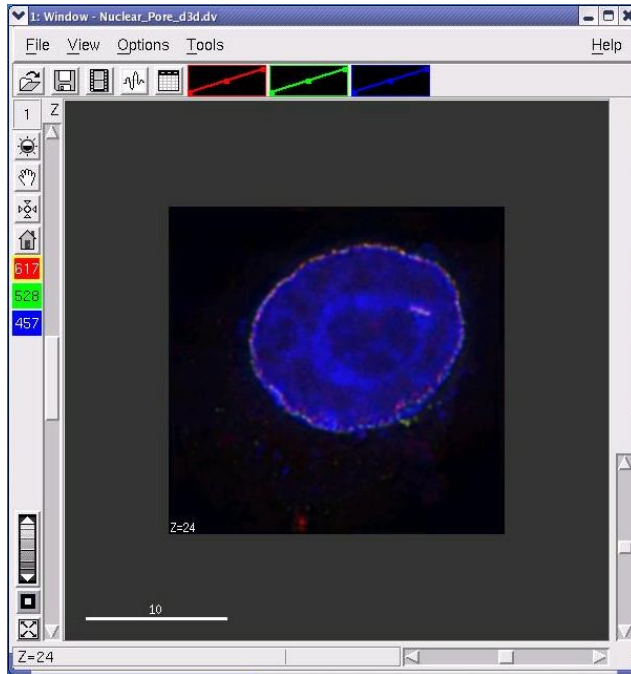


The image opens in the Image window

 **Tip** You can also open an image by clicking Data Folder and then double-clicking on the .dv file in the data1 directory.

The Image Window

DeltaVision images are displayed in the Image window.

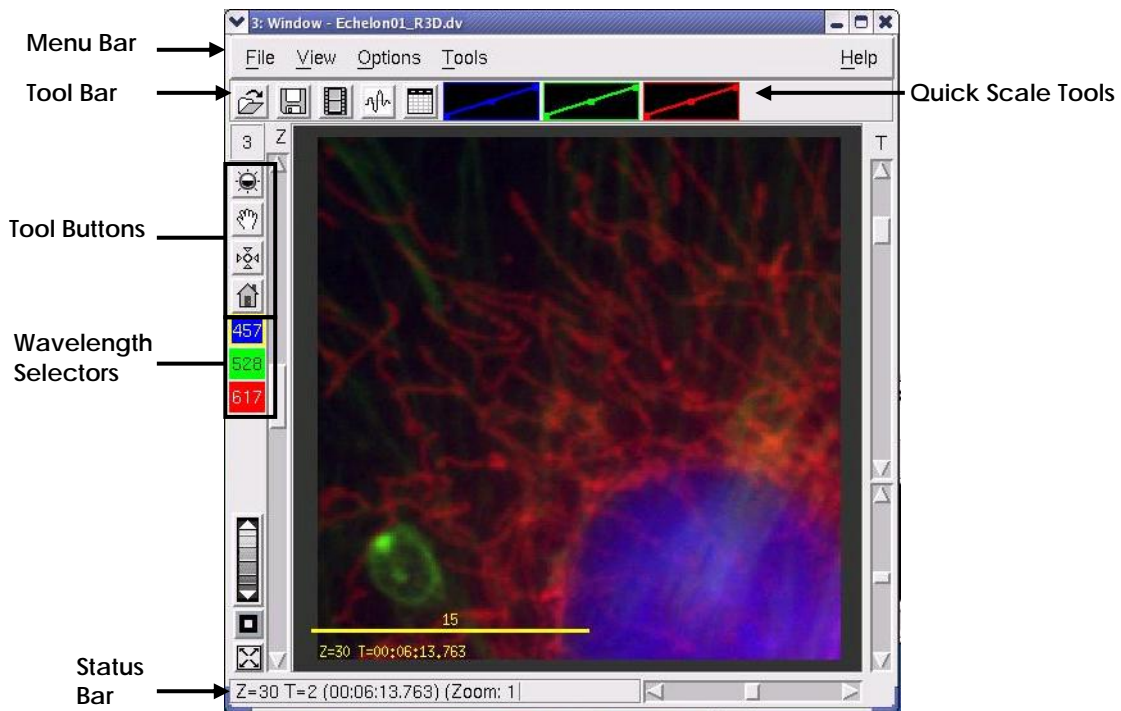

The Image Window

The Image window provides controls and tools that you can use to view and analyze image data:

- The **Menu Bar** allows you to open or save images, control the display of the data, and open tools to analyze intensity data.
- The **Toolbar** provides buttons as alternatives to the menu bar above. The buttons can be used to open or save image data files, play movies, view intensity line profiles, or inspect intensity data.
- The **Tool Buttons** are used to change the image view. These controls allow you to scroll through Z sections and time points, zoom, pan the image vertically or horizontally, select which channels to view, and scale intensity to adjust brightness and contrast.
- The **Wavelength Selectors** show or hide the wavelengths (channels) in the Image window. When the Image is displayed in color, each button has the same color as the wavelength that it controls. The number on each control button indicates its wavelength. When the data is displayed as grayscale, use these controls to choose which wavelength is displayed. The buttons are white with black numbers when on and black with white numbers when off.
- The **Status Bar** shows which Z section or time interval is displayed. It also displays the intensity of the pixel currently under the mouse pointer.
- The **Quick Scale Tools** provide a convenient method by which you can adjust the intensity scale of each wavelength in the image.

Viewing 5D Images

After opening an image file, use the *softWoRx* Image window controls for:

- Viewing Z sections and time points in the data.
- Viewing different areas of a section.
- Zooming in on a selected area.
- Displaying or hiding channels.


Viewing Z Sections and Time Points

Use the Z and T sliders on the left side of the Image window to display different Z sections and time points. If images contain both Z and T data, you can use the Z

slider to show all of the sections at a time point or use the T slider to show how a section changes with time.

To navigate through Z sections and time points:

- ▶ Move the Z slider down to show deeper Z sections (the top section is displayed when the slider is at the top). Move the T slider down to show data acquired at later time points (delta time increases as the slider is moved down).

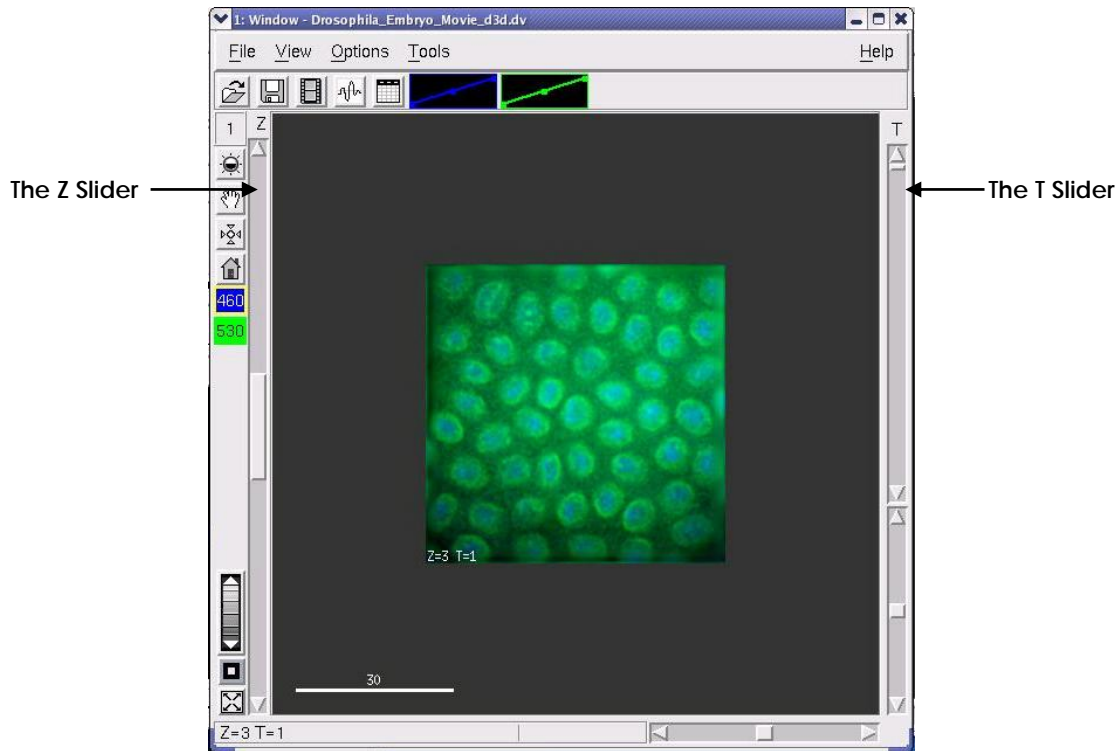

Move the Z and T sliders to display Z sections and time points


Tip You can also use the right and middle mouse buttons to scroll through Z sections. To scroll through time points, hold down the CTRL key and press the right or middle mouse buttons.

Viewing Z Sections in Several Image Windows

The Section slider allows you to scroll through Z sections of all open windows simultaneously. (Only the windows that are open at the time that you activate the Section slider are affected.)

To scroll Z sections simultaneously:

1. Choose **View | Slider** from the *softWoRx* main menu to open the Section Slider tool.

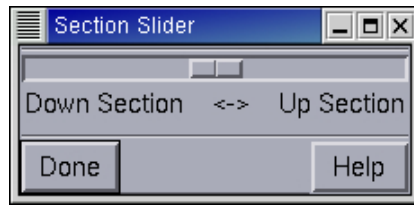


The Section Slider

2. Move the slider to the left to decrease the current Z Section or to the right to increase it.





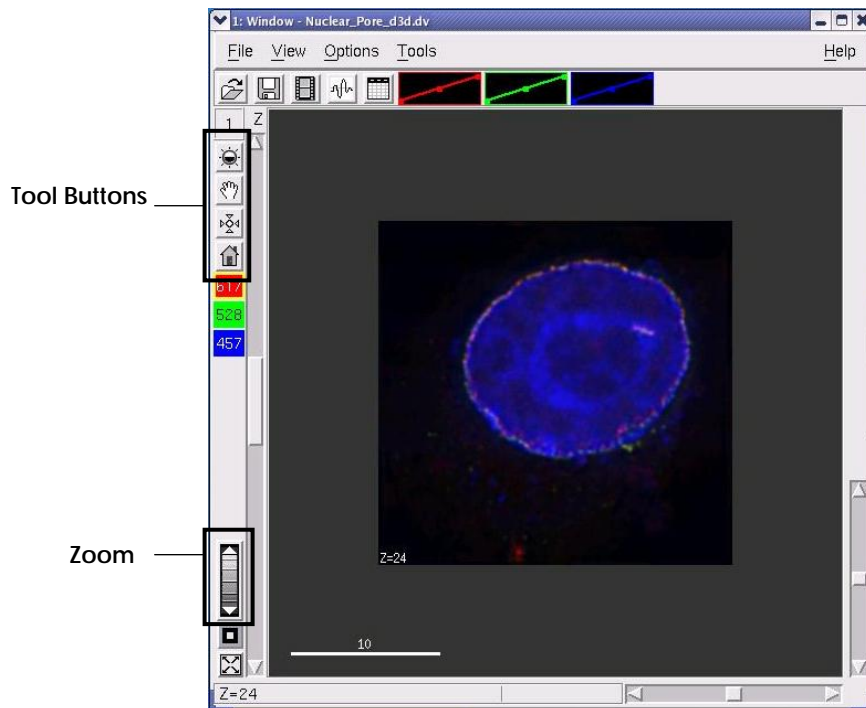


Viewing Areas

You can view different areas of a section by sliding the vertical and horizontal scroll bars and the tool buttons to reposition the image.



Use tool buttons, scroll bars, and zoom to view different image areas.

You can also use the following tool buttons to change display characteristics or reposition the image.


Use this tool	To...
	Adjust the intensity scale of the image.
	Pan across the image.
	Center the image on a point. Select the tool and click the point that you want to center.
	Position the image on its original center.
	Zoom in or out. Move the Zoom wheel up to zoom out or down to zoom in.
	Return to a 1:1 zoom level.
	Zoom to fit image to window.


 **Tip** From the Image window, you can also use **Alt + Mouse Wheel** to zoom the view in and out.

Zooming In or Out on Specific Points


One common way to use the tool buttons is to position the image and zoom in on a specific point.

To zoom in on a specific point:


1. On the Image window, click the **Choose New Window Center**  button and then click the point in the image on which you want to zoom. The image is centered on the selected point.

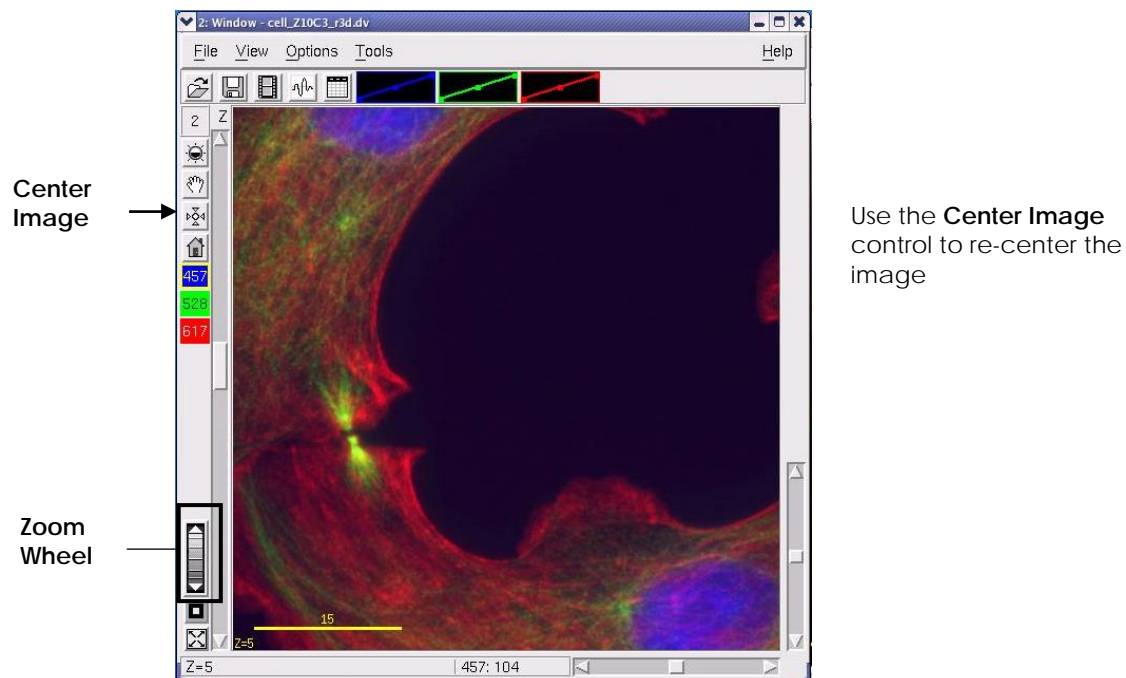


- Use the Zoom wheel to zoom in on the point that you selected. The new zoom level is displayed in the status bar.

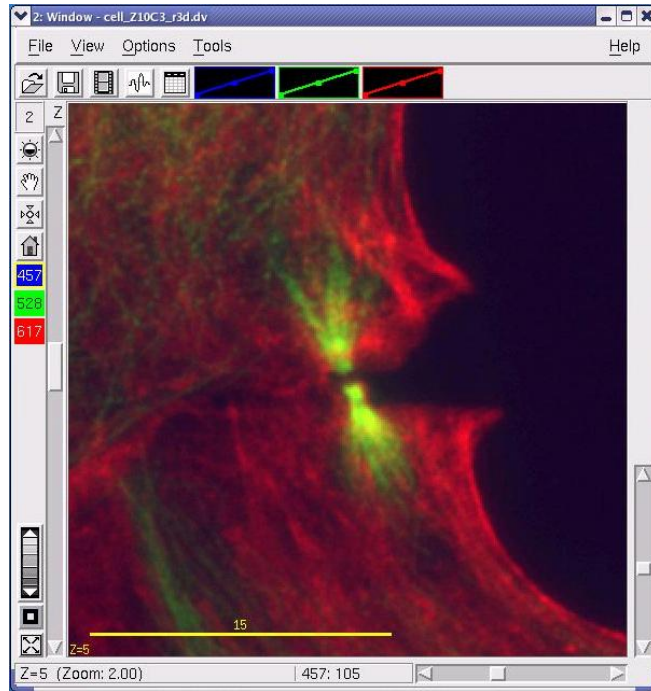


Use the Zoom wheel to zoom on the point that is selected as the center point

- To return the zoom to a 1:1 display, click the **Reset Zoom to 1:1**  button.

 **Tip** You can specify to interpolate (smooth) images when the zoom level is greater than 1. Interpolated images provide better quality results but take longer to display. To interpolate, choose **Options | Display** and select the Interpolate zoom option.

Displaying or Hiding Channels

You can display or hide channels with the Wavelength Selectors on the Image window. When the channel is displayed, the channel displays the color that is assigned to it. When the channel is not displayed, the selector is black.

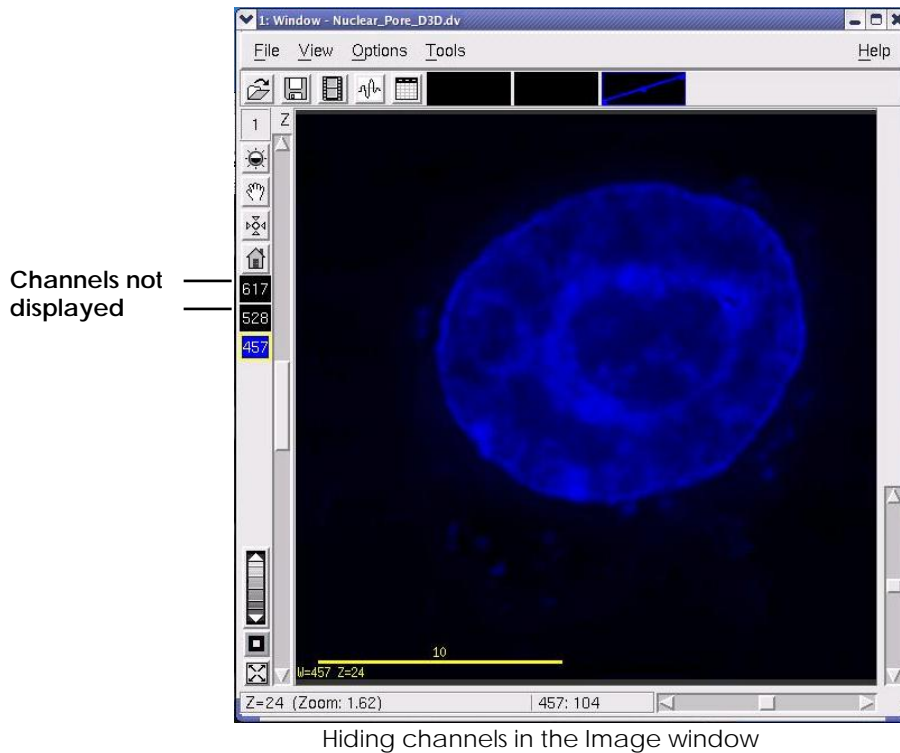

To hide a channel:

- ▶ Double-click the Wavelength Selector of a displayed channel. The Wavelength Selector turns black to indicate that the channel is not displayed.

To display a channel:

- ▶ Click the Wavelength Selector of an undisplayed channel. The Wavelength Selector displays the color of the channel and the channel is displayed.

Adjusting Brightness and Contrast

You can improve the contrast of selected data in a channel by changing the channel's intensity scale.

softWoRx uses shades of the color selected for a channel to display an intensity scale. The darkest shade is mapped to the lowest (dimmiest) intensity value in the wavelength and the lightest shade is mapped to the highest (brightest) intensity value. The remaining shades are mapped to values between the lowest and highest values.


Color shades can be mapped to create linear or nonlinear intensity scales. In linear scales, the color shades are mapped to values that are distributed evenly from the minimum to the maximum intensity values. In nonlinear scales, the shades are mapped to values that are distributed unevenly throughout the range.

You can adjust the intensity scale for an individual image window or for all open image windows simultaneously.

Note Changing intensity scale values only adjusts the display of the data. It does not alter the image data.


To change the intensity scale for the current image window:

1. On the Image window, click the **Scale Image Intensities**  button to open the Image Scaling dialog box. This dialog box shows the image intensity scale.

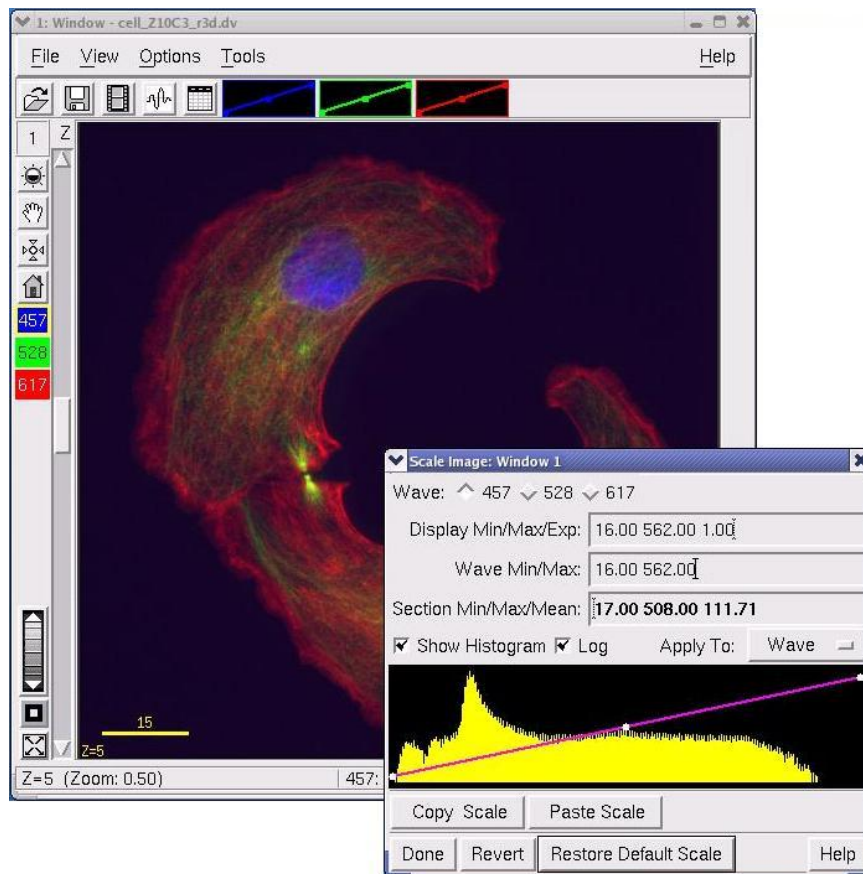

The Scale Image dialog box is used to adjust the intensity scale. The histogram on this dialog box is a frequency plot that shows the distribution of pixel intensities in the image file. The Y-axis shows the number of pixels for a given intensity.

2. In the **Wave** field, select which channel to scale.
3. To change the minimum or maximum scale value, click and drag on the Left or Right handles.

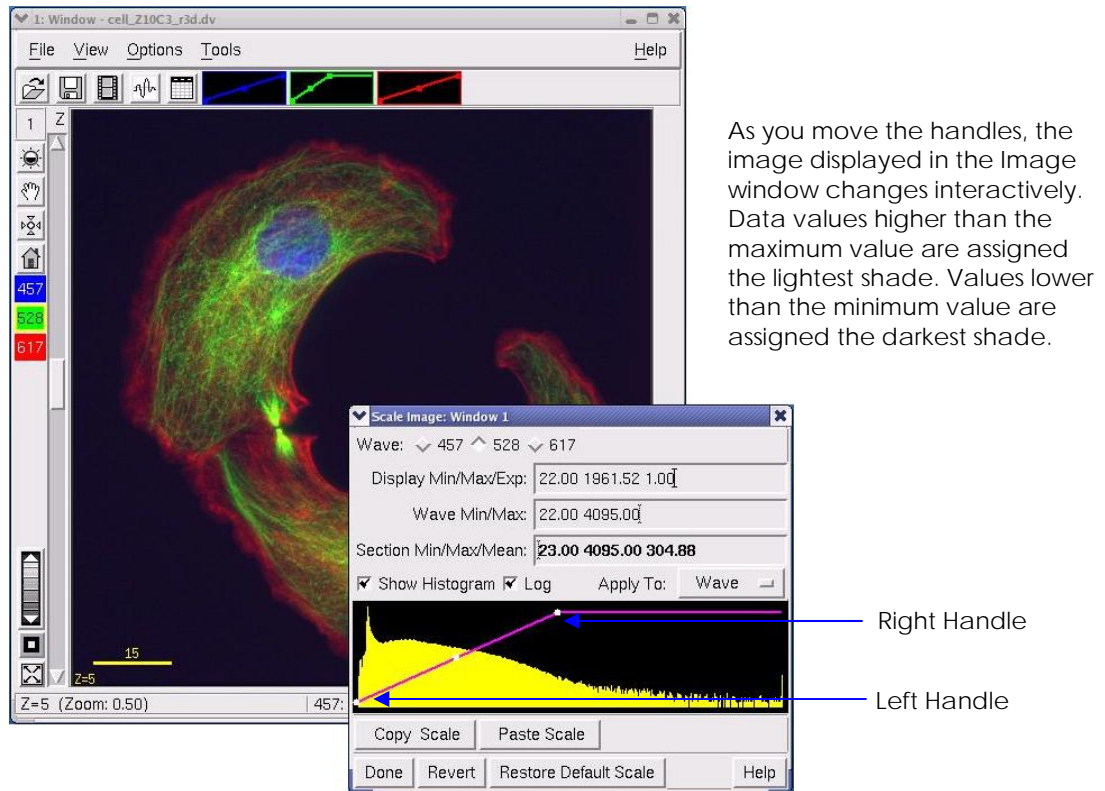


As you move the handles, the image displayed in the Image window changes interactively. Data values higher than the maximum value are assigned the lightest shade. Values lower than the minimum value are assigned the darkest shade.

Right Handle

Left Handle

The Left and Right handles change the scale range.


4. To slide the range to the left or right, click and drag on the Center handle.

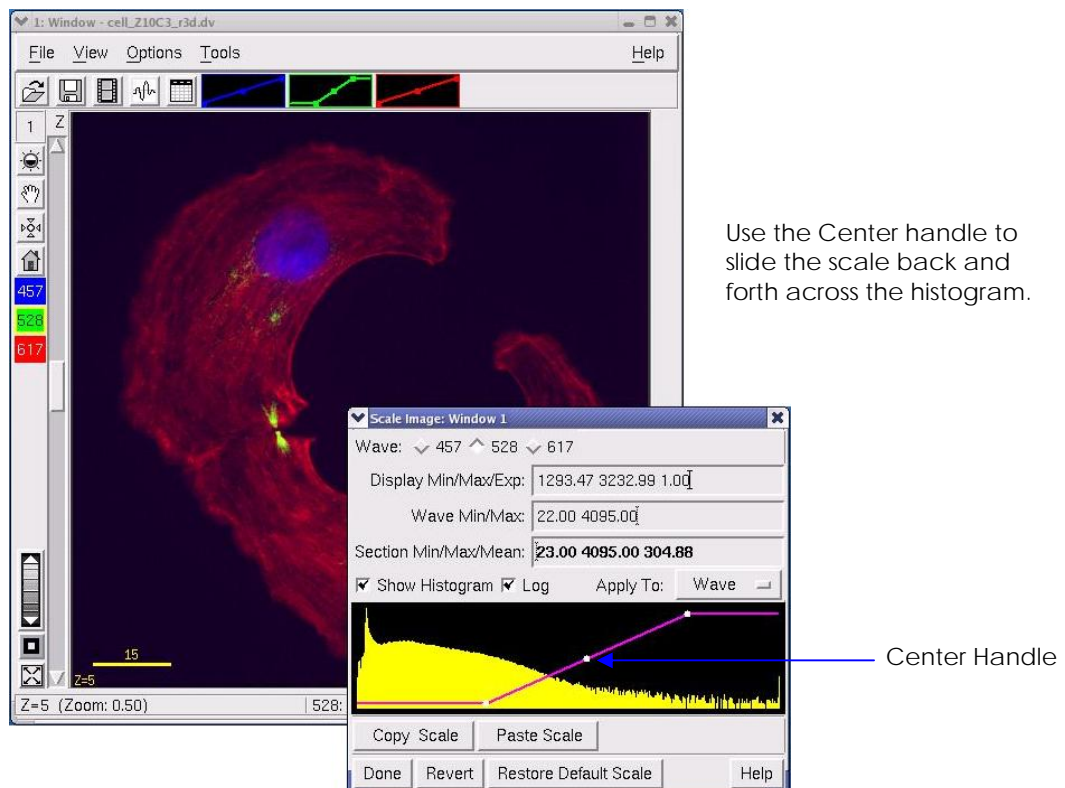

Use the Center handle to slide the scale back and forth across the histogram.

Center Handle

- To change the intensity scale distribution, click anywhere in the histogram (except on the handles) and drag the mouse up or down.

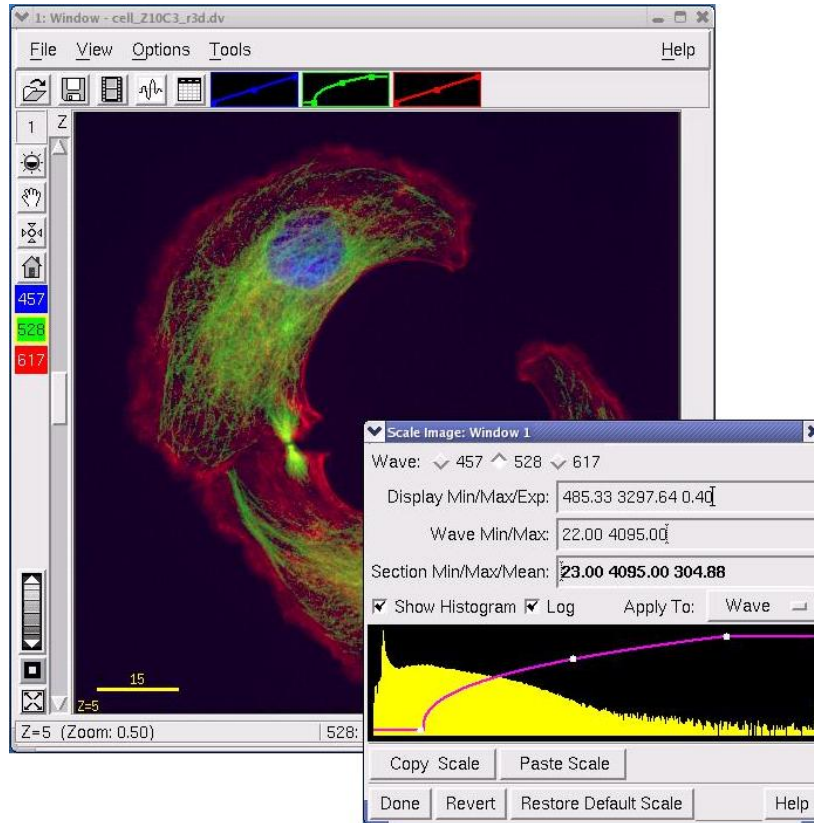


Dragging the mouse up or down in the histogram changes the scale distribution.

Tips

- #1 You can improve contrast at the low end of the intensity range by reducing the gamma value. To improve contrast at the high end of the range, increase the gamma value.
- #2 Another way to scale the image is to enter values into the Min/Max/Exp fields.
- #3 You can restore all of the default values by clicking **Restore Default Scale**.

To change the intensity scale for all open image windows:

- Select **View | Scale All Windows** from the softWoRx main menu. The Scale All Windows dialog box is displayed.

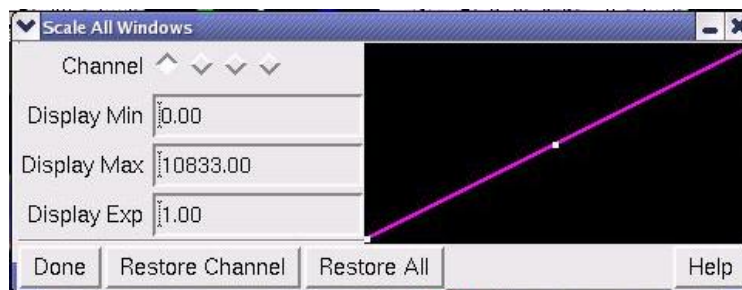

2. Similar to the Image Scaling dialog box, the right side of the Scale All Windows dialog box includes a line with three nodes (or handles) that graphically represent the intensity scale. Use the mouse to click and drag the white handles on the line as follows:
 - To change the minimum or maximum scale value, click and drag on the left or right handle.
 - To slide the scale range to the left or right, click and drag on the center handle.

As you move the handles, the images displayed in all of the currently open Image windows change accordingly. Changing the minimum and maximum intensity values changes brightness and contrast by mapping all 256 color shade values to a larger or smaller range of data. Data values higher than the maximum value are assigned the brightest shade. Data values lower than the minimum value are assigned the dimmest shade.

3. To change the intensity scale distribution, change the shape of the curve by clicking anywhere on the histogram (except on one of the handles) and dragging the mouse up or down.
 - To improve contrast at the low end of the intensity range, increase the slope of the curve at the left side of the graph.
 - To improve contrast at the high end of the range, increase the slope of the curve at the right side of the graph.

Changing the scale distribution increases the contrast at one end of the data range and decreases it at the other end.

4. Use the **Restore Channel** button to restore the selected channel to its original intensity scale.
5. Use the **Restore All** button to restore all channels to their original intensity scales.
6. Press **Done** when finished with the Scale All Windows dialog box.

Assigning Colors or Grayscale to Channels

You can view image data in grayscale or in two different color modes:

Grayscale Mode is useful for studying detail in a single wavelength. Because of the way the eye reacts to colors, you may be able to see more detail in Grayscale than in a Color mode. You can only view one channel at a time in Grayscale mode.

Color Mode can be used to visually compare intensities of two or three wavelengths. It also allows you to use the Volume Viewer with RGB opacity,

improve speed of volume rendered images, or save multi-channel *DeltaVision* images as a series of TIFF images.

Blended Color Mode allows you to overlay nonfluorescent data such as Differential Interference Contrast (DIC) data sets onto fluorescent data sets or to visually compare intensities of more than three channels simultaneously.


Grayscale Mode

You can switch between Color and Grayscale modes. Because of the way the eye reacts to colors, you may be able to see more detail in Grayscale than in Color. In general:

- Use Grayscale when you want to see more detail in a single wavelength of an image.
- Use Color mode when you want to visually compare intensities of two or more wavelengths.

To switch between Grayscale and Color mode:

- ▶ On the Image window, choose **View | Color** to toggle between Color and Grayscale.

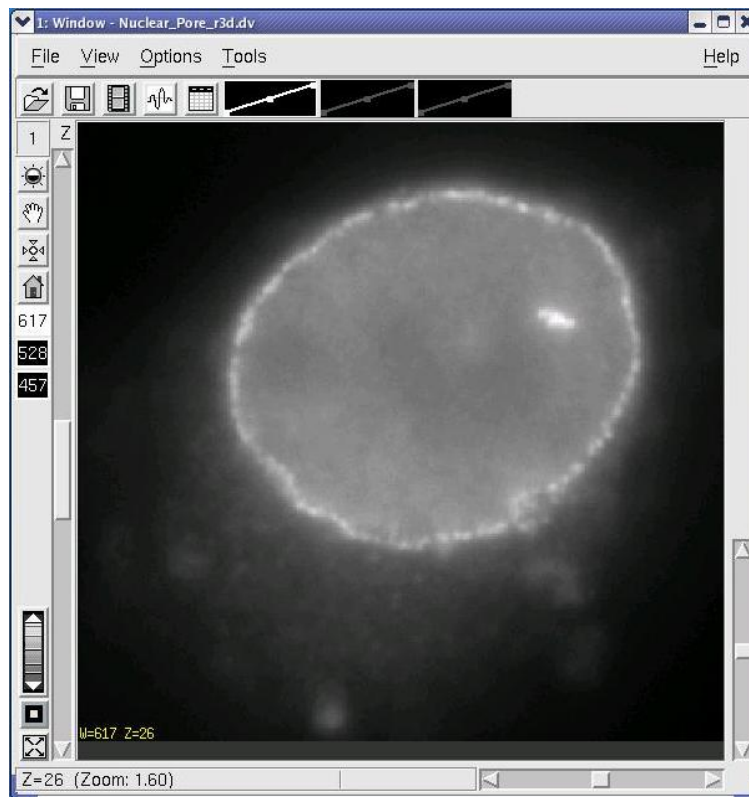


Grayscale shows more detail in a single wavelength

Color Mode

If you choose a basic color for each channel (red, green, or blue), you can display up to three channels in this mode. If you select any other colors (e.g., cyan, magenta, or yellow), the two colors used to create these mixtures are disabled for the other channels and they are turned off (Black). Colors such as Cyan, Magenta, or Yellow can be viewed in combination with only one other color.

To assign a basic color to a channel:

1. Choose **View | Select Image Colors** on the Image window menu to open the Select Image Colors dialog box.

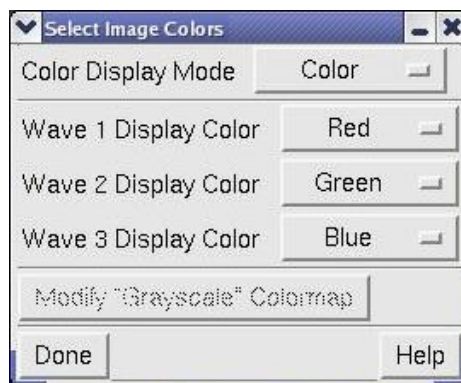

Set channel color options in Select Image Colors

2. Select the wavelength for each channel in the **Display Color** option lists. The new colors are displayed in the Image window as they are selected.
3. Click **Done** to set the colors that you selected.

Note You can assign basic red, blue and green colors to as many as three channels. You can assign other colors to two channels.


Blended Color Mode

In Blended Color Mode, you can assign any color to each channel. You can view up to five channels as separate colors. You can also assign grayscale as a color (this is useful for DIC data).


You can select an arbitrary color for each wavelength or you can specify to use the true color that is normally associated with each wavelength in the color spectrum.

To set Colors in Blended Color Mode:

1. On the Image window menu, choose **View | Blended Color** to set Blended Color mode.
2. Choose **View | Select Blended Colors** to open the Blended Colors dialog box.

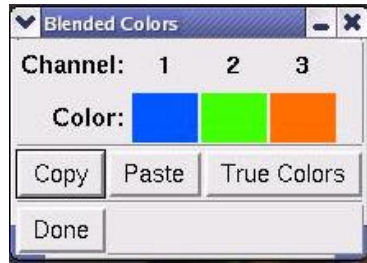

3. To specify the colors normally associated with a wavelength, click **True Colors**.
4. To specify a custom color for a channel, click the color under the channel to open the Choose Color for... dialog box and use the **Red**, **Green**, and **Blue** sliders to choose the color.

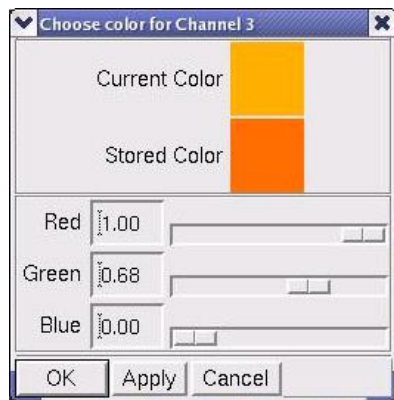

Choose Color for the selected channel

To assign grayscale to a channel in Blended Color Mode:

1. On the Image window menu, choose **View | Blended Color** to set Blended Color mode.
2. Choose **View | Select Blended Colors** to open the Blended Colors dialog box.
3. Click the color under the channel you want to modify and move all three color sliders all the way to the right to assign White as the Current Color, then press **OK**.


Controlling the Image Window Display

You can hide the Image window Display controls, toolbar, and scroll bars. You can also display and set a scale on your images. This is useful for preparing images for presentations.

Hiding or Displaying Image Window Border Tools

The border tools are the icons and controls on the left of the Image window. The toolbar is the set of icons above the Image window. These tools are displayed by default. You can hide them to capture a JPEG of the image.

Hiding border tools and the toolbar can focus the screenshot on your data.

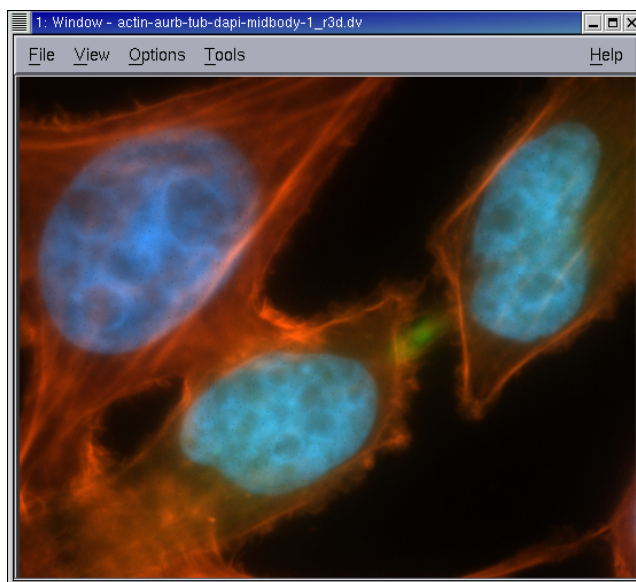

The Image window with the border tools and toolbar hidden

To switch border tools on and off:

- ▶ Choose **Options** on the Image window menu and display or hide the border tools as follows:

To display the tools, select the **Show Border Tools** toggle on the **Options** menu.

To hide the tools, clear the **Show Border Tools** toggle on the **Options** menu.

To switch the toolbar on and off:


- ▶ Choose **Options** on the Image window menu and display or hide the toolbar as follows:

To display the toolbar icons, select the **Show Toolbar** toggle on the **Options** menu.

To hide the tools, clear the **Show Toolbar** toggle on the **Options** menu.


The Image Window Scale Bar

You can display or hide a scale bar to show the scale of the image. You can also set options to move the scale bar (to show the scale of a point of interest in the image) or to control how the scale bar is displayed. The scale bar is displayed by default.

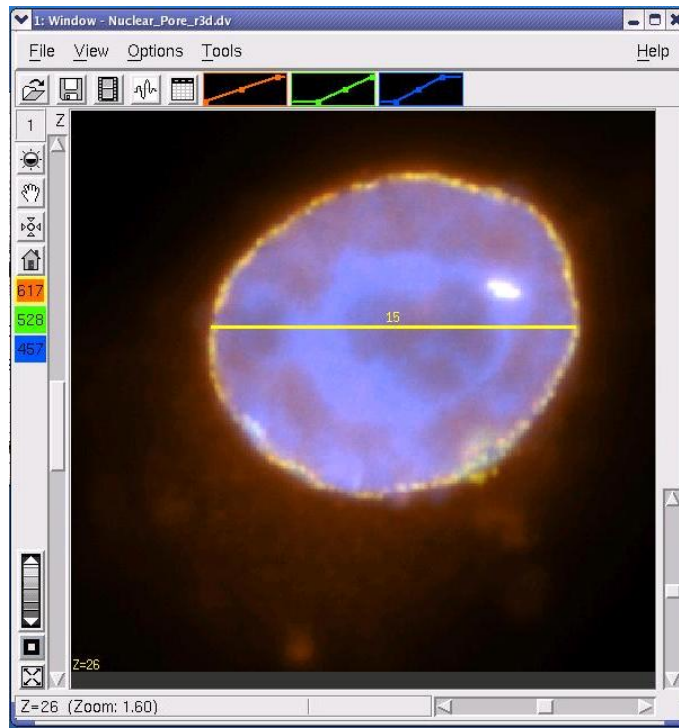

To change the image scale bar:

1. From the Image window menu, choose **Options | Display** to open the Display Attributes dialog box.

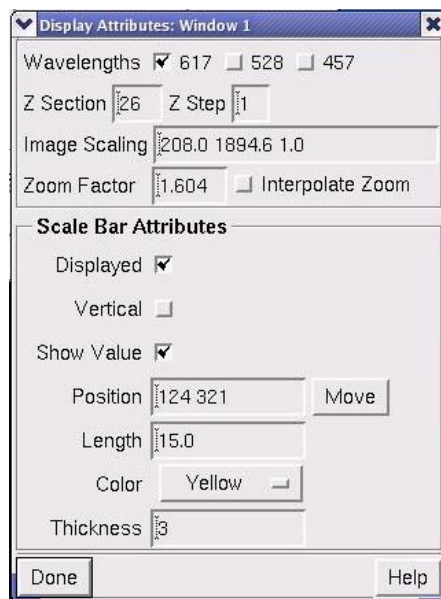

Set Scale Bar Attributes

2. Choose whether to display or hide the scale bar.

To display the scale bar, select the **Displayed** option.

To hide the scale bar, clear the **Displayed** option.

3. To display the scale as a vertical bar, select **Vertical** (the default is horizontal).
4. Use the **Show Value** option to select whether or not you want the scale bar value displayed.
5. Adjust the **Position**, **Length**, **Color**, and **Thickness** of the scale bar. The scale bar changes interactively as you set these properties.
6. Select the **Move** button and use the mouse to drag the scale bar to any position within the Image window.

Resizing or Reorienting an Image


You can use the Resample2D tool to resize an image or to reorient an image in X, Y, and Z directions.

Resizing an Image

When Resample2D magnifies an image, it interpolates values to add pixels to the image. When it reduces an image, it combines or eliminates pixels to create a subset of the original pixel data.


To resize an image:

1. Choose **View | Resample2D** from the *softWoRx* main menu to open the Resample2D dialog box.

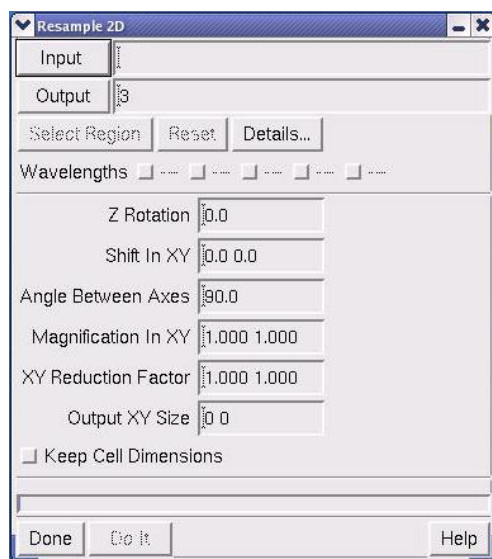


2. Enter the window number in the **Input** field.

- To resize part of the image, click **Select Region** and then drag the mouse across the area in the Image window.

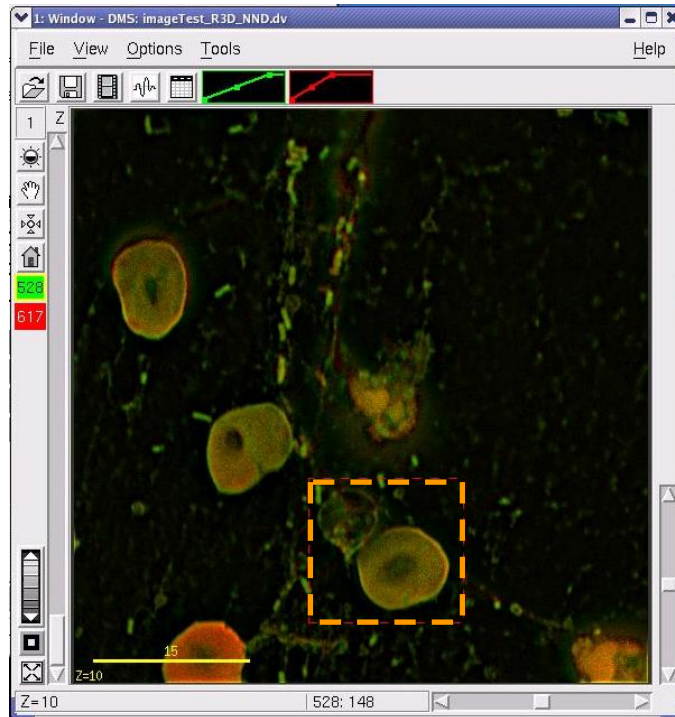


Select a region in the Image window

- To magnify the image, enter a magnification factor in the **Magnification in XY** field.
- To reduce an image, enter a reduction factor in the **XY Reduction Factor** field.
- Select **Keep Cell Dimensions** to keep the X and Y pixel size constant. (If unselected, a new size is calculated, based on the current settings for magnification.)

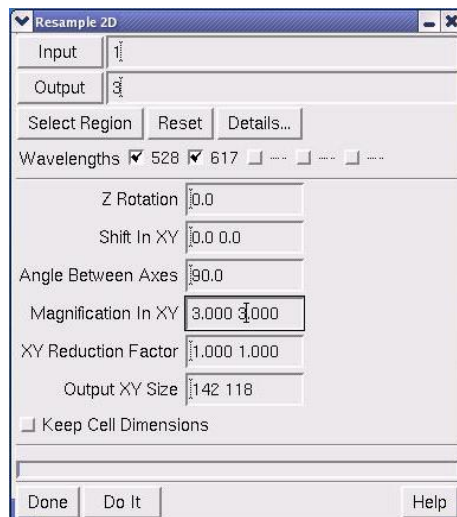

7. Click Do It.

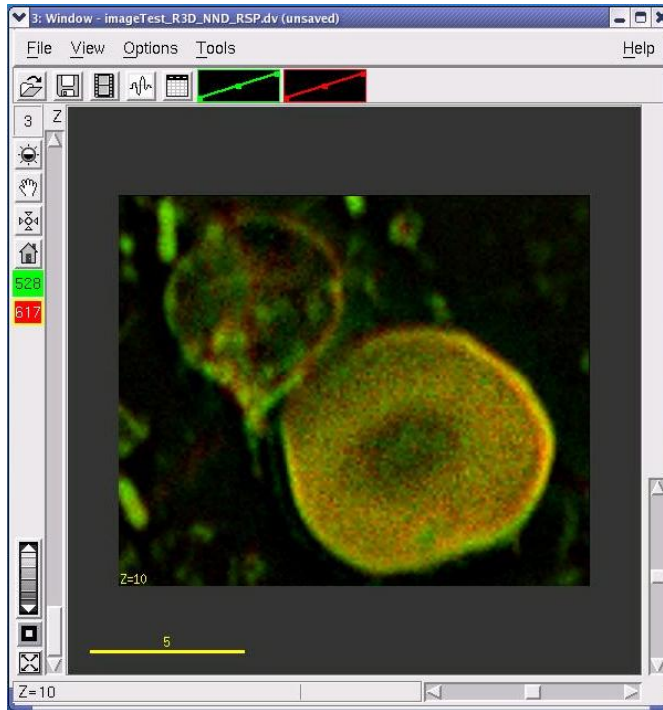

Rotating an Image

To rotate an image:


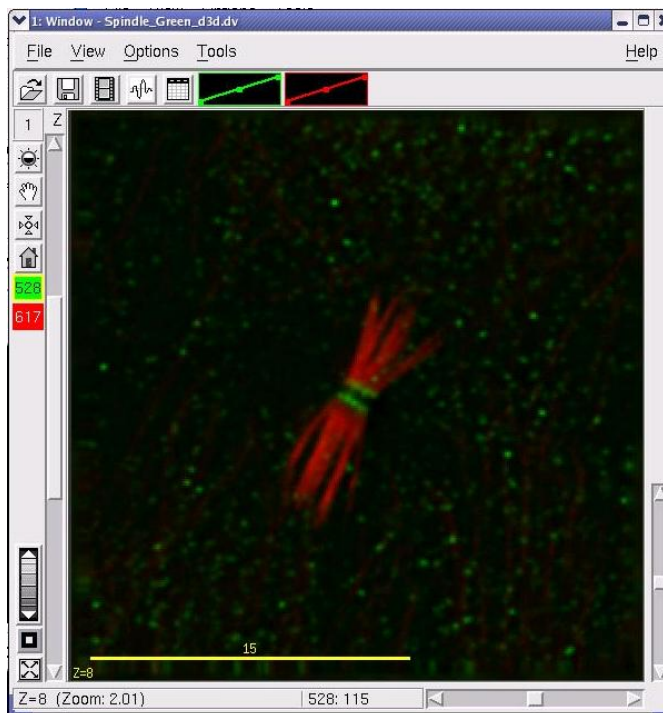


Image before rotation

1. Choose **View | Resample2D** from the *softWoRx* main menu to open the Resample2D dialog box.
2. Enter the window number in the **Input** box.
3. If you want to reorient part of the image, click **Select Region** and then drag the mouse across the area in the Image window.
4. Select the wavelengths that you are interested in.
5. Enter the angle to rotate the image in the Z direction in the **Z Rotation** field.
6. Enter the distance to shift the data in the **Shift in XY** field.
7. Enter the angle to rotate the data on the XY plane in the **Angle Between Axes** field.
8. If you want to maintain the same relative pixel size in the rotated image, activate the **Keep Cell Dimensions** checkbox.

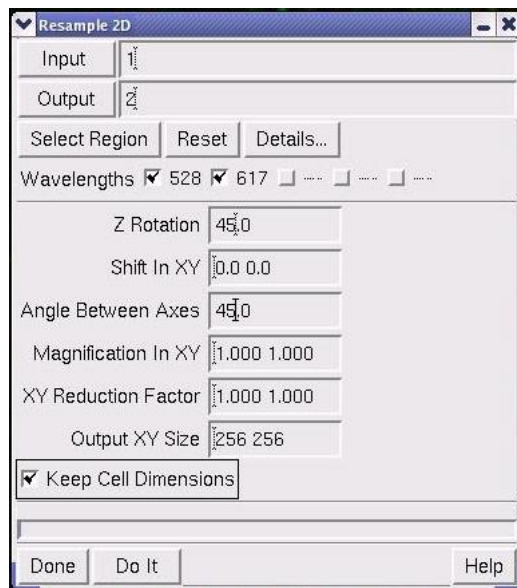

Enter the angle to rotate the axes

9. Click **Do It**.

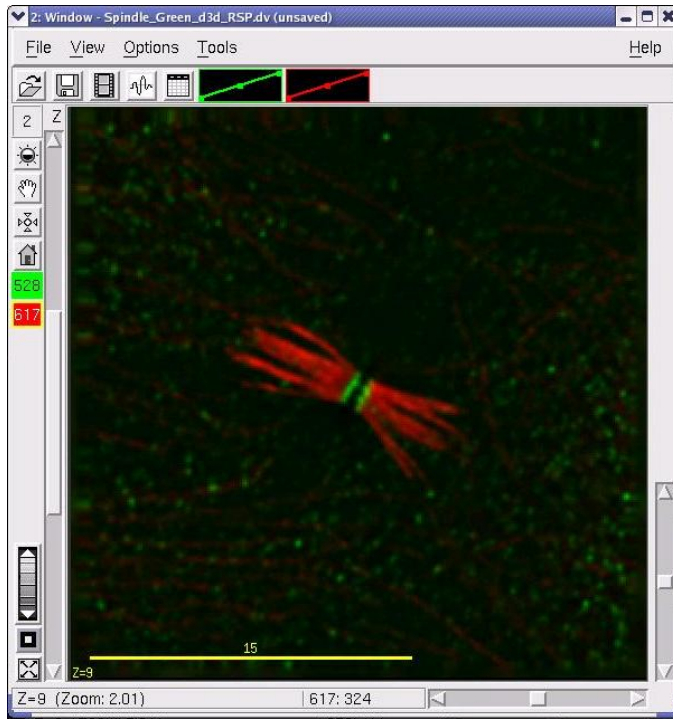


Image after rotation

To reorient images:



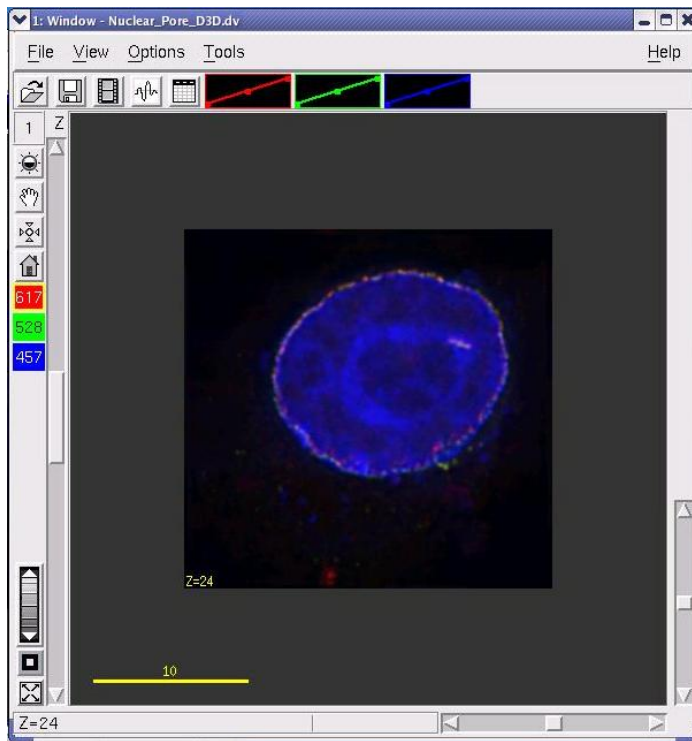


Image before reorientation

1. Open the Rotate3D dialog box by choosing **View | Rotate3D** on the main *softWoRx* menu.

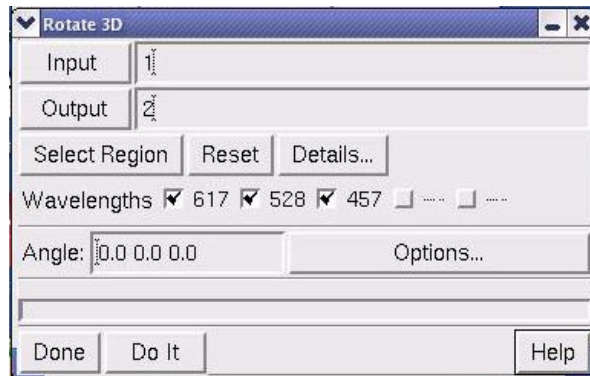

2. Enter the Image window number in the **Input** field.
3. Specify the X, Y, and Z rotation angles to apply to the data in the **Angle** field. When viewing an image in a window, a positive X rotation rotates the top of the image towards you, and the bottom away from you. A positive Y rotation rotates the right side of the image towards you and the left away from you. A positive Z rotation indicates counter-clockwise rotation.
4. To modify the output size that *softWoRx* creates from the rotation angle and the dimensions of the input image, select **Options** from the Rotate 3D dialog box. The Rotate Options dialog box is displayed. Enter the new dimensions in the **Output Dimensions** field.

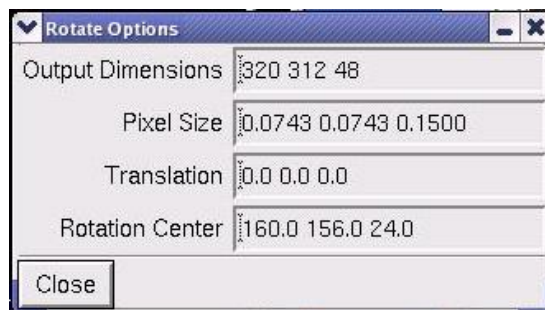

5. Specify an X,Y,Z vector for translating the image (pixels) in the **Translation** field. The size of each pixel in real-world coordinates (usually microns) is displayed in the **Pixel Size** field.
6. Specify a center point about which the rotation occurs in the **Rotation Center** field. The **Rotation Center** is by default the center of the image. You can specify a different center point (in pixel coordinates).
7. Click **Do It** to generate the image.

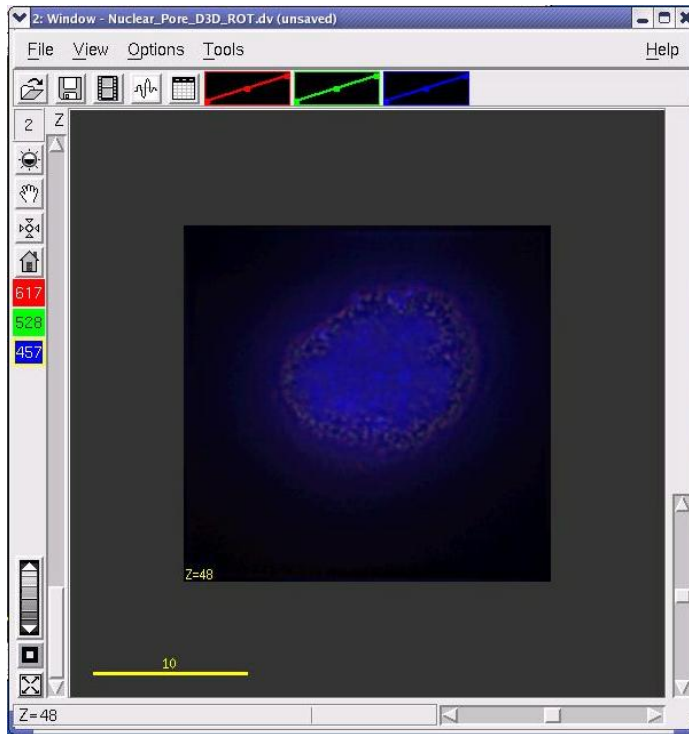



Image after reorientation

Viewing Cross Sections

You can view cross sections of the data by creating orthogonal projections. The Orthogonal Viewer allows you to interactively view YZ and XZ plane cross sections.

To view YZ or XZ cross sections:

1. Open the image in the desired Image window.
2. Open the Orthogonal Viewer by choosing **Tools** | **Orthogonal Viewer** from the Image window. The orthogonal projection is displayed. The new Image window displays the original image. It also displays an XZ projection (at the bottom of the window) and a YZ projection (on the right side of the window). Projection lines show the areas of the image that are displayed on the cross sections.

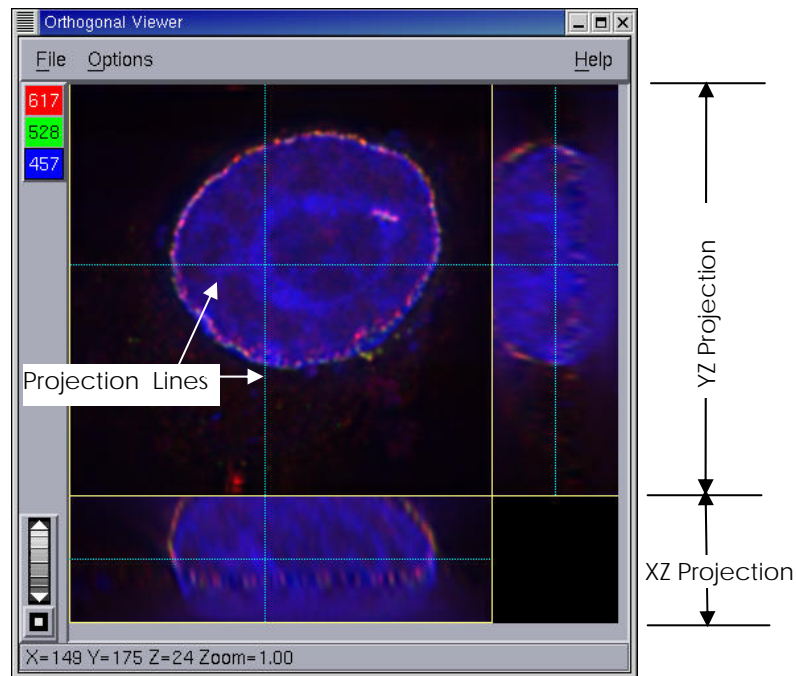

3. To orient the image in "real world" coordinates, select **Options** on the Orthogonal Viewer, and then select the **Cover Slip at Bottom/Right** toggle. This orients the image so that the display in the window represents the orientation of the sample when the data was collected (e.g., on the XZ projection, the cover slip is down).
4. To change the cross sections that are displayed on the projections, use the mouse to drag the projection lines across the Image window.

8. Viewing Movies

Movies greatly enhance the analysis of certain types of image data. When used on a volume rendering, a movie shows the relationships between objects in 3-D space. When used with a time-lapse data file, a movie allows you to visualize the course of events captured in the study. You can also use movies to trace particles in time-lapse data.

In This Chapter

Viewing Volumetric or Time-Lapse Movies	113
Tracking Particle Movement with Trails Movies.....	114

Viewing Volumetric or Time-Lapse Movies

At least one Image window must be open in order to apply the movie function. There are two ways to access the movie function: from the *softWoRx* main toolbar and from the Image window. Accessing the movie function from the *softWoRx* toolbar applies the movie function to all open windows. Opening the movie function from an Image window applies it only to that Image window.

To view one image data set as a movie:

1. Open the file that you want to view as a movie in an Image window.
2. Click **View**.
3. Click **Movie**.
4. Unselect the **Paused** check box.

To view two or more image data sets as movies:

1. Open the files that you want to view as movies in Image windows.
2. Click **View** on the *softWoRx* main toolbar.
3. Click **Movie**.
4. Unselect the **Paused** check box.

To view time-lapse Z series data:


1. Open the file that you want to view as a movie in an Image window.
2. Click **View**.
3. Click **Movie**.
4. Unselect the **Paused** check box.
5. Select the **Time Lapse** check box.
6. Adjust the **Manual Z Control for Time Data** slider.

Tracking Particle Movement with Trails Movies


You can trace the movement of particles in time-lapse data with the Trails Movie tool.

To trace particle movement:

1. Open an image that contains time-lapse data in the Image window.

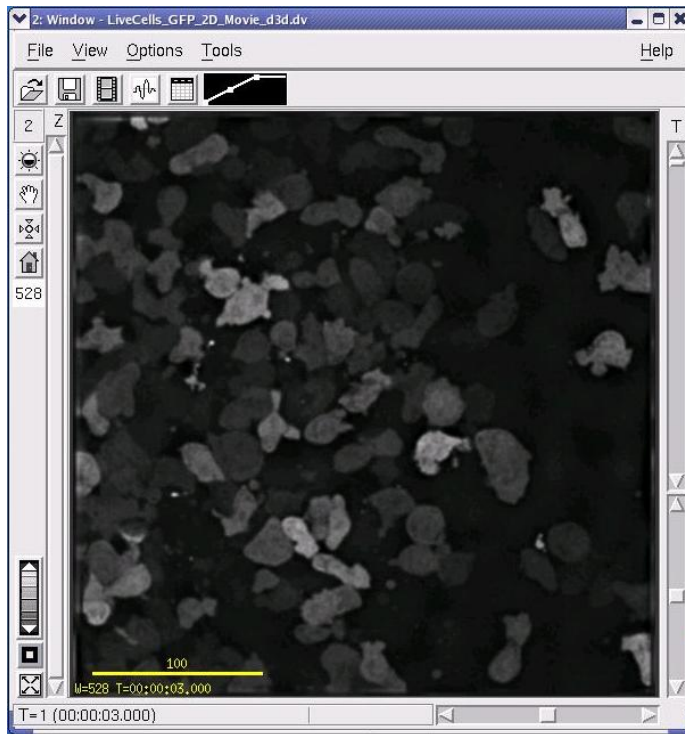


- From the main *softWoRx* window, choose **View | Trails Movie**. The Trails Movie dialog box is displayed.

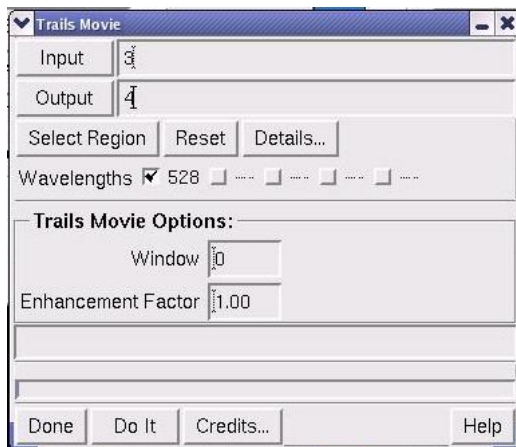


- Drag the Image window number into the **Input** field.
- Click **Select Region** to select a region of interest. Then select which wavelengths to include in the movie.
- Enter the length of the trace history in the **Window** field (e.g., for a value of 5, the trace includes the previous 4 time frames and the current time frame).
- To emphasize the display of the current point, enter a value greater than 1.0 in the **Enhancement Factor** field (typical values for this field are 1.0 - 2.0).

Note The displayed intensity value of each point is a weighted average of the corresponding points in the previous time frames. The previous points all have a weight of 1.0. The Enhancement factor is assigned as the weight for the intensity of the current value.

- Click **Do It** to create the trails display.

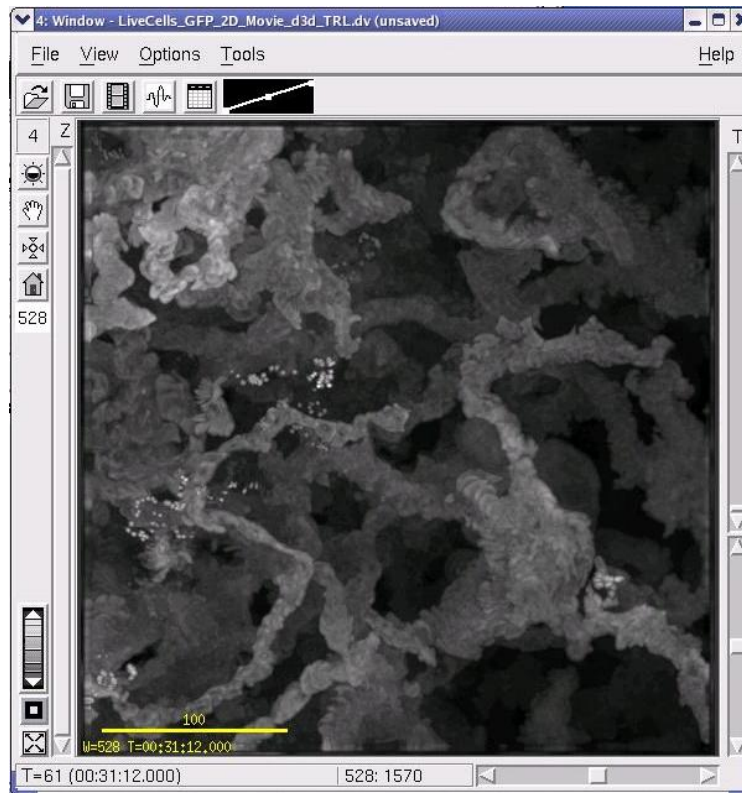

- To create a Trails movie, choose **File | Save As Movie** on the Image window and save the movie. The Save As Movie dialog box is displayed.

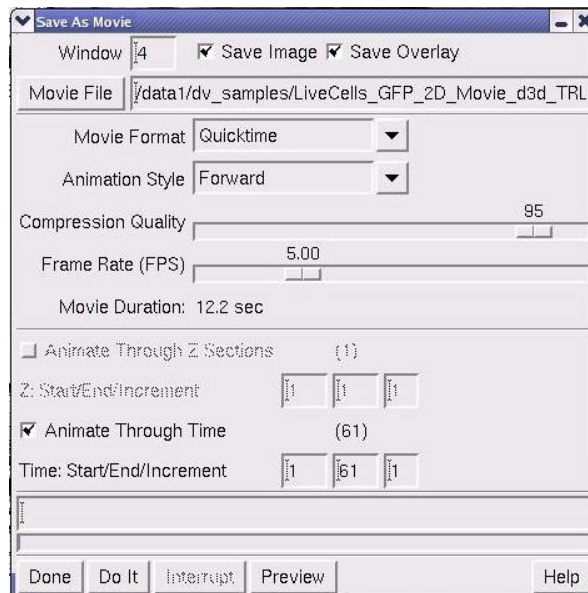

9. From this dialog box, you can use the appropriate fields to select the movie format (**Quicktime**, **AVI**, or **MPEG**), the animation style (**Forward**, **Backward**, or **Forward and Back**), the compression quality, the frame rate, and the time increment to use. You can also select whether to animate the movie through Z sections or time.

9. Viewing Projections and Volumes

You can create two types of data projections of multiple Z sections:

Two-dimensional projections can help you to visualize how the data are oriented in the XY direction. These projections allow you to view the paths of individual fibers, chromosomes, or other types of linear data.


Volume projections can help you to understand the three dimensional nature of the data.

In This Chapter

Creating 2D Projections.....	119
Creating Volume Projections.....	122


Creating 2D Projections

Use the Quick Projection tool to quickly combine information from multiple Z Sections. Averaging all of the sections into one provides an approximation of a volume rendering of the image looking directly down the Z axis.

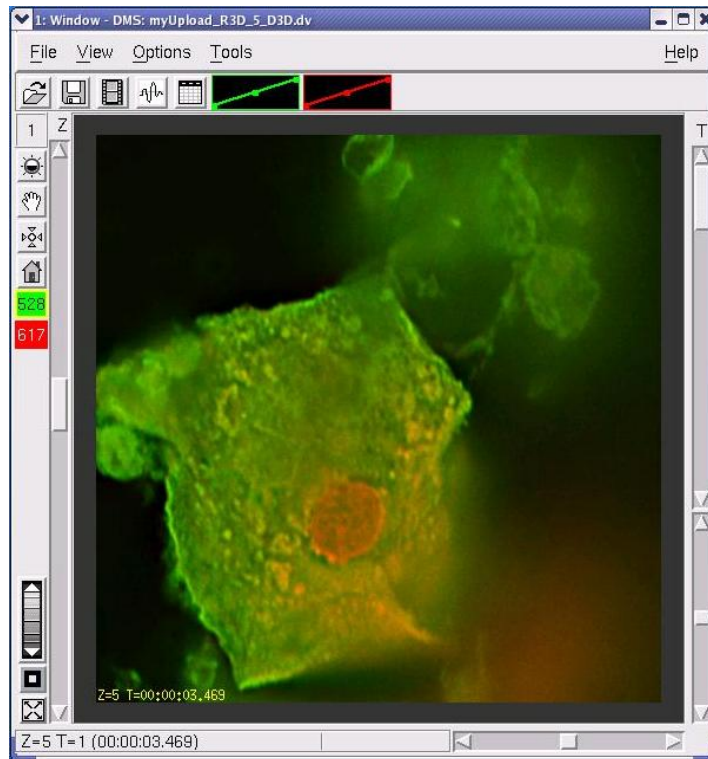


A multiple Z section image before projection


1. Choose **View | Quick Projection** from the main *softWoRx* menu to open the Quick Projection dialog box.

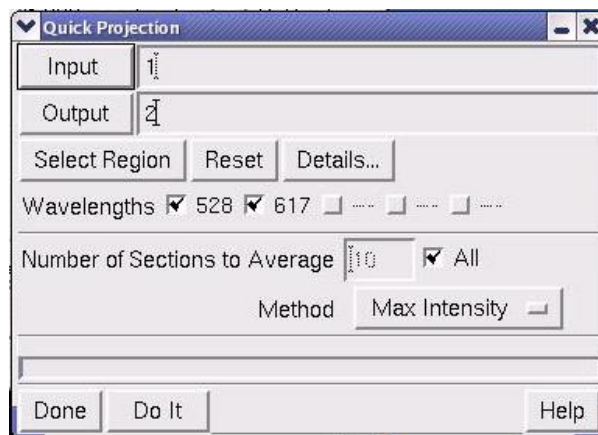

2. Enter an image file name or window number in the **Input** field.
3. If you want to include only selected data, click **Details** to open the Region Details dialog box. Then specify the ranges of data that you want to include in the **Output Options** fields.

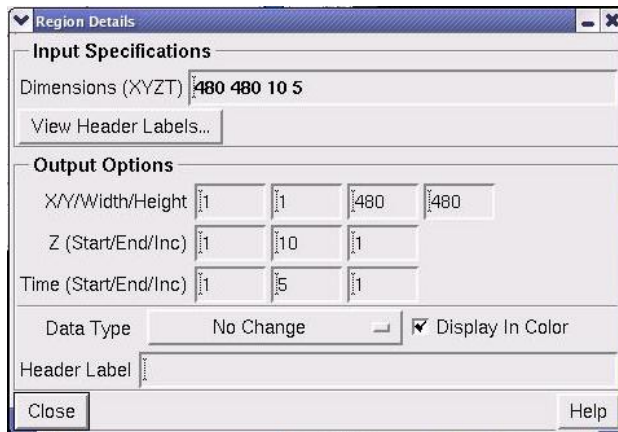

4. In the Quick Projection dialog box, select which wavelengths to include.
5. In the Quick Projection dialog box, choose how to group sections as follows:
 - To group sequential sets of sections into output sections, specify how many input sections to average for each output section in the **Number of Sections to Average** field.
 - To group all of the sections into one section, select **All**.
6. Choose one of the following ways to combine the sections in the **Method** list:
 - To add the intensity of each pixel to create the output values, choose **Sum**. (Be careful when using this option. If the output intensity values are too large for the output data type specified, the output image will appear to be saturated.)
 - To average the input data values to create the output image, choose **Average**.
 - To use the largest intensity value of all the input intensities to create an output value, choose **Max Intensity**. (This method may give you the most realistic representation of a volume rendered image, especially when combining all of the images in the input data set.)
7. Click **Do It**. The projected image is displayed in a new Image window.

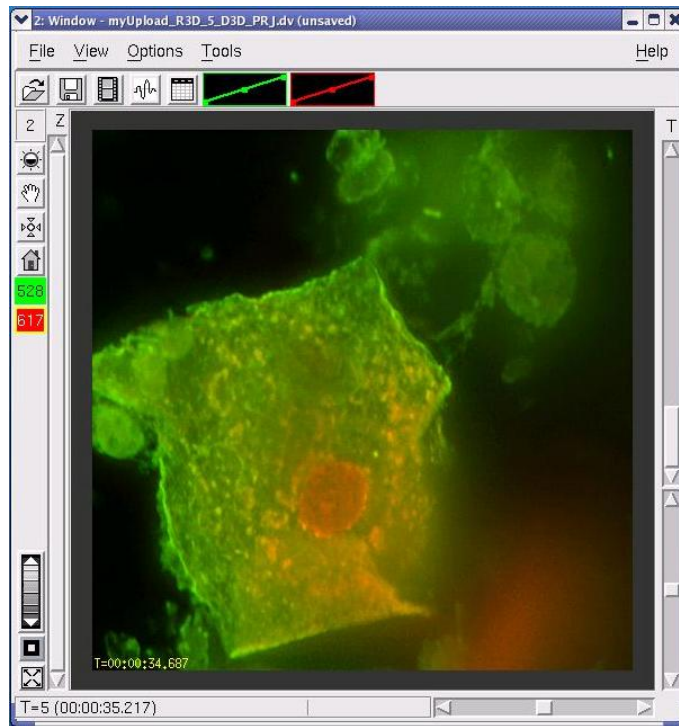

A single Z section image after projection


Creating Volume Projections

The Volume Viewer provides you with the ability to view the image data in 3-D. This tool allows greater visual understanding of the image data and comparison of features within the image data. It also allows quantitative assessment of structures throughout the entire data set on a single image.

Volume Rendering

A brief understanding of how *softWoRx* creates a volume rendering will help you utilize this tool for various image data sets. Theoretically, a set of parallel rays is sent through the data set at various angles to analyze the data in those paths and collect new data from that perspective. Each time a set of rays is passed through the data, a projection is created from the resulting data. These projections constitute the volume rendering.

Figure 4. Volume Rendering Projections

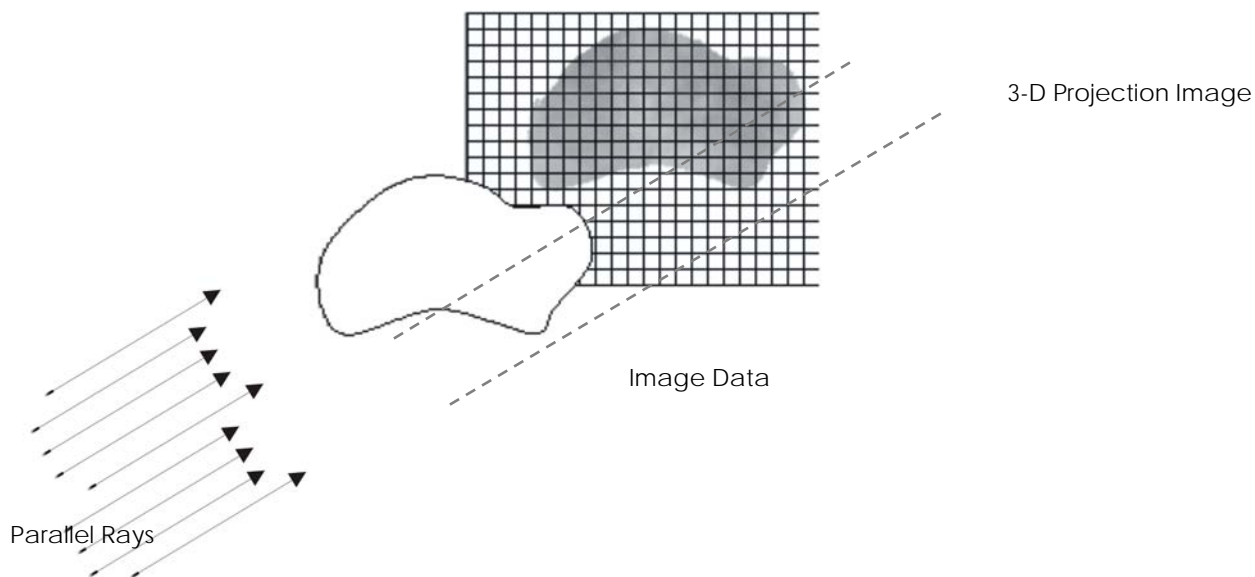


Axes of Rotation

Volume Viewer enables you to create a movie of the data rotating around an axis. This axis of rotation can be any of the three common axes, X, Y, or Z.

Given a Z series of data, the coordinates are defined as shown in the following figure.

Figure 5. Axes of Rotation

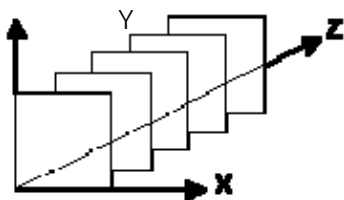

Rotations about the axes are as shown in the following figures.

Figure 6. Rotation About the X Axis

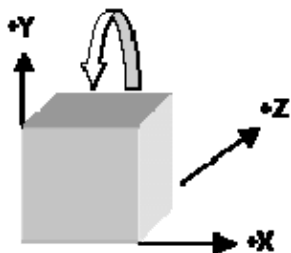


Figure 7. Rotation About the Y Axis


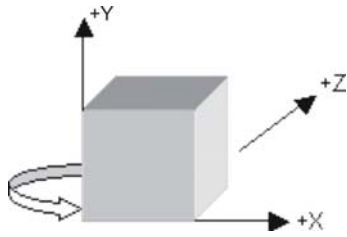


Figure 8. Rotation About the Z Axis

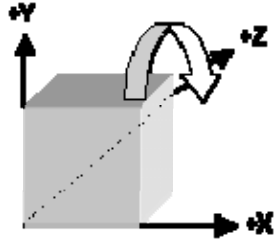






According to the parameter settings, *softWoRx* enables you to see the desired portion of the Z section, in the desired wavelength, using a variety of methods. Options in Volume Viewer help you to perform a volume rendering.


The most important options are **Select Region**, **Details**, and **Method**. By limiting the size of the data set using **Select Region** and **Details**, the time needed to create projection images can be drastically reduced.

For detailed information about each of Volume Viewer options, refer to the online Help.

Methods for Projecting Volumes


The following table summarizes the six methods of data collection.

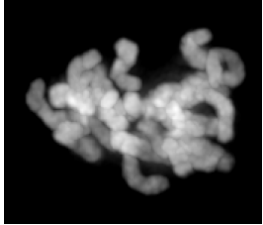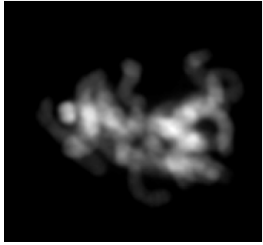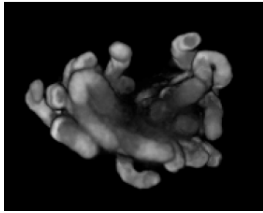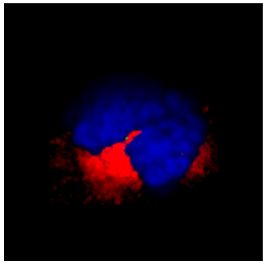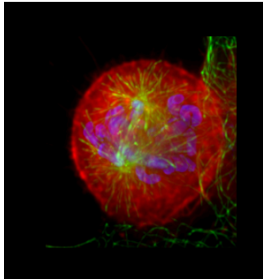
Method	Functionality	Application
Maximum Intensity 	Each ray collects the data from the voxel with maximum intensity.	Best choice for showing internal detail of a translucent image.
Additive 	Each ray collects and sums data from all the voxels in its path and scales it down to an appropriate intensity between 0-255.	Generates quantitative projections. This data can be used for comparison of intensity in various structures within the image data.
Progressive 	Each ray collects the data in the voxel that is closest to the front of the image.	Clearly displays opaque features within an image. Works well for objects whose internal details are not needed, such as metaphase chromosomes.
RGB/Opacity 	If the image data has been processed using Blend Colors, with the Opacity option toggled on, then this data can be used to form a volume rendering in multiple wavelengths. The resulting image captures the positional information and relationship among all the wavelength data sets.	Realistically renders a volume in a multi-wavelength image. Clearly relates the data points of the various wavelengths.
Mixed 	Different methods (maximum intensity, additive, and progressive) can be assigned to different wavelengths.	One wavelength contains diffuse, cloudy features and is set to the maximum intensity method and another wavelength has opaque features and the method is set to progressive.

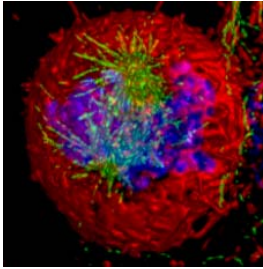Method	Functionality	Application
VolPack 	Generates images that use lighting techniques to highlight surfaces in the 3D rendered image.	Substantially faster than the other methods supported, although the method may not be optimal for all image types.

Note VolPack uses libraries obtained from the Stanford Computer Graphics Laboratory. It is an implementation of the shear-warp volume rendering algorithm as described in *Lacroute, P. and Levoy, M., Fast Volume Rendering Using a Shear-Warp Factorization of the Viewing Transformation, Proc. SIGGRAPH 1994, ACM.*

To render a volume:


1. Open an image file. Click  to open the intensity scaling dialog box and subtract background from the image by setting the left control point at the middle of the first intensity peak.
2. Choose **View | Volume Viewer** in the *softWoRx* main menu to open the Volume Viewer dialog box.

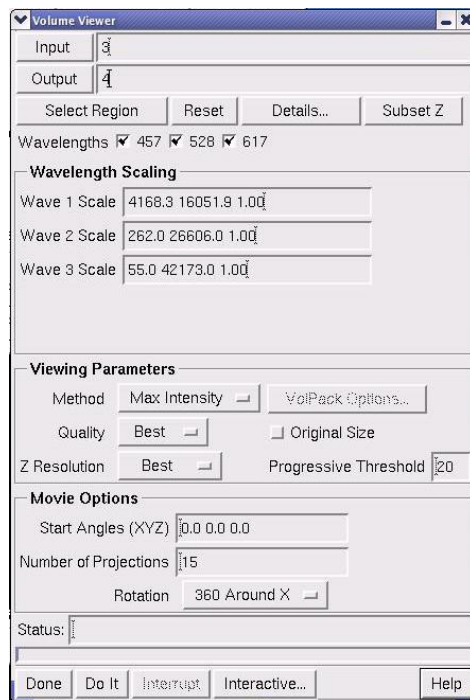


Use Volume Viewer to render volumes

3. Drag the Image window number into the **Input** field.

4. Click **Select Region** and drag the mouse across the region of interest to select a rectangular region. Scroll through the Z sections to insure the region of interest includes all the image data that you want to include in the data set.

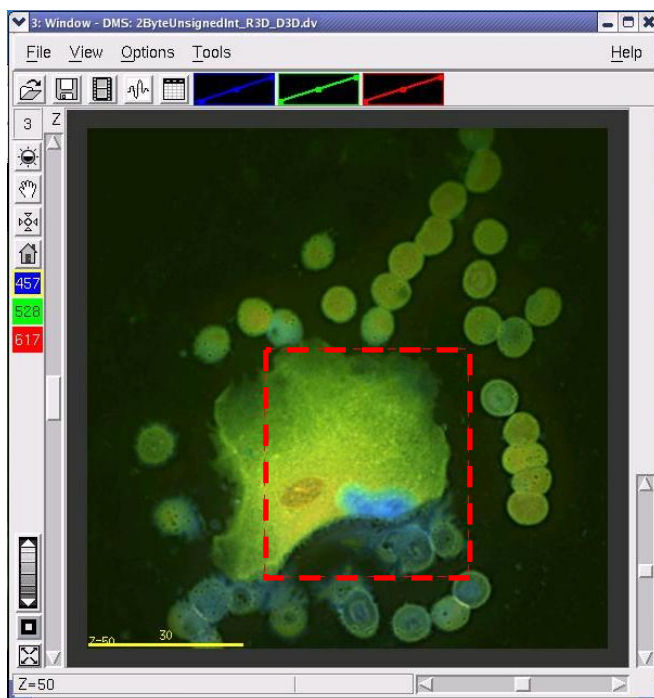

5. Specify the Viewing Parameters and Movie Options.
6. Click **Do It** to process the image data.

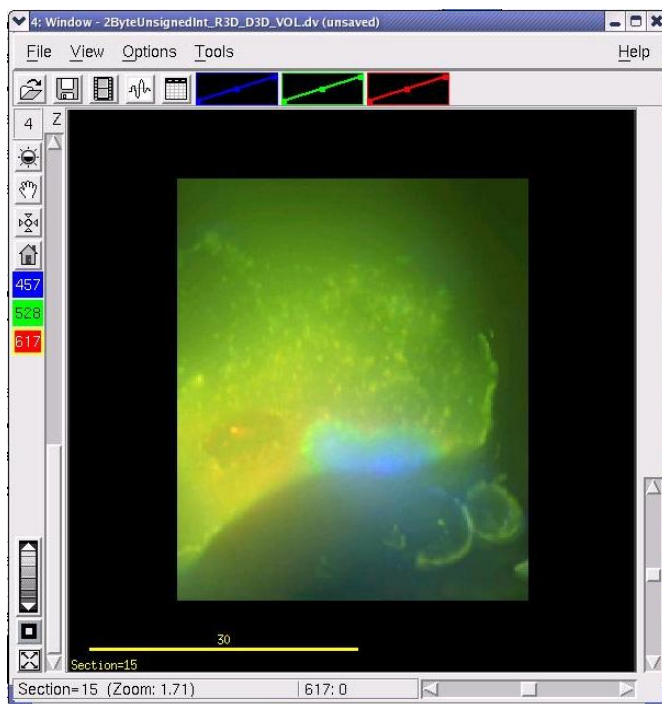


Volume created with the Max Intensity method

7. Move the Z slider up and down to view the projections or use the Movie tool.

To rotate the image using interactive image rotation:

1. Open the image file and choose **View | Volume Viewer** in the *SoftWoRx* main menu to open the Volume Viewer dialog box.

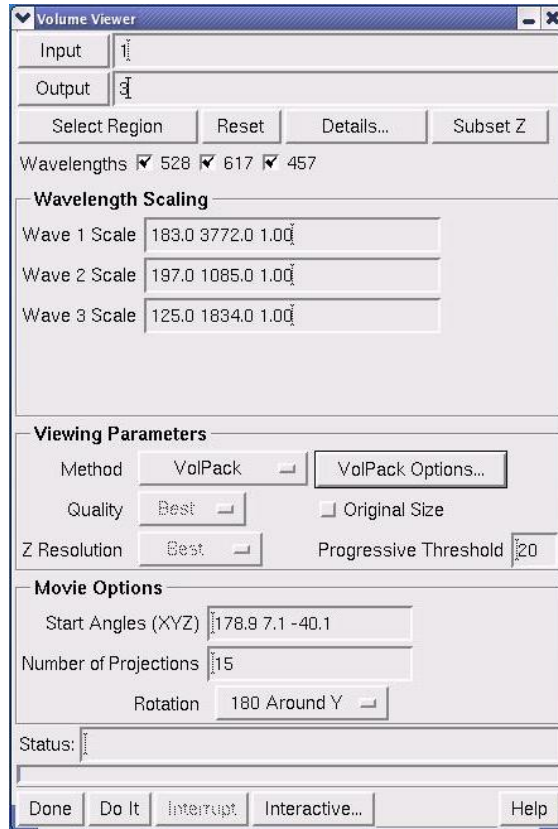

2. Drag the Image window number into the **Input** field.



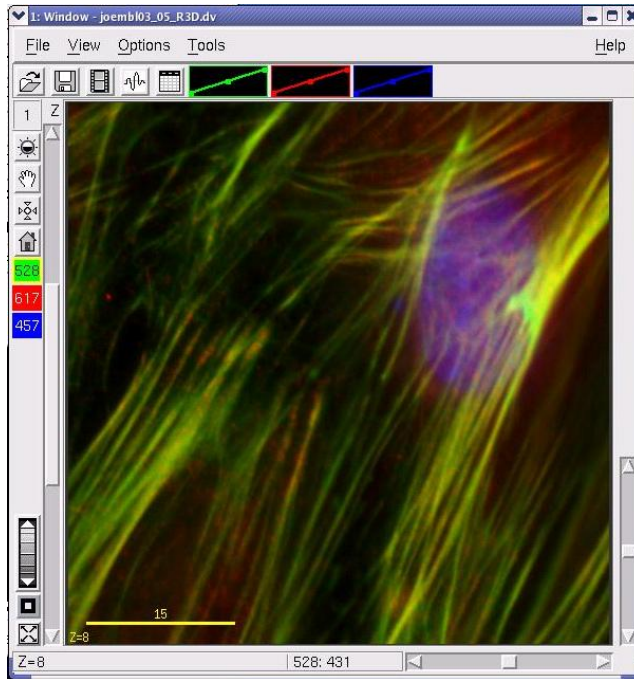


Image before interactive rotation


3. Click **Select Region** and drag the mouse across the region of interest.

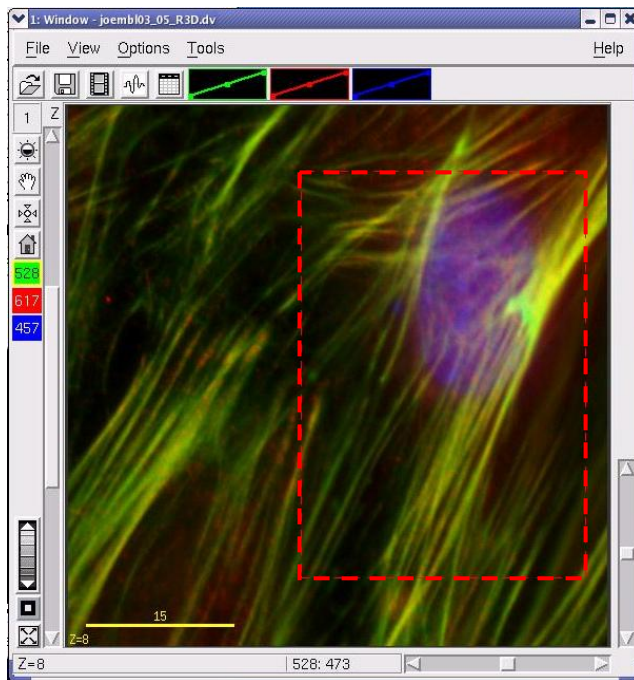


4. Click **Interactive** to open the Interactive Volume Viewer Parameters dialog box.

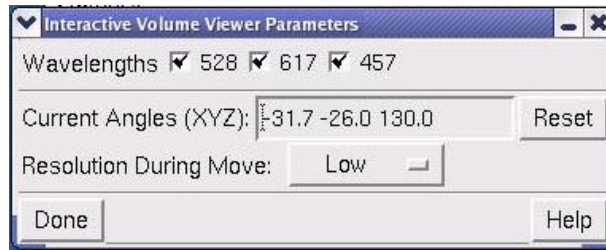

5. Choose which wavelengths to display in the window in the **Wavelengths** options.

 **Tip** VolPack is the preferred method for interactive viewing. With VolPack, you can view up to three wavelengths at a time. To use VolPack, select this option under Viewing Parameters in the Volume Viewer dialog box.

6. Select **Low** in the **Resolution During Move** option.
7. Drag the cursor on the image to rotate the image to the desired orientation.

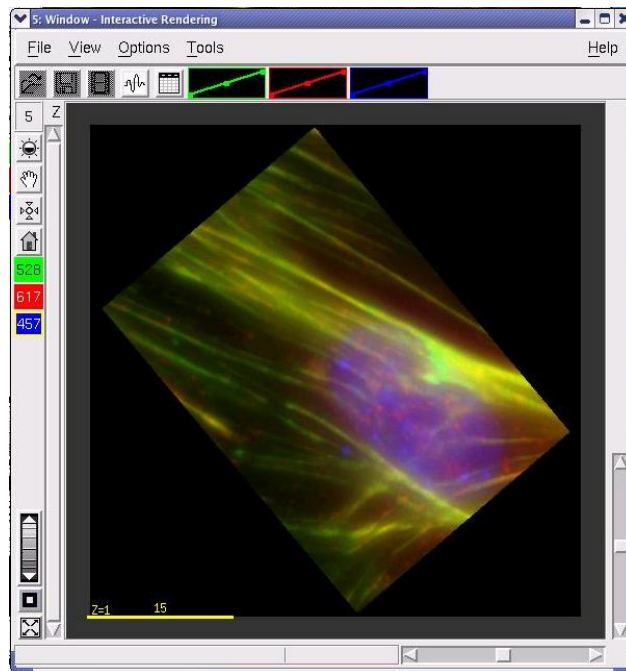

8. Click **Done** in the Interactive Volume Viewer Parameters dialog box.
9. Click **Do It** in the Volume Viewer to render the volume with the new orientation.


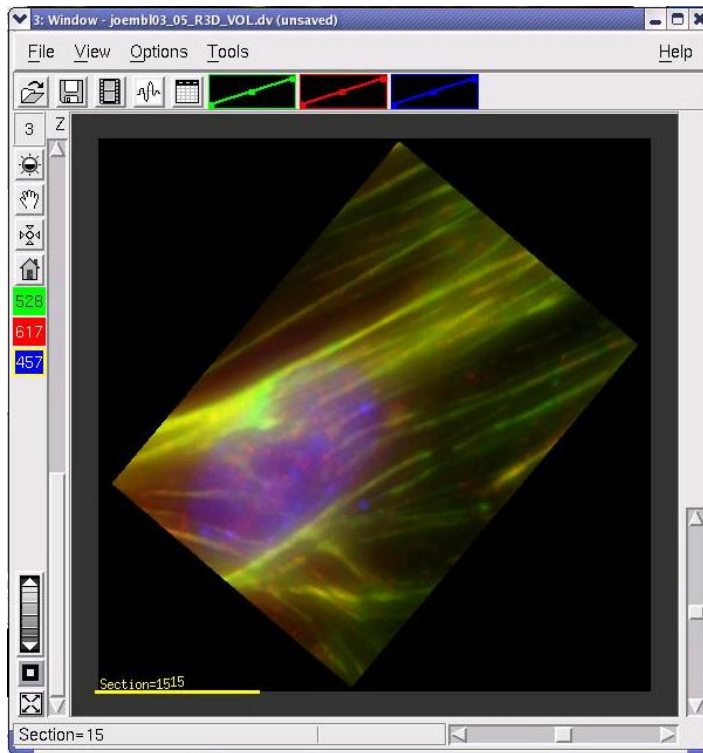



Image after interactive rotation

To render a volume using the RGB Opacity Method:


1. Open the desired image file.

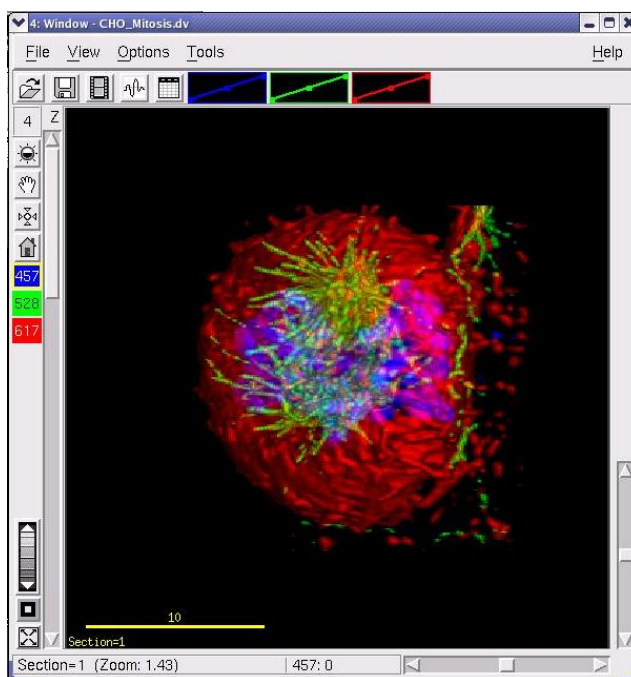


- Choose **View | Blend Colors** in the *softWoRx* main menu. The Blend Colors dialog box is displayed.

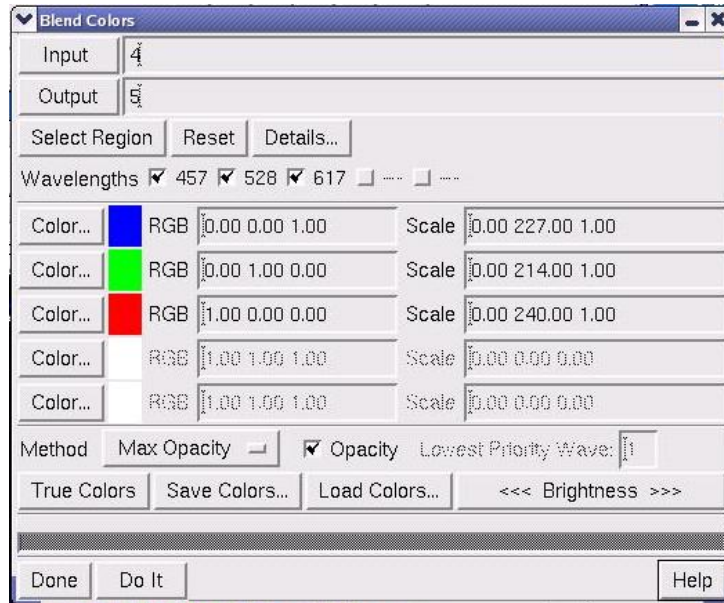


- Drag the Image window number to the **Input** field.
- Select **Max Opacity** in the **Method** option and select the **Opacity** option.
- Click **Do It** to create a blended color image.

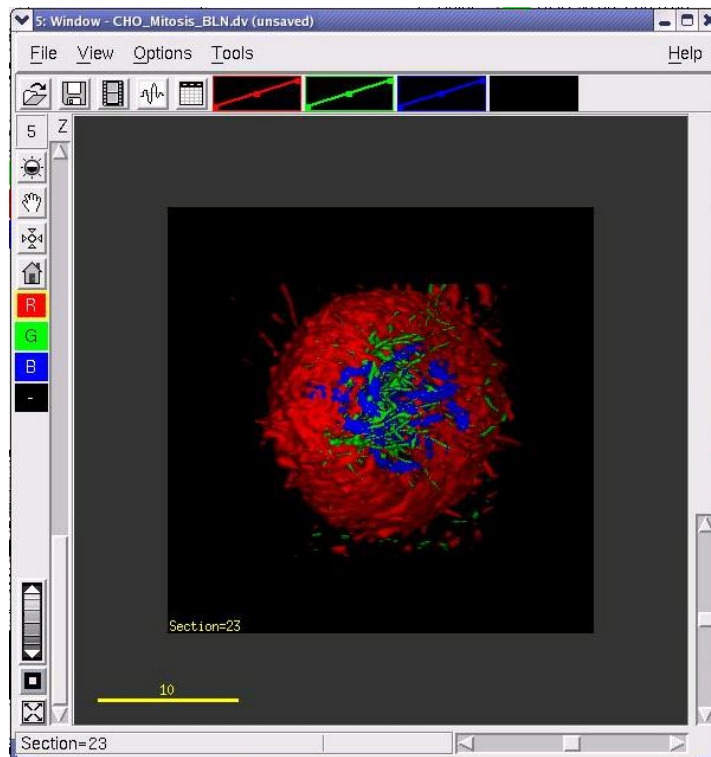


- In the *softWoRx* main menu, select **View | Volume Viewer** to open the Volume Viewer dialog box.

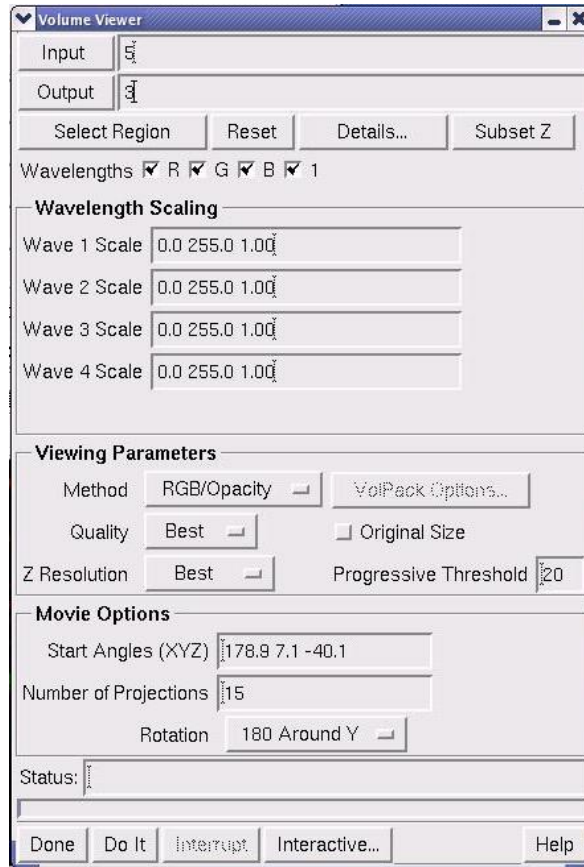

- Drag the Image window number of the blended image to the Input field in the Volume Viewer dialog box.
- To select which data to view, click **Select Region** and drag the mouse across the Image window. (The area inside of the rectangle that is displayed as you drag the mouse is the region of interest.) Scroll through the Z sections to insure the region of interest includes all the image information you want to include in the data set.

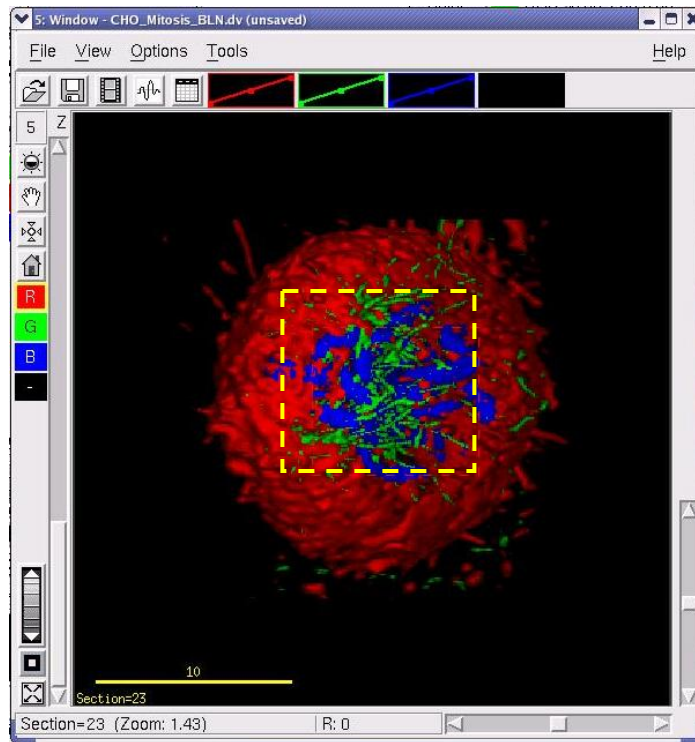

9. In the Viewing Parameters section of the Volume Viewer dialog box, select **RGB/Opacity** as the **Method** option and adjust the Movie Options.
10. Click **Do It** to create the RGB opacity volume.

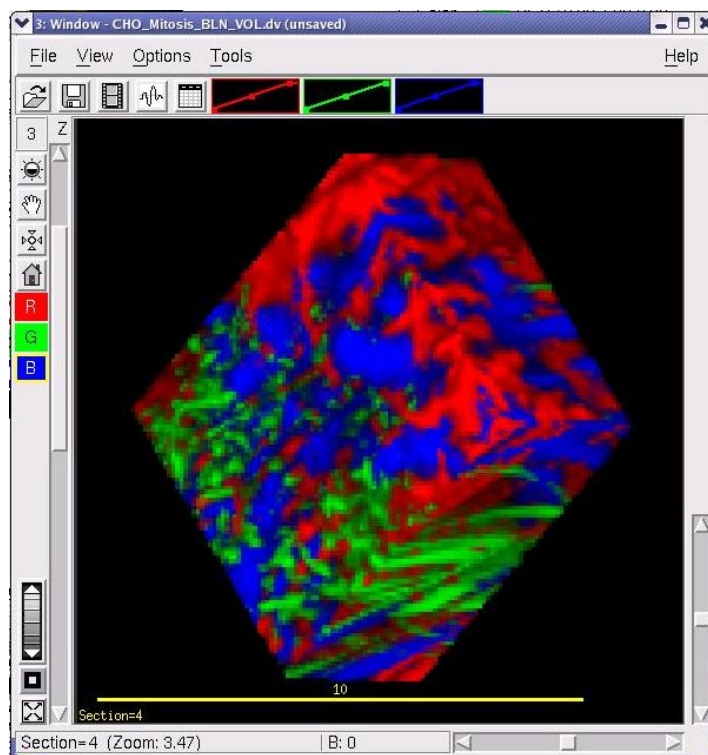

11. To view the volume, move the vertical Z scroll bar up and down or use the Movie tool.

To save the volume rendered images:

1. Click **File | Save** in the Image window menu to open the Save to File window.
2. Examine the **Output** field to ensure that the desired name is used. The *softWoRx* software automatically adds a `_VOL` tag near the end of the filename.
3. Click **Do It** to save the image.

10. Filtering Image Data

Use the *softWoRx* filters to prepare data for modeling and other types of analysis.

In This Chapter

About <i>softWoRx</i> Filters.....	138
Using Convolution Filters	138
Enhancing Object Boundaries.....	139
Using 2D Statistical Filters	141
Using Image Arithmetic	142
Scaling Pixel Intensity to Enhance Local Contrast.....	143
Setting an Intensity Threshold.....	144


About *softWoRx* Filters

softWoRx provides the following types of filters:

- Convolution filters perform high pass, low pass, and other digital filtering.
- The Edge Enhancement filter enhances object boundaries.
- 2D Filter limits noise-like intensity.
- The Image Arithmetic filter scales images, combines information from images, or subtracts images to isolate features.
- The Local Contrast Enhancement filter enhances local contrast around pixels.
- The Threshold filter removes data below an intensity threshold.

Using Convolution Filters


Use the Convolution³ tool to perform basic digital filtering, such as high-pass and low-pass filtering. When you select a filter, the fields in this dialog box are updated to indicate the kernel values.

 **Tip** You may add or change convolution kernels to the list by modifying the CONVOLUTION_FILTERS file in the *softWoRx* configuration directory. Use the Revise Convolution Kernels menu item within the *softWoRx* Utilities menu to access this file.

To use Convolution filters:

1. Choose **Filter | Convolution** from the *softWoRx* main menu to open the Convolution dialog box.

³ For more about Convolution filters, see *Digital Image Processing*, by Kenneth R. Castleman, Prentice Hall, 1995.

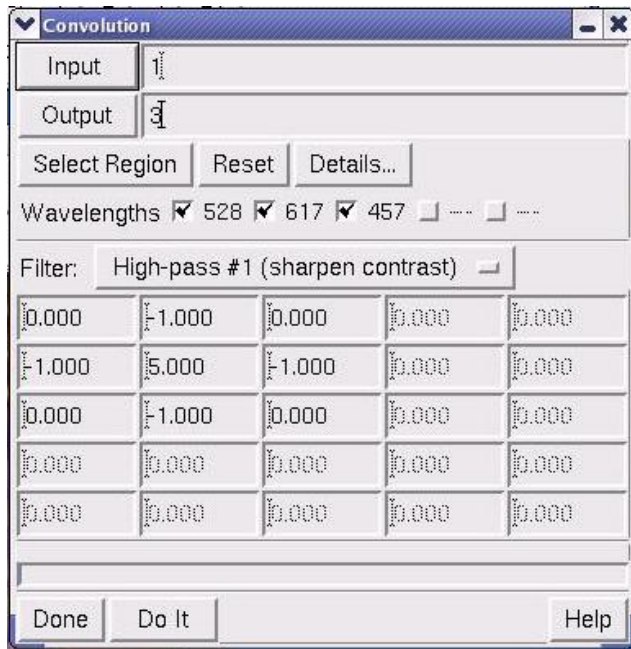

2. Enter an image file name or Window number in the **Input** field.
3. If you want to include only selected data, use the **Select Region** button and drag the mouse across the portion of the image you want to include. Alternatively, you can click **Details** to open the **Region Details** dialog box. Then specify the ranges of data that you want to include in the **Output Options** fields.
4. Select which wavelengths to filter.
5. Select which filter to use in the **Filter** option list.
6. To customize the filter, change the values in the numbered grid.
7. Click **Do It** to run the filter.

Enhancing Object Boundaries


Use Edge Enhance to enhance object boundaries. This tool uses the image intensity gradient to calculate boundaries. The result is calculated from the following expression:

$$\text{Result} = \text{Input} * \{ \text{Fraction} + (1 - \text{Fraction}) * \text{gradient}[F(\text{Input})] \}$$

where *Fraction* is a number between 0 and 1, and *F* is either a linear or an arc tangent function.

To enhance object boundaries:

1. Choose **Filter | Edge Enhance** from the *softWoRx* main menu to open the Edge Enhancement dialog box.

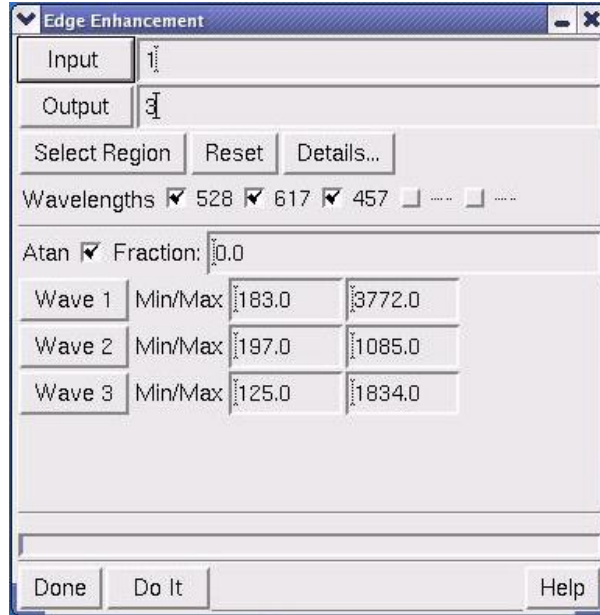


2. Enter an image file name or window number in the **Input** field.
3. If you want to include only selected data, use the **Select Region** button and drag the mouse across the portion of the image you want to include. Alternatively, you can click **Details** to open the Region Details dialog box. Then specify the ranges of data that you want to include in the **Output Options** fields.
4. Select which wavelengths to filter.
5. If you want to use the arc tangent method to calculate the gradient, select **Atan**. To use a linear method, unselect this option.
6. Specify the relative contribution of the original image and the gradient image by entering a value between 0 and 1 in the **Fraction** field. Increasing the **Fraction** value increases the relative contribution of the original image.
7. Click **Do It** to apply the filter.

Using 2D Statistical Filters

Use 2D filters to remove pixels with noise-like intensity or to achieve other effects.

To use statistical filters:

1. Choose **Filter** | **Filter2D** from the *softWoRx* main menu to open the 2D Filter dialog box.

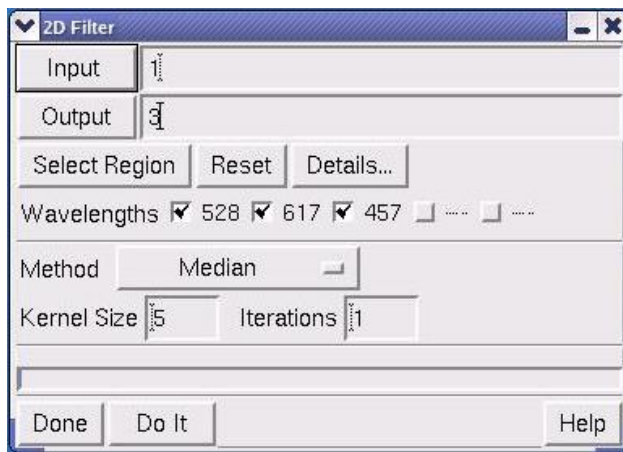

2. Enter an image file name or window number in the **Input** field.
3. If you want to include only selected data, use the **Select Region** button and drag the mouse across the portion of the image you want to include. Alternatively, you can click **Details** to open the Region Details dialog box. Then specify the ranges of data that you want to include in the **Output Options** fields.
4. Select which wavelengths to filter.
5. Select one of the following filters in the **Method** option list:
 - **Median** uses the image intensities within a box-shaped region around each pixel and selects the median value for the resulting image. (This is useful for removing noise.)
 - **Mean** is similar to the median filter except that the mean value is used instead of the median.
 - **Variance** calculates the statistical variance of the image about the mean intensity within the local region.
 - **Weighted Mean** is similar to the mean filter, except that mean is weighted by the variance.
6. Specify the size (pixels) of the square box used to calculate the filtered images in the **Kernel Size** field. (Kernel sizes of 3 and 5 are most useful.)
7. Specify the number of times to run the filter in the **Iterations** field.

- Click **Do It**. The filtered image is displayed in a new Image window.

Using Image Arithmetic


Use Image Arithmetic to scale images, combine information from multiple images, and subtract images in order to isolate features.

Image Arithmetic calculates an output image from one or two input images by applying one of several operations. The input images may come from one or two windows or files. To create the image, specify the input images (files or windows), the wavelength number of Image 1, the wavelength number of Image 2, the arithmetic operation, and any necessary coefficients.

Note In order for the calculation to be processed, the input images must have the same X, Y, and Z dimensions.

To perform an image arithmetic calculation:

- Click **Filter** | **Image Arithmetic** in the *softWoRx* main menu to open the Image Arithmetic dialog box.

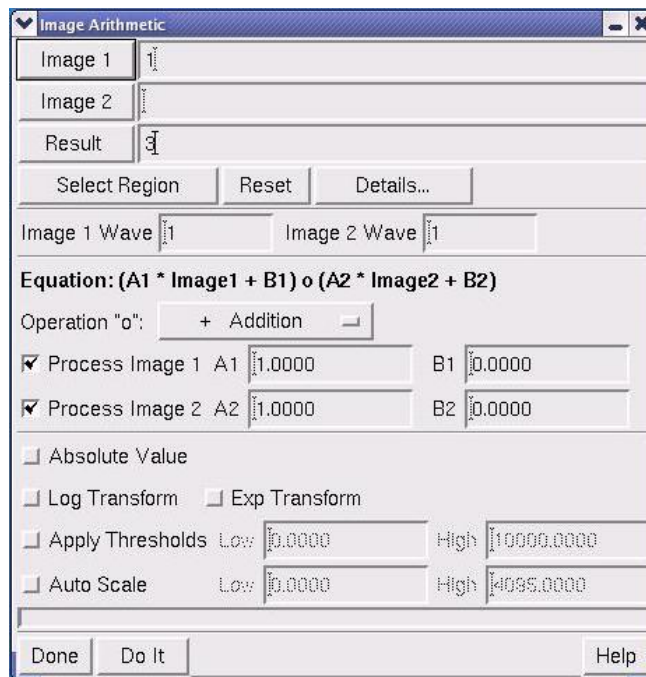

- Drag a window number from an Image window into the **Image 1** field, or click **Image 1** and browse to an image file.
- If desired, repeat Step 2 for the **Image 2** field.
- Define the destination file or window for the calculated result. (The default is **Window 1**.)


5. Enter the appropriate input wavelength from **Image 1** into the **Image 1 Wave** field.
6. If you are using a second image, enter the appropriate wavelength from **Image 2** to use as input into the **Image 2 Wave** field.
7. Select the appropriate operation from the **Operation "o"** list.
8. Set the **Process Image 1** and **Process Image 2** options to enable or disable the processing of the appropriate images.
9. Enable or disable the other options in the dialog box as necessary for your calculation.
10. Click **Do It**. The results of the calculation are displayed in the file or window defined in the **Results** field.

Scaling Pixel Intensity to Enhance Local Contrast

Use Local Contrast Enhancement to scale the intensity of each pixel in the image based on the **Local Mean** and **Local Contrast** of the pixel. (The term "local" refers to the fact that the mean and contrast are calculated from the pixel elements that form a box around the pixel of interest.)

To use nonlinear local contrast enhancement:

1. Choose **Filter | Enhance Contrast** from the *softWoRx* menu to open the Local Contrast Enhancement dialog box.

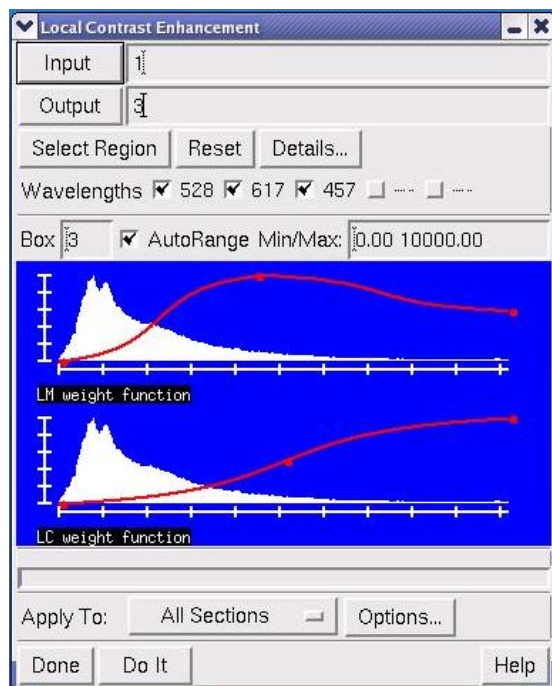


2. Enter an image file name or window number in the **Input** field.
3. If you want to include only selected data, use the **Select Region** button and drag the mouse across the portion of the image you want to include. Alternatively, you can click **Details** to open the Region Details dialog box. Then specify the ranges of data that you want to include in the **Output Options** fields.
4. Select which wavelengths to filter.
5. Specify the box size (in pixels) that determines the size of the region used for the local mean and contrast calculations. Only odd values are valid.
6. To reposition the intensity scaling curves based on the minimum/maximum values of the local mean and contrast, select **AutoRange**.
7. Specify the minimum and maximum values of the intensity scaling curve in the **Min/Max** fields.
8. To modify the intensity scaling curves displayed on the **LM weight function** and **LC weight function** histograms, select one of the three points on the curves with the mouse and move the point to the left or right.
9. Choose one of the following options for applying the contrast enhancement in the **Apply To** option list:
 - **All Sections** creates a regular output image file or window.
 - **Current Section** stores the result in a temporary ("scratch") window.
10. Click **Do It** to create the filtered output.

Setting an Intensity Threshold

Use the Threshold dialog box to remove all data in an image that is below a certain intensity cutoff value. You can choose to output either the image values above the threshold or a binary mask, where all pixels below the cutoff are 0 and all above are set to 1.

To remove data below an intensity threshold:

1. Choose **Filter | Threshold** from the *softWoRx* main menu to open the Threshold dialog box.

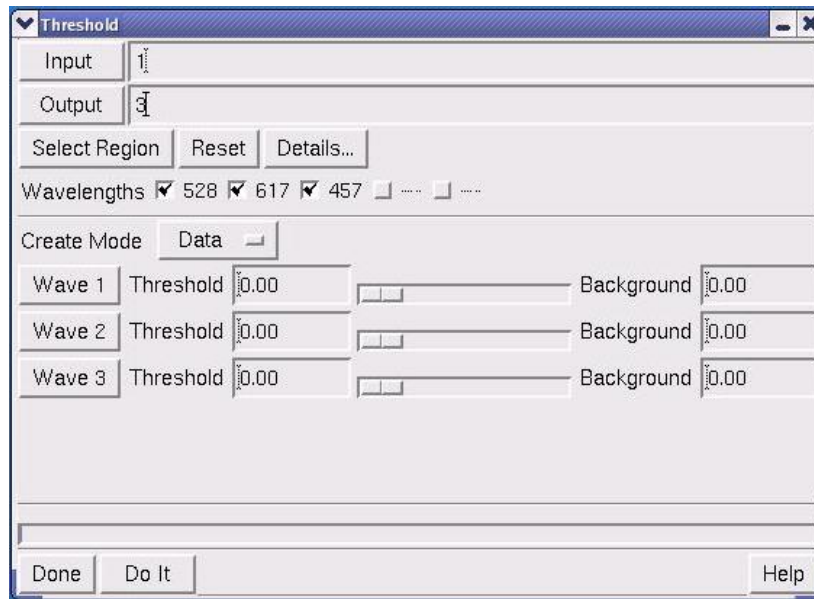

2. Enter an image file name or window number in the **Input** field.
3. If you want to include only selected data, use the **Select Region** button and drag the mouse across the portion of the image you want to include. Alternatively, you can click **Details** to open the Region Details dialog box. Then specify the ranges of data that you want to include in the **Output Options** fields.
4. Select which wavelengths to filter.
5. Select one of the following options in the **Create Mode** list:
 - Choose **Data** to create an output image of all intensities equal to or above the threshold (the cutoff intensity value). If a pixel has an intensity value below the threshold, the pixel in the output image is set to this value (usually 0).
 - Choose **Mask** to put zeros in all pixels where the intensity is below the threshold and ones for all pixels equal to or above the threshold.
6. Specify a **Threshold** and **Background** value for each wavelength.
7. Click **Do It**

11. Saving, Exporting, and Printing

softWoRx provides several options for saving, exporting, and printing image data. You can save complete *DeltaVision* files or save only selected image data (e.g., selected wavelengths or an XYZ region). You can also export *DeltaVision* files to TIFF, Movie, or PhotoShop formats and capture screenshots of images.

After you save images or data, you can archive them to a CD. You can also print TIFF or *DeltaVision* files as images directly from the Image window.

In This Chapter

Image Data Files and Image Graphic Files	148
Saving <i>DeltaVision</i> Files	148
Exporting <i>DeltaVision</i> Files	151
Capturing Screen Shots	156
Archiving Files to CD/DVD	157
Printing Images.....	159

Image Data Files and Image Graphic Files

softWoRx distinguishes between image data files and image graphic files. Image data files contain more information than image graphic files; image data files can be used for analysis while image graphic files are used as graphics.


- 5D Image data files include spatial and possibly temporal and spectral data.
- Image graphic files that are created when you save or copy Image windows include only the 2D pixel data that is displayed inside the window.

Saving *DeltaVision* Files

You can save complete *DeltaVision* files or you can choose to save only specific data. *softWoRx* allows you to select an area and to specify which Z sections and time points to include. You can also choose which channels to save.

To save a .dv file:


1. From the Image window, choose **File | Save** to open the Save File dialog box.

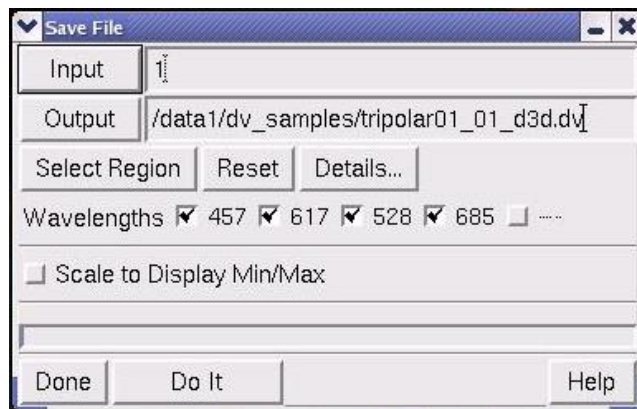

2. Enter an input file or window number in the **Input** field.

Tip You can also specify an existing window by dragging an Image window button from the main *softWoRx* menu to the Input field, or by dragging the window number indicator in the upper left hand corner of the Image window to the Input field.


3. Enter an output file or window in the **Output** field. If the window or file exists, you will need to choose whether to overwrite the data or to append the new data as new channels.
4. To save a region of a window, choose **Select Region** and select an area in the Image window by dragging the mouse across the area. Adjust the rectangle you've created until it contains the desired area. Then click outside the Image window with the mouse.

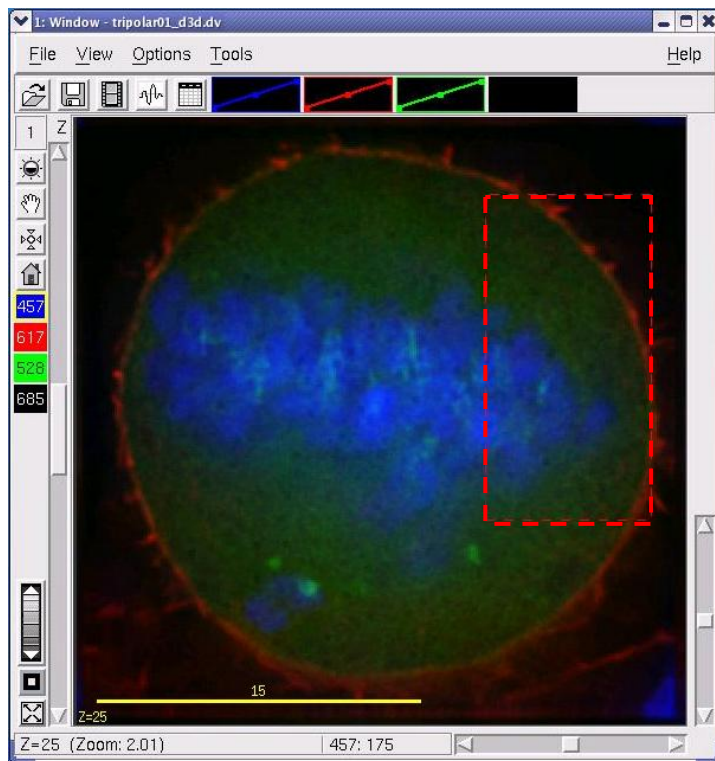

The Region of Interest

Note Selecting a region is optional and you can do this only when your input is a window.

- Click **Details** to open the Region Details dialog box. Specify the ranges of the X, Y, Z, and time data to save in the selected region. Then set any other output details.

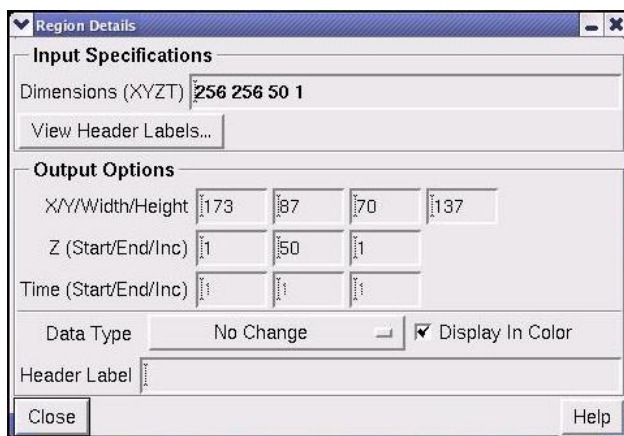


- On the Save File dialog box, choose which wavelengths of the input data to process and include in the output data set.

7. If you want to scale the output data according to the current minimum and maximum intensity scale factors, click **Scale to Display Min/Max**. (Scaling only works when the output data type is 16-Bit integer or 8-Bit).
8. Click **Do It** to save the file.

Saving to the Data Management System (DMS)


You can save *DeltaVision* image files from your local file system to the DMS database if you have this feature set up. See Chapter 6, “DMS Integration” for details.

4. From the Image Window, select **File | Save to DMS**. The Save To Data Management System menu is displayed as shown.

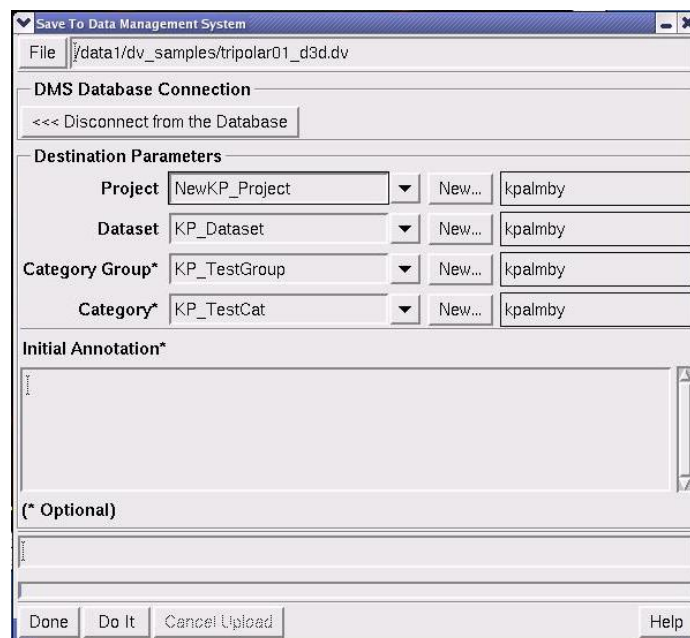

5. From this menu, select the Project and Dataset to which you want to save the data on the DMS database. You can optionally select a Category Group and Category for the saved image data and add initial annotation to the file for future reference. When you are satisfied with the image data to be uploaded, click **Do It**.

When the upload is complete, “Finished” is displayed on the status line as shown.

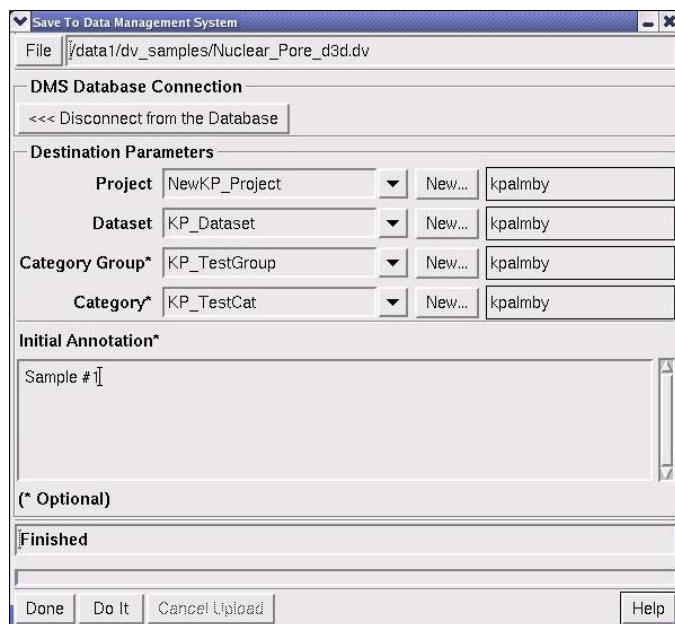

6. Click **Done** to exit the Save To Data Management System menu.

Exporting DeltaVision Files


You can export *DeltaVision* files to TIFF, PhotoShop® or MPEG Movie files. Exporting to PhotoShop and Movie files saves only the pixel image data (the intensity data is not preserved).

Exporting to TIFF Files

Exporting to TIFF files allows you to share data sets in a nonproprietary format.

To export a *DeltaVision* file to a TIFF file:

1. Open the image that you want to export in a *softWoRx* Image window.
2. From the Image window menu, select **File | Save As TIFF** to open the Save as TIFF dialog box.

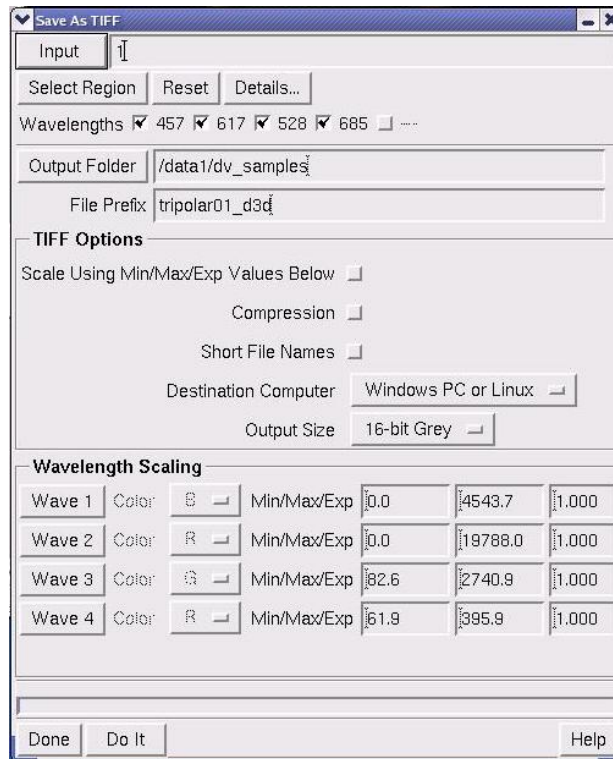

Use Save As TIFF to specify export options

3. If you want to specify a particular time point or a single Z section, click **Details** and set the applicable options in the Region Details dialog box.
4. Enter the directory and file name in the **Output** directory. The default is the directory in which the input file is located.
5. Enter the prefix of the file name in the **File Prefix** field.
6. Select the TIFF options from the lists on the dialog box. (The default settings should work well for most applications.)
7. Under **Output Size**, select one of the three possible TIFF output formats:
 - **8-bit Grey scale** – generates compressed data, with each channel separate.
 - **16-bit Grey scale** – generates uncompressed data. Use this option for quantitation.
 - **24-bit RGB** – generates compressed data; a 3-channel color TIFF with 8-bits per channel.
8. Click **Do It** to export the data to a TIFF file.

Exporting to PhotoShop Files


You can save the visible contents of an Image window as an Adobe PhotoShop 24-bit, RGB color, image file. You can choose to:

- Save the image with the overlay graphics (scale markers, etc) merged into a single image.
- Save the image without the overlay graphics.
- Save only the overlay graphics.

Exporting to PhotoShop saves only the pixel image data (the original intensity data is not preserved).

To export to a PhotoShop file:

1. Open the folder containing the image file that you want to export. Then double-click the image file name to automatically open that image in a *softWoRx* Image window.
2. From the Image window **File** menu, choose **Save As Photoshop** to open the Save Photoshop dialog box.

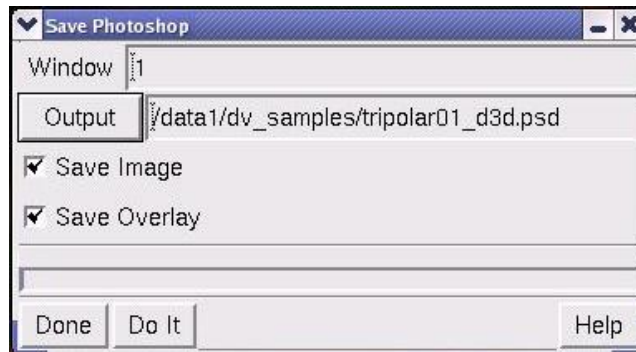

3. Enter the directory and file name in the **Output** directory. The default is the directory in which the input file is located.
4. Select the one or both of the following options for saving the file:
 - To save the image, select **Save Image**.
 - To save the overlay graphics, select **Save Overlay**.
5. Click **Do It** to complete the export.

Exporting to Movie Files


You can save the contents of an Image window as an MPEG movie in 24-bit, RGB color. You can choose to:

- Save the image with the overlay graphics (e.g, scale markers) merged into a single image, save it without overlay graphics, or save only the overlay graphics.
- Save movies that include Z sections, time-lapse data, or both types of data.
- Specify a range of data to include in the movie file.

You can play the MPEG movie format on the QuickTime viewer, the Windows Media Player, or a variety of Linux or Macintosh movie players. You can also import movies into PowerPoint. Double clicking on an MPEG file in a file browser opens the PlayMPEG viewer, which you can use to vary the speed of the movie.

To save a movie:


1. From the Image window, choose **File | Save As Movie** to open the Save Movie dialog box.

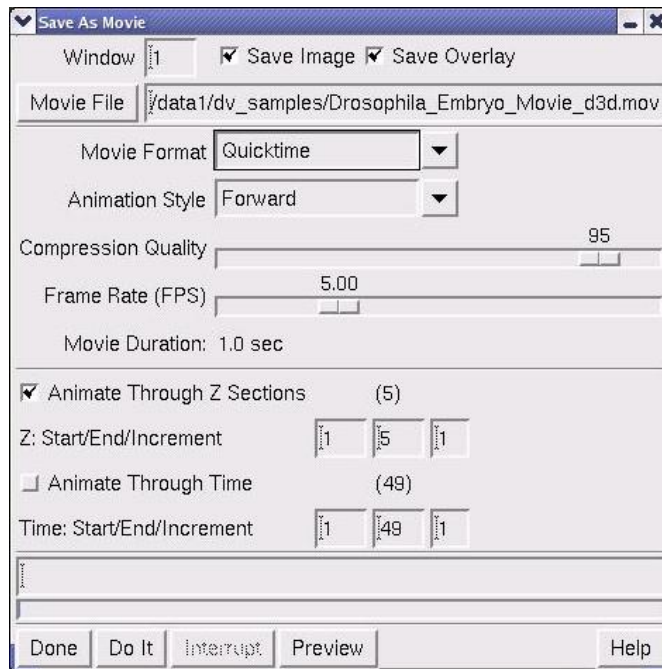

2. Drag the Image window number into the Movie File field, or click Movie File and select a name and path for the movie.
3. At the top of the window, select one or both of the following options:
 - To save the image, select **Save Image**.
 - To save the overlay graphics, select **Save Overlay**.
4. In the **Movie Format** list, select from **Quicktime**, **AVI**, or **MPEG** movie formats.

5. In the **Animation Style** list, set the movie looping mode by selecting one of the following options:
 - **Forward & Back** records a movie that starts on the first frame selected, plays to the last frame, and then back again to the last frame.
 - **Forward** records a movie that starts on the last frame, plays to the first frame, and stops.
 - **Backwards** records a movie that starts on the last frame selected, plays to the first frame, and stops.
6. Adjust the **Compression Quality** slider to specify image quality and file size. Moving the slider to the left produces better quality and bigger file sizes. Moving the slider to the right produces lower quality and smaller file sizes.
7. Adjust the **Frame Rate** slider to specify the playback speed.
8. Select whether to animate through Z sections. Then select the range of data to include in the movie and the increment between frames.
9. Select whether to animate through Time. Then select the range of data to include in the movie and the increment between frames.
10. Click **Do It** to save the movie.

 **Tip** Select the **Preview** button to preview the speed and other settings before saving.

 **Note** *softWoRx* Task Builder provides the capability of exporting Image window contents to AVI and QuickTime movies. For details on using the Task Builder, see Page 54.


Exporting to JPEG Files

You can save the contents of any window that includes an image as a JPEG file. This includes:

- The Image window
- The 3DModel window
- The Statistics window

To save a JPEG file:

1. In a window that contains graphical data, choose **File | Save JPEG Snapshot** to open the JPEG Snapshot dialog box.

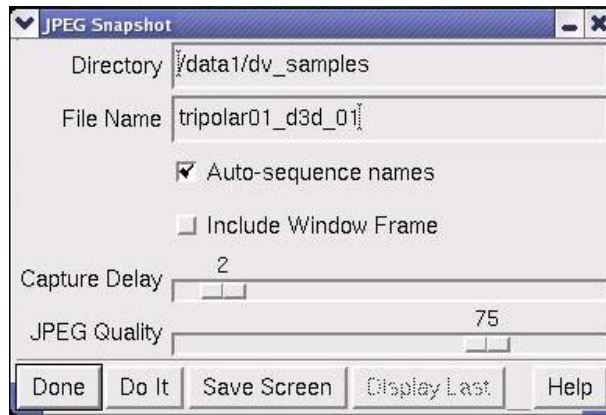


2. Enter a directory and file name and click **Do It**.

Capturing Screen Shots

You can capture a screen shot of any window and save it as a JPEG image. This is a useful way to save images for presentations.


To capture a screen shot:

1. From the main window, choose **Utilities | Image Snapshot** to open the JPEG Snapshot window.

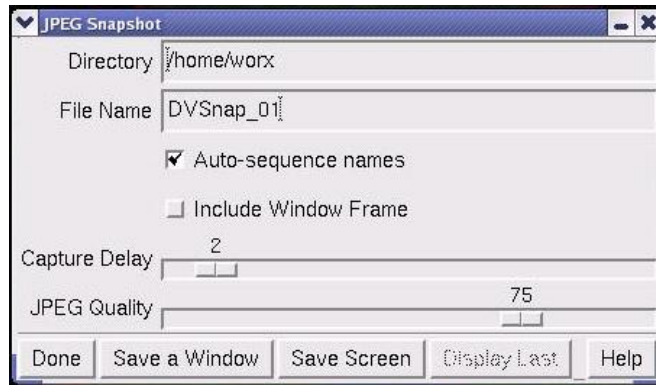

2. Select the directory and file name for the file.
3. To use the same prefix for a series of screen shots, click **Auto sequence names**. (A sequential number is added to the prefix to create a unique name for each file.)
4. To set the delay between when the image is selected and captured, move the **Capture Delay** slider.
5. Set the JPEG slider to the right to increase the quality of the screen shot. (Higher quality results in larger file sizes.)
6. Choose whether to save a window or the entire screen.

7. Click the window (or screen) to capture the image.
8. To view the screen shot, click **Display Last**.

 **Tip** You can also open Snapshot from the Image window, the 3DModel window, the Chromatic Correction dialog box, and the Orthogonal Viewer window.

Archiving Files to CD/DVD


You can open the Linux K3b “CD Kreator” tool directly from the *softWoRx* main menu to archive your *softWoRx* files to CDROM or DVD discs. This tool supports 650 or 700 MB capacity CD-RW discs and 4.3 GB DVD discs.


To copy files to a CD using K3b:

1. Place a blank CD into the CD drive.
2. On the *softWoRx* main menu, choose **Utilities | Archive Data to CD** to open the K3b CD/DVD creation window.




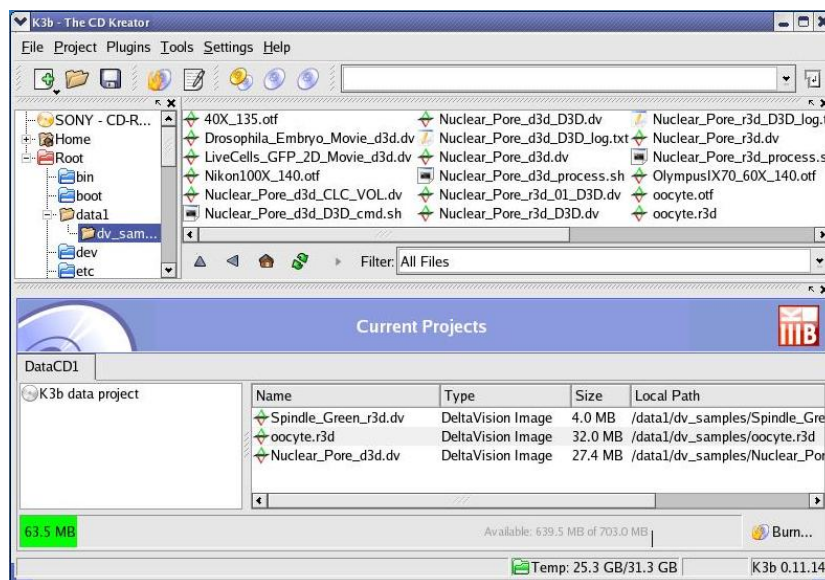
3. Click on the **New Data CD Project** icon. The Current Projects window opens.

 **Note** The process for archiving files to a DVD is the same, except you choose **New Data DVD Project** in this step.

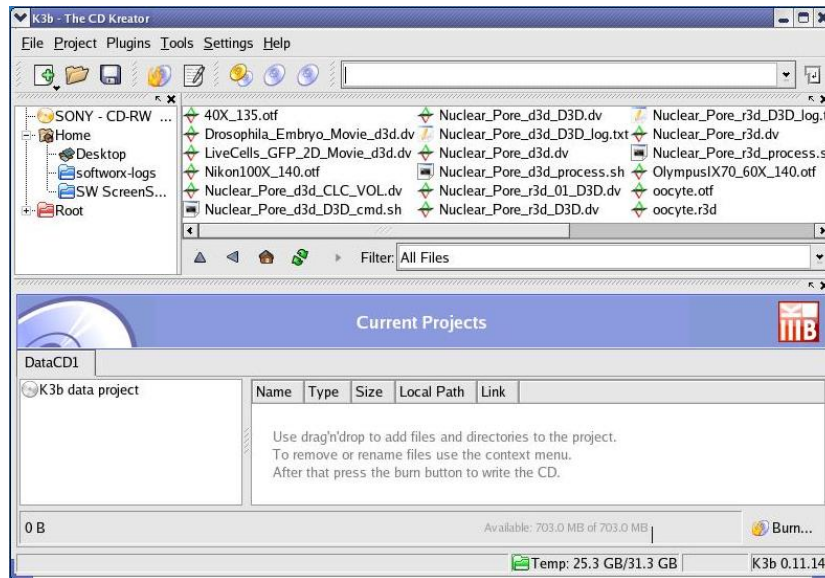


4. Navigate to the files you want to archive and double-click on each file you want to copy to the CD, or drag and drop the files into the Current Projects window.

 **Tip** You can select multiple files in the K3b window by holding down the CTRL key while selecting files.



5. When you are satisfied with your selections, click the **Burn...** button in the bottom right corner of the window. The Data Project – K3b dialog box is displayed.

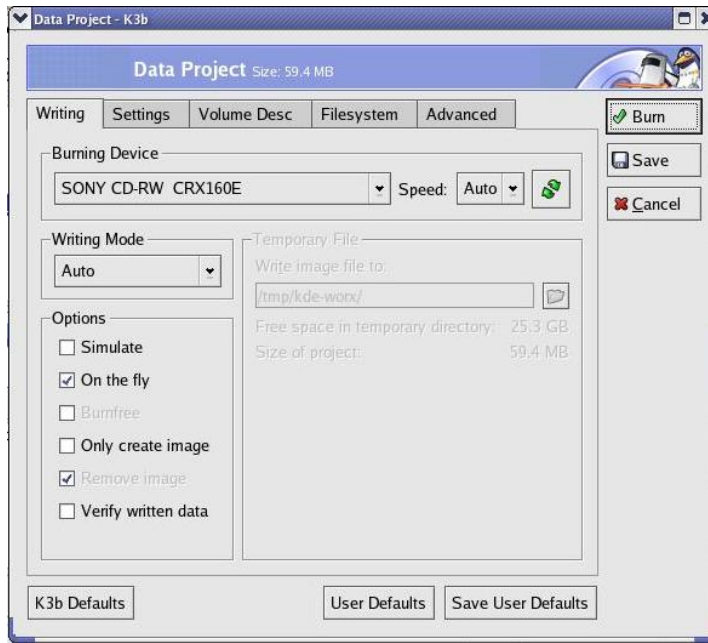


6. Click the **Burn** button at the top right corner of the window to begin the CD creation process. For more complete information on using the K3b “CD Kreator” tool, select **Help | K3b Handbook** on the K3b main window to view the entire user’s manual for K3b.

Printing Images

softWoRx allows you to print *DeltaVision* images from an Image window.

To print a *DeltaVision* image:

1. Open the .dv image that you want to print in the Image window.
2. From the Image window menu, choose **File | Print**.

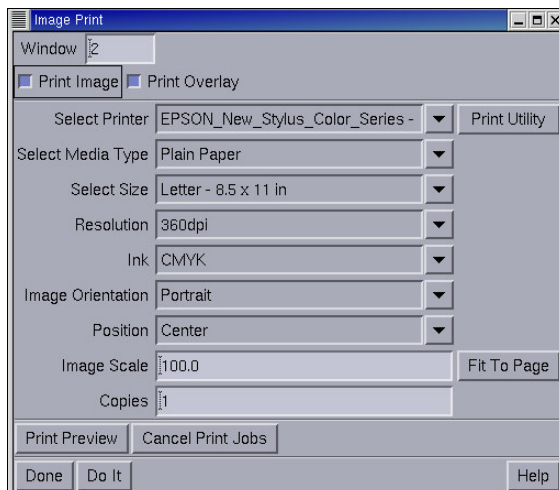

3. In the Image Print dialog box, enter the window number of the image in the **Window** field.
4. Click **Do It** to print the image.

Part Three

ANALYZING RESULTS

Part Three shows how to use *softWoRx* tools to perform quantitative analysis.

In Part Three

Chapter 12. Examining Intensity Data162

Chapter

13. Measuring Distance and Velocity..... 175

Chapter

14. Volume Modeling 181

Chapter

15. Detecting and Analyzing Colocalization.....193

Chapter

16. Other Applications203

12. Examining Intensity Data

This chapter introduces the tools used to examine intensity data.

In This Chapter:

Examining Point Values	162
Examining Intensity Data With Data Inspector	163
Viewing Intensity Line Profiles	166
Calculating Statistics	169


Examining Point Values

You can examine intensity values for individual pixels in the Image window. The wavelength and intensity value of the point under the mouse are displayed at the bottom of the Image window.

To view point values:

- ▶ From the main *softWoRx* toolbar, choose **Measure | Point Values**. Select a channel (**Wave**) and move the mouse across the image to view individual


point values for that channel. The point coordinates and intensity of the point under the mouse are displayed in the Point Values dialog box.

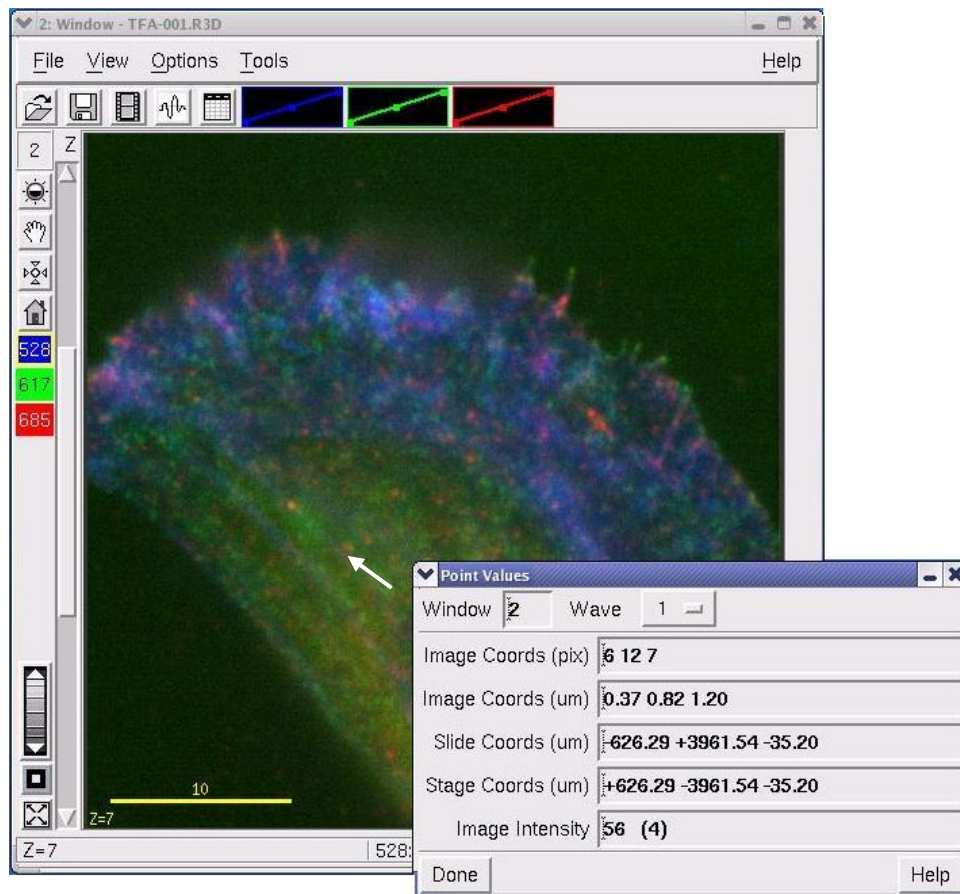

Tip You can also view the intensities of individual points for the selected channel on the status bar (e.g., 457:1216 in this example).


Examining Intensity Data with Data Inspector

The Data Inspector includes several tools for examining the intensity data in the image file. You can simultaneously view a graphical image, a table of intensity values, a 3-D graph, and a histogram of intensity values. With these views open, you can select various regions of interest (ROIs) to explore the data. As you select an ROI, the data in each view is updated for that ROI.

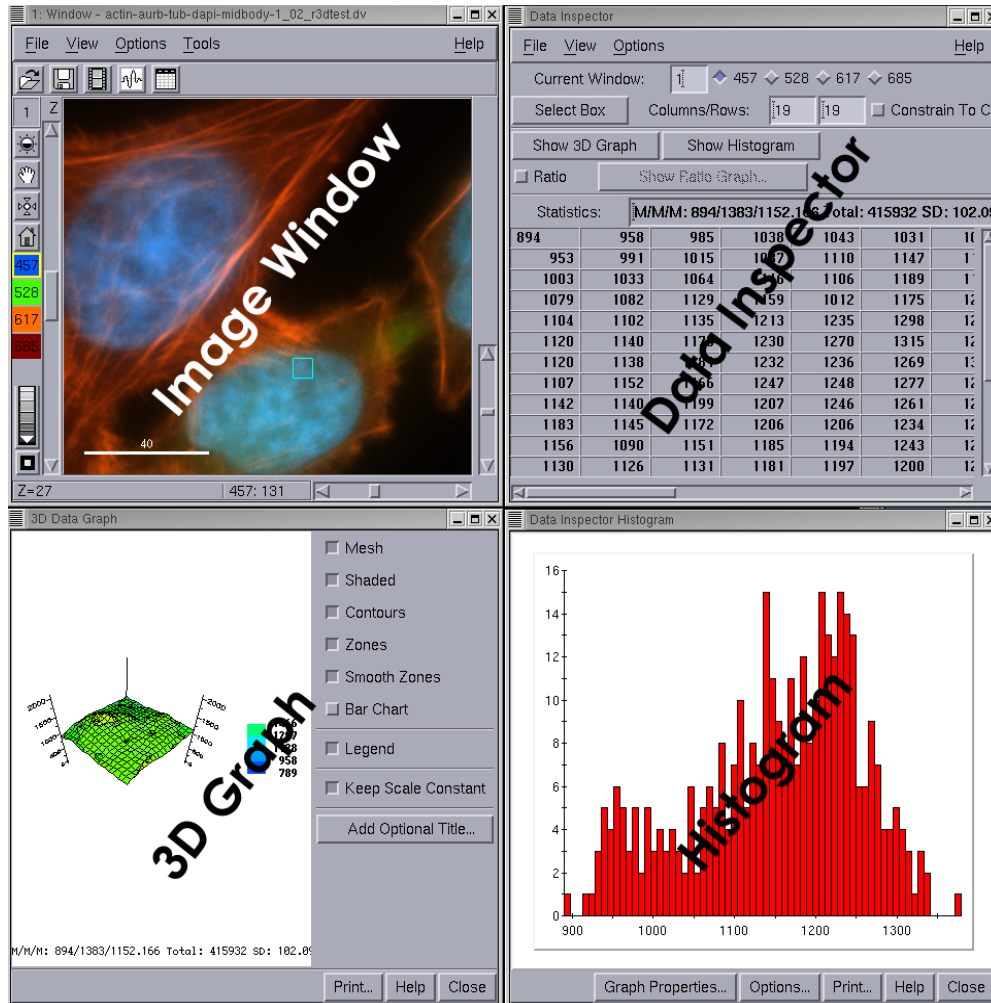

To open the Data Inspector tools:


1. Open an image file in the Image window.
2. Choose **Tools** | **Data Inspector** on the Image window to open the Data Inspector window.

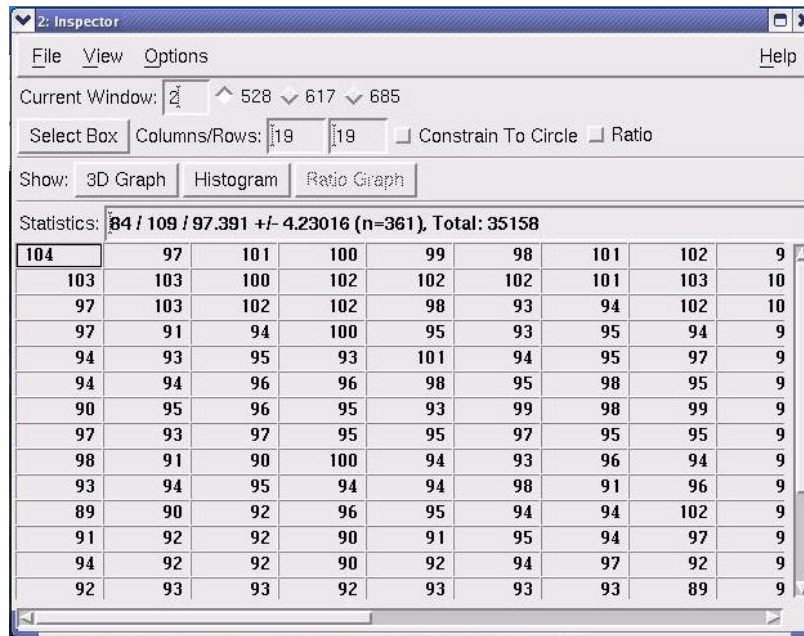

3. In Data Inspector, click **Show 3D Graph**.
4. Click **Show Histogram**.
5. Arrange the windows on the screen.

Selecting a Region of Interest

The region of interest, or ROI, is a region of the image that you can resize and move to visually examine image details. You can select a region of interest (ROI) in the Image window to display its intensity values in the 3D graph, the Histogram, and the table in Data Inspector. The display in each window changes automatically as you change the ROI in the Image window.

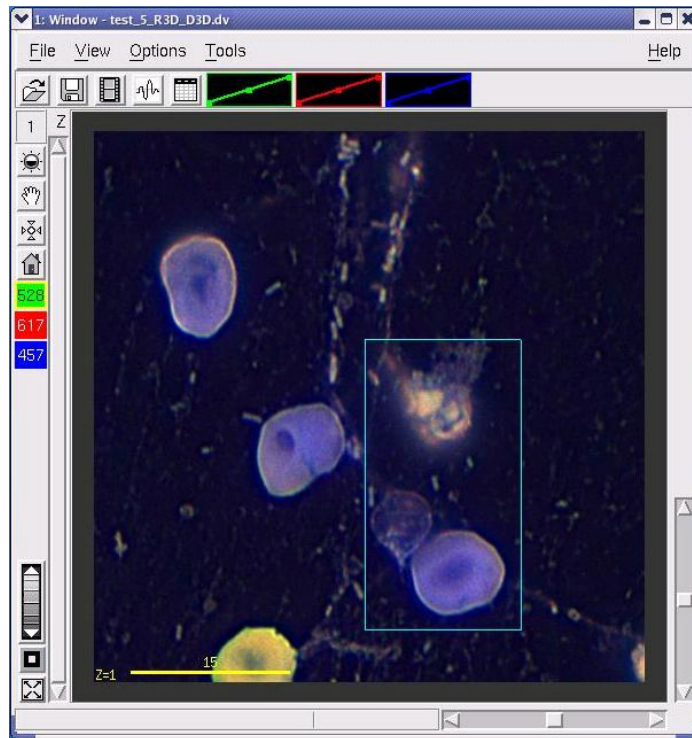

The ROI is the area within the box

To define an ROI:

1. Choose **Tools | Data Inspector** on the Image window to open the Data Inspector window.
2. On the Data Inspector window, click **Select Box**.
3. In the Image window, click the top left corner of the ROI and drag the mouse to enclose a region of interest with the ROI box.
4. To make the ROI a circle instead of a square, select **Constrain to Circle**.
5. To change the position of the ROI, click on any point in the Image window. (The ROI is centered on the point that you click.)

Viewing Intensity Line Profiles


A line profile is a plot of intensity values for pixels along a straight line. Two types of line profile tools are available:

- **Line Profile** displays a plot of intensity values for pixels in a row or column of the Image window. This profile is overlain on the image. You can view a profile of a line of pixels or a band of pixels. This data can be saved as a text file.


- **Arbitrary Profile** displays a plot of intensity values along a line segment that is oriented at any angle in the Image window. This profile is displayed in a separate window. You can interactively change the orientation of the line. This data can be saved in an .s1k spreadsheet compatible file.



Viewing the Line Intensity of a Row or Column

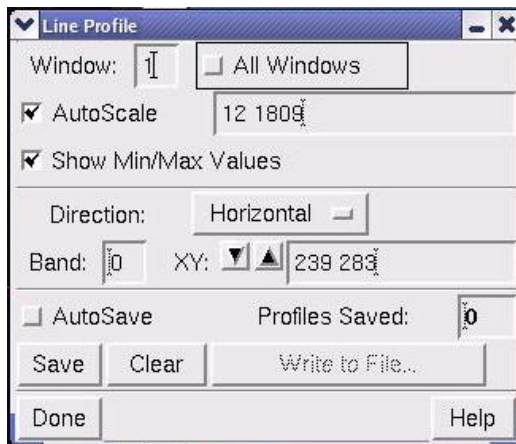1. Open an image in the Image window and choose **Tools | Line Profile** on the Image window to view the Line Profile dialog box.

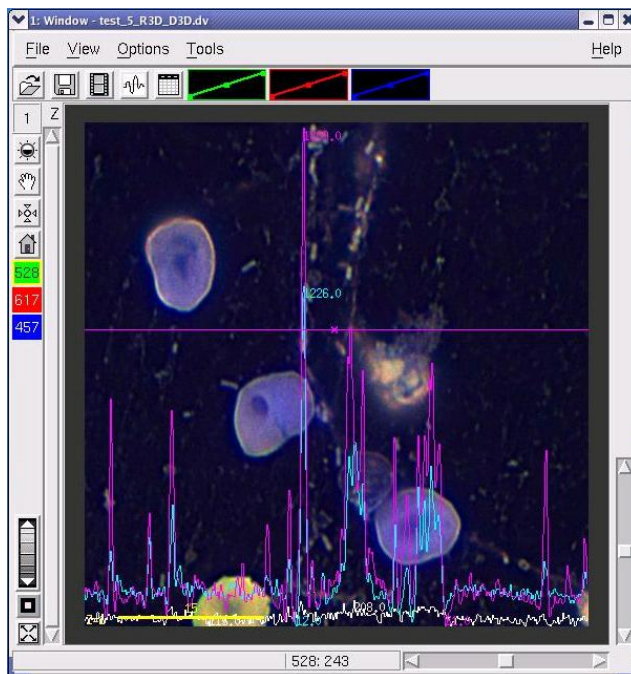

2. Drag the window number from the Image window to the **Window** field in the Line Profile dialog box.
3. Click on the image to display a horizontal line profile.



4. To get the average line profile of several rows and columns of pixels, enter the width in the **Band** field.
5. To display a vertical line profile, choose **Vertical** in the **Direction** list.
6. To change the position of the profile, click another point on the image or use the  and  buttons next to the XY field.
7. Create a group of profiles to save in a file as follows:
 - To save only selected profiles, unselect **AutoSave**. Then click **Save** after each profile that you want to save is displayed.
 - To save all profiles, select **AutoSave** before you create the profiles.
8. To save a group of profiles, click **Write to File**, enter the file name and other options in the Profile Output Options dialog box, and click **Do It**.




Tip You can change the Line Profile colors that are associated with different wavelengths. From the Image window, choose **View > Select Graphics Colors**.

Viewing the Line intensity in Any Direction

To display an arbitrary line profile:

1. Open an image in the Image window and choose **Tools | Arbitrary Profile** to open the Arbitrary Line Profile dialog box.
2. Click and drag on the image to display a line profile.

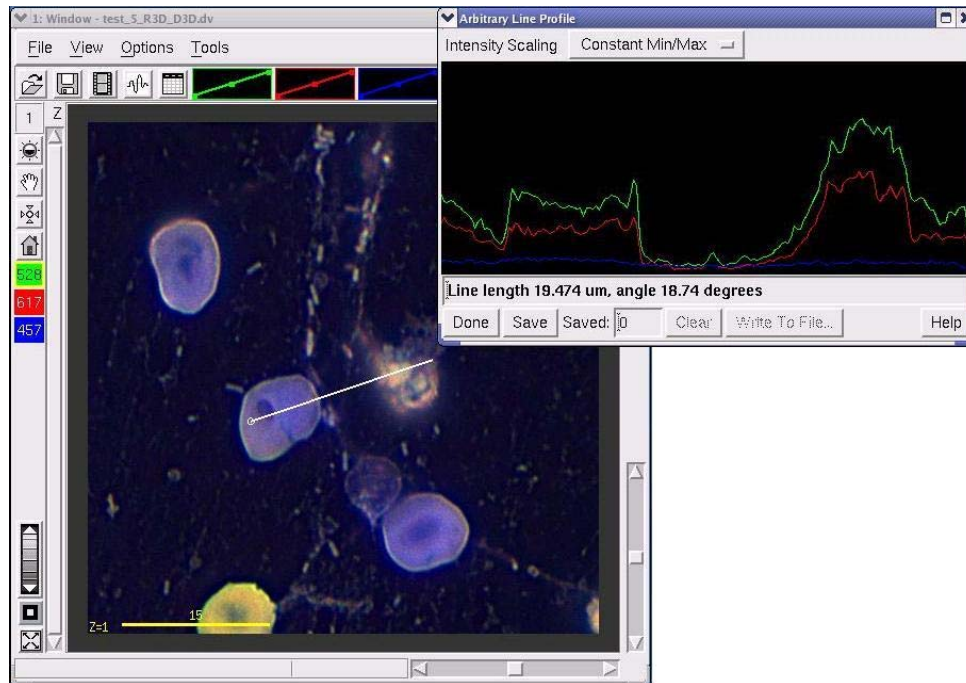

3. Choose how to scale the profile in the **Intensity Scaling** list.
 - **Constant Min/Max** sets the scale range to the minimum and maximum intensity values in the image.
 - **Autoscale** sets the range to the minimum and maximum intensity values in the profile.
4. To save a group of profiles in a spreadsheet-compatible text file, click **Save** after each profile that you want to save is displayed. When you are finished collecting profiles, click **Write to File**, enter the file name and other options in the **Save as SYLK spreadsheet** dialog box, and click **OK**.

Calculating Statistics

Calculating Statistics for Selected Areas

Use Data Inspector to calculate data for selected areas.

To open Data Inspector:


- ▶ Select **Tools | Data Inspector** on the Image window to open the Data Inspector window.

The Data Inspector window displays a grid of intensity data with the following statistics:

Statistics: M/M/M: 1092/1250/1181.515 Total: 426527 SD: 95.14202								
1163	1171	1145	1163	1172	1210	1201	1198	1174
1147	1160	1174	1149	1165	1167	1197	1188	1170
1149	1163	1161	1133	1173	1176	1168	1192	1176
1143	1161	1156	1185	1188	1169	1189	1154	1185
1195	1198	1196	1186	1179	1175	1191	1173	1160
1167	1181	1164	1195	1167	1177	1182	1181	1166
1176	1181	1175	1175	1168	1175	1173	1168	1177
1220	1198	1190	1172	1205	1179	1180	1188	1194
1198	1195	1186	1193	1193	1187	1185	1176	1180
1209	1208	1197	1199	1216	1192	1185	1178	1194
1184	1211	1213	1244	1218	1197	1194	1191	1181
1225	1211	1220	1216	1230	1200	1188	1192	1181
1221	1222	1217	1223	1212	1190	1167	1194	1207
1205	1227	1231	1220	1223	1218	1193	1213	1209

To gather statistics:

1. Click on a point of interest in the Image window and press the space bar. (The statistics for a rectangular area around that point are displayed in the Data Inspector **Statistics** field.)

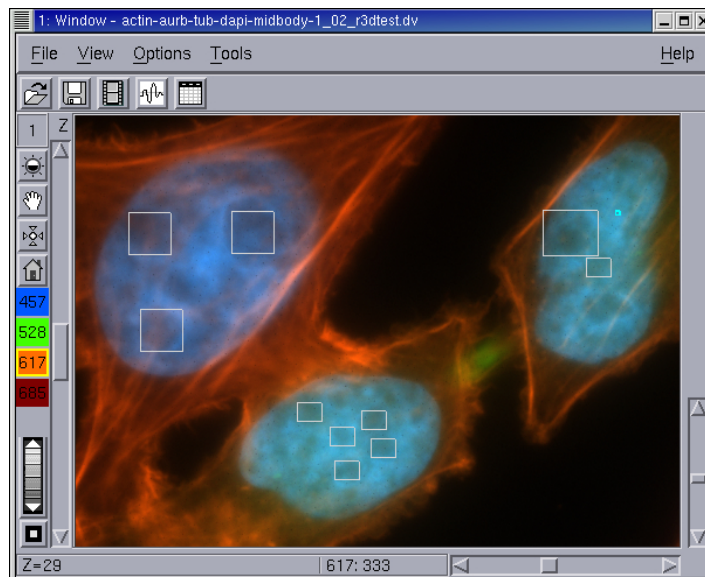

2. From the Data Inspector window, choose **File | Save Statistics Record** to open the Statistics Record File dialog box.
3. To save the statistics file, enter a file name in the Statistics Record File dialog box and click **OK**. The file is saved as a text file similar to the following.

File Generated By softWoRx DataInspector
/data1/statisticsRecord_sample2 Tue Feb 17 16:55:01 2004


wave	X	Y	spot diameter	box width	incl_points	min	max	mean	total intensity	SD
685	-4196.678	1215.97	39	39	1070	190	262	215.49	230575	10.77
685	-4193.034	1188.47	39	39	1070	188	273	233.24	249563	16.83
685	-4167.192	1215.63	39	39	1070	180	232	201.01	215080	7.95
685	-4128.099	1157.66	24	24	408	175	206	190.15	77580	5.39
685	-4136.05	1150.04	24	24	408	180	229	199.39	81353	7.39
685	-4129.755	1166.27	24	24	408	187	222	202.22	82505	5.73
685	-4066.146	1195.09	36	36	960	171	270	199.19	191227	16.18
685	-4075.422	1215.30	36	36	960	169	218	186.62	179153	7.57
617	-4191.709	1189.13	36	36	960	592	1199	945.76	907931	134.62
617	-4195.022	1216.63	36	36	960	612	1057	811.14	778695	86.70
617	-4165.867	1216.96	36	36	960	557	858	672.46	645557	52.15
617	-4136.713	1152.69	16	16	176	443	631	535.04	94167	41.04
617	-4126.442	1158.98	16	16	176	447	571	510.66	89876	28.96
617	-4129.093	1166.93	16	16	176	540	719	626.86	110328	38.53
617	-4138.038	1162.30	16	16	176	472	558	520.51	91609	15.31
617	-4147.315	1169.25	16	16	176	496	755	628.27	110575	46.45
617	-4065.483	1210.33	16	16	176	441	612	523.47	92131	35.21
617	-4077.742	1216.30	39	39	1070	399	745	508.66	544263	46.83

Calculating Statistics for Irregular Areas

The Edit Polygon tool can be used to calculate statistical values for the area of data that is inside of a polygon. You can draw multiple polygons on different Z sections and view the statistics for each. You can also save statistical reports to text files or to SYLK spreadsheet compatible files.


To select polygon areas:

1. Choose **Model | Edit Polygon** on the main *softWoRx* menu.

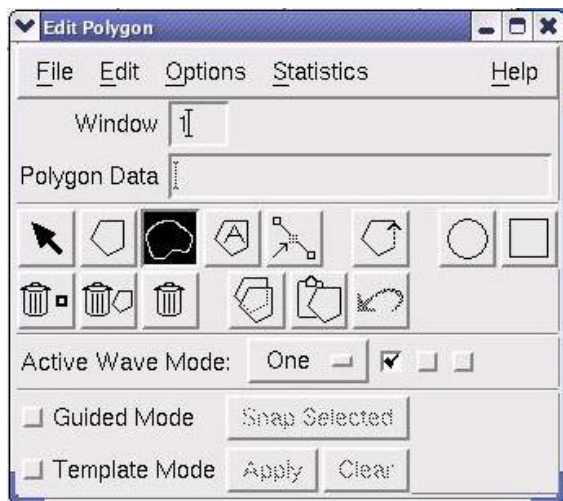


2. Click a polygon option (e.g., ) on the Edit Polygon window.


3. To automatically find the borders around objects (based on changes in intensity values) select the **Guided Mode** option.
4. Drag the mouse to draw the polygons on the Image window. You can draw sets of polygons on the same Z section and on different Z sections.

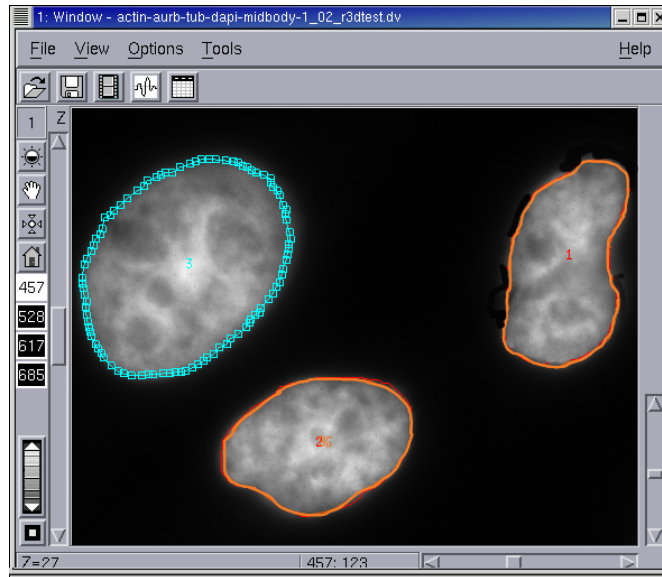

5. To copy a polygon to other sections or wavelengths, use the arrow tool in the Edit Polygon window to select the polygon. Then choose **Edit | Propagate Polygons**.

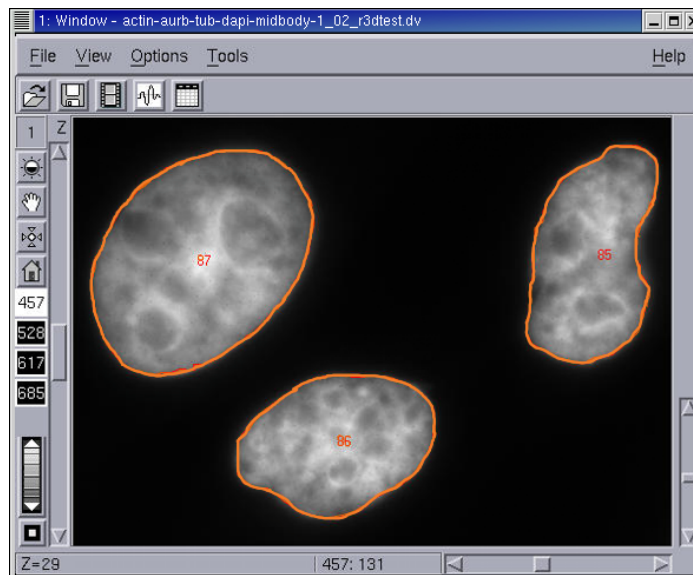


6. Choose **Statistics | Table** to display a table that shows statistical values for each polygon.

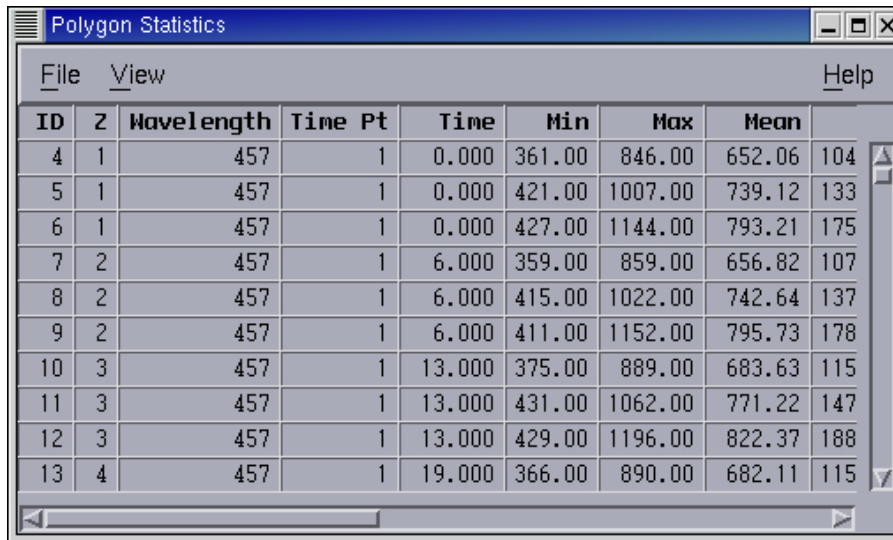

ID	Z	Wavelength	Time Pt	Time	Min	Max	Mean	
4	1	457	1	0.000	361.00	846.00	652.06	104
5	1	457	1	0.000	421.00	1007.00	739.12	133
6	1	457	1	0.000	427.00	1144.00	793.21	175
7	2	457	1	6.000	359.00	859.00	656.82	107
8	2	457	1	6.000	415.00	1022.00	742.64	137
9	2	457	1	6.000	411.00	1152.00	795.73	178
10	3	457	1	13.000	375.00	889.00	683.63	115
11	3	457	1	13.000	431.00	1062.00	771.22	147
12	3	457	1	13.000	429.00	1196.00	822.37	188
13	4	457	1	19.000	366.00	890.00	682.11	115

- Choose **Statistics | Graph** to open the Plot Polygon Graph dialog box.

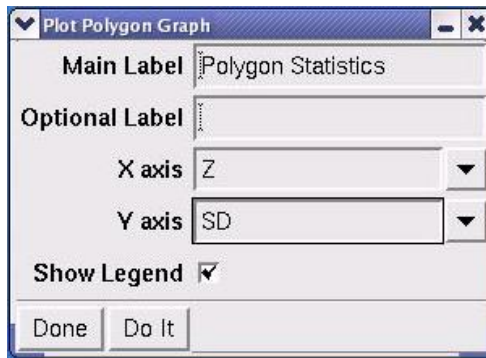

Plot Polygon Graph

Main Label: Polygon Statistics

Optional Label:


X axis: Z

Y axis: SD

Show Legend:

Done Do It

- In the Y axis field, choose which type of statistical parameter (e.g., SD) to plot.
- Choose other options in the Plot Polygon Graph dialog box and click **Do It** to display a graph of the statistics values.

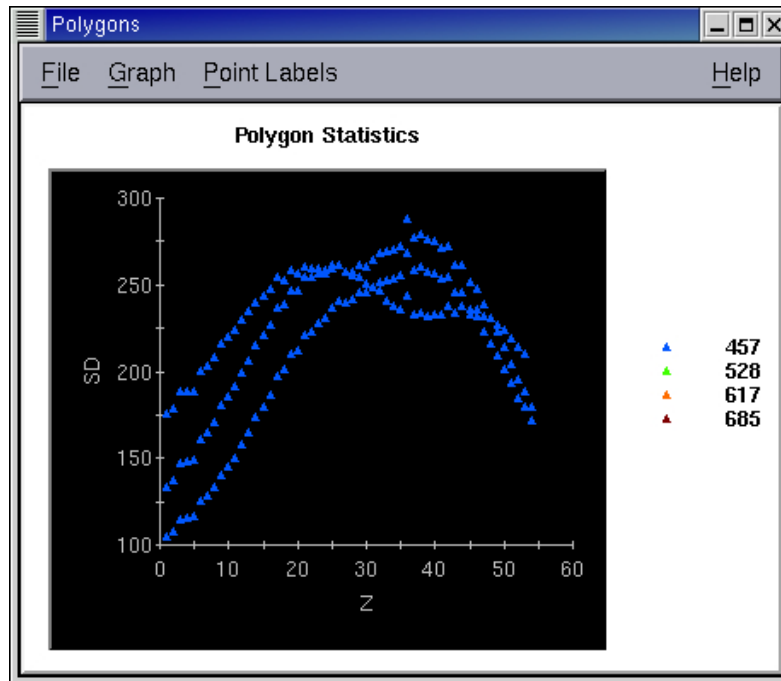

10. From the Polygons window **File** menu, choose **Save As SYLK** to save the file as a spreadsheet compatible file.

13. Measuring Distance and Velocity

This chapter shows how to measure distance and velocity.


In This Chapter

Measuring Distances	175
Measuring Velocity	177

Measuring Distances

The Distance tool allows you to measure distances between points in one Z plane or between points in many different Z planes. The measurement data can be saved to a file for off-line analysis.

The Measure Distances dialog box contains the options for measuring distances. This section briefly describes these options and contains step-by-step instructions for completing the most common procedures.

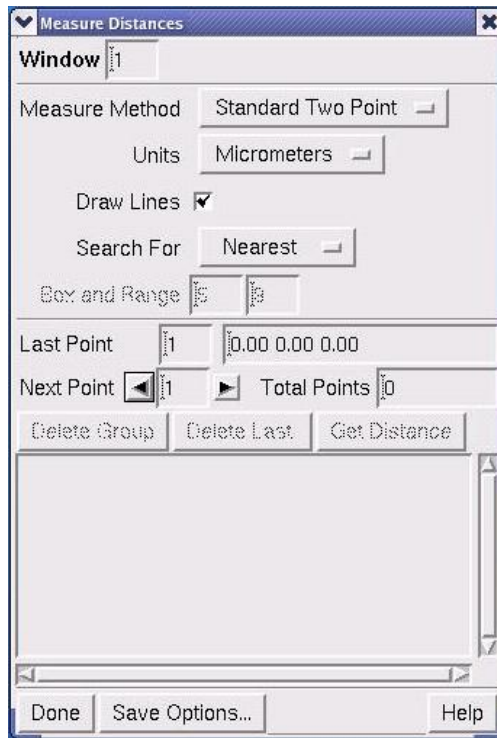

Options in Measure Distances are summarized below and are described in further detail in the *softWoRx* online Help.

Select	To...
Standard Two Point	Measure the distance between two points.
Single Reference	Measure the distance from a single reference point to other points.
Leap Frog	Measure the distance between two consecutive selected points.
Multiple Segment	Measure the sum of the distance between consecutive selected points.

The following procedures describe the most common measurement tasks performed using *softWoRx*.

To measure distance using the Standard Two Point method:

1. Open an image in the Image window.
2. On the Image window menu, click **Tools | Measure Distances** to open the Measure Distances dialog box.
3. Enter the desired window number in the **Window** field.
4. In the **Measure Method** list, select **Standard Two Point**.
5. Select the appropriate unit of measurement from the **Units** option.

6. Select two points using the mouse. The distance will be displayed in the gray information window.

To measure distance using the Multiple Segment method:


1. Open an Image in the Image window.
2. On the Image window menu, click **Tools | Measure Distances** to open the Measure Distances dialog box.
3. Enter the desired window number in the **Window** field.
4. In the **Measure Method** list, select **Multiple Segment**.
5. Select the appropriate unit of measurement from the **Units** option.
6. Drag the mouse across the image to select consecutive points.
7. Click **Get Distance** to display the coordinates of all the selected points and the sum of the distance in the gray information window.

 **Note** If an error occurs, click **Delete Last**.


Measuring Velocity

You can use the distance tool to measure the velocity of objects in time-lapse data. Use the Measure Distance tool to measure the velocity of particle movement.

1. Open an image in the Image window.
2. On the Image window menu, click **Tools | Measure Distances** to open the Measure Distances dialog box.

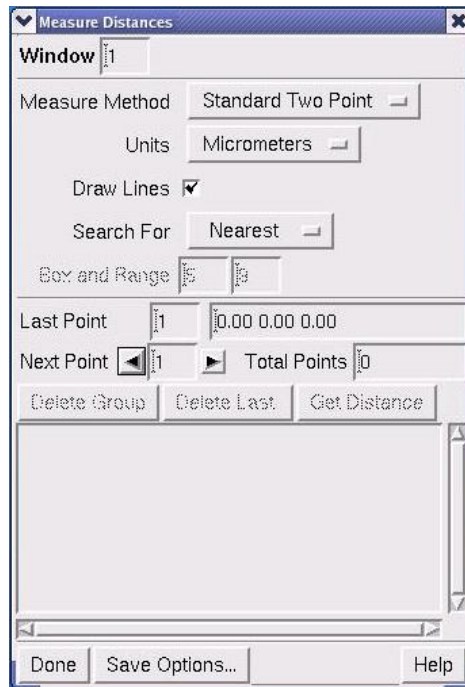


3. In the **Measure Method** list, select the **Standard Two Point** method.
4. In the **Units** list, select the appropriate unit of measurement.
5. In the Image window, click on the particle that you want to measure.

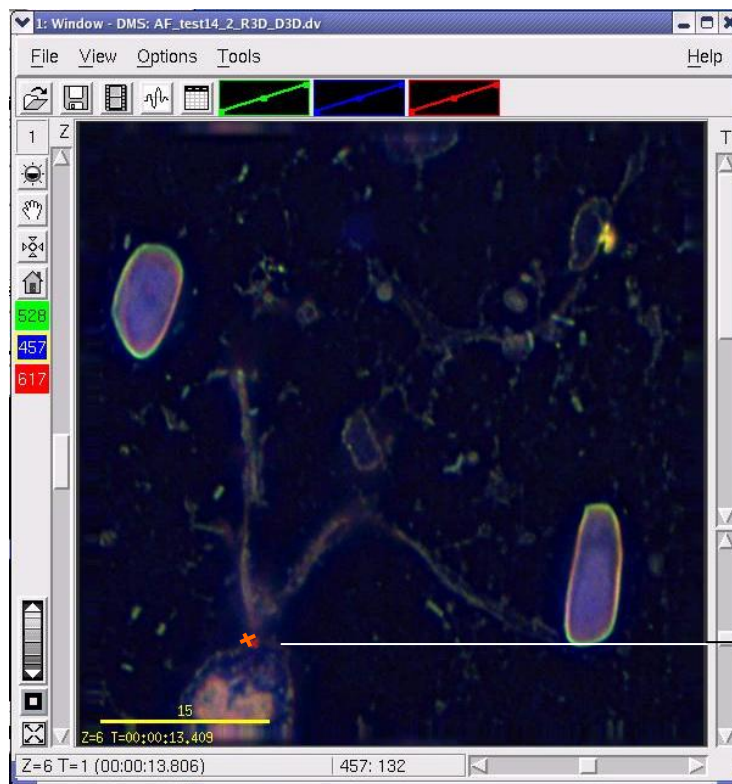


Original Time Point

6. Move the T (time) slider on the Image window to display the particle at a different time point.
7. Click the same particle that you selected at the original time point.

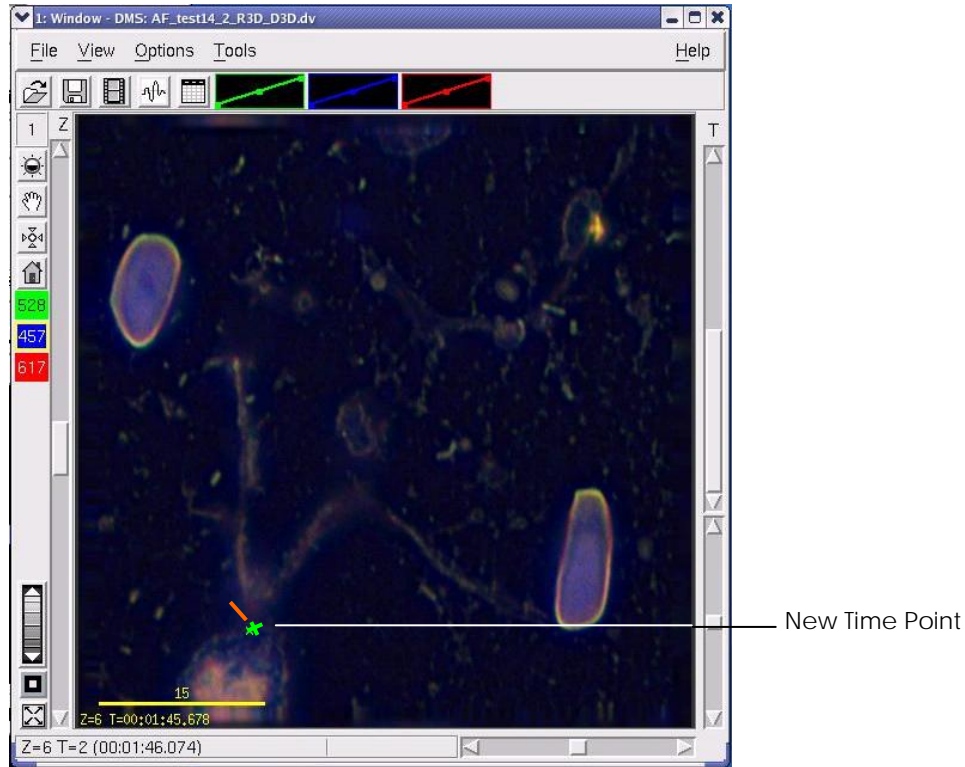


The distance and velocity of the particle movement are displayed at the bottom of the Measure Distances dialog box.

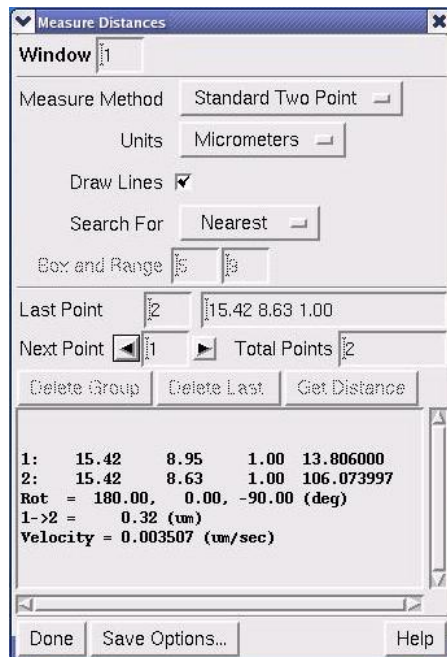

14. Volume Modeling

Modeling features of your image data can help you to understand the three-dimensional nature of your data. Volume models have been used to study objects such as nuclei and cell boundaries.

 **Note** *softWoRx* also provides Line modeling tools that have proven to be useful for studying chromosomes, neurons, and other complex three-dimensional structures. For information about Line modeling, see the online Help.


In this Chapter:

About Volume Modeling	182
Edit Polygon Dialog Box	182
2D Polygon Finder	185
3D Object Builder	187
Volume Modeling Example	190

About Volume Modeling

The 2D Polygon Finder and the 3D Object Builder are used to create three-dimensional models of features within the image data and obtain quantitative information. Initially, the 2D Polygon Finder is used to specify the object of interest. The 3D Object Builder is then used to create a three-dimensional model from the two-dimensional polygons in each Z section. Finally, the real space coordinates are saved to an ASCII file (which can be viewed from a table of measurements) and the 3D Model Display can be viewed.


In summary, this is the process:

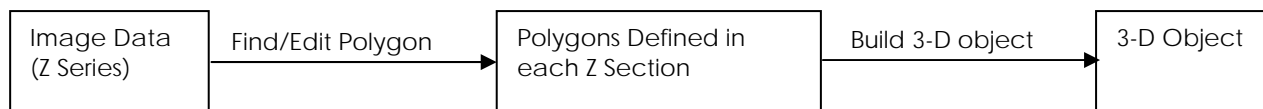

3D Modeling Flow

Edit Polygon Dialog Box

When a satisfactory polygon is not obtained using the 2D Polygon Finder, you can use the Polygon Editor to define a polygon.

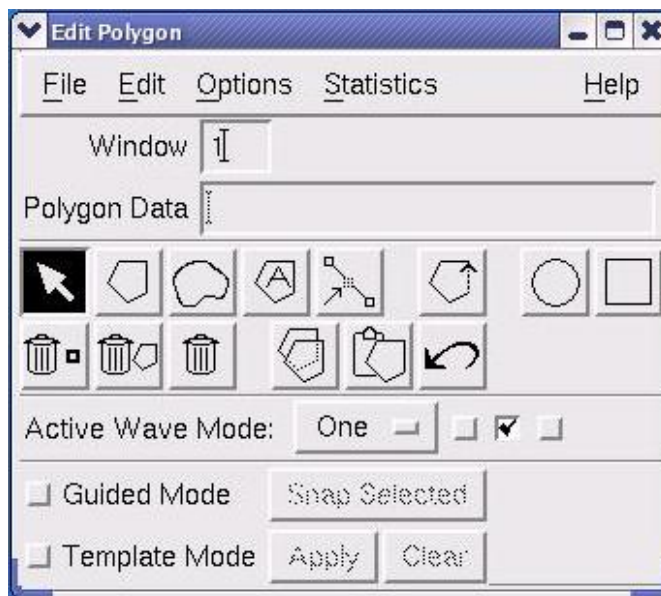

Several controls enable the Polygon Editor to define the intended feature. Descriptions of the most useful controls follow. For additional information, refer to the online Help.

Window

Sets the Image window number of the image data to be processed.

Polygon Data

Defines the Image window to be used by a file name.

Snap Selected

Determines the boundaries of the polygon after you have manually selected an approximation of the polygon. Polygon Editor finds the closest pixel to your manually selected polygon that matches the criteria of the Guide Options.

Options | Guide Mode Options

Defines the parameters for the Polygon Editor to evaluate when you use **Snap Selected**. For additional information, refer to the online Help.

Done








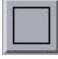


Closes the Edit Polygon dialog box.





File | Save

Allows you to specify a name for the polygon file and save the new data.



The following table provides a brief description for each of the buttons on the Edit Polygon dialog box.


Edit Polygon Tools

Button	Name	Functionality
	Select polygon	Selects, moves, or modifies existing polygons or points within a polygon.
	Add polygon	Connects points which you select using the mouse.
	Add polygon (freehand)	Allows you to draw the polygons freehand. It is especially useful when used with Guided mode.
	Automated polygon creation tools	Allows you to define parameters for the automatic detection of polygons.
	Insert point into polygon	Adds a point to a line segment in a selected polygon. Drag and drop this point to change the shape of the polygon.
	Close current polygon	Connects the last point selected with the first point in order to close the polygon.
	Add circle	Allows you to define a circular area.
	Add box	Allows you to define a rectangular area.
	Delete selected polygon point	Deletes the selected or most recently added point.
	Delete selected	Deletes the selected polygon.

polygon		
	Delete all polygons	Deletes all polygons.
	Copy polygon for pasting	Allows you to select a polygon to copy.
	Paste polygon	Pastes the polygon that is selected for copying.
	Undo last action	Restores the last deleted object or moves the object to its former position.

To create polygons using Snap Selected in Polygon Editor:

1. Open the desired image data file.
2. Click **Model | Edit Polygon** in the *softWoRx* main menu.
3. Drag the Image window number to the **Window** field.
4. Click the **Add Polygon**  button.
5. Click points to form the polygon. To connect the last point with the first, click the **Close Polygon**  button.
6. Select **Guided Mode**.
7. Select **Options | Guided Mode Options** on the Edit Polygon dialog box. The Guided Polygon Select Parameters dialog box is displayed.

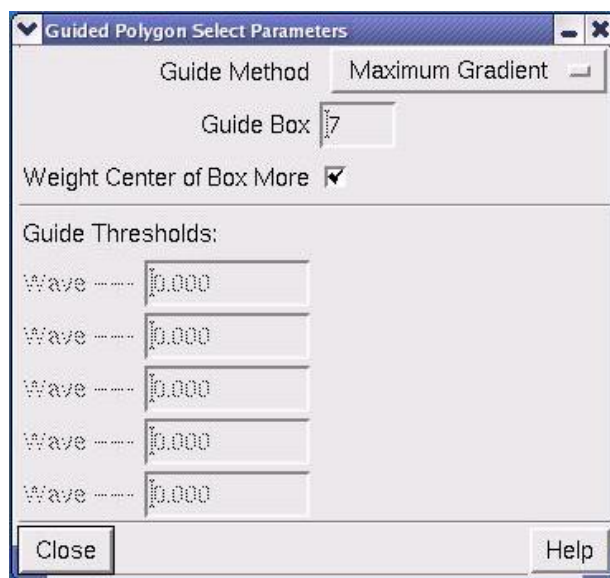


8. Adjust the values in the Guided Polygon Select Parameters dialog box. For more information on this topic, consult the online Help.
9. Click **Close**.
10. In the Edit Polygon dialog box, click **Snap Selected**.
11. Repeat this process for each Z section that will be used in the 3-D model.
12. In the Edit Polygon dialog box, select **File | Save**.
13. Enter a name for the polygon file and click **OK**.
14. Use the polygon file to create a 3-D model using the 3D Object Builder.

2D Polygon Finder

The first task that must be performed is to isolate the object that you want to study. This is accomplished by finding the 2-D representation in each Z section and combining them into a 3-D object. Setting a threshold value in the wavelength intensity and allowing 2D Polygon Finder to create a polygon in each Z section can often isolate two-dimensional features in a Z series. A simple adjustment of the threshold value allows you to modify the polygon created. The *softWoRx* Polygon Editor includes additional tools for more complex adjustments.

The 2D Polygon Finder Dialog Box

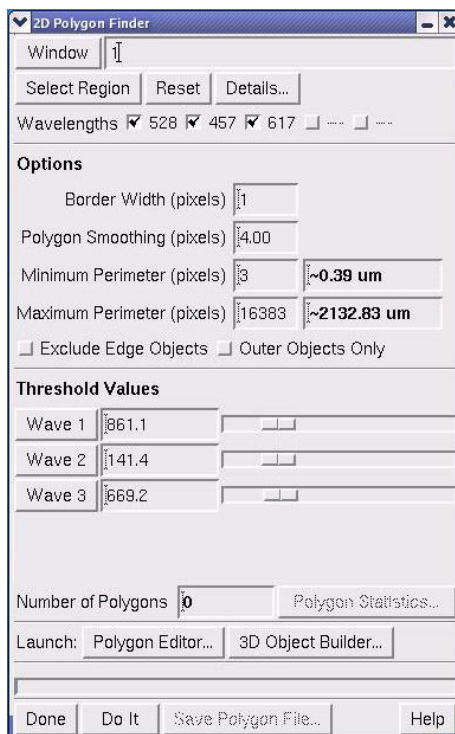

The main options in the 2D Polygon Finder dialog box are summarized below. For information regarding the other options and for additional details on these, refer to the online Help.

Window

Sets the Image window number of the image data to be processed.

Select Region

Defines the region of interest.

Reset

Erases the selected region.

Minimum Perimeter

Sets a value for the minimum number of points that can define a polygon. Polygons with less than this number of points will be discarded.

Maximum Perimeter

Sets a value for the maximum number of points that can define a polygon. Polygons with more than this number of points will be discarded.

Polygon Smoothing

Sets the desired number of pixels between points on the polygon. The 2D Polygon Finder will use greater detail when able.

Exclude Edge Objects

Specifies not to use those polygons that touch the edge of the image.

Outer Objects Only

When enabled, specifies that 2D Object Finder will only find the outermost continuous polygons. When disabled, specifies that polygons that are fully contained within others will be created if detected.

Wave # Threshold

Defines the minimum intensity to be included in the polygon. (Default Value: Minimum Value for the Wavelength + 20% of the Dynamic Range).

Launch

Opens either the Polygon Editor, to enable more sophisticated selection of a polygon, or 3D Object Builder, to continue creating the three-dimensional model.

Done

Closes the 2D Polygon Finder.

Do It

Begins the process of creating polygons.

Save Polygon File

Allows you to specify a name for the polygon file and save the new data.


To create polygons using 2D Polygon Finder:

1. Open the desired image data file.
2. Select **Model | 2D Polygon Finder** in the *softWoRx* main menu. The 2D Polygon Finder dialog box is displayed.
3. Drag the Image window number to the **Window** field.
4. Click **Select Region**.
5. Using your mouse, draw a box around the region of interest. It may be helpful to scroll through the Z sections to ensure that all of the desired areas are included.
6. Select the desired wavelengths in the **Wavelengths** check boxes.
7. Enter the desired values in the **Threshold** fields. In order to estimate an initial threshold value, use **Point Values** (located in the **Tools** menu of the Image window) to view the image intensity at various points.
8. Click **Do It**.
9. Click **Save Polygon File**. Then enter a name for the polygon file and click **OK**. By default, *softWoRx* uses the previous file name and replaces the file extension with POL.
10. Create a 3-D model using the 3D Model Builder or use Polygon Editor to create new polygons or modify existing ones.

3D Object Builder

The 3D Object Builder joins the 2-D polygons together to create a three-dimensional model. This is helpful for quantitative analysis as well as visual understanding of the image data.

The 3D Object Builder Dialog box

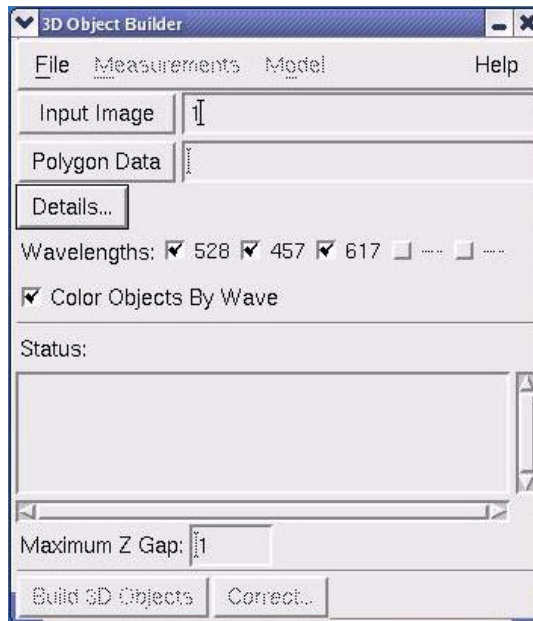

Information loaded into 3D Object Builder must include the polygon data. The two methods of accomplishing this are described in the following procedures.

To load data from an Image window:

- ▶ Select **Model | 3D Object Builder** from the *softWoRx* main menu and drag the desired Image window number into the **Input Image** field.

To load data from a saved polygon file:

1. Load the image file which corresponds to the polygon file into the **Input Image** field by dragging and dropping the Image window number.
2. Click **Polygon Data** and select the appropriate polygon file. (You may move up a directory level by clicking on the path bar above the desired directory.)
3. Click **OK**.

Creating and Viewing the 3-D Object

Once the polygon file is loaded into the 3D Object Builder, you are ready to create the 3-D object. Before a 3-D object can be viewed, it must be saved as a solid model. Before the measurements can be viewed, it must be saved as a measurement file. These measurement files are saved as tab delimited text files that can easily be exported to spreadsheet programs such as Microsoft Excel.


To build and save a 3-D object:

1. Select **Model | 3D Object Builder** on the *softWoRx* main menu.

2. Load the data into the 3D Object Builder as described in the previous section.
3. Choose the desired wavelengths to be modeled.
4. Click **Build 3D Objects**.
5. Click **Model** in the 3D Object menu.
6. Click **Save Solid Model**.
7. Enter the desired name in the field. By default, *softWoRx* will use the previous file name and replace the file extension with SOL.
8. Click **OK**.

To view a 3-D object (not available without optional 3D Model):

1. Click **Model** in the 3D Object Builder menu.
2. Click **View Model**.

 **Note** You must have the original Image window open in order to view the 3-D model.

To measure the area of an object:

1. Click **Measurements** in the 3D Object Builder menu.
2. Click **Table of 2D Measurements**. The Save Measurements File dialog box is displayed.
3. Type the desired name for the 2-D measurement file.
4. Click **OK**.
5. Open the folder containing the saved measurement file.
6. Double-click the icon of the desired measurement file to view a text file similar to the following.

wave	time	sec	poly	pixs	integr_int	area	centroid	center_mass
wave 1 time 1								
1	1	1	1	31	18150	0.17	21.75 9.61	21.75 9.61
1	2	1	2	57	33576	0.31	21.34 19.35	21.34 19.35
2	1	2	1	27	15996	0.15	21.63 9.41	21.63 9.41
2	2	2	2	43	25218	0.24	21.42 19.50	21.42 19.50
3	1	3	1	26	15340	0.14	21.52 9.20	21.52 9.20
3	2	3	2	38	22233	0.21	21.64 19.26	21.64 19.26
4	1	4	1	18	10615	0.10	21.56 9.36	21.55 9.36
4	2	4	2	24	14149	0.13	21.64 19.22	21.64 19.22
5	1	5	1	36	21250	0.20	21.45 9.30	21.45 9.30
5	2	5	2	42	24669	0.23	21.53 19.26	21.53 19.26
6	1	6	1	29	17030	0.16	21.58 9.67	21.57 9.66
6	2	6	2	24	14156	0.13	21.38 19.39	21.38 19.39
7	1	7	1	29	17069	0.16	21.47 9.57	21.46 9.57
7	2	7	2	38	22388	0.21	21.40 19.28	21.40 19.28
8	1	8	1	29	16953	0.16	21.62 9.87	21.62 9.87
8	2	8	2	25	14748	0.14	21.59 19.10	21.58 19.10
9	1	9	1	41	24188	0.23	21.40 9.57	21.40 9.57
9	2	9	2	47	27705	0.26	21.43 19.18	21.42 19.18
wave 2 time 1								
1	1	1	1	263	173347	1.45	19.61 14.99	19.61 14.98
1	2	1	2	6	4014	0.03	13.49 20.04	13.49 20.04
2	1	2	1	261	171848	1.44	19.88 14.32	19.88 14.32
3	1	3	1	215	141459	1.19	19.94 14.68	19.94 14.68
4	1	4	1	217	142986	1.20	19.91 15.05	19.91 15.05
5	1	5	1	185	122070	1.02	20.18 13.84	20.18 13.84
6	1	6	1	221	145557	1.22	20.07 14.66	20.06 14.67
7	1	7	1	184	121316	1.02	20.30 14.09	20.30 14.09
8	1	8	1	211	139238	1.16	20.09 14.97	20.09 14.96
9	1	9	1	202	133238	1.12	20.23 14.62	20.23 14.63
wave 3 time 1								
sec	poly	pixs			integr_int	area	centroid	center_mass

To measure the volume of an object:

1. Select **Measurements** in the 3D Object Builder menu.
2. Select **Table of 3D Measurements**. The Save Measurements File dialog box is displayed.
3. Enter the desired name for the 3-D measurement file.
4. Click **OK**.
5. Open the folder containing the saved measurement file.
6. Double-click the icon of the desired measurement file to view a text file containing the data.

Volume Modeling Example

The following steps show how to use the Polygon Finder and 3D Object Builder in a specific set of image data. The image data file `Nuclear_Pore_D3D` (included with your system) is used in this tutorial.

1. Open Nuclear_Pore_D3D.
2. Select **Model | 2D Polygon Finder** in the *softWoRx* main menu.
3. Drag the Image window number to the **Window** field.
4. Click **Select Region**.
5. Using your mouse, draw a rectangle around the region of interest. It may be helpful to scroll through the Z sections to ensure that all of the desired areas are included.
6. Select the wavelength 457 in the **Wavelengths** check box.
7. Type **750** in the **Wave 3: Threshold** field. In order to estimate an initial threshold value, use **Point Values** in the **Tools** menu in the Image window menu to view the image intensity at various points.
8. Click **Do It**.
9. Click **Save Polygon File**.
10. Type a name for the polygon file and click **OK**.
11. In the **Launch:** field, click **3D Object Builder**. The 3D Object Builder dialog box is displayed. Notice that *softWoRx* automatically loads the polygon data from the open Image window which was used to make the polygon.
12. Select the wavelength 457 in the **Wavelengths** check box.
13. Click **Build 3D Objects**.
14. Click **Model | Save Solid Model** in the 3D Object Builder menu.
15. Type the desired name in the **File Name** field and click **OK**.
16. Select **Model | View Solid Model** to view the model.
17. Use the center mouse button to rotate the model for viewing from other angles.

15. Detecting and Analyzing Colocalization

softWoRx provides two tools that you can use together to detect and analyze colocalization.

First, use the Colocalization tool to examine the entire image and identify areas that appear to have colocalized data. You can use the data generated by this tool to create volume views and graphically examine them to find structures or specific areas that appear to be colocalized.

Then use the ROI Colocalization tool to examine the specific structures or areas that you have identified. The data selection features of this tool can be used to select many types of areas.

In This Chapter


Examining the Entire Image	194
Identifying Potential Colocalized Areas	197
Detecting Colocalization with ROIs.....	201

Examining the Entire Image

Use the Colocalization tool to identify possible areas of colocalization throughout the data set. This tool generates a product image of two channels after subtracting a threshold value for each. Then a scatter plot of the results is created and the Pearson Coefficient of Correction is measured.


To use Image Colocalization:

1. Open the image to analyze in the Image window.

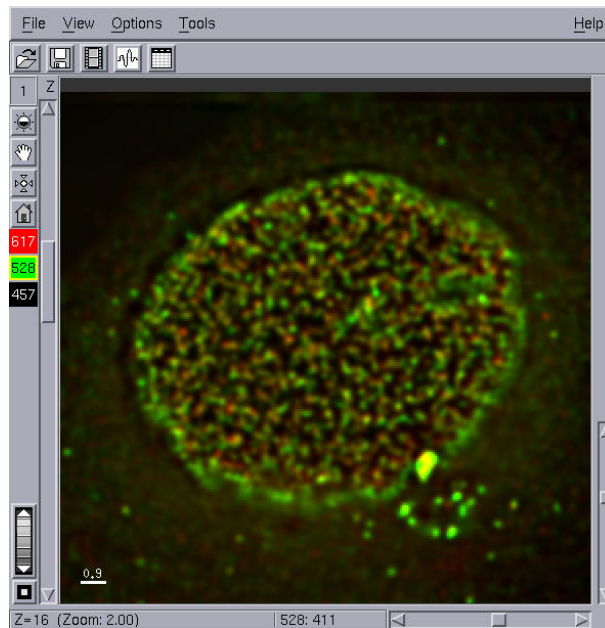


The Nuclear Pore image displays two proteins: channel 528 is tagged to a protein that regulates the gateway to the cell. Channel 617 is VOM, an HIV protein.

2. Choose **Measure | Colocalization** from the *softWoRx* main menu to open the Colocalization dialog box and enter the number of the Image window that you want to analyze in the Input field.

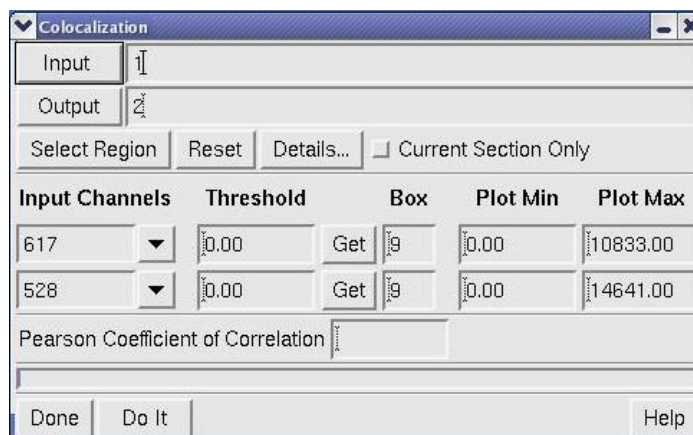

- To analyze a region of a window, click **Select Region** and select an area in the Image window by dragging the mouse across the area. Adjust the rectangle you've created until it contains the desired area. Then click outside the Image window with the mouse.

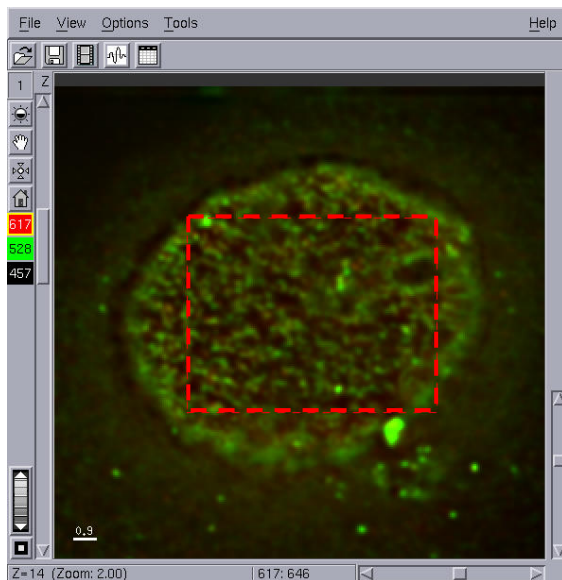


- Click **Details** to open the Region Details dialog box. Specify the ranges of the X, Y, Z, and time data to analyze in the selected region.




For example, in this case we are interested in colocalization that is occurring below the surface of the cell.

- Select which channels to analyze in the **Input Channels** lists.
- Select a background threshold for each channel by clicking **Get** and then clicking on a background area of the image. The background is an averaged value within a box of the size specified. Click **Get** again when you are finished selecting the background.

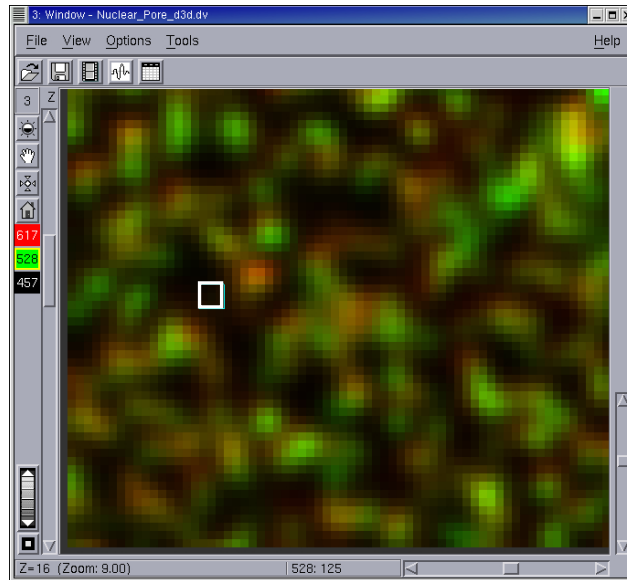


Tip You can change the size of the box.

- Click **Do It** to run the colocalization analysis for the selected data.

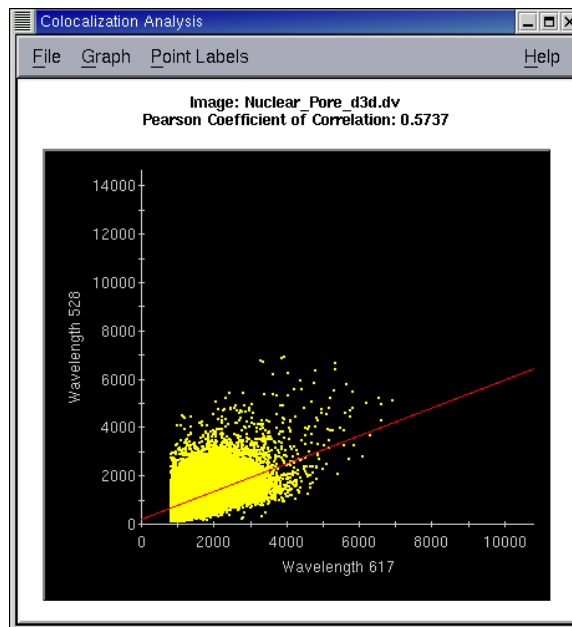

The Colocalization Graph and a new window with the colocalization data are displayed

The Colocalization graph is a plot of the two intensities on a pixel-by-pixel basis (each spot is a pixel). The Pearson Coefficient of Correlation indicates how closely the two intensities are colocalized (full colocalization is 1.0) and calculates the Pearson Coefficient of Correlation. This value is displayed in the Colocalization dialog box.

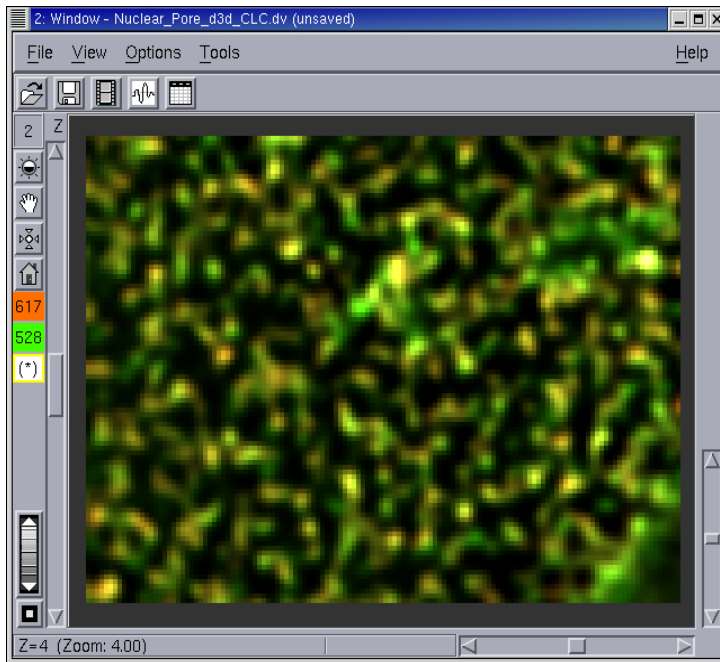


The new window contains a third channel (*) that is the product of the two intensities at each data pixel. This channel indicates possible areas of colocalization. (If the intensities of both channels are high for a given pixel, the product of the intensities is high. If one of the intensities is low or zero, the product is much lower.)

Identifying Potential Colocalized Areas


To visually identify areas that may be colocalized, you can select specific points or groups of points to display them on the three channel image. You can also render a volume projection of the three channel image that includes the selected product channel in white.

To identify colocalized structures in Image Colocalization Tool data:


1. Select the points on the colocalization graph that have higher intensities by dragging the mouse across the graph as shown below.

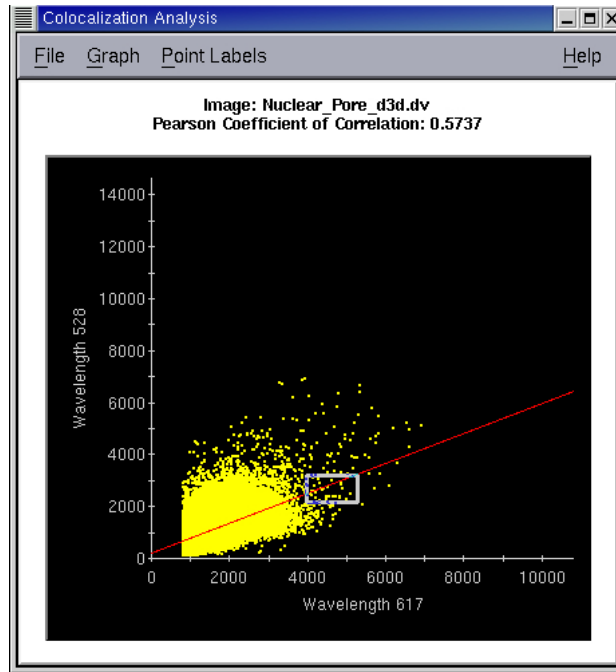


2. Select additional points by holding the CTRL button as you drag the mouse across additional areas.

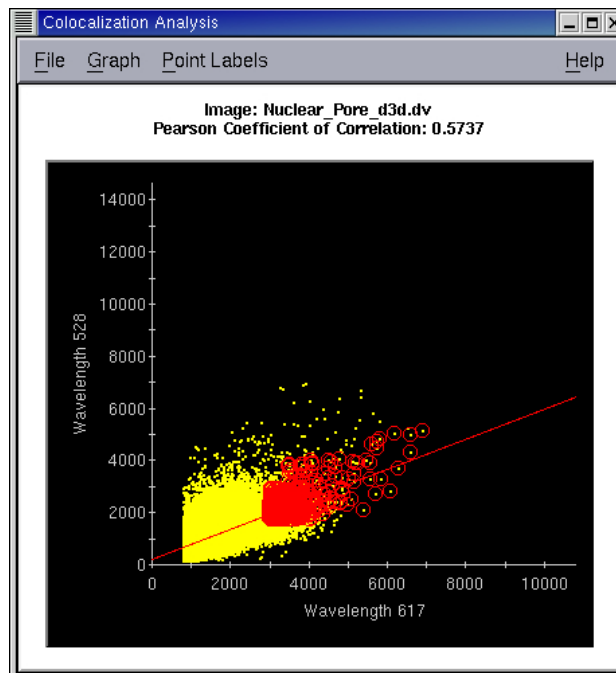


As you select points, they are highlighted in red on the image graph and displayed in white on the three channel image.

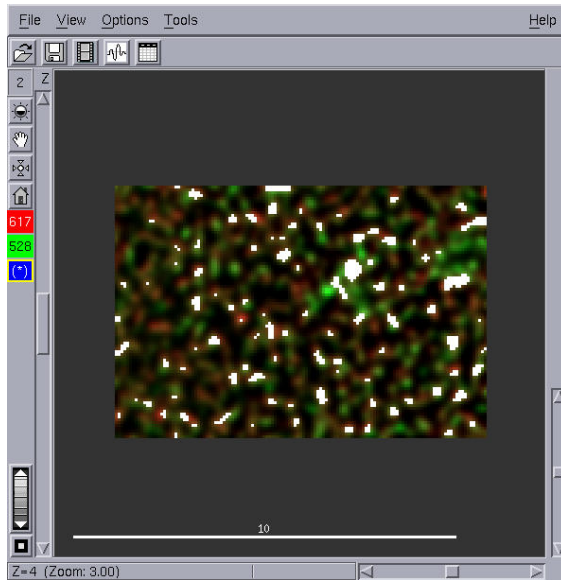

3. Choose **View | Volume Viewer** on the *softWoRx* main menu to open the Volume Viewer window and drag the output window number (post-colocalization) into the **Input** field.
4. Select Volume Viewer parameters and click **Do It** to render the volume.

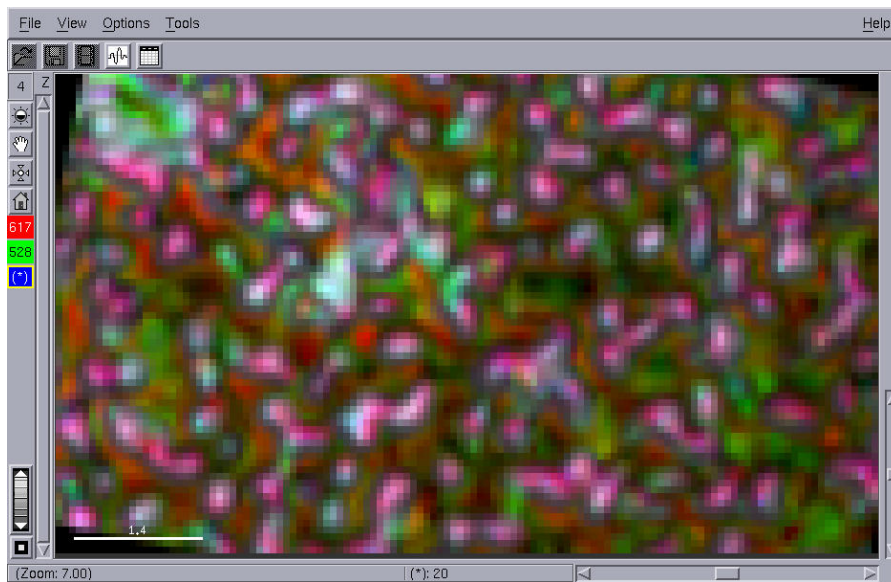

5. Click **Interactive** on the Volume Viewer and examine the image from several angles to find intense white areas that indicate potential colocalization.

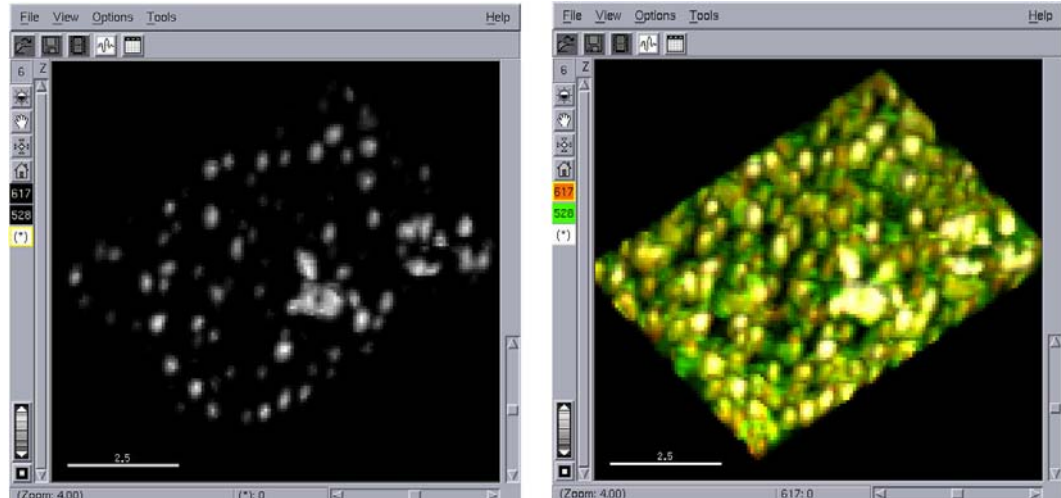


Interactive Viewer shows product channel and original data

6. To view only the colocalized channel, select the (*) channel.


Alternatively, you can view the colocalized channel by changing the grayscale color map for the (*) channel to a rainbow or cold to hot color map so the bright intensity is displayed in a different color.

To change the grayscale color map (from the previous example):


1. While viewing only the (*) channel, choose **View | Select Image Colors** on the output window menu. The Select Image Colors dialog box is displayed.

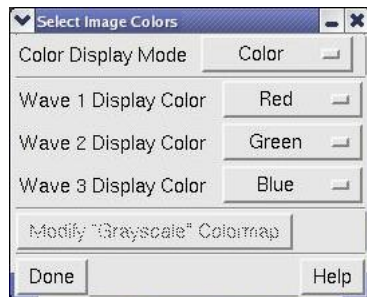

2. In the Color Display Mode field, select **Grayscale**. The **Modify \"Grayscale\" Colormap** button is activated.

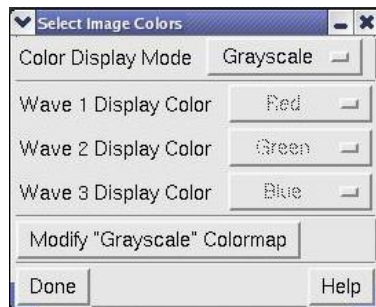

- Click the **Modify “Grayscale” Colormap** button. The Select Image Colormap dialog box is displayed.

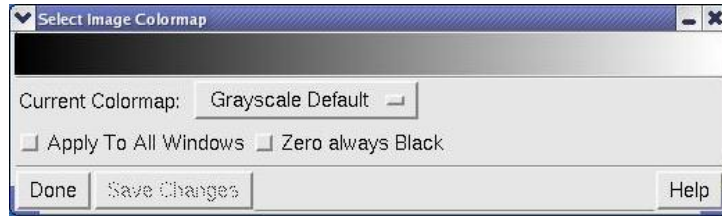

- In the Current Colormap field, change the color map from **Grayscale Default** to one of the rainbow or temperature color maps (for example, **Rainbow 1** or **Cold To Hot**).


Detecting Colocalization with ROIs

After you use the Colocalization tool to identify possible areas of colocalization, you can use Region-Of-Interest (ROI) Colocalization to selectively analyze those areas.


ROI Colocalization allows you to create ROI polygons from which a scatter plot and Pearson Coefficient of Correlation are derived.

To use ROI colocalization:


- Choose **Measure | Colocalization (ROI)** from the *softWoRx* main menu to open the Colocalization Analysis dialog box. Enter the number of the Image window that you want to analyze in the **Input Window** field.

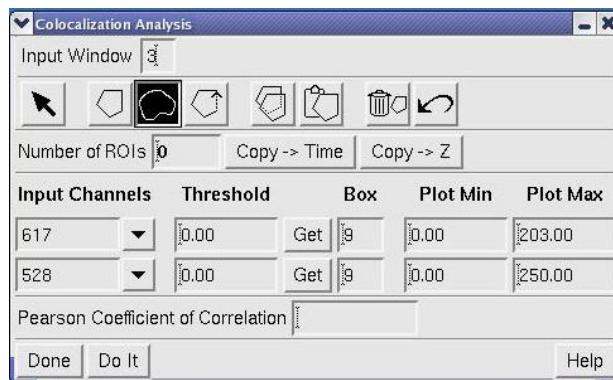

- Select which channels to analyze in the **Input Channels** lists.
- To select specific areas of the image, use the **Create Freehand ROI Polygon** button to create polygons on each section that you want to analyze. If the position and shape of the structure are consistent through time and Z intervals, you can use the **Copy Selected Polygon** and the **Paste Polygon** buttons to copy and paste the polygons through time points or Z sections.

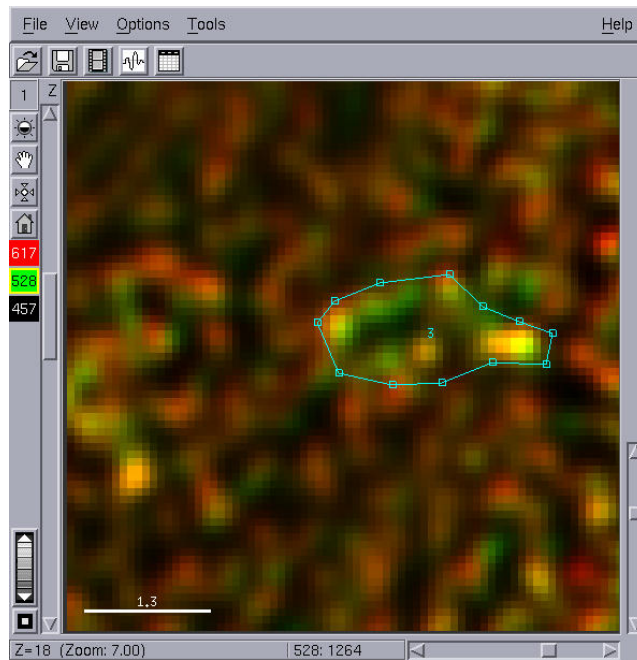

4. Select a background threshold for each channel by clicking **Get** and then clicking on a background area of the image. The background is an averaged value within a box of the size specified. Click **Get** again when you are finished selecting the background.
5. Click **Do It** to plot the Colocalization graph and calculate the Pearson Coefficient of Correlation value for the selected region (this value is displayed in the Colocalization Analysis dialog box).

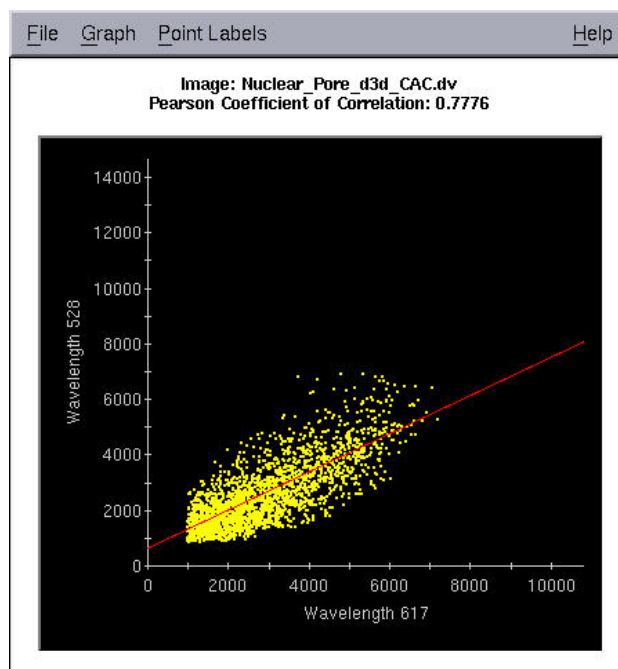

16. Other Applications

This chapter shows how to analyze experiments that are performed with the optional Quantifiable Laser Module (QLM).

Contents in this Chapter

About Photokinetics.....	203
Analyzing Fluorescence Recovery After Photobleaching	204
Analyzing Fluorescence Resonance Energy Transfer	208

About Photokinetics

Photokinetics refers to the reactivity of fluorescent molecules while they are in the excited state. Photokinetic reactions can be used to study the interactions of molecules within living cells.

Photo-bleaching, FRET, and photo-activation are examples of photokinetic reactions.

The following table shows photokinetic experiment methods and which biological applications can be studied with those methods.


Photokinetic Methods and Applications

Photokinetic Methods	Biological Applications
FRAP Single-Point and Multi-Point	Affinity, Biomolecular Cycling, Biomolecular Environment, Structural Kinetics
Pattern Bleaching	Compartmental Analysis, Biomolecular Cycling, Transport
FLIP	Compartmental Analysis, Biomolecular Cycling, Transport, Structural Visualization
Background Reduction	Structural Visualization
Combinations -FRAP/FRET -Repeat during cell cycle -Rapid Repeat	Compartmental Analysis, Biomolecular Cycling, Transport Affinity, Biomolecular Cycling, Biomolecular Environment
FRET -Sensitized emission -Donor Photo-bleaching -Acceptor Depletion	Affinity, Biomolecular Environment
Photo Activation	Compartmental Analysis, Affinity, Biomolecular Cycling, Biomolecular Environment, Transport, Cell Fate, Structural Kinetics, Structural Visualization

Analyzing Fluorescence Recovery After Photo-bleaching

The Fluorescence Recovery After Photo-bleaching (FRAP) experiment method consists of photo-bleaching a point (or points) of interest and then observing the recovery of fluorescence in the bleached area. For detailed instructions, see the *Fluorescence Recovery After Photo Bleaching* Product Note at: www.appliedprecision.com.

An example of a Single-Point FRAP experiment is shown below.

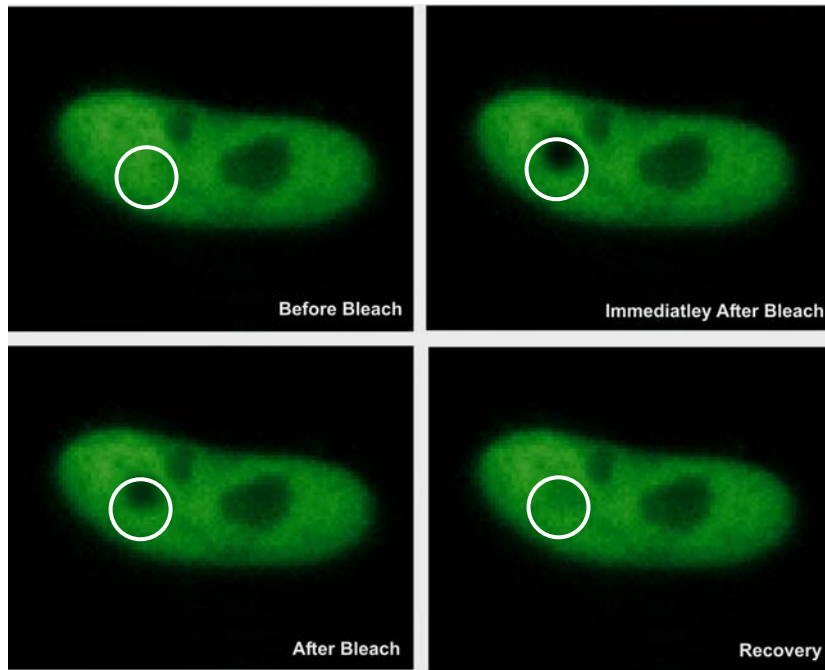

A point of interest is photo bleached and monitored

About FRAP Experiments

There are two types of FRAP data:


- Single-Point FRAP data is collected in experiments that monitor a single location or monitor several locations in a sequential fashion.
- Multi-Point FRAP data is collected in experiments that monitor several locations in the sample at the same time.


Single and Multi-point FRAP experiments can be used for the following types of studies:

- Affinity Studies
- Biomolecular Cycling
- Biomolecular Studies
- Environment Studies
- Structural Kinetics

To analyze FRAP data:


1. Open an image in the Image window. From the *softWoRx* main menu, choose **Measure | PK Analysis**. The Photokinetic Data Analysis dialog box is displayed.

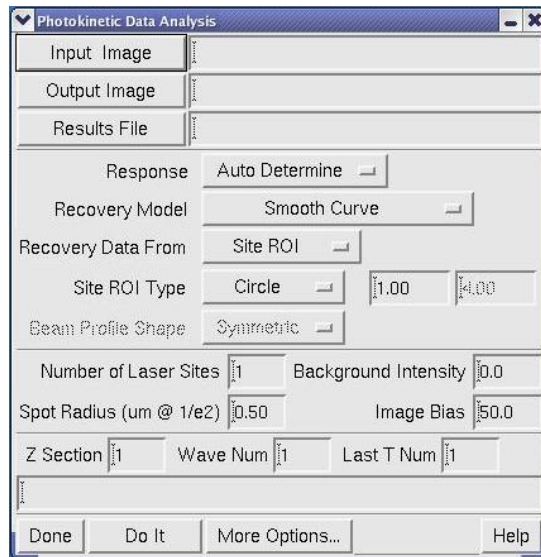

 **Tip** You can use the Photokinetic Data Analysis dialog box to specify a recovery model, type of ROI, beam profile shape, and number of sites. You can also use it to remove background intensity, select which Z sections and wavelengths to include in the analysis, and specify other options.


2. Drag the window number into the Input Image field.
3. Select the desired Response type, Recovery Model, and Data Recovery options.
4. If you are analyzing a Multi-point FRAP data set, enter the number of laser sites in the **Number of Laser Sites** field.
5. If you are using background subtraction, enter a background value in the **Background Intensity** field.
6. To determine a number to enter in the **Spot Radius** field, use the Measure Distances tool to make an approximate measurement of the bleach spot.
7. Click **Do It** to run the analysis. The following files are generated:

The three channel output image file includes the original time lapse image, the ratio of the current time point to average pre-bleach time points, and the ratio data at the location used for analysis.

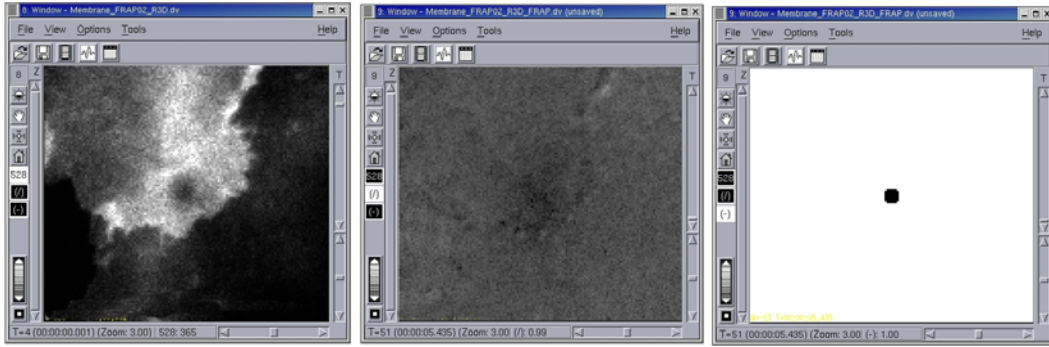

The three channel image generated by the Photokinetic Data Analysis tool

A JPEG file contains a recovery graph that is a plot of the fluorescence intensity before and after the event.

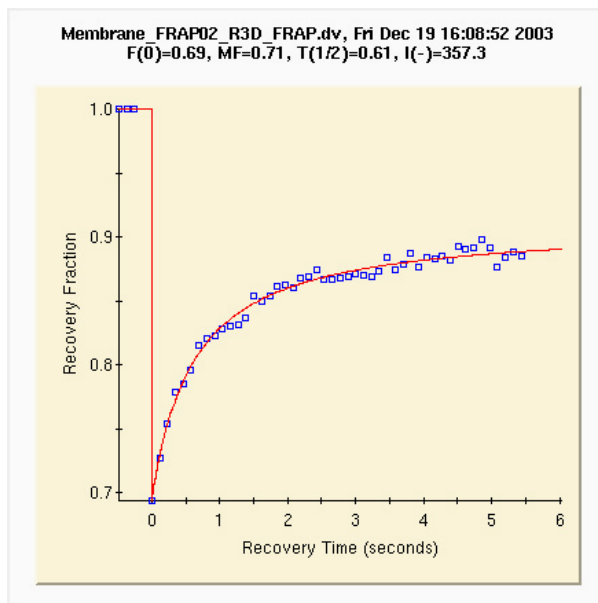

The Fluorescence Recovery Graph

A log file contains the analysis results.

```

Membrane_FRAP02_R3D_FRAP.log
-----
FRAP Results:
Fri Dec 19 16:08:46 2003
-----
Input Image: /data1/InSitu2003/Membrane_FRAP02_R3D_FRAP.dv
Result Image: 9.dv
Result File: /home/worx/Membrane_FRAP02_R3D_FRAP.log

FRAP Best-Fit Results - Single Component:
-----
Bleach Point: 1
Location (X,Y): 16.3, 16.3 um
Post Bleach Fraction: 69.388 % (Bleached Frac= 30.61%, K=0.781, B=1.076)
Mobile Fraction: 70.677 %
Final Fraction: 91.023 %
Beam Radii (X,Y): 0.500, 0.500 um
Time Constant: 0.565 secs
Half-time: 0.608 secs

Chi-squared: 2301.473
Avg Error: 0.022 +/- 1.757 counts
Avg Abs Error: 1.435 +/- 0.991 counts

FRAP Measurement Information:
First Measurement: 69.435 % at 0.0010 secs (0.002 half-times)
Last Measurement: 88.465 % at 5.4350 secs ( 8.94 half-times)
Experiment Duration: 5.434 secs
Num Pre-bleach Pnts: 3
Num Post-bleach Pnts: 48
Time Zero Offset: 0.000 secs
Image Bias: 50.000 counts
Background Intensity: 0.000 counts
Photosensor Normalization: Off
Mean Inten. Normalization: On
Pre-bleach Statistics: 1.000000 +/- 0.000141 (fractional)
Pre-bleach Statistics: 357.273 +/- 0.051 (counts)

Recovery Measurements:
Tn Ts ModelF ModelI RecoveryF RecoveryI
-3 -0.482 1.000 357.273 1.000 357.216
-2 -0.365 1.000 357.273 1.000 357.292
-1 -0.252 1.000 357.273 1.000 357.312

```


The FRAP Analysis log file

Analyzing Fluorescence Resonance Energy Transfer

FRET (Fluorescence Resonance Energy Transfer) is a method for determining whether two types of molecules are in close proximity. FRET occurs when there is a quantum physical exchange of energy between dipoles. The presence of FRET indicates that the molecules are within 60 \AA (6 nm).

FRET is orientation specific. Negative FRET does not mean the molecules are not interacting. Positive FRET means that they are close, but does not necessarily mean that the molecules are interacting.

There are many ways to measure FRET. Experiments are simple, but controls are essential.

 **Tip** Before you conduct a FRET experiment, consider using the Colocalization tool to determine whether the molecules that you are studying are close to each other. (Colocalization experiments provide more approximate results than FRET, but they are much simpler to prepare and analyze.)

Using the FRET Analysis Tool

The FRET analysis tool provided assists in analysis of sensitized emission FRET experiments. Acceptor photo-bleaching can also be done with Deltavision. Use the Ratio Analysis tool or Polygon Editor to analysis acceptor photo-bleaching experiments.

Before You Begin

Be sure to follow the conventions described in the online Help topic *Acquiring and Preparing FRET Data*.

Before you begin, make sure that you have the following data:

You must have the following three image files:


- A Donor Control data file that contains data from a slide prepared for the Donor probe only.
- An Acceptor Control data file that contains data from a slide prepared for the Acceptor probe only.
- The FRET Experiment file that contains data from a slide prepared for both Donor and Acceptor probes.

Each image must have the following three channels:

Channel	Description
1	Donor (Donor Excitation, Donor Emission)
2	Acceptor (Acceptor Excitation, Acceptor Emission)
3	FRET (Donor Excitation, Acceptor Emission)

Analyzing FRET Data

Analysis of Direct FRET data consists of determination of Donor and Acceptor crosstalk factors, calculation of Net FRET and FRET efficiency, and analyzing the Net FRET and FRET Efficiency statistics. Use the FRET analysis tool and the following process to analyze the FRET data:

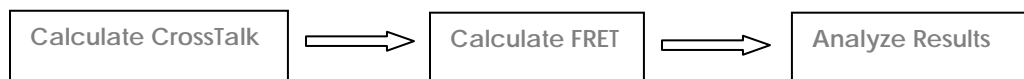


Calculating Crosstalk

Calculation of Donor and Acceptor crosstalk factors involves creating Region of Interest (ROI) polygons on Donor and Acceptor Control images and determining representative background values to be subtracted from intensities in ROIs.

To calculate crosstalk:

1. From the *softWoRx* main menu, choose **Measure | FRET Analysis**. The FRET Analysis dialog box is displayed. Make sure Calculate Crosstalk is the active tab.

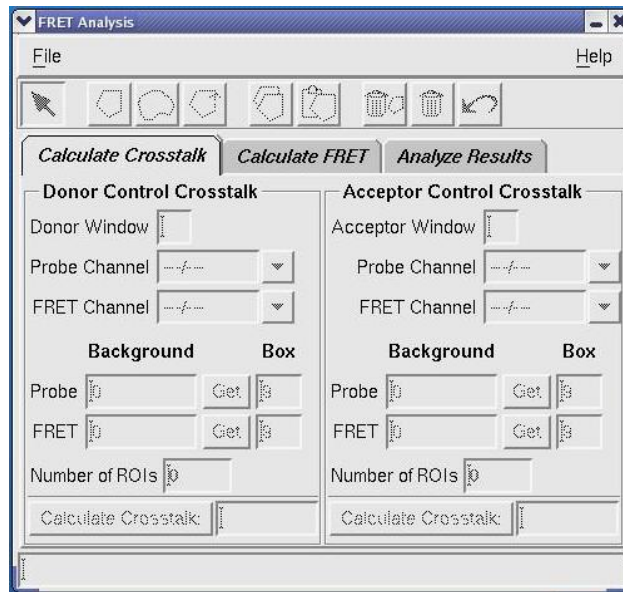

2. Open the Donor and Acceptor images in Image windows.
3. Drag the Image window number icon from each of the Donor and Acceptor Image windows to the appropriate Donor and Acceptor **Window** fields in the FRET Analysis dialog box to connect to these windows.
4. Validate that the Probe Channel for the Donor Control Crosstalk is set to the correct (Donor) channel and the Acceptor Channel for the Acceptor Control Crosstalk is set to the correct (Acceptor) channel.
5. Validate that the **FRET Channel** is set to the correct channel for each control image.
6. Specify the background for the **Probe** and **FRET** channels of each image using the **Get** buttons. The "Get" functions average intensity values in a box (specified in the **Box** fields) while you drag the cursor around in the Image window with the left mouse button held down. Click the **Get** button a second time to disable the "Get" function.
7. For each image, define ROIs using the tools on the toolbar at the top of the FRET Analysis dialog box to define representative areas where FRET would occur if these were Experiment images.
8. Click **Calculate Crosstalk** to analyze the background and defined ROIs to generate a crosstalk factor for both the Donor and Acceptor control images. Typical Donor Crosstalk factors are 50-70%; typical Acceptor crosstalk factors are 15-30%, if using CFP and YFP and the FRET pair.


The Donor and Acceptor crosstalk factors are loaded in to the appropriate fields for FRET calculation.

Calculating FRET

Calculation of Net FRET and FRET Efficiency uses a FRET Experiment image as input along with Donor and Acceptor Crosstalk factors and background values for each channel. It generates an image with 2 channels: Net FRET and FRET Efficiency (%E).

To calculate FRET:

1. Choose the Calculate FRET tab on the FRET Analysis dialog box.

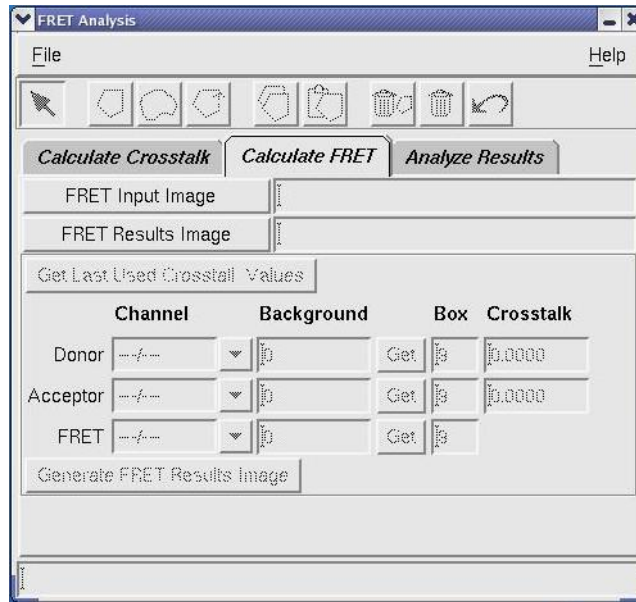


2. Open the FRET Experiment image you would like to analyze in an Image window.
3. Drag the Image window number icon from the FRET Experiment Image window to the **FRET Input Image** field in the FRET Analysis dialog box. An output FRET Results image file name is automatically generated. (The FRET Results image is saved to the disk and an Image window containing the saved image is displayed after the calculation is completed).
4. Validate that the correct channels have been assigned to the Donor, Acceptor and FRET channels of the input image.
5. Validate that the crosstalk factors are reasonable (These factors were calculated when you calculated crosstalk).
6. Use the "Get" function to specify a background value for each of the 3 channels (see Step 6 in *To calculate crosstalk* on page 209).
7. Once you are satisfied that all parameters are set up correctly, click **Generate FRET Results Image**. When this process is finished, the image opens.

Analyzing Results

Analyzing Net FRET and FRET Efficiency statistics involves specifying a FRET Results Image and one or more ROI polygons to generate a table and graph of statistics.

To analyze FRET results:

1. Choose the Analyze Results tab of the FRET Analysis dialog box.

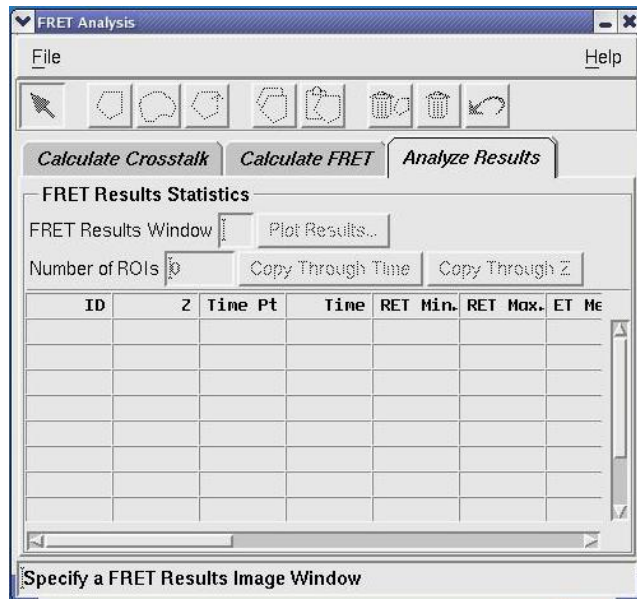

2. If it isn't already being viewed, open the FRET Results image that you would like to analyze in an Image window.
3. If it isn't already specified, drag the window number of FRET Results Image window to the appropriate **Input** field in the FRET Results Statistics area.
4. Use the ROI creation tool icons at the top of the FRET Analysis dialog box to specify one or more regions of interest.
5. If you have an image with multiple time points or multiple Z sections, you may want to propagate ROI polygons through time or Z. To do this, create a polygon and make sure it is selected. Then, choose **Copy Through Time** or **Copy Through Z** to propagate the polygons. (As changes to ROI polygons are made, the table of statistics is updated to reflect the statistics of the chosen ROI set.)
6. To export the table of numbers in a form that can be used in a spreadsheet, choose **Save Results As SYLK** (Symbolic Link format) or **Save Results As CSV** (Comma-Separated Values) from the **File** menu on the FRET Analysis dialog box.

7. To view the results in a graph, choose **Plot Results** to generate an X/Y plot of parameters that you choose. If the X axis is time-related, the software associates ROIs from time point to time point and plots the values as connected sets on the graph.

 **Tips**


#1 You can modify the details of how the graph is displayed after it is created with the Graph-Properties tool.

#2 You can optionally save the FRET Results graph by selecting Save As JPEG from the FRET Results graph's File menu.

Appendix A: Image Quality

Because many restoration problems can be attributed to problems with data collection, you'll find it worthwhile to become familiar with the common problems and their solutions.

All of these methods are simple to perform on a regular basis.

 **Note** Although the *softWoRx* restoration algorithms contain many refinements that improve their ability to handle experimentally obtained optical sections, there are certain types of data that simply cannot be properly restored.

In This Appendix

Using Deconvolution Residuals	216
Visually Evaluating Images	217

Using Deconvolution Residuals

What is a Residual?

The deconvolution residual is a measure of the difference between the measured image and the solution convolved with the point spread function (PSF). In mathematical terms:

$$\text{Residual} = (\text{Measured Image}) - (\text{Deconvolved Image} * \text{PSF})$$

where * represents convolution.

In principle, the two quantities on the right side of the equation are equal, so that the residual should be zero. Not surprisingly, however, the use of experimental data prevents perfectly precise results and the residuals are not zero.

For the purpose of digital deconvolution, the residual is calculated from the average of the residuals measured at each point in the three-dimensional image. As a general rule, a small residual is better than a large residual.

The most useful form of the residual is the "Average Counts Residual," which is the average difference between the measured image and the result convolved with the PSF.

The Standard Residual

The "standard residual" is the sum of all residuals divided by the sum total intensity of the image. Images with a large total intensity may therefore yield an unrealistically small residual. For this reason, the standard residual may not be as useful as the average count residual. Use the standard residual to compare deconvolution performance results with deconvolutions prior to version 2.10 of the *DeltaVision softWoRx* software.

Refer to the following table when assessing the "standard residual."

Value	Quality	Suggested Action
> 0.1	Poor deconvolution quality	Check experiment conditions and data quality.
0.1 -0.05	Marginal Deconvolution	May not be appropriate for this data.
0.05-0.01	Reasonable	Typical for data with a low signal-to-noise ratio or a large amount of spherical aberration.
< 0.01	Good	

Normalized Residual

As a means of watching the deconvolution progress, the "Normalized Residual" is also displayed. By definition, the normalized residual equals 1 after the first

iteration. Subsequent residuals are then scaled in the same way as the first iteration to yield numbers between 0 and 1. A residual larger than 1 indicates that the deconvolution algorithm has encountered serious difficulties. The table below gives a guideline for assessing normalized residuals.

Value	Quality	Suggested Action
> 1.0	Worse than before the first iteration.	Review experimental conditions.
1.0	No improvement.	
0.5	Reasonable improvement.	
0.25	Substantial improvement.	
0.10	Excellent improvement.	

Most deconvolutions yield residuals somewhere between 0.1 and 0.3. Consistent results between 0.3 and 1.0 should prompt a review of experimental conditions and the deconvolution problems listed in *Visually Evaluating Images*. In particular, you should check pixel sizes, wavelength, and PSF selection.

Visually Evaluating Images

The second method of assessing a deconvolution is to simply study the resulting images. Comparison of the measured and deconvolved images is an excellent way of verifying that structures present in the results are a valid representation of the actual object. With the aid of the deconvolution image, it is usually possible to understand the structures present in the measured image.

It is expected that certain deconvolutions will be less successful than others, due to the dependence upon experimental data. It is not always possible to meet the conditions required for deconvolution microscopy. Fortunately, there are only a few characteristic problems.

Weak Convergence of the Residual

A common cause of poor convergence is that the optical sections were measured in the presence of spherical aberration. As a consequence, the standard PSF is not appropriate for deconvolution. Flip the image on its side (using the "Flip" or "Rotate" program) and study the quality of the image along the optical axis ("Z"). Asymmetric blurring and greatly elongated points indicate spherical aberration.

Use of very low intensity images (with a corresponding low signal-to-noise ratio) can also limit convergence.

Dark Halo Around Bright Structures

The halo problem is often caused by the presence of excessive spherical aberration. As always, ensure that the image properties are correct before beginning an exhaustive study of this sort of problem.

Use the Line Profile tool to assess the relative magnitude of the dark intensity region. Although very noticeable, the intensity of the halo is often quite small (e.g., about 5% below the adjacent background).

Another source of halos is refractive index changes.

Bright Spot Problem

In situations where the object contains concentrated areas of fluorescent probe, it is natural for the deconvolution process to yield even brighter spots in the resulting image. For example, a region with 2000 counts of fluorescence could deconvolve to a brightness of 8000 counts. The intensity of such areas can be so great that low intensity structures present in the 16-bit image are not visible on a standard 8-bit computer screen. Although these low intensity data exist, these data are simply not visible next to the bright spots. To view low intensity areas, adjust the image contrast, brightness, and intensity scaling factor.

Pebble Grain Texture

Low intensity texture patterns are often visible in low intensity images where the signal-to-noise ratio has dropped below about 10 to 1. Increasing the image intensity is the obvious method of avoiding this problem.

This can also be caused by problems with the CCD camera elements.

Deconvolution Holes

A bright ringed hole in background areas of the image is probably a result of subtracting too much background intensity during deconvolution. The "deconvolution hole" is the edge of the region that has reached the minimum possible intensity (typically 0). Deconvolve the image again using a background subtraction of 0, rather than the default value.

Increasing the "Z Transform Size" can also help control deconvolution holes. In some situations, the measured data may actually have a low intensity threshold caused by the loss of detector response. In this case, it is not possible to simply add intensity to the image. The edge of the low intensity threshold is indistinguishable from the edges being sought by the deconvolution process.

Poor Z Resolution

Elongated blurring in the Z directions is characteristic of spherical aberration. Also, note that the football shape along the Z-axis is normal for an optical microscope. The conical extensions above and below an object represent the uncertainty involved with measuring light through a lens with less than a 90 degree cone angle. At present, the best available lenses have a cone angle of approximately 68 degrees. Information between 68 and 90 degrees is not measured by the lens.

Section Intensity Fluctuation

Since each optical section is measured separately, it is common for the image intensity to be slightly different. To view this experimental situation, flip the optical sections on their side and look at the sections in the XZ plane. Striations perpendicular to the Z-axis indicate that the relative intensity of the sections varies. The striations are typically caused by arc lamp fluctuations during data acquisition.

Use the image correction program before deconvolution to compensate for arc lamp fluctuation. In some cases, the striations are visible after the image correction process. This is a serious problem that needs to be addressed! Such images will yield disastrous deconvolutions since the striations will be enhanced by the image processing.

Invalid Optical Transfer Function

There are at least two ways that an optical transfer function (OTF) can be invalid.

- The corresponding PSF may not have been well measured.
- Incorrect OTF may have been applied for deconvolution. Check that the lens ID number of the measured image corresponds to the lens ID of the OTF.

Index

- 2D Polygon Finder, 182–85
 - Dialog options, 186
 - Excluding edge objects, 186
 - introduction, 182
 - Minimum perimeter, 186
 - Smoothing pixel distance, 186
 - Specifying outer object only, 186
- 3D graph, 163–65
- 3D Object Builder, 187–90
 - creating the object, 188
 - example, 190
 - introduction, 182
 - loading the input image, 188
 - measuring the area of an object, 188
 - measuring the volume of an object, 188
 - viewing the object, 188
- 3D, viewing data in. *See* Volume Viewer
- 5D image data files, 148
- Acquisition
 - data, 2
 - workstation, 2
- Additive method of data collection, 125
- Adjacent images, 21–22
- Adjusting thumbnail images, 88
- Aligning adjacent images, 18
- Analysis workstation, 2
- Analyzing data, 4
- Applied Precision, LLC, contacting, iii
- Arbitrary profile, 166
- Arbitrary profiles, 168–69
- Archiving, 157–59
- Area of object, measuring, 188
- ASCII file, saving coordinates to, 182
- Assigning colors, 101
- Auto-Polygon Creation icon tool, 182
- Axes of rotation, 123–24
- Background value, 145
- Bad pixel file, 18, 20–21
- BioRad MRC-600 Pic, 35
- Blended color mode, 100, 101–2
- Border rolloff, 27
- Bright spot problem, 218
- Brightness, 95–99
- Burning discs, 157–59
- Calibration, 20–21
- Capturing screens, 156–57
- CCD defects, 17
- CD Kreator tool, 157–59
- CDROM discs, 157–59
- Channels, assigning colors to, 101
- Chromatic aberration, 21–22
- Chromatic Aberration Corrector, 18
- Close Polygon icon tool, 182
- Colocalization
 - graph, 196
 - identifying potential areas of, 197–201
 - Pearson Coefficient of Correction, 194
 - ROIs, 201–2
 - tool, 194–97
- Color
 - assigning to channels, 99–102
 - blended, 101–2
 - true, 101
- Color mode, 100
- Colormap, 95–99
 - changing, 100
- Combining data, 119–22
 - with image arithmetic, 142–43
- Combining data files, 52–54
- Combining multiple tasks, 54–57
- Connecting to a DMS database, 74
- Constrained iterative algorithm, 16
- Contact Applied Precision, iii
- Contrast, 95–99
- Conventions, document, ii
- Converting file formats, 31
- Convolution filters, 138–39
- Correct Image tool, 17, 18–19
- Correcting images, 17–25
 - chromatic aberration, 18, 21–22
 - inconsistent illumination intensity, 17
 - motion artifacts, 18
 - photo-bleaching, 17
- Cover slip, 112
- Create Polygon icon tool, 182
- Cropping regions, 44–50

- Cross sections, 111–12
- Crosstalk, calculating, 209
- Customer Service Hotline, iii
- Cut Mask, 49
- Dark halo, 218
- Data
 - Presenting, 85
 - Visualizing, 85
- Data collection, methods of, 125
 - additive, 125
 - maximum intensity, 125
 - mixed, 125
 - progressive, 125
 - RGB/opacity, 125
 - VolPack, 125
- Data files, 148
 - combining, 52–54
 - defined, 148
 - filtering, 137–45
 - opening, 88–89
- Data Inspector, 163–65
 - 3D graphing with, 165
 - and statistics, 169–74
- Database integration, 73–84
- Deconvolution, 9–16
 - about, 10
 - constrained iterative algorithm, 16
 - Deconvolve dialog box, 12
 - holes, 218
 - nearest neighbor method, 10, 16
 - options, 14–16
 - residuals, 216–17
 - run options, 13
 - single image, 11–13
 - standard method, 10, 16
 - using a queue, 13–14
 - visually assessing, 217
- Delete Selected Point icon tool, 182
- Delete Selected Polygon icon tool, 182
- DeltaVision
 - archiving files from, 157–59
 - exporting files from, 151–56
 - saving files for, 148–51
- Display border tools, 103
- Display channel, 95
- Distance, measuring, 175–77
 - options, 176
 - Tutorial, 176
- DMS
 - about, 73
 - connecting to database, 74
 - integration, 73–84
 - uploading images, 74–81
- Document conventions, ii
- DVD discs, 157–59
- Edge enhance, 139–40
- Edge objects, excluding in 2D Polygon Finder, 186
- Edit polygon tools, 182–85
- Edit Polygon tools
 - for gathering statistics, 41–44
- Enhance contrast, 143–44
- Equalize intensities. *See* Equalize time points
- Equalize time points, 18, 19–20
- Experiments
 - FRAP, 204–8
 - FRET, 208–13
- Explorer, 5
- Exporting
 - DeltaVision files, 151–56
 - to JPEG files, 155
 - to Movie files, 153–55
 - to PhotoShop, 153
 - to TIFF, 151
- Files
 - archiving, 157–59
 - data, 148
 - exporting DeltaVision, 151–56
 - formats of, 31
 - graphic, 148
 - saving as JPEG, 155
 - saving as movie, 153–55
 - saving as TIFF, 151
 - saving DeltaVision, 148–51
- Filter2D, 140–42
- Filtering images, 137–45
 - convolution, 138–39
 - enhancing boundaries, 139–40
 - removing noise, 140–42
 - softWoRx filters, 138
- Filters
 - activating multiplexed sets, 63–67
- Fluctuation of intensity, 219
- Fluorescence recovery after photo-bleaching.
See FRAP
- Fluorescence resonance energy transfer. *See* FRET
- Fluorescence, punctate, 15
- FRAP, 204–8
- FRET, 203, 208–13
 - analysis tool, 208–13
 - analysis workflow, 209
 - analyzing results, 212

- calculating, 211
- calculating crosstalk, 209
- efficiency statistics, 212
- prerequisites, 209
- Graph, ratio, 62
- Grayscale, 100
 - about, 99
 - assigning from blended color mode, 102
 - colormap, 100
- Hide border tools, 103
- Hide channel, 95
- Histogram, 163–65
- Image arithmetic, 142–43
- Image fusion, 52–54
- Image graphic files, 148
- Image viewer
 - cross sections, 111–12
- Image window, 102–7
 - about, 89–90
 - border tools, 103
 - display controls, 92
 - menu, 90
 - scale bar, 103–5
 - status bar, 90
 - toolbar, 90
 - zoom wheel, 94
- Images
 - adjacent, 21–22
 - analyzing quality, 215–19
 - and colocalization, 194–97
 - and colocalized structures, 197–201
 - and intensities, 162–74
 - and line intensity, 167–68
 - archiving, 157–59
 - arithmetic, 142–43
 - calibrate, 20–21
 - calibrating, 18
 - combining data, 52–54
 - contrast and brightness, 95–99
 - correction, 17–25
 - cropping, 27, 44–50
 - cross sections, 111–12
 - deconvolve single, 11–13
 - deconvolving, 9–16
 - enhancing boundaries, 139–40
 - filtering, 137–45
 - graphic files, 148
 - Inovision ISee, 33
 - intensity threshold, 144–45
 - interpolation, 93, 94
 - local contrast enhancement, 143–44
 - opening, 88–89
 - ratio, 57–62
 - reorienting, 105–11
 - repositioning, 92
 - resizing, 105–7
 - RGB opacity method, 131–35
 - rotating, 107–9
 - saving as JPEG, 155
 - saving as movie, 153–55
 - saving as PhotoShop files, 153
 - saving as TIFF files, 151
 - saving volume rendered, 135
 - scaling, 95–99
 - scaling pixel intensity, 143–44
 - selecting, 40–44
 - stitching, 25–29
 - time-lapse, 19–20
 - uploading to DMS, 74–81
 - viewing as movies, 107–12
 - viewing data, 87–112
 - visually evaluating, 217–19
 - window, 102–7
 - with bright spots, 218
 - with dark halo issues, 218
 - with deconvolution holes, 218
 - with invalid OTF, 219
 - with multiple Z sections, 27–29
 - with pebble grain texture, 218
 - with poor Z resolution, 218
- Importing
 - BioRad MRC-600 Pic files, 35
 - file formats, 31
 - Inovision ISee files, 33
 - TIFF images, 32
- Inconsistent illumination intensity, 17
- Inovision ISee™, 33
- Insert Point icon tool, 182
- Intensity data, 162–74
 - and viewing rows or columns, 167–68
- Intensity fluctuation, 219
- Intensity scale, 95–99
 - restore defaults, 98
- Intensity threshold, 144–45
- Intensity values, plotting, 166
- Interactive image rotation, 128–31
- Internet Address, Applied Precision, iii
- Interpolated images, 93, 94
- Invalid OTF, 219
- Irregular data regions
 - cropping, 47–50
 - selecting, 41–44
- JPEG files, 155
- K3b, 157–59

- Large images, 25–29
- Line intensity, 167–69
 - viewing in any direction, 168–69
- Line Profile tool, 167
- Line profiles, 167–68
- Linear intensity scales, 95–99
- Live specimen tracking, 114–18
- Local contrast enhancement, 143–44
 - dialog box, 143
- Local mean, 143
- Maximum intensity, 125
- Maximum perimeter, 186
- Measuring
 - area of an object, 188
 - chromatic correction, 21–22
 - colocalization, 194–97
 - colocalized ROI, 201–2
 - distance. *See* Distance, measuring
 - FRAP data, 204–8
 - FRET data, 208–13
 - residuals, 216–17
 - velocity of particle movement, 177–79
 - volume of an object, 188
- Methods of data collection. *See* Data collection, methods of
- Minimum perimeter, 186
- Mixed method of data collection, 125
- Modeling, 181–92
 - 2D Polygon Finder, 182–85
 - 3D Object Builder, 187–90
 - Edit Polygon, 182–85
 - Edit Polygon tool, 182–85
 - example, 190
 - summary of process, 182
- Modes
 - blended color, 101–2
 - color, 101
 - color and grayscale, 99–102
 - switching between, 100
- Motion artifacts, 18, 21
- Movie files, 153–55
- Movie players, 154
- Movies
 - time-lapse, 113–14
 - trails, 114–18
 - viewing, 107–12
 - volumetric, 113–14
- Multi-layered images, 27–29
- Multiple Segment method, 177
- Multiplexed wavelength
 - experiment design, 68–71
 - filter set activation, 63–67
 - option, 62–71
 - overview, 63
- Multi-Point FRAP, 205
- Navigating Z sections, 90–92
- Nearest neighbor, 16
 - deconvolution, 10
- Nonlinear intensity scales, 95–99
- Nonlinear local contrast enhancement. *See* Local contrast enhancement
- Normalize time points, 18, 19–20
- Normalized residual, 217
- Opening files, 88–89
- Optical File Transfer function, 11
- Optional components, 5
- Orthogonal viewer, 111–12
- OTF, 11
 - invalid, 219
- Panel collection feature, 25
- Panels, 25–29
- Particle movement, 114–18
 - measuring velocity of, 177–79
- Pearson Coefficient of Correction, 194
- Pebble grain texture, 218
- Photo-activation, 203
- Photo-bleaching, 203
 - correcting, 17
 - for FRAP, 204
- Photokinetics, 3, 203–4
- PhotoShop files, 153
- Point values, 162
- Polygon Editor, 182–85
 - maximum points, 186
 - minimum points, 186
- Polygon smoothing, 186
- Polygons. *See* 2D Polygon Finder, 3D Object Builder, Modeling
- Poor Z resolution, 218
- Presenting data, 85
- Primary light path, 63
- Printing, 159
 - screens, 156–57
- Process chains, 54–57
- Processing data, i, 3, 7
- Progressive method of data collection, 125
- Projections, 119–35
 - quick, 119–22
 - volume, 122–35
- Punctate fluorescence, 15
- QLM. *See* Quantifiable Laser Module

- Quality of images, 215–19
- Quantifiable Laser Module, 3, 203
- Queue
 - for deconvolution, 13–14
 - manager dialog box, 14
- Queue manager, 56, 79
- Quick projection, 119–22
- Ratio graph, 62
- Ratio imaging, 57–62
- Rectangular data regions
 - cropping, 45–47
 - selecting, 40–41
- Reducing time data, 50–52
- Region of interest, 165
- Removing noise, 140–42
- Rendering with volume viewer, 122–24
- Reorienting images, 105–11
- Repositioning images, 92
- Resample2D, 105–7
- Residuals, 216–17
 - weak convergence of, 217
- Resizing images, 105–7
- Resolve3D, 2
- RGB opacity method, 131–35
- RGB/opacity method of data collection, 125
- ROI. *See* Region of interest
- ROI colocalization, 201–2
- Rotating images, 107–9, 128–31
- Rotation angles, 107–9
- Rotation axes, 123–24
- Rotation center, 110
- Save as SYLK spreadsheet dialog box, 169
- Saving
 - as JPEG, 155
 - as Movie, 153–55
 - as PhotoShop, 153
 - as TIFF, 151
 - DeltaVision files, 148–51
 - statistics records, 170
 - volume rendered images, 135
 - with overlay graphics, 153–55
- Scale bar, 103–5
- Scaling pixel intensity, 143–44
- Screen shots, 156–57
- Scrolling Z sections, 91
- Secondary light path, 63
- Section intensity fluctuation, 219
- Select Polygon icon tool, 182
- Selecting
 - irregular regions, 41–44
 - rectangular regions, 40–41
- Selecting data, 40–44
- Show channel, 95
- Single-Point FRAP, 205
- Smoothing pixel distance in 2D Polygon Finder, 186
- Snapshots, 156–57
- Spherical aberration, 218
- Standard residual, 216
- Standard Two Point method, 176
- Statistics
 - calculating for selected areas, 169–74
 - using Data Inspector for, 170
 - with Edit Polygon tools, 41–44
- Status bar, 90
- Stitching, 25–29
 - border rolloff, 27
 - example, 26
 - images with multiple Z sections, 27–29
- Table of 3-D measurements, 190
- Task Builder, 54–57
- Task chains, 54–57
- Technical support, iii
- Threshold
 - default value for 2D Polygon Finder, 186
- Thumbnails, adjusting, 88
- TIFF output formats, 152
- TIFF, importing, 32
- Time points
 - reducing, 50–52
 - viewing, 90–92
- Time-lapse image data, 19–20, 59
- Time-lapse movies, 107–12, 113–14
- Tools
 - Align Image, 18, 21–22
 - calibrate, 18
 - calibration, 20–21
 - chromatic aberration corrector, 18, 21–22
 - colocalization, 194–97
 - convolution, 138–39
 - copy region, 53
 - Correct Image, 17, 18–19
 - correction, 17–25
 - Cut Mask, 49
 - Data Inspector, 163–65
 - deconvolve, 9–16
 - edge enhance, 139–40
 - edit header, 53
 - Edit Polygon, 41–44
 - equalize time points, 18, 19–20
 - Filter2D, 140–42

- for calculating statistics, 169–74
 - for identifying colocalized structures, 197–201
 - for measuring velocity, 177–79
 - image arithmetic, 142–43
 - image fusion, 52–54
 - image window, 90, 102–7
 - intensity threshold, 144–45
 - interactive image rotation, 128–31
 - Line Profile, 167
 - local contrast enhancement, 143–44
 - measuring distances, 175–77
 - nearest neighbor, 10
 - orthogonal viewer, 111–12
 - Point values, 162
 - Polygon Editor, 182–85
 - projections, 119–35
 - queue manager, 56, 79
 - quick projection, 119–22
 - removing data below a threshold, 144–45
 - reorienting, 105–11
 - Resample2D, 105–7
 - resizing, 105–7
 - ROI colocalization, 193, 201–2
 - rotating, 107–9
 - scale bar, 103–5
 - scale image, 95–99
 - switching border on/off, 103
 - time-lapse movies, 113–14
 - trails movie, 114–18
 - viewing movies, 107–12
 - volume rendering, 122–24
 - volume viewer, 122–35
 - volumes, 119–35
- Tracking particle movement, 114–18
- Trimming time data, 50–52
- True color, 101
- Tutorials
- Measuring distances, 176
 - modeling, 190
- UIC MetaMorph STK, 36
- Undo Last icon tool, 182
- Uploading images to DMS, 74–81
- URL, Applied Precision, iii
- Velocity, 177–79
- measuring, 175
- Viewing data in 3D. *See* Volume Viewer
- Viewing image data, 87–112
- Viewing movies. *See* Movies
- Viewing time points, 90–92
- Viewing Z sections, 90–92
- Visualizing data, 3
- VolPack, 126
- Volume models, 181–92
- Volume of object, measuring, 188
- Volume projections, 122–35
- Volume rendering, 122–24
- Volume viewer, 122–35
- axes of rotation, 123–24
 - interactive image rotation, 128–31
 - limiting size of data set, 124
 - rendering, 122–24
 - RGB opacity method, 131–35
 - using with colocalization, 199
- Volumes, 119–35
- Volumetric movies, 113–14
- Voxels, 27
- Weak convergence, 217
- Web site, Applied Precision, iii
- Z sections
- and blurring, 218
 - data projections, 119–35
 - measuring points in, 175–77
 - slider, 91
 - viewing, 90–92
 - viewing as movie, 107–12
- Zooming, 93
- Z-stack image, 10